XKM 3200 WL LMD

da	Brugsanvisning Kommunikationsmodul
de	Gebrauchsanweisung Kommunikationsmodul
en	Operating instructions Communication module
fi	Käyttöohje Tiedonsiirtomoduuli
fr	Mode d'emploi Module de communication
it	Istruzioni d'uso Modulo di comunicazione
nl	Gebruiksaanwijzing Communicatiemodule
no	Bruksanvisning kommunikasjonsmodul
sv	Bruksanvisning kommunikationsmodul

da	5
de	25
en	45
fi	65
fr	85
it	105
nl	125
no	145
sv	165

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Notes about these instructions	45
Queries and technical problems	45
Definition of terms	45
Module	45
Intended use	46
Items supplied	47
Scope of delivery	47
Technical data	47
EU Conformity declaration.....	47
Guide to the appliance	48
Module overview	48
Network connection	48
Switching on and off	49
LED indicator.....	49
Button	49
WPS button.....	49
Reset button.....	49
Data protection and data security	50
Factory default settings for network configuration	50
Copyright and licences	50
Warnings and safety notes	51
Fitting and installation	52
Getting ready to use WiFi	52
Range and availability of WiFi signal.....	52
LAN cable	52
WiFi password	52
Installation.....	53
Module slot.....	53
Inserting the module	54
Laying data cables	54
Switching the cleaning machine on	54
Network connection	55
Establishing the WiFi connection	55
LAN cable	55
WPS login	55
WiFi access point	55
Configuring the interface.....	55
Opening the menu	56
Checking the connection status.....	56
Selecting settings	57
Activating the configuration	58
Fault 559.....	58
Web interface	59
Password	59
Configuring the module	59
Firmware update.....	60
PIN for Cloud Service	61

en - Contents

Caring for the environment **62**
Disposal of the packing material 62

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

- Select an option using the arrow buttons and save your choice with *OK*.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings  menu.

Queries and technical problems

In the event of queries or technical problems, please contact Miele. Contact details can be found at the end of the operating instructions for your cleaning machine or at www.miele.com/professional.

Definition of terms

Machine

In these operating instructions, the laboratory washer is referred to as “the machine”.

Module

The communication module is generally referred to as the module in these operating instructions.

The XKM 3200 WL LMD communication module supplied enables an interface to be set up between a Miele washer-disinfector and a terminal device or local network. The data connection is a standard Ethernet or WiFi connection. The interface serves a number of purposes, including process documentation. Further information on process documentation can be obtained from Miele.

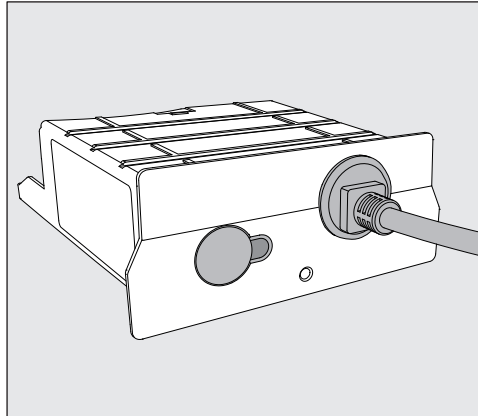
Only terminal devices (PCs, tablets, printers, etc.) compliant with IEC/EN 62368 should be used.

Only specialists are permitted to configure the interface.

The communication module can be used in the following washer-disinfectors:

- PG 8562
- PG 8581
- PG 8582
- PG 8582 CD
- PG 8583
- PG 8583 CD
- PG 8591
- PG 8592
- PG 8593

Scope of delivery

- XKM 3200 WL LMD communication module

Technical data

Voltage, rated load, fuse rating	See data plate
Mains cable	Approx. 1.5 m
Ethernet baud rate	10/100 Mbit/s
WiFi standard	802.11 b/g/n
WiFi frequency band	2400-2483.5 MHz
Maximum WiFi transmission power	< 100 mW
Temperature during operation	Up to max. 80 °C
CE marking	2014/53/EU Radio Equipment Directive

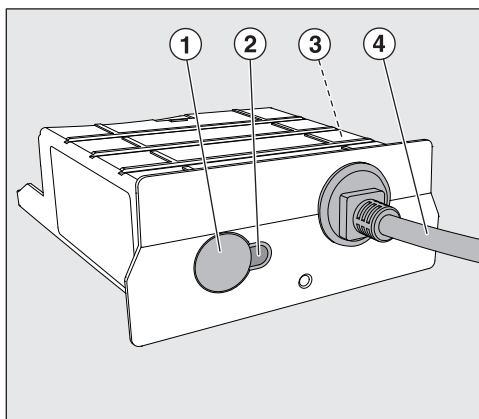
EU Conformity declaration

Miele hereby declares that the XKM 3200 WL LMD communication module complies with Directive 2014/53/EU.

The complete text of the EU declaration of conformity can be downloaded from the following website by quoting the product name or serial number:

- www.miele.com/professional
 - > Products > Download
- or
- > Service & Support > Operating instructions

Module overview

- ① Reset/WPS button
- ② LED indicator
Indicates operational readiness after installation and active communication when in operation
- ③ Data plate (located on the side)
- ④ Data cable

Network connection

Integrating the cleaning machine into the on-site network requires the following:

- Ethernet network or WPA2-secured 2.4 GHz WiFi 4 network
- DHCP server or specifications for static IPv4 settings
- WiFi: network name (SSID) and password
- Network printer (compatible with Internet Printing Protocol and PostScript 3): printer URL and port

Switching on and off

The module does not have a separate On/Off switch. It is switched on and off together with the cleaning machine.

The module requires approx. 90 s for initialisation after switch-on. Until this has happened data cannot be transferred and interface settings cannot be changed.

LED indicator

The LED indicator (light emitting diode) provides status information about the communication module.

LED	Status
Lit up GREEN	Connection to the network was successful. The module is ready for operation.
Lit up YELLOW	Network is configured but there is no connection to the network.
Lit up RED	The module is not configured.
Off	Power is not being supplied to the module. Check that the module is in the correct position in the module slot.

Button

The button on the module has several functions.

WPS button

The WPS button can be used to establish a WiFi connection between the module and your router or repeater (WPS, WiFi, Protected Setup). WPS-enabled routers or repeaters must be used.

Reset button

The button on the module can be used to reset the module to the factory default settings.

- Press the button for approx. 5–9 seconds to restore the factory default settings.

Data protection and data security

When you activate the networking function and connect your machine to the Internet, your machine sends the following data to the Miele Cloud:

- Machine serial number
- Machine model and technical features
- Machine status
- Information about the software version of your machine

Initially, this data cannot be assigned to a specific user and is not saved permanently. Data cannot be saved permanently or assigned to a specific user until after you have linked your machine to a user. Data transmission and processing are governed by Miele's strict security standards.

Factory default settings for network configuration

You can reset all of the settings on the communication module or your integrated WiFi module to the factory default settings. The network configuration should be reset whenever an appliance is being disposed of or sold, or if a used appliance is being put into operation. This is the only way to ensure that all personal data has been removed and the previous owner will no longer be able to access the appliance.

Copyright and licences

For the purpose of operating and controlling the communication module, Miele uses proprietary or third-party software that is not covered by open source licensing terms. This software/these software components are protected by copyright. The copyrights held by Miele and third parties must be respected.

Furthermore, this communication module contains software components which are distributed under open source licensing terms. The open source components contained in the machine along with the corresponding copyright notices, copies of the licensing terms valid at the time and any further information can be accessed locally by IP using a web browser (*<https://<IP address>/Licenses>*). The liability and warranty arrangements for the open source licences displayed in this location only apply in relation to the respective rights holders.

To avoid the risk of accidents and damage to this communication module please read these instructions carefully before using it for the first time.

Keep these instructions in a safe place.

 Please also read the operating instructions for your washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ This communication module must only be used for applications specified in the operating instructions. Changes or modifications to it and incorrect use of it are not permitted as this can be dangerous.
- ▶ The communication module must only be installed, commissioned and serviced by Miele service or by a suitably qualified person. Repairs are only to be undertaken by Miele Service. Repair work by unqualified persons could be dangerous.
- ▶ Check the communication module carefully before installation for any visible damage. A damaged module could be dangerous and must not be installed.
- ▶ Before installing the communication module compare the connection data (voltage and frequency) on the data plate with that on the data plate of your washer-disinfector. If the data does not match, the module could sustain damage. If in any doubt, consult an electrician.
- ▶ The communication module housing must not be opened on any account.
Touching live electrical components or altering the electrical or mechanical components is dangerous and could cause the module to malfunction.
- ▶ Make sure that the data cable for the communication module is assembled and routed correctly. The cable must not be kinked or trapped, for example. It must not be routed around sharp edges or be placed under tension.
- ▶ The WiFi technology supported by the module can be temporarily or completely disrupted e.g. by structural conditions or the set-up location. For this reason availability of featured functions cannot always be guaranteed.

Getting ready to use WiFi

If you wish to use the WiFi connection on the module, please note the following prior to installation.

Range and availability of WiFi signal

The WiFi range or signal strength is dependent on the structural environment. Reinforced concrete walls and ceilings, metal surfaces, metal-vaporised films, glass screens and fire protection constructions can muffle or divert the signal, therefore reducing the range.

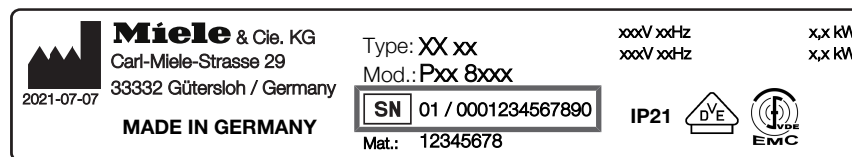
If the cleaning machine is located under a metal worktop or is surrounded by metal cabinets, for example, the signal will be almost completely blocked.

LAN cable

Do not connect the LAN cable to the network. Plugging in the LAN cable switches off the WiFi signal automatically. The technology does not support the simultaneous use of WiFi and LAN.

WiFi password

- Take a note of the serial number of your cleaning machine. This is required as the password for establishing the WiFi connection. You will find the serial number **SN** on the data plate on the cleaning machine. The position of the data plate is indicated in the machine description for the cleaning machine.
Example:

The serial number of the module, without the initial zeroes, acts as the password. Example:

SN 01 / **0001234567890**

Password: **1234567890**

Installation

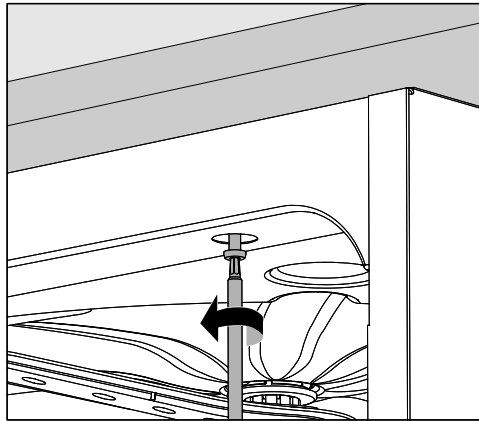
⚠ Before fitting the communication module compare the connection data (voltage and frequency) on the data plate with that on the data plate of your cleaning machine. If the data does not match, the module could sustain damage. If in any doubt, consult a qualified electrician.

Module slot

- Disconnect the cleaning machine from the mains electricity supply.

The module is inserted into the module slot on the cleaning machine. The module slot is located on the back of the cleaning machine.

- If the cleaning machine is installed in a niche:
 - Pull the cleaning machine forwards out of the niche.
- If the cleaning machine is installed under a worktop:
 - Open the door.

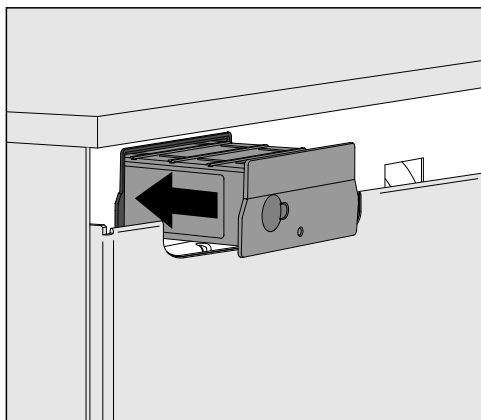
- Unscrew the fixing screws.
- Pull the cleaning machine out from under the worktop. Take care not to damage the hoses and cables when moving the machine.

en - Fitting and installation

Inserting the module

- Remove any stickers or protective caps as necessary from the slot aperture.

- Push the module into the module slot. Gently push the module until it clicks into place.

Laying data cables

The technology does not support the simultaneous use of WiFi and LAN.

WiFi

If you wish to use the WiFi connection, the LAN cable must not be connected to a network or terminal device. The WiFi signal will only be available if the LAN cable is not plugged in.

- We recommend that you roll the LAN cable up and secure it somewhere at the back of the machine. This will stop it from getting in the way when the cleaning machine is being installed or undergoing maintenance work.

LAN

If you wish to use the LAN cable for data transfer instead of the WiFi connection, connect the LAN cable to your terminal device. Plugging in the LAN cable switches off the WiFi signal automatically.

- Take care to ensure that the cable is laid correctly; it should not be kinked or trapped, for example. Keep the cable away from sharp edges and do not place it under tension.

Only terminal devices (PCs, printers, etc.) which comply with EN/IEC 62368 should be used.

Switching the cleaning machine on

- Connect the cleaning machine to the electricity supply.
The module is supplied with mains voltage by the cleaning machine. No additional power connection is required.

After around 90 s, the module is ready for operation and the interface can be configured.

Establishing the WiFi connection

LAN cable

Do not connect the LAN cable to the network. Plugging in the LAN cable switches off the WiFi signal automatically. The technology does not support the simultaneous use of WiFi and LAN.

WPS login

WPS login requires the use of a WPS-enabled WiFi router or repeater.

- Switch the cleaning machine on.
After around 90 s, the module is ready for operation.
- Briefly press the button on the module (1–4 seconds).
Never press the button for more than 4 seconds, as this will reset the module to the factory default settings; see Reset button.
- Activate the WPS function on your WiFi router or repeater within 2 minutes.

The LED on the module will light up green if the WPS login is successful.

WiFi access point

If you do not wish to set up a WPS connection and the module has not yet been configured for a WiFi network, a WiFi access point will open automatically when the cleaning machine is switched on. You can connect the access point to your PC.

- To do this, switch the cleaning machine on.
After around 90 s, the module is ready for operation.
- Configure the interface of your cleaning machine. To do this, follow the instructions in the “Configuring the interface” chapter.
- Open the module’s web interface and configure the interface under “WiFi Settings”; see “Configuring the interface”.

Configuring the interface

Unauthorised access poses a risk.

Settings in the cleaning machine, e.g. parameters for disinfection or the dispensing of process chemicals, may be changed as a result of unauthorised access via the network.

Under no circumstances should it be possible to access the machine via the Internet or other public or unsecured networks, either directly or indirectly (e.g. using port forwarding).

Only specialists are permitted to configure the interface.

en - Network connection

Opening the menu ■ Switch the cleaning machine on using the  button.

You must wait approx. 90 seconds before starting to configure the interface.

■ Open the menu as follows:

▸ button

▸ Further settings

▸ Interface

■ Select the Ethernet interface and confirm your choice with *OK*.

The following options are available:

- Module status

Displays the connection status.

- DHCP

The Ethernet interface can either be implemented via a Dynamic Host Configuration Protocol (DHCP) or by setting the individual parameters.

Checking the connection status

■ Select the Module status option and confirm your choice with *OK*.

The following messages may appear in the display:

- Active

The module is ready for operation.

- Initialising, module not yet active

The module has not finished initialising. This process takes approx. 90 s.

- Inactive

No module has been detected.

Check the installation of the module. Switch the cleaning machine off and then back on again.

If the message continues to appear, contact the Miele Customer Service Department.

Selecting settings

Document all settings in case technical support is required at a later date.

DHCP must be deactivated prior to manual intervention.

- Select the DHCP option and confirm your choice with *OK*.

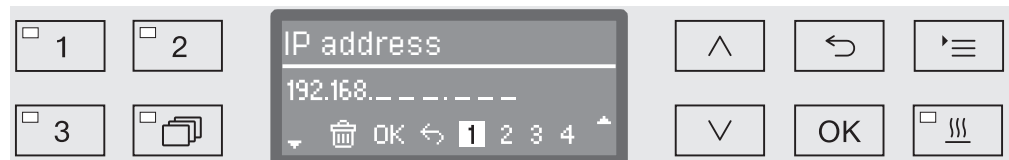
- Select the Off option and confirm your choice with *OK*.

The following settings are then available for selection:

- IP address
- Subnet mask
- Standard gateway
- DNS server automatic
- DNS server 1
- DNS server 2
- Port type (not relevant)
- Port (not relevant)

- Select an option and confirm your choice with *OK*. You can then make the following settings according to the information displayed on the screen.

To enter the IP address, Subnet mask etc., the display contains the following information:

The current setting is shown on the middle line of the display. It can be changed using the options shown in the bottom line. The top line shows which option has been selected from the bottom line. The following data entry options are available:

- Numbers from 0 to 9.
- Decimal point “.”.
- Use the  symbol to delete the last position.
- Select the *OK* symbol on the display to save the setting. The display will then revert to the main menu.
- Select the  symbol on the display or use the  button to exit the process without saving the setting(s) made. The display reverts to the main menu.

The IP address, Subnet mask, etc. must always be entered in full. If the IP address is 192.168.1.23, for example, enter 192.168.001.023.

- Use the arrow buttons \wedge (right) and \vee (left) to move the cursor to the option you require.
- Confirm each entry with *OK*.
- To save your entries, select the *OK* symbol on the display and confirm with the *OK* button.

Activating the configuration

The interface configuration is only effective after the cleaning machine has been restarted.

- Wait approx. 10 seconds before switching the cleaning machine off for a restart.
- Switch the cleaning machine off and then back on again with the  button.
- Wait approx. 90 seconds for the module to initialise.
- Then check if the interface is working.

Fault 559

If the fault message Fault 559 is displayed during commissioning or a programme is interrupted with this message, the RS232 interface is activated and must be deactivated.

- Open the menu for configuring the RS232 interface via Further settings / Interface.
- Select the Ethernet option and confirm your choice with *OK*.

The RS232 interface is deactivated. After around 90 s, the module is ready for operation.

Web interface

The XKM 3200 WL LMD module has a dedicated web interface via which the network settings can be configured, printers can be set up or the password can be changed.

The interface can be accessed from any workstation on your local network. A configured interface in the controller of the cleaning machine, an up-to-date web browser and an existing network connection to the cleaning machine are required.

The web interface is divided into websites. The output language is English.

- Enter one of the following addresses in the address bar of your browser (where *<IP-address>* corresponds to the IP address of your cleaning machine).

Password

<https://<IP-address>/password.html>

When the web interface is opened for the first time, the “Admin” user must set up a password for their profile (once only). A module reset will reset the module settings to the factory default settings; this will include resetting the password for the “Admin” user.

- Change Password

Changing the module password

The “Admin” user and the password must always be set at the factory as a first step before further settings can be made. When setting the password for the first time, the field for the old password must be left blank.

Configuring the module

<https://<IP-address>/index.html>

To change the settings, you must enter a user name and a password. You cannot access the settings without doing this.

- WiFi Settings

Selecting a WiFi network

- IP Configuration

Modifying the interface configuration

- Network Printer

Specifying a network printer for printing reports

- Reset

Restart the module or restore the factory default settings (this will also delete user profiles and module passwords).

Firmware update *<https://<IP-address>/update.html>*

You can use this option to update the firmware for your XKM module. The software version of your cleaning machine will not be affected. There is a different process for updating cleaning machines.

Updates should be carried out when the interface is not being used to query process data, for example. Updates must not be carried out while a cleaning programme is in progress.

For firmware updates, the cleaning machine must be connected to the Internet.

- NewerAvailable
 - true
An update is available.
 - false
The installed firmware version is up to date.
- CurrentVersion
Designation of the firmware version that is currently installed.
- AvailableVersion
Designation of the available firmware version if an update is available (NewerAvailable: true).
- Type
Model identifier of the XKM module in use.
- UpdateInProgress
 - true
An update is being installed.
 - false
An update is available but installation has not yet started.
- *Check for Update* button
This button can be used to check whether an update is available.
- *Start Update* button
Starts installing the update if an update is available (NewerAvailable: true). The process may take several minutes. The button is greyed out if no update is available (NewerAvailable: false).

PIN for Cloud Service

<https://<IP-address>/tan.html>

This page allows you to view the PIN for accessing the Miele Cloud Service and to request a new PIN if necessary. Access to this page is protected with the module administrator password.

You must register your cleaning machine with the Miele Cloud in order to use the Miele Cloud Service. Once cleaning machines have been registered, they will log into the Cloud automatically each time they are switched on, provided they have an active Internet connection.

- Serial number

 - Serial number of the cleaning machine (see data plate).

- PIN

 - PIN for unique identification with the Cloud Service (pairing).

- *Refresh* button

 - To request a new pairing PIN. The PIN is automatically generated by the Cloud.

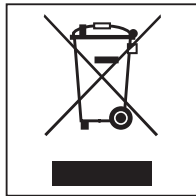
Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

Disposing of your old appliance

Electronic and electrical appliances contain many valuable materials. They also contain certain materials, compounds and components which were essential for their correct functioning and safety. These could be hazardous to your health and to the environment if disposed of with general waste or if handled incorrectly. Please do not, therefore, dispose of your old appliance with general waste.

Instead, please make use of officially designated collection and disposal points to dispose of and recycle electrical and electronic appliances in your local community, with your dealer, with Miele or with the manufacturer. By law, you are solely responsible for deleting any personal data from the appliance prior to disposal. You are legally obliged to remove any old batteries which are not securely enclosed by the appliance and to remove any lamps without destroying them, where this is possible. These must be taken to a suitable collection point where they can be handed in free of charge. Please ensure that your old appliance poses no risk to children while being stored prior to disposal.

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