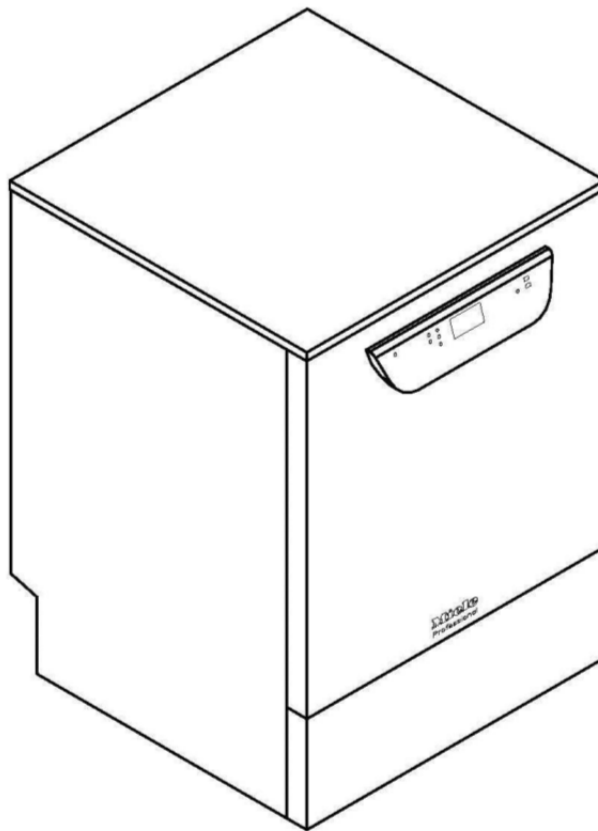
Miele

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PG 8591 - 8593

Legend:

Circled, bold-type abbreviations:
 Connection required

Abbreviations surrounded by broken circle:
 Connection optional or required, depending on model

(EL)	Electrical connection PG 8591 PG 8592	Voltage (default) Rated load Fuse rating Minimum cross-section Length of electrical cord (UL 921) with plug Electrical connection	2 AC 208V 60 Hz 6.2 kW 2 x 30 A 3 x AWG 10 approx. 5'9" NEMA L6-30	2 AC 208V 60 Hz 6.2 kW 2 x 30 A 3 x 5.2 mm ² approx. 1.8 m NEMA L6-30
(EL)	Electrical connection PG 8591 PG 8592	Voltage (convertible) Rated load Fuse rating Minimum cross-section Length of electrical cord (UL 921) with plug Electrical connection	3 AC 208V 60 Hz 6.2 kW 3 x 20 A 4 x AWG 12 approx. 5'9" NEMA L15-20	3 AC 208V 60 Hz 6.2 kW 3 x 20 A 4 x 3.3 mm ² approx. 1.8 m NEMA L15-20
(EL)	Electrical connection (240V Option PG8591)	Voltage (default) Rated load Fuse rating Minimum cross-section Length of electrical cord (UL 921) with plug Electrical connection	2 AC 240V 60 Hz 6.2 kW 2 x 30 A 3 x AWG 10 approx. 5'9" NEMA L6-30	2 AC 240V 60 Hz 6.2 kW 2 x 30 A 3 x 5.2 mm ² approx. 1.8 m NEMA L6-30
(EL)	Electrical connection (240V Option PG8591)	Voltage (convertible) Rated load Fuse rating Minimum cross-section Length of electrical cord (UL 921) with plug Electrical connection	3 AC 240V 60 Hz 6.2 kW 3 x 20 A 4 x AWG 12 approx. 5'9" NEMA L15-20	3 AC 240V 60 Hz 6.2 kW 3 x 20 A 4 x 3.3 mm ² approx. 1.8 m NEMA L15-20
(EL)	Electrical connection (PG 8593)	Voltage (default) Rated load Fuse rating Minimum cross-section Length of electrical cord (UL 921) with plug Electrical connection	3 AC 208V 60 Hz 6.2 kW 3 x 20 A 4 x AWG 12 approx. 5'9" NEMA L15-20	3 AC 208V 60 Hz 6.2 kW 3 x 20 A 4 x 3.3 mm ² approx. 1.8 m NEMA L15-20
(EL)	Electrical connection (PG 8593)	Voltage (convertible) Rated load Fuse rating Minimum cross-section Length of electrical cord (UL 921) with plug Electrical connection	2 AC 208V 60 Hz 6.2 kW 2 x 30 A 3 x AWG 10 approx. 5'9" NEMA L6-30	2 AC 208V 60 Hz 6.2 kW 2 x 30 A 3 x 5.2 mm ² approx. 1.8 m NEMA L6-30

The machine is supplied with electrical cord and plug. The receptacle must be accessible after machine installation. For hard-wired machines, connection should be made via a suitable mains switch with all-pole isolation. The contact opening between all open contacts must be at least 3 mm wide and the mains switch must be lockable in the open position. If necessary, an equipotential bond with good contact connection must be provided in accordance with all appropriate national and local regulations.

Splash protection class IP 21

	Equipotential bonding and earthing	<p>There is a screw connection point at the back of the machine, to which equipotential bonding can be connected if required.</p> <p>Connect equipotential bonding and protective earth conductor!</p>																													
	Network/printer connection (optional)	<p>Optional machine interface for transmission or process data output</p> <p>Communication module (not supplied)</p> <p>Ethernet module or RS232 module (install electrical sockets near machine if required)</p> <p><i>Connections/installation must comply with IEC 60950.</i></p>																													
	Cold Water	<p>Two cold water connections are required for machines with steam condenser (DK). Use supplied Y-adaptor if only one cold water connection is available.</p> <table border="1" data-bbox="456 645 1402 1120"> <tr> <td>Max. temperature</td> <td>68 °F</td> <td>20 °C</td> </tr> <tr> <td>Max. water hardness</td> <td>73 gpg</td> <td>70 °dH</td> </tr> <tr> <td>Recommended flow pressure</td> <td>29 psi</td> <td>200 kPa</td> </tr> <tr> <td>Min. flow pressure</td> <td>14.5 psi</td> <td>100 kPa</td> </tr> <tr> <td>Max. pressure</td> <td>145 psi</td> <td>1,000 kPa</td> </tr> <tr> <td>Throughput</td> <td>2 gal/min</td> <td>7.5 l/min</td> </tr> <tr> <td>Connection thread - on site (flat sealing)</td> <td>3/4" Male Garden Hose Thread</td> <td>3/4" Male Garden Hose Thread</td> </tr> <tr> <td>Length of connection hose (supplied)</td> <td>approx. 5' 7"</td> <td>approx. 1.7 m</td> </tr> <tr> <td>Length of the steam condenser hose</td> <td>approx. 5' 7"</td> <td>approx. 1.7 m</td> </tr> </table>			Max. temperature	68 °F	20 °C	Max. water hardness	73 gpg	70 °dH	Recommended flow pressure	29 psi	200 kPa	Min. flow pressure	14.5 psi	100 kPa	Max. pressure	145 psi	1,000 kPa	Throughput	2 gal/min	7.5 l/min	Connection thread - on site (flat sealing)	3/4" Male Garden Hose Thread	3/4" Male Garden Hose Thread	Length of connection hose (supplied)	approx. 5' 7"	approx. 1.7 m	Length of the steam condenser hose	approx. 5' 7"	approx. 1.7 m
Max. temperature	68 °F	20 °C																													
Max. water hardness	73 gpg	70 °dH																													
Recommended flow pressure	29 psi	200 kPa																													
Min. flow pressure	14.5 psi	100 kPa																													
Max. pressure	145 psi	1,000 kPa																													
Throughput	2 gal/min	7.5 l/min																													
Connection thread - on site (flat sealing)	3/4" Male Garden Hose Thread	3/4" Male Garden Hose Thread																													
Length of connection hose (supplied)	approx. 5' 7"	approx. 1.7 m																													
Length of the steam condenser hose	approx. 5' 7"	approx. 1.7 m																													
	Hot Water Connection	<table border="1" data-bbox="456 1120 1402 1554"> <tr> <td>Max. temperature</td> <td>150 °F</td> <td>65 °C</td> </tr> <tr> <td>Max. water hardness</td> <td>73 gpg</td> <td>70 °dH</td> </tr> <tr> <td>Recommended flow pressure</td> <td>29 psi</td> <td>200 kPa</td> </tr> <tr> <td>Min. flow pressure</td> <td>5.8 psi</td> <td>40 kPa</td> </tr> <tr> <td>Max. pressure</td> <td>145 psi</td> <td>1,000 kPa</td> </tr> <tr> <td>Throughput</td> <td>2 gal/min</td> <td>7.5 l/min</td> </tr> <tr> <td>Connection thread - on site (flat sealing)</td> <td>3/4" Male Garden Hose Thread</td> <td>3/4" Male Garden Hose Thread</td> </tr> <tr> <td>Length of connection hose (supplied)</td> <td>approx. 5' 7"</td> <td>approx. 1.7 m</td> </tr> </table>			Max. temperature	150 °F	65 °C	Max. water hardness	73 gpg	70 °dH	Recommended flow pressure	29 psi	200 kPa	Min. flow pressure	5.8 psi	40 kPa	Max. pressure	145 psi	1,000 kPa	Throughput	2 gal/min	7.5 l/min	Connection thread - on site (flat sealing)	3/4" Male Garden Hose Thread	3/4" Male Garden Hose Thread	Length of connection hose (supplied)	approx. 5' 7"	approx. 1.7 m			
Max. temperature	150 °F	65 °C																													
Max. water hardness	73 gpg	70 °dH																													
Recommended flow pressure	29 psi	200 kPa																													
Min. flow pressure	5.8 psi	40 kPa																													
Max. pressure	145 psi	1,000 kPa																													
Throughput	2 gal/min	7.5 l/min																													
Connection thread - on site (flat sealing)	3/4" Male Garden Hose Thread	3/4" Male Garden Hose Thread																													
Length of connection hose (supplied)	approx. 5' 7"	approx. 1.7 m																													
	Demineralised water (DI)	<table border="1" data-bbox="456 1554 1402 1980"> <tr> <td>Max. temperature</td> <td>150 °F</td> <td>65 °C</td> </tr> <tr> <td>Recommended flow pressure</td> <td>29 psi</td> <td>200 kPa</td> </tr> <tr> <td>Min. flow pressure</td> <td>4.4 psi</td> <td>30 kPa</td> </tr> <tr> <td>Max. pressure</td> <td>145 psi</td> <td>1,000 kPa</td> </tr> <tr> <td>Connection thread - on site (flat sealing)</td> <td>3/4" Male Garden Hose Thread</td> <td>3/4" Male Garden Hose Thread</td> </tr> <tr> <td>Throughput</td> <td>2 gal/min</td> <td>7.5 l/min</td> </tr> <tr> <td>Length of connection hose (supplied)</td> <td>approx. 5' 7"</td> <td>approx. 1.7 m</td> </tr> </table> <p>Hose from supply container to machine to be fitted on site.</p>			Max. temperature	150 °F	65 °C	Recommended flow pressure	29 psi	200 kPa	Min. flow pressure	4.4 psi	30 kPa	Max. pressure	145 psi	1,000 kPa	Connection thread - on site (flat sealing)	3/4" Male Garden Hose Thread	3/4" Male Garden Hose Thread	Throughput	2 gal/min	7.5 l/min	Length of connection hose (supplied)	approx. 5' 7"	approx. 1.7 m						
Max. temperature	150 °F	65 °C																													
Recommended flow pressure	29 psi	200 kPa																													
Min. flow pressure	4.4 psi	30 kPa																													
Max. pressure	145 psi	1,000 kPa																													
Connection thread - on site (flat sealing)	3/4" Male Garden Hose Thread	3/4" Male Garden Hose Thread																													
Throughput	2 gal/min	7.5 l/min																													
Length of connection hose (supplied)	approx. 5' 7"	approx. 1.7 m																													

(AW)	Drain Water	Drainage temperature	200 °F	93 °C
		Length of drain hose	approx. 4' 6"	approx. 1.4 m
		Drain hose, max. drainage length	approx. 13' 1"	approx. 4.0 m
		Drain pump head height from floor level, max.	approx. 3' 3"	approx. 1 m
		Drain hose delivery, max.	4.2 gal/min	16 l/min
		On-site hose connector (external diameter x length)	7/8" x 1 3/16"	22 mm x 30 mm
		It is advisable to lay the drain hose in a curve min. 600 mm above floor level to the drainage point.		
(DS)	External dispensing (optional) (note machine model)	Connection for up to two dispensing units for liquid agents on the back panel.		
	Machine feet	Height-adjustable	0-2 3/8"	0 – 60 mm
		Foot diameter	1 3/8"	35 mm
		Entry point into housing	8 M	8 M
	Heat Dissipation to Room	Heat transfer by emission	1,194 btu/h	0.35 kWh/h
		Load on removal	1,365 btu/h	0.40 kWh/h
	Machine data	Height	32 7/8"	835 mm
		Height without lid	32 1/4"	820 mm
		Width	23 5/8"	598 mm
		Depth	23 5/8"	598 mm
		Door height	24 1/2"	622 mm
		Net weight	172 lbs	78 kg
		Dynamic floor load	1,200 N	1,200 N
		Width including transport pallet	26 3/8"	670 mm
	Depth including transport pallet	29 3/8"	740 mm	
	Height including transport pallet	36 1/4"	920 mm	
	Set-up Conditions	Ambient temperature	40-105 °F	5-40 °C
		Max. humidity	80%	80%
		Max. altitude above sea level	4,921'	1,500 m
<p>It is essential to observe the installation instructions when installing the machine.</p> <p>In order to reduce the risk of steam damage, the area around the machine should be limited to furniture and fittings that are designed for use in commercial environments.</p> <p>Installation should only be performed by Miele Service or authorized installers in accordance with valid regulations, relevant standards and health and safety codes!</p> <p>Subject to change!</p>				

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
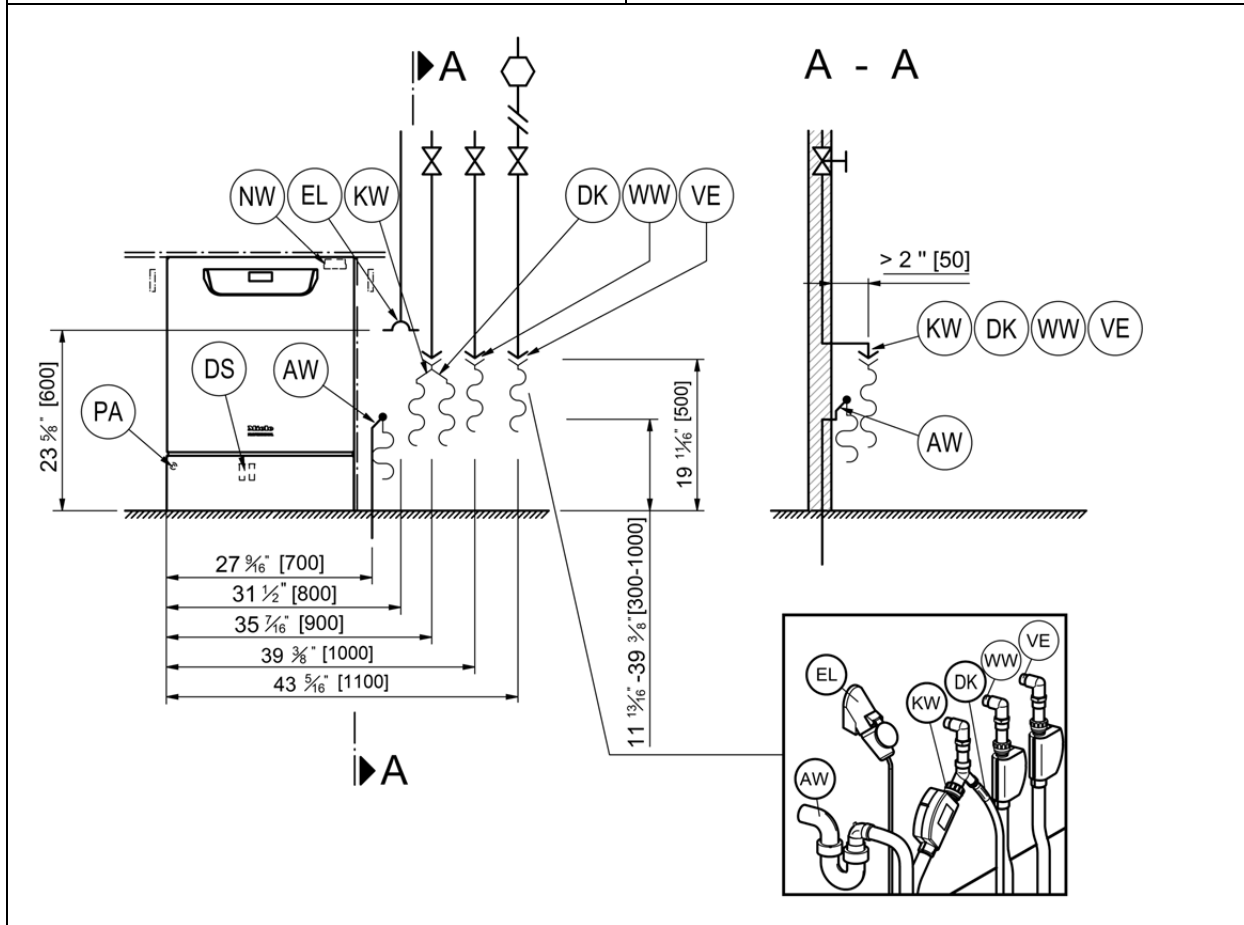
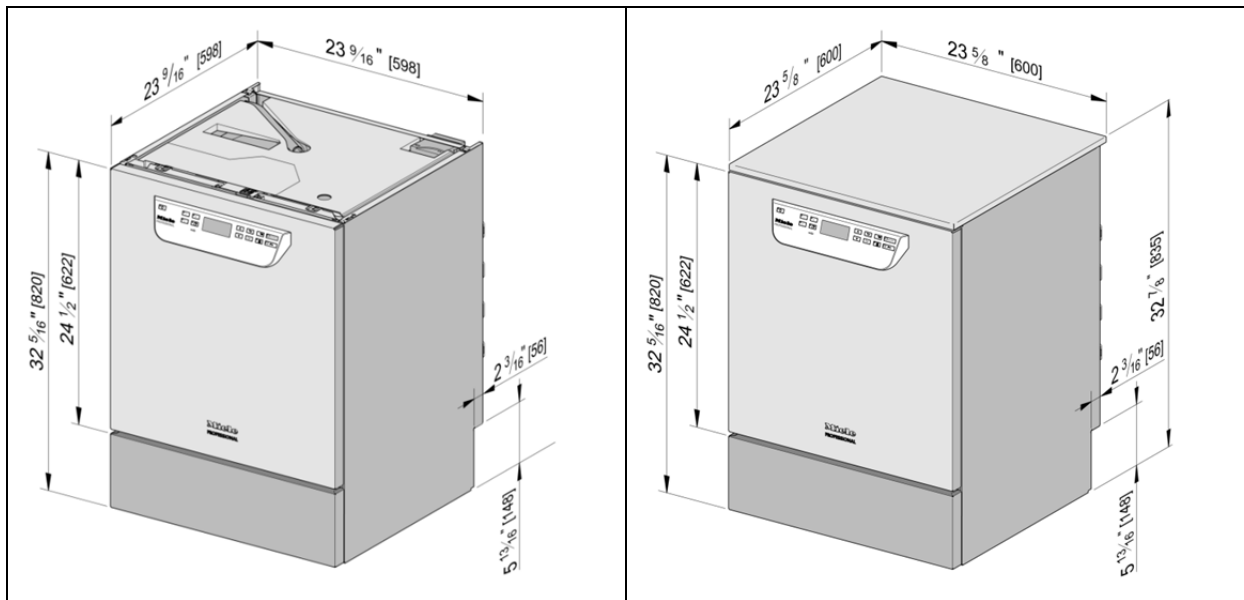
AW	Drain water	NW	Network and printer connection
DK	Cold water connection, steam condenser	PA	Equipotential bonding
DS	External dispensing system connection	VE	Demineralized water connection
EL	Electrical connection	WW	Hot water connection
KW	Cold water connection		

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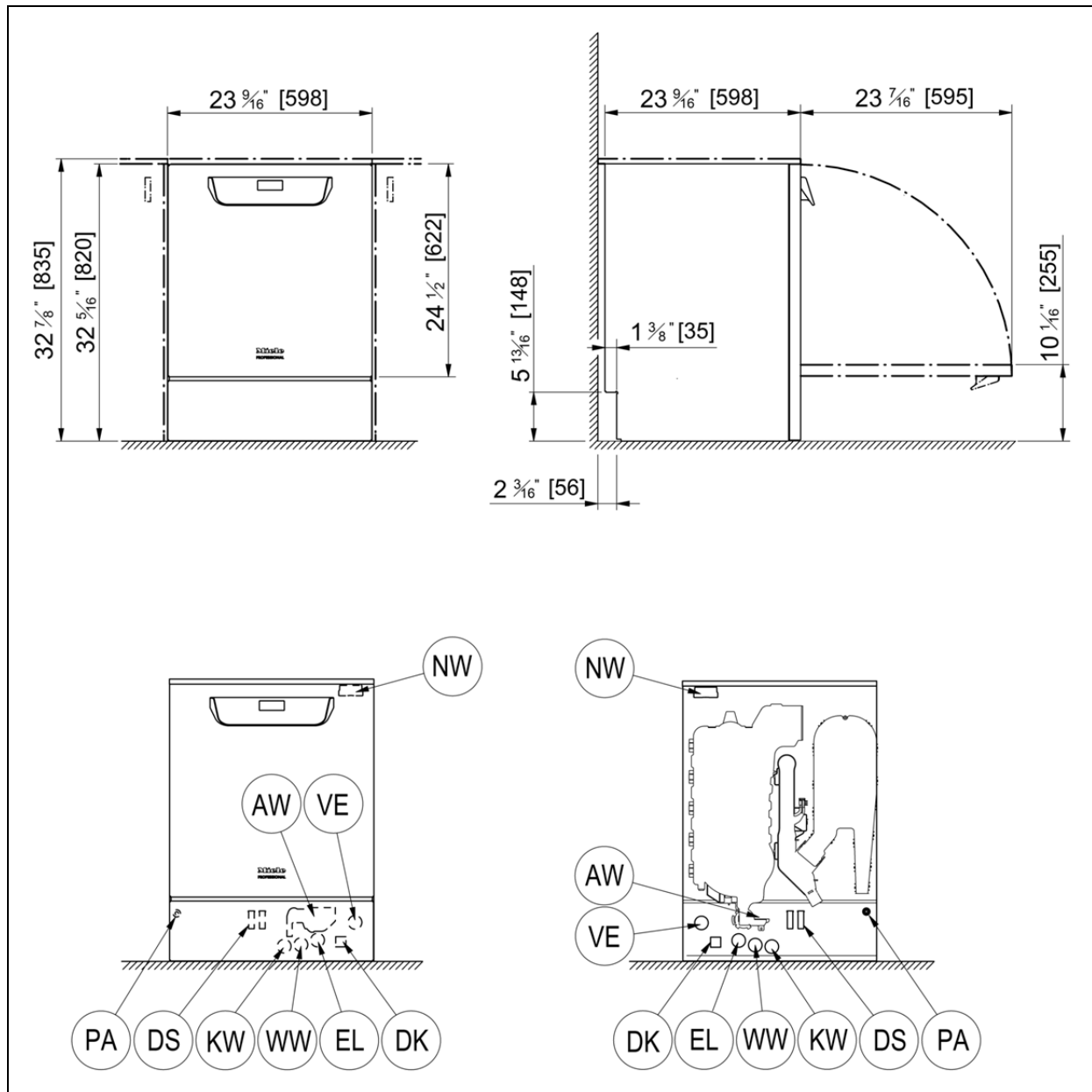AW	Drain water	NW	Network and printer connection
DK	Cold water connection, steam condenser	PA	Equipotential bonding
DS	External dosing system connection	VE	Demineralized water connection
EL	Electrical connection	WW	Hot water connection
KW	Cold water connection		

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