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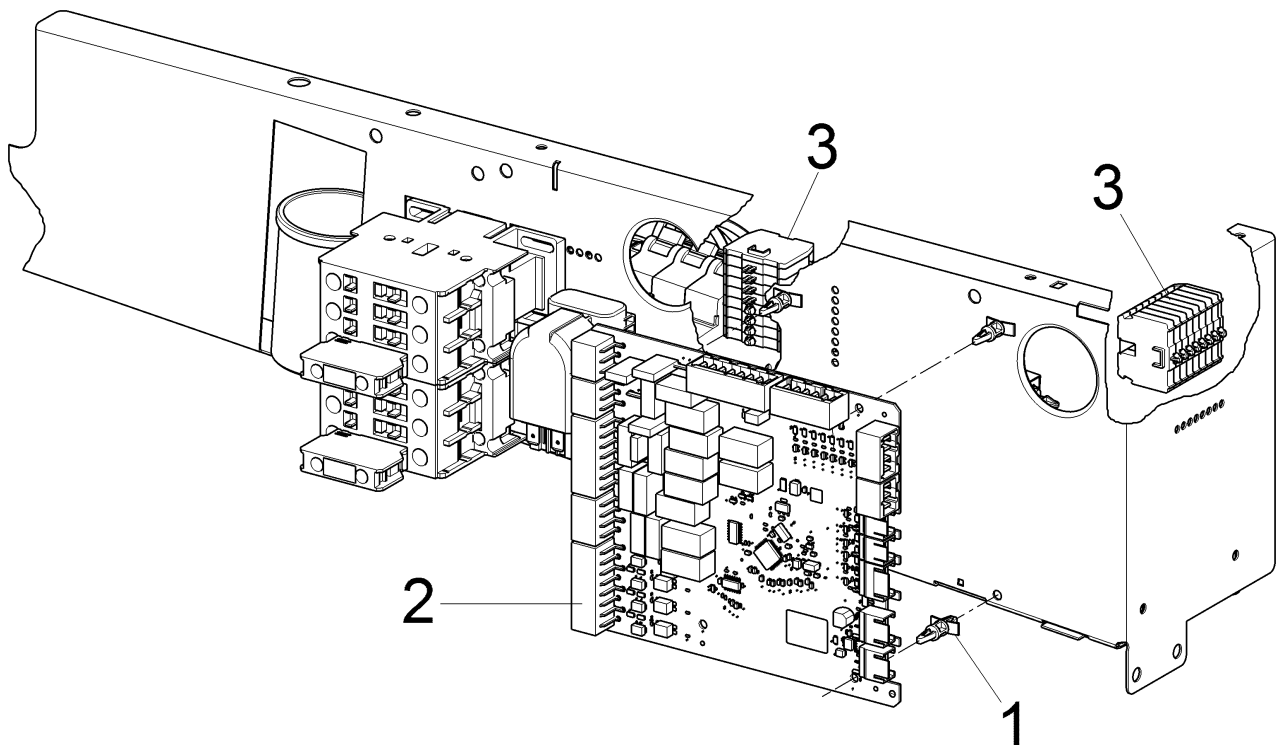
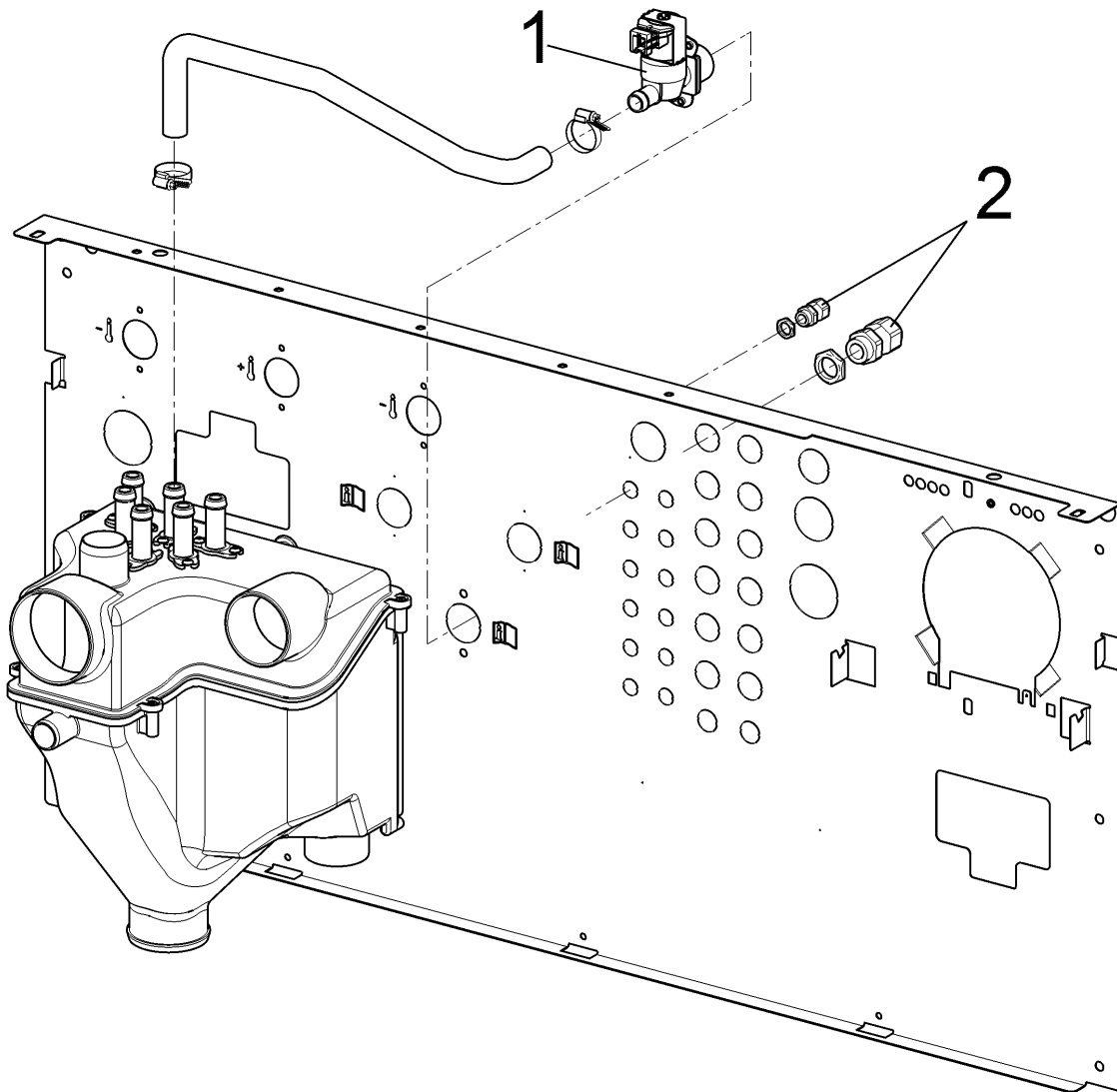
M.-Nr. 10717781

1 von 35


<b>de</b>	Montageanweisung Modul 5 für Flüssigdosierung 1-3 und Wasserrückgewinnung	<b>en</b>	Installation instructions - Module 5, liquid dispensing 1 - 3 and water recycling (WRG)	<b>cs</b>	Montážní návod modulu 5 pro dávkování tekutých prostředků 1-3 a rekuperaci vody
<b>da</b>	Monteringsanvisning modul 5 til dosering af flydende midler 1-3 og vandgenvinding	<b>el</b>	Οδηγίες τοποθέτησης μονάδας 5 για τη δοσομέτρηση υγρού απορρυπαντικού 1-3 και την ανακύκλωση νερού	<b>en-US</b>	Installation instructions - Module 5, liquid dispensing 1 - 3 and water recycling (WRG)
<b>es</b>	Instrucciones de montaje del módulo 5 para la dosificación líquida 1-3 y la recuperación de agua	<b>fi</b>	Asennusohje – Moduuli 5, nestemäisten aineiden annostus 1-3 ja vedenkierrätyslaite	<b>fr</b>	Notice de montage module 5 dosage liquide 1-3 et récupération d'eau
<b>it</b>	Indicazioni di montaggio modulo 5 per dosaggio liquido 1-3 e modulo recupero acqua	<b>nl</b>	Montage-instructie module 5 voor dosering 1-3 van vloeibaar reinigingsmiddel en waterterugwinning	<b>no</b>	Monteringsveiledning for Modul 5 til flytende dosering 1-3 og vanngjenvinningssystem
<b>pt</b>	Instruções de montagem do módulo 5 para dosagem de líquido 1-3 e recuperação de água	<b>ru</b>	Монтажная инструкция / модуль 5 для жидкого дозирования 1-3 и повторного использования воды	<b>sv</b>	Monteringsanvisning modul 5 för flytande dosering 1-3 och vattenåtervinning
<b>zh-CN</b>	安装说明——模块 5，液体药剂配给 1-3 及水循环 (WRG)				

1

2

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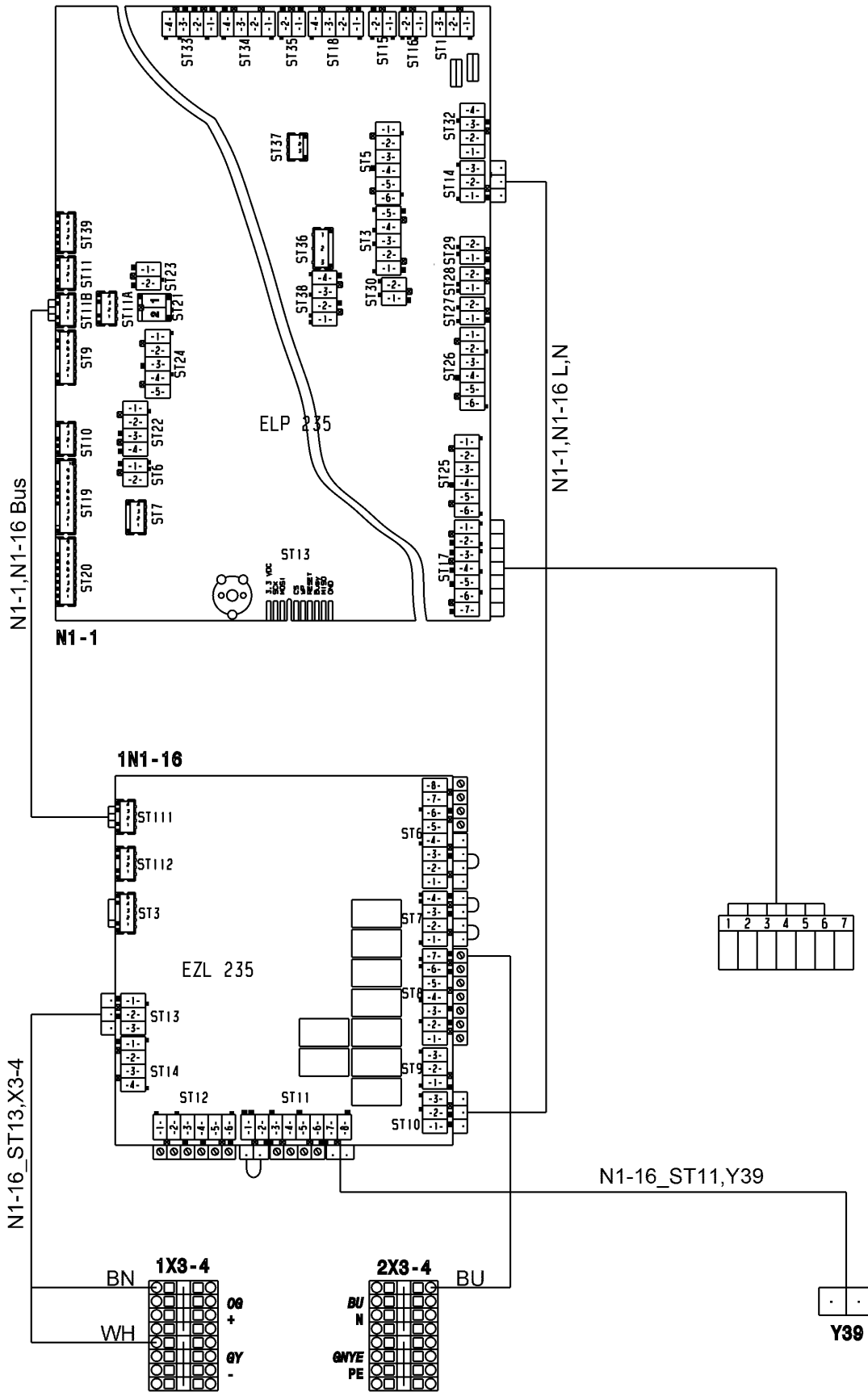**Reason:** the module is required for connecting dispenser pumps for liquid dispensing and water recycling to the washing machine.

#### Included parts

No.	Mat. no.	Designation
1		EZL 235 electronic unit, module 5
4		Electronic board holder
1		Coding plug MFM 1-2-5
1		Coding plug WRG 1
1		Coding plug MFM 53X
1		Plug N1Q_ST11A
1		Wiring harness N1-1, N1-16 L, N
1		Wiring harness N1-1_ST13, X3-4
1		Wiring harness N1-16_ST11, Y39
1		Wiring harness N1-1, N1-16 Bus
1		Connection wire, blue 0.75 x 280
2		Raised-head screws, M4 x 10
1		Intake valve 3/4"
1		Hose
2		Hose clips, 8 - 16 mm
1		Screw connector strip, 4-pole, neutral
1		Screw connector strip, 4-pole, red
1		Screw connector strip, 6-pole, blue
1		Screw connector strip, 7-pole, green
1		Terminal strip 1X3-4, orange / grey
1		Terminal strip 2X3-4, blue / green / yellow
1		Liquid dispensing accessory
1		Wiring diagram

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##### **WARNING**

Danger due to incorrectly carried out service and repair work

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✂ Service and repair work should only be carried out by a suitably qualified electrician with specialist training, knowledge and experience.

✂ All appropriate safety regulations must be taken into account.

✂ First read the Technical Service Documentation (TSD), then commence work.

#### Risk of injury due to sharp edges

##### **CAUTION**

There may be a risk of injury due to sharp edges.

Risk of injury due to sharp edges!

✂ Protective gloves should be worn and the edge protection, Mat. no. 05057680, should be used.

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- ✚ Remove the machine lid, see the model-specific TSD.
- ✚ Secure module (1) with holders (2) onto the cross support, see Fig. 1.
- ✚ Fit terminal strips (3) in the pre-punched holes of the cross support, see Fig. 1.
- ✚ Fix inlet valve (1) to the top of the back panel. To do this push the relevant embossed hole out of the top of the back panel, see Fig. 2.
- ✚ Fix cable screw connectors (2) to the top of the back panel. To do this push the relevant embossed holes out of the top of the back panel, see Fig. 2.
- ✚ Connect screw connector strip 4-pole, neutral to ST6 on the module, see Fig. 3.
- ✚ Connect screw connector strip 7-pole, green to ST8 on the module, see Fig. 3.
- ✚ Connect screw connector strip 4-pole, red to ST11 on the module, see Fig. 3.
- ✚ Connect screw connector strip 8-pole, blue to ST12 on the module, see Fig. 3.
- ✚ Connect coding plug WRG1 to ST6 pin 2-3 on the module, see Fig. 3.
- ✚ Connect coding plug MFM 53X to ST7 on the module, see Fig. 3.
- ✚ Connect plug N1Q\_ST11A to ST11 Pin 1-2 on the module, see Fig. 3.
- ✚ Connect coding plug 1-2-5 to ST3 on the module, see Fig. 3.

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- ✚ Remove any wiring harness connected to ST17 on the ELP 235 electronic unit, see Fig. 3.
- ✚ Lay wiring harness N1-16\_ST11, Y39 through the opening in the cross support beside the module, along the valve wiring harness and connect it, see Fig. 3.
- ✚
- ✚ Fasten wiring harness N1-16\_ST11, Y39 to the valve wiring harness with cable ties.
- ✚ Lay and connect wiring harness N1-16\_ST13, X3-4, see Fig. 3. To do this, feed the wiring harness through the opening in the cross support beside the module.
- ✚ Lay and connect wiring harness N1-1, N1-16 bus, see Fig. 3.
- ✚ Lay and connect wiring harness N1-1, N1-16 L, N, see Fig. 3.
- ✚ Lay and connect blue connection wire (bu), see Fig. 3.
- ✚ Connect the dispenser pumps in accordance with the wiring diagram.
- ✚ Fit the machine lid.
- ✚ Log the module on, see Service mode in the model-specific TSD.
- ✚ Calibrate the dispenser pumps, see Supervisor level in the Operating instructions for the machine.

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el

## en-US

### Installation instructions - Module 5, liquid dispensing 1 - 3 and water recycling (WRG)

**Reason:** The module is required for connecting dispenser pumps for liquid dispensing and water recycling to the washing machine.

#### Parts included

Quantity	Mat. no.	Designation
1		EZL 235 electronic, module 5
4		Standoff
1		MFM 1-2-5 coding plug
1		WRG 1 coding plug
1		MFM 53X coding plug
1		Plug N1Q_ST11A
1		"N1-1, N1-16 L, N" wiring harness
1		"N1-1_ST13, X3-4" wiring harness
1		"N1-16_ST11, Y39" wiring harness
1		"N1-1, N1-16 bus" wiring harness
1		Connection wire, blue 0.75 x 280
2		Raised-head screws, M4 x 10
1		Inlet valve, 3/4"
1		Hose
2		Hose clips, 8 - 16 mm
1		Beige 4-pin screw-terminal strip
1		Red 4-pin screw-terminal strip
1		Blue 6-pin screw-terminal strip
1		Green 7-pin screw-terminal strip
1		Orange/gray terminal strip 1X3-4
1		Blue/green/yellow terminal strip 2X3-4
1		Liquid-dispensing accessory
1		Wiring diagram

#### Danger due to incorrectly carried out service and repair work

##### WARNING

Danger due to incorrectly carried out service and repair work.

There is a risk of serious damage, injury and even death if service and repair work is carried out incorrectly.

- ✂ Service and repair work should only be carried out by a suitably qualified electrician with specialist training, knowledge and experience.
- ✂ All appropriate safety regulations must be taken into account.
- ✂ Read the technical service documentation (TSD) before starting service work.

#### Risk of cuts

##### CAUTION

Components are sharp.

Risk of cuts.

- ✂ Protective gloves should be worn and the edge protection, mat. no. 05057680, should be used.

## Installing module 5 for liquid dispensing 1-3 and water recycling

- ✚ Remove the machine lid; see the model-specific TSD.
- ✚ Secure the module (1) with standoffs (2) onto the crossbar; see Fig. 1.
- ✚ Install the terminal strips (3) in the pre-punched holes on the crossbar. See Fig. 1.
- ✚ Secure the inlet valve (1) to the top of the rear panel. To do this, push the appropriate pre-punched hole out of the top of the rear panel; see Fig. 2.
- ✚ Secure cable screw connectors (2) to the top of the rear panel. To do this, push the appropriate pre-punched holes out of the top of the rear panel; see Fig. 2.
- ✚ Connect the beige 4-pin screw-connector strip to plug ST6 on the module; see Fig. 3.
- ✚ Connect the green 7-pin screw-connector strip to plug ST8 on the module; see Fig. 3.
- ✚ Connect the red 4-pin screw-connector strip to plug ST11 on the module; see Fig. 3.
- ✚ Connect the blue 8-pin screw-connector strip to plug ST12 on the module; see Fig. 3.
- ✚ Connect the WRG1 coding plug to ST6 pin 2-3 on the module; see Fig. 3.
- ✚ Connect the MFM 53X coding plug to plug ST7 on the module; see Fig. 3.
- ✚ Connect the N1Q\_ST11A plug to ST11 pin 1-2 on the module; see Fig. 3.
- ✚ Connect the MFM 1-2-5 coding plug to plug ST3 on the module. See Fig. 3.

The power (L, N) and control (bus) wiring harnesses must be laid separately in order to avoid interference in the control circuit (electromagnetic compatibility).

Excess cable should be gathered so that current flows in opposite directions and tied to avoid any antenna effect.

- ✚ Remove any wiring harness connected to ST17 on the ELP 235 electronic; see Fig. 3.
- ✚ Lay the "N1-16\_ST11, Y39" wiring harness through the opening in the crossbar beside the module, along the valve wiring harness and connect it; see Fig. 3.
- ✚
- ✚ Fasten the "N1-16\_ST11, Y39" wiring harness to the valve wiring harness with cable ties.
- ✚ Lay and connect the "N1-16\_ST13, X3-4" wiring harness. See Fig. 3. To do this, feed the wiring harness through the opening in the crossbar beside the module.
- ✚ Lay and connect the "N1-1, N1-16 bus" wiring harness; see Fig. 3.
- ✚ Lay and connect the "N1-1, N1-16 L, N" wiring harness; see Fig. 3.
- ✚ Lay and connect the blue connection wire (bu); see Fig. 3.
- ✚ Connect the dispenser pumps in accordance with the wiring diagram.
- ✚ Reinstall the machine lid.
- ✚ Log the module on; see the "service mode" section in the model-specific TSD.
- ✚ Calibrate the dispenser pumps; see "Supervisor level" in the operating instructions for the machine.

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