A 303

de	Gebrauchsanweisung Injektormodul
da	Brugsanvisning Injektormodul
en	Operating instructions Injector module
es	Instrucciones de manejo Módulo inyector
fi	Käyttöohje Suorasuihkumoduuli
fr	Mode d'emploi Module à injection
hr	Upute za uporabu Injektorski modul
it	Istruzioni d'uso Modulo iniettore
nl	Gebruiksaanwijzing injectormodule
no	Bruksanvisning injektormodul
pt	Instruções de utilização Módulo injetor
sl	Navodila za uporabo Injektorski modul
sv	Bruksanvisning injektormodul

de	4
da	11
en	18
es	25
fi	32
fr	39
hr	46
it	53
nl	60
no	67
pt	74
sl	81
sv	88

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Notes about these instructions	18
Appropriate use	19
Queries and technical problems	19
Items supplied	20
Disposal of the packing material	20
Safety instructions and warnings	21
Areas of application.....	22
Injector modules.....	22
Inserting an injector module	22
Removing an injector module	22
To check before loading the machine and before starting a programme	23
Loading pipettes	23
Volumetric pipettes	23

Important warnings

⚠ Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property.
Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

- Select an option using the arrow buttons and save your choice with *OK*.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings | menu.

This module can be used to reprocess machine-reprocessible laboratory glassware and utensils in a Miele cleaning machine for laboratory glassware and utensils. Follow the operating instructions for your cleaning machine as well as the instructions of the glassware and utensil manufacturer on how to reprocess their items by machine.

The A 303 injector module is designed for reprocessing pipettes.

The module can hold up to 98 pipettes. The maximum length of the pipettes is 450 mm in the A 150 and 605 mm in the APLW 106. During reprocessing, the bodies of volumetric pipettes must not touch each other. This reduces the capacity of the module.

Use the Pipettes programme for reprocessing.

To ensure sufficient wash pressure, only one A 303 module may be inserted in the A 150 or APLW 106 lower basket. The module can only be used on the right-hand side. The second module connection can be used for another module.

Upper baskets cannot be used with this module.


This module is suitable for use in the following:

- Lower basket A 150
- APLW 106 basket

The cleaning machine is generally referred to as “the machine” in these operating instructions. Reprocessible laboratory glassware and utensils are referred to as “items” if they are not more closely defined.

Queries and technical problems

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your dishwasher or the washer-disinfector or go to www.miele-professional.com.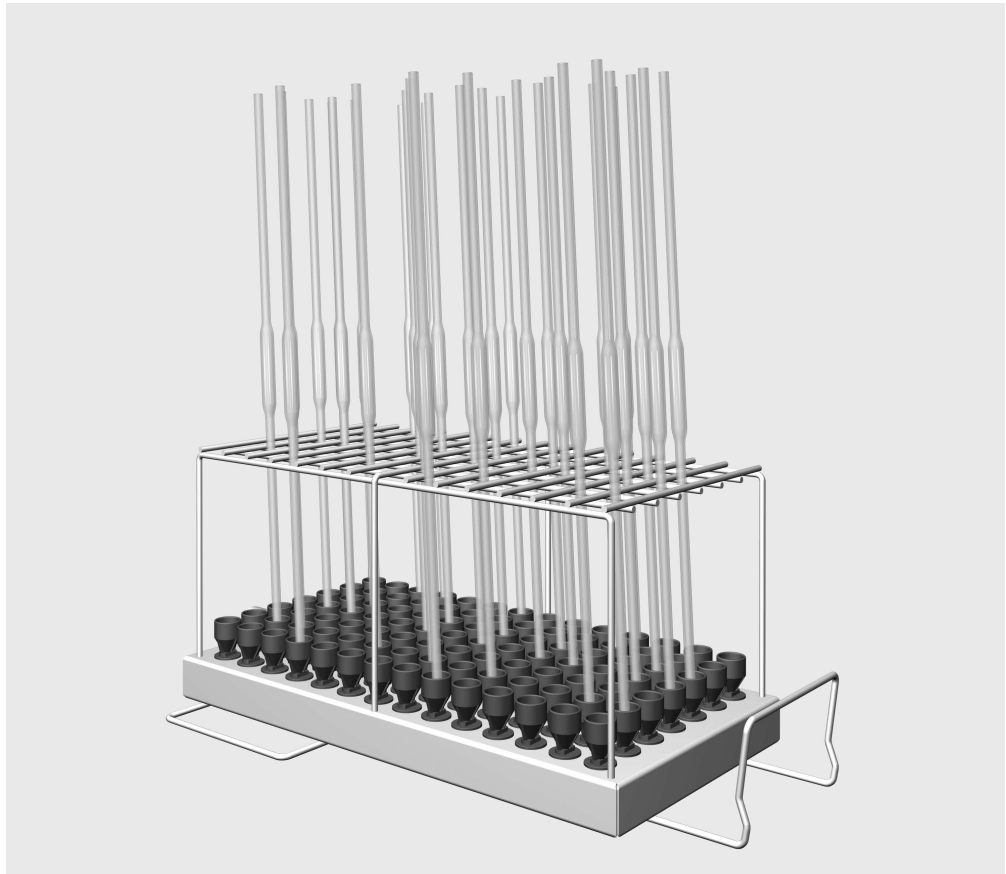

- Injector module A 303, height 185 mm, width 225 mm, depth 471 mm.

Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

To avoid the risk of accidents and damage to this module please read these instructions carefully before using it for the first time. Keep these instructions in a safe place where they are accessible to users at all times.

⚠ Please also read the operating instructions for your dishwasher or the machine and pay particular attention to the Warning and Safety instructions.

- ▶ The module is approved solely for the applications specified in the “Application technology” chapter of these operating instructions. Components such as nozzles may only be replaced with Miele accessories or genuine Miele original spare parts. Conversions, modifications and any other use are not permitted.
- ▶ Inspect all mobile units, baskets, modules and inserts daily as described in the “Maintenance” section in the Operating instructions for the washer-disinfector.
- ▶ Only items which have been declared by their manufacturer as suitable for machine reprocessing may be processed. The manufacturer's specific reprocessing instructions must be observed.
- ▶ Broken glass can result in serious injury when loading or unloading. Broken glass items must not be processed in the machine.
- ▶ Only place empty, unloaded modules in baskets. Check that they are correctly engaged before loading them. Modules must be completely emptied before taking them out. Placing loaded modules in baskets or removing them can cause damage to the wash load and cause injury should glassware be broken.
- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

Miele cannot be held liable for damage or injury caused by non-compliance with the warning and safety notes.

en - Areas of application


Before placing pipettes in the injector unit, please make sure that all hollow sections are free of blockages.

Injector modules

⚠ Only place empty, unloaded modules in baskets. Check that they are correctly engaged before loading them. Modules must be completely emptied before taking them out. Placing loaded modules in baskets or removing them can cause damage to the wash load and cause injury should glassware be broken.

Inserting an injector module

This injector module is only suitable for use in lower basket A 150. The injector nozzles must point upwards.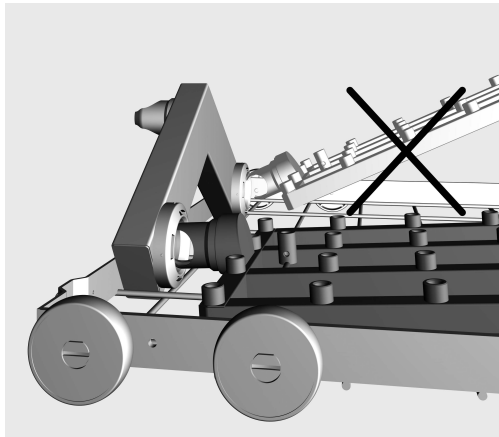

- Carefully slide the injector module in and connect it to the water connection point.
- Then press down on the module to secure it on the rails in the basket.

Removing an injector module

⚠ Do not hold the basket by the injector module when removing it from the cleaning machine. The module could come loose or the water connection could be damaged.


- Release the module from the locking mechanism and pull it out of the water connection at a flat angle. If the module is removed at too steep an angle, the lever action may damage the connection point and the connector.

To check before loading the machine and before starting a programme

- Check that the baskets and modules are correctly connected to the water supply.

Loading pipettes


The module can hold up to 98 pipettes. Please note the maximum length of 450 mm (in the A 150) or 605 mm (in the APLW 106). During reprocessing, the bodies of volumetric pipettes must not touch each other. This reduces the capacity of the module.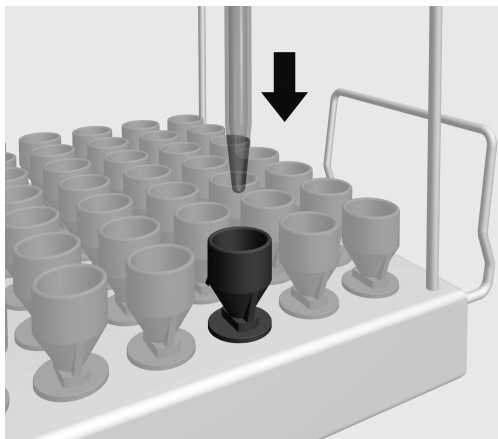

- Insert the pipettes through the frame and into the funnels with the pointed end facing downwards.

Volumetric pipettes

The bodies of volumetric pipettes must not touch each other.

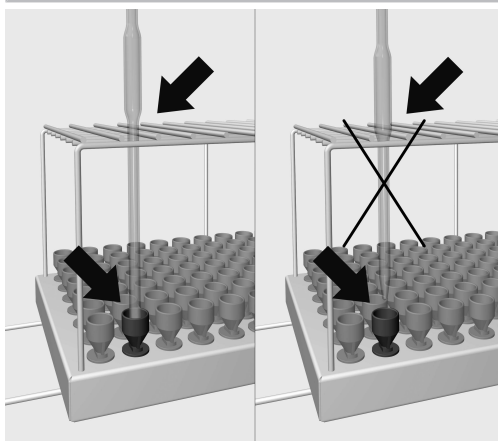- If the body of a pipette does not fit through the frame, the outflow tube will need to be long enough for the pipette to reach the funnel.

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