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

M.-Nr. 10950153

1 von 43


de	Ablaufpumpe (DP)	en	Drain pump (DP)	cs	Vypouštěcí čerpadlo (DP)
da	Afløbspumpe (DP)	el	Αντλία αποχέτευσης (DP)	en-US	Drain pump (DP)
es	Bomba de desagüe (DP)	fr	Pompe de vidange (DP)	it	Pompa di scarico (DP)
nl	Afvoerpomp (DP)	pl	Pompa odpływowa (DP)	pt	Bomba de esgoto (DP)
ru	Сливной насос (DP)	sl	Odtočna črpalka (DP)	sv	Avloppspump (DP)
zh-CN	排水泵 (DP)				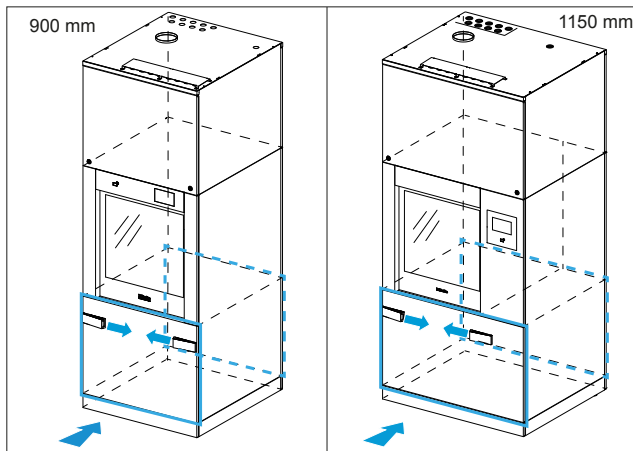


1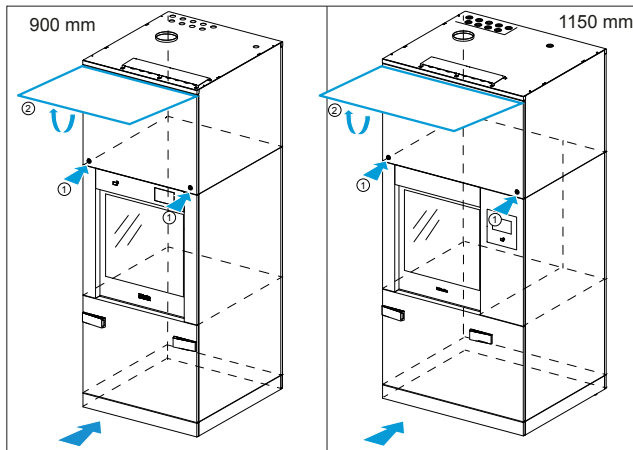


2

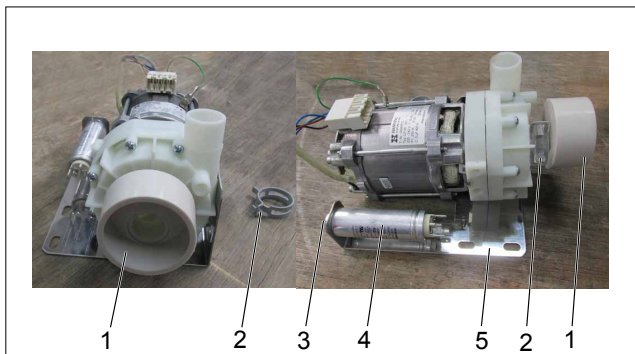
3

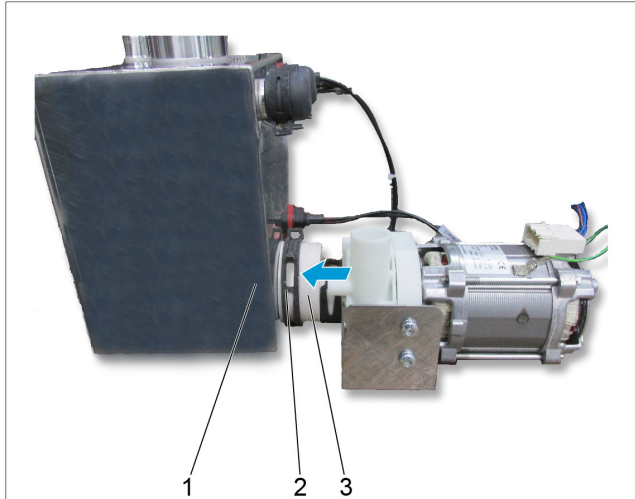
4

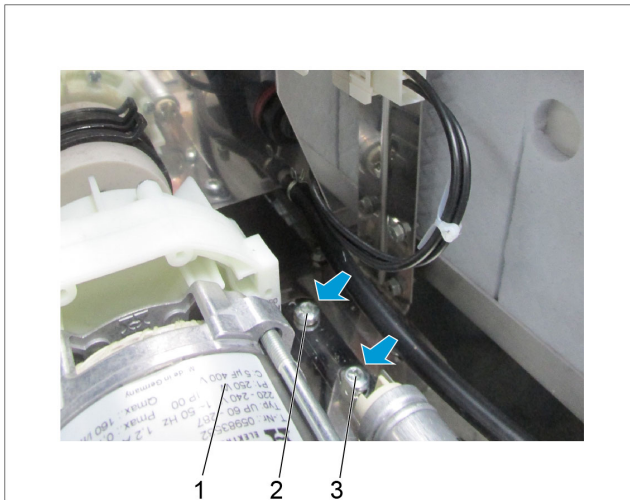
5

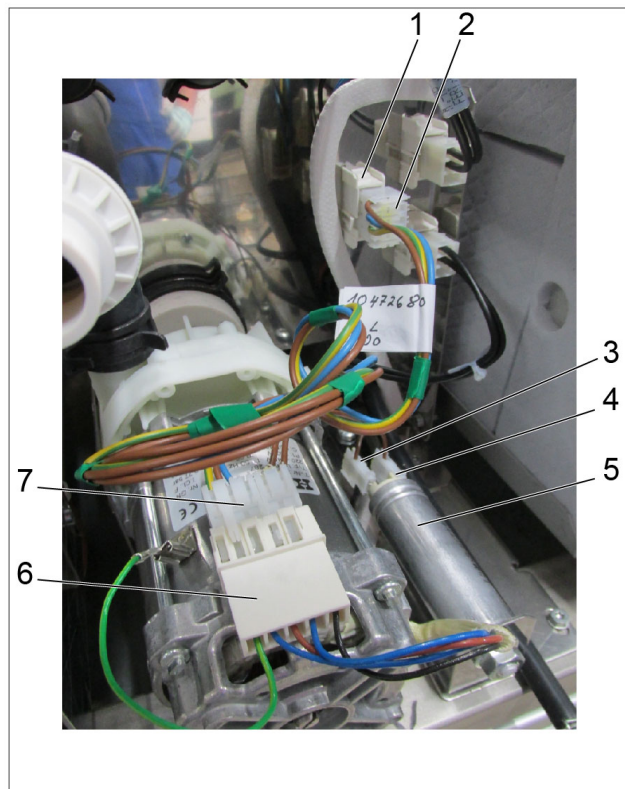
6

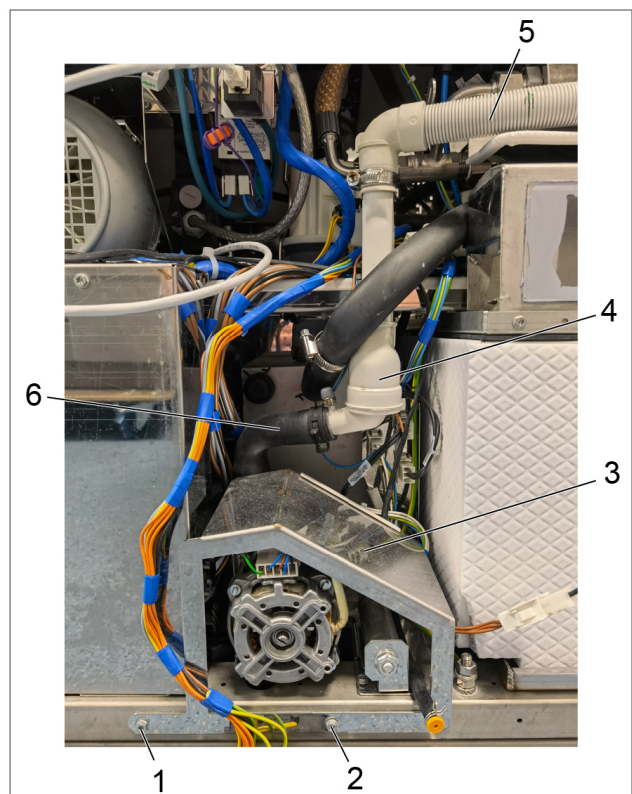
7

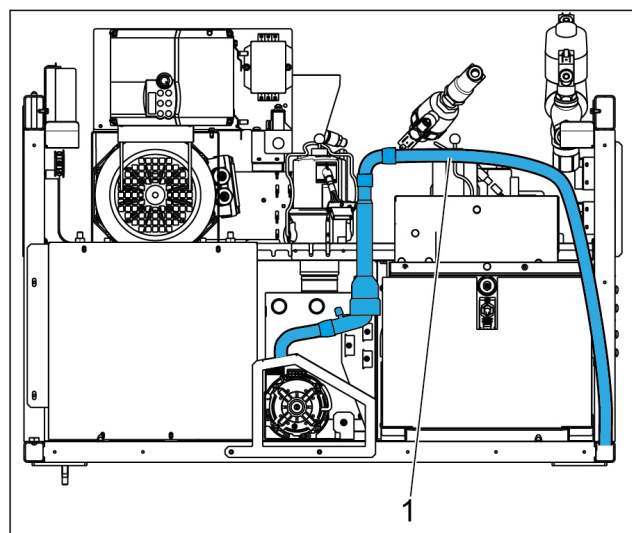8

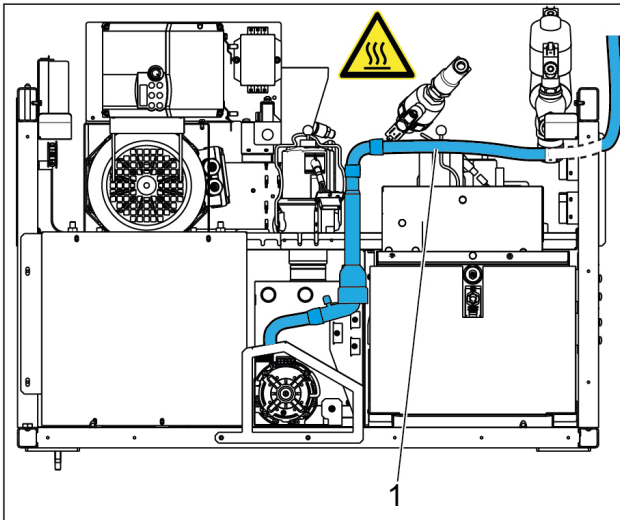9

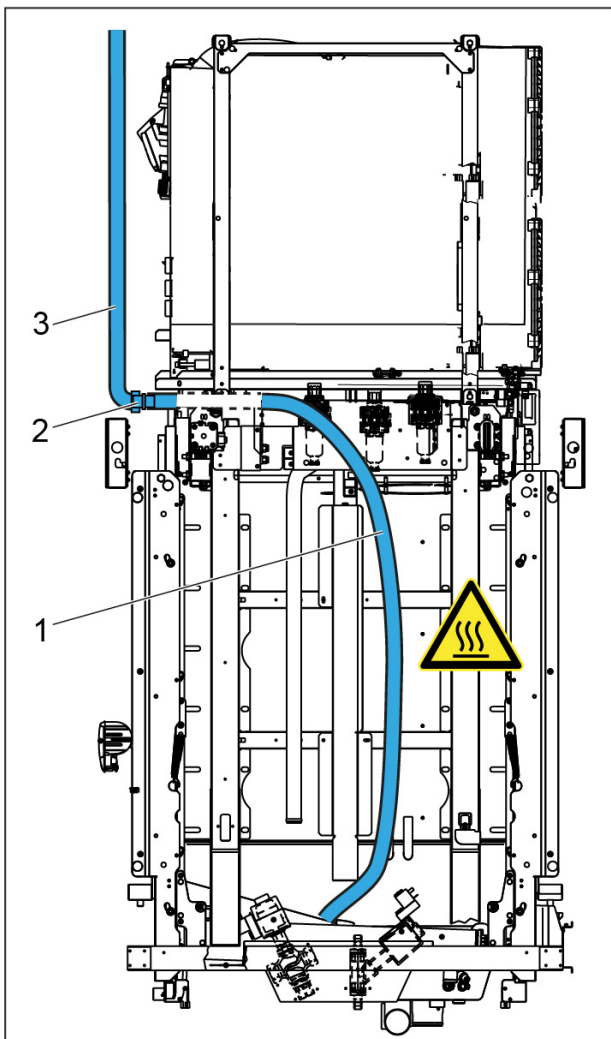10

de

Ablaufpumpe (DP)

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en

Drain pump (DP)

Reason: this conversion kit enables a drain pump (DP) to be installed in machine models PLW 8615, PLW 8616, PLW 8617, PWD 8626 or PWD 8628.

50 Hz drain pump:

Parts required

No.	Mat. no.	Designation
1	10931330	Drain pump (DP) PLW/PWD 86xx

Alternative 60 Hz drain pump:

Parts required

No.	Mat. no.	Designation
1	11516100	DP60 Hz drain pump PLW/PWD 86xx

The “Miele Service Documentation (MSD)” service application is required when carrying out maintenance work on the appliance

The safety instructions and warnings in the Miele Service Documentation (MSD) for the corresponding model must be observed.

Work on the appliance must be carried out in accordance with the work instructions.

Disconnect the machine from the mains and ensure power cannot be switched on again in error.

Wear protective gloves and goggles.

Do not lay cables over sharp edges or hot surfaces.

Lay the hoses without kinking or straining them.

Opening the service panel lower part

The service panel lower part is secured with two bars, which each clip into the frame at the sides of the machine.

✂ Slide the service panel retaining bolts towards the centre of the machine at the same time; see Fig. 1.

✂ Open the service panel slightly at the top and remove it.

Opening the top-box panelling service flap

✂ Open the 2 locks on the right and left of the top-box panelling service flap using the appropriate key (1), see Fig. 2.

✂ Fold up the top-box panelling service flap (2) and secure to prevent it from falling, see Fig. 2.

Fitting the drain pump

Preparing the drain pump for installation

✂ Fit the spring clip (2) on the small side of the adapter (1), see Fig. 3.

✂ Slide the adapter (1) onto the drain pump connector and secure it with the spring clip (2), see Fig. 3.

✂ Fit the interference suppression capacitor (4) with the M8 nut (3) to the holding bracket (5), see Fig. 3.

✂ Fit the drain pump on the holding bracket (5), see Fig. 3.

Fitting the drain pump in the machine

✂ Fit the large side of the adapter (3) onto the drain connection of the intermediate container (1) and secure it with a hose clip (2), see Fig. 4.

✂ Fit the drain pump (1) with the fixing screws (2) and (3), see Fig. 5.

Fitting the drain pump minor wiring harness

✂ Insert the connection plug (7) into the connection socket (6) of the drain pump, see Fig. 6.

✂ Connect the connection plugs (3) and (4) to the connections of the interference suppression capacitor (5), see Fig. 6.

✂ Connect the connection plug (2) to the connection socket (1) of the machine wiring harness, see Fig. 6.

✂ Fit edge protection (15 mm) to the holding plate for the plugs to protect the cables.

Connecting the waste water line to a drain pump

Installing the non-return valve

✂ Attach the hose elbow bend (6) to the drain pump connector and secure it with a hose clip, see Fig. 7.

✂ Push the short end of the non-return valve (4) into the hose elbow bend (6) and secure it with a hose clip, see Fig. 7.

✂ Slide the drain hose (5) onto the long end of the non-return valve (4) and secure it with a hose clip, see Fig. 7

Installing the splash protection plate

✂ Fit the splash protection plate (3) with the fixing screws (1) and (2) to the machine frame, see Fig. 7.

✂ Fit edge protection (30 mm) to the splash protection plate to protect the cables.

Routing the drain hose to a floor drain or side drain

If the drain hose is routed to a floor drain or side drain, care must be taken to ensure that waste water cannot flow back. To do this, the drain hose must be routed with a bend.

✂ Connect the drain hose to the non-return valve (1) of the drain pump and secure using a hose clip, see Fig. 8.

- ✂ Lay the drain hose (1) to the drain in such a way that there is a loop in the machine to prevent the waste water from flowing back, see Fig. 8.

Routing the drain hose upwards

If the drain hose is routed upwards, part of the drain hose is already laid in the machine ex works.

NOTE

Risk of damage due to hot pipes and sharp edges.

The drain hose can become damaged and leak.

- ✂ Do not route the drain hose over hot parts.
- ✂ Do not route the drain hose over sharp edges.
- ✂ Cover sharp edges with edge protection.

- ✂ Connect the pre-laid drain hose (1) to the non-return valve of the drain pump and secure using a hose clip, see Fig. 9.

The pre-laid drain hose runs to the side of the machine and then upwards. Below the top-box panelling, the drain hose is routed out of the machine towards the outfeed side. Here, the enclosed drain hose can be connected together with the hose connector and routed upwards out of the top-box panelling.

- ✂ Insert the hose connector (2) into the pre-laid drain hose (1) at the top and secure using a hose clip, see Fig. 10.
- ✂ Push the angled end of the enclosed drain hose (3) onto the hose connector (2) and secure using a hose clip, see Fig. 10.
- ✂ Route the drain hose (3) upwards out of the opening in the top-box panelling, see Fig. 10.
- ✂ When feeding the drain hose through one of the openings on the top-box panelling, remove the edge protection from the opening first.
- ✂ Pull the edge protection over the drain hose.
- ✂ Route the drain hose through the opening.
- ✂ Refit the edge protection in the opening.
- ✂ Connect the drain hose to the on-site drain and secure using a hose clip.

Drain pump activation

- ✂ Connect the service computer to the machine.
- ✂ Log in to the Web User Interface (WebUI) with a service technician with administrator rights.
- ✂ Request write rights in the WebUI.
- ✂ Select the System menu.
- ✂ Select the Machine configuration menu.
- ✂ Select the Drainage submenu.
- ✂ Select the Drain pump option parameter.

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en-US

Drain pump (DP)

Reason: This conversion kit enables a drain pump (DP) to be installed in models PLW 8615, PLW 8616, PLW 8617, PWD 8626, or PWD 8628.

50 Hz drain pump:

Parts required

Quantity	Mat. no.	Designation
1	10931330	Drain pump (DP) PLW/PWD 86xx

Alternative 60 Hz drain pump:

Parts required

Quantity	Mat. no.	Designation
1	11516100	60 Hz drain pump (DP) PLW/PWD 86xx

The "Miele Service Documentation" (MSD) service application is required for all service work

The warning and safety instructions given in the model-specific technical service documentation must be observed.

Service work must be done following the model-specific work instructions.

Disconnect the machine from power and ensure that power cannot be switched on again in error.

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Do not lay cables or wires over sharp edges or hot surfaces.

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Opening the bottom service panel

The bottom service panel is secured with 2 bolts which each clip into the frame at the sides of the machine.

- ✂ Slide the service panel retaining bolts towards the center of the machine at the same time; see Fig. 1.
- ✂ Open the service panel slightly at the top and remove it.

Opening the top service panel

- ✂ Open the 2 locks on the right and left of the top service panel using an appropriate key (1); see Fig. 2.
- ✂ Fold the top service panel up (2) and secure it to keep it from falling down; see Fig. 2.

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- ✂ Install the spring clip (2) on the small side of the adapter (1); see Fig. 3.
- ✂ Slide the adapter (1) onto the drain pump connection and secure it with the spring clip (2); see Fig. 3.
- ✂ Secure the interference suppression capacitor (4) to the mounting bracket (5) with the M8 nut (3); see Fig. 3.
- ✂ Install the drain pump on the mounting bracket (5); see Fig. 3.

Installing the drain pump in the machine

- ✂ Install the large side of the adapter (3) onto the drain connection of the intermediate container (1) and secure it with a hose clip (2); see Fig. 4.
- ✂ Install the drain pump (1) with the retaining screws (2) and (3); see Fig. 5.

Installing the drain pump minor wiring harness

- ✂ Insert the connection plug (7) into the connection socket (6) of the drain pump; see Fig. 6.
- ✂ Connect the connection plugs (3) and (4) to the connections of the interference suppression capacitor (5); see Fig. 6.
- ✂ Insert the connection plug (2) into the connection socket (1) of the machine wiring harness; see Fig. 6.
- ✂ To protect the wiring, install 15 mm edge protection on the plug's mounting bracket.

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Installing the non-return valve

- ✂ Attach the hose elbow bend (6) to the drain pump connector and secure it with a hose clip; see Fig. 7.
- ✂ Push the short end of the non-return valve (4) into the hose elbow bend (6) and secure it with a hose clip; see Fig. 7.
- ✂ Slide the drain hose (5) onto the long end of the non-return valve (4) and secure it with a hose clip; see Fig. 7

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- ✂ Install the splash protection plate (3) with the fixing screws (1) and (2) on the machine frame; see Fig. 7.
- ✂ Install edge protection (30 mm) on the splash protection plate to protect the wiring.

Installing the drain hose in a floor drain or side drain

If the drain hose is routed to a floor drain or side drain, care must be taken to ensure that waste water cannot flow back. The drain hose must have a bend in it.

- ✂ Connect the drain hose to the non-return valve (1) of the drain pump and secure with a hose clip; see Fig. 8.
- ✂ Lay the drain hose (1) to the drain in such a way that there is a loop in the machine to prevent the waste water from flowing back; see Fig. 8.

Feeding the drain hose upward to a drain

If the drain hose is routed upward, part of the drain hose is already laid in the machine ex works.

NOTE

Risk of burns and cuts from hot pipes and sharp edges.

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- ✂ Do not lay the drain hose over hot components.
- ✂ Do not lay the drain hose over sharp edges.
- ✂ Cover sharp edges with the edge protection.

- ✂ Connect the pre-existing drain hose (1) to the non-return valve of the drain pump and secure with a hose clip; see Fig. 9.

The pre-laid drain hose runs to the side of the machine and then upward. Below the top-box paneling, the drain hose is routed out of the machine toward the outfeed side. From here, the supplied drain hose can be connected together with the hose connector and fed up and out of the top-box paneling.

- ✂ Push the hose connector (2) into the top of the pre-existing drain hose (1) and secure with a hose clip; see Fig. 10.
- ✂ Push the angled edge of the supplied drain hose (3) onto the hose connector (2) and secure with a hose clip; see Fig. 10.
- ✂ Feed the drain hose (3) up and out of the opening on the top-box paneling; see Fig. 10.
- ✂ When feeding the drain hose through one of the openings on the top-box paneling, remove the edge protection from the opening first.
- ✂ Pull the edge protection over the drain hose.
- ✂ Feed the drain hose through the opening.
- ✂ Reinstall the edge protection in the opening.
- ✂ Connect the drain hose to the on-site drain and secure with a hose clip.

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