
de	Sockel fest (FP)	en	Fixed plinth (FP)	cs	Pevný sokl (FP)
el	Σταθερή Βάση (FP)	en-US	Permanent stand (FP)	es	Zócalo fijo (FP)
fr	Socle fixe (FP)	it	Zoccolo fisso (FP)	nl	Vaste sokkel (FP)
no	Fast sokkel (FP)	pt	Base fixa (FP)	ru	Неподвижное основание (FP)
sv	Fast sockel (FP)				

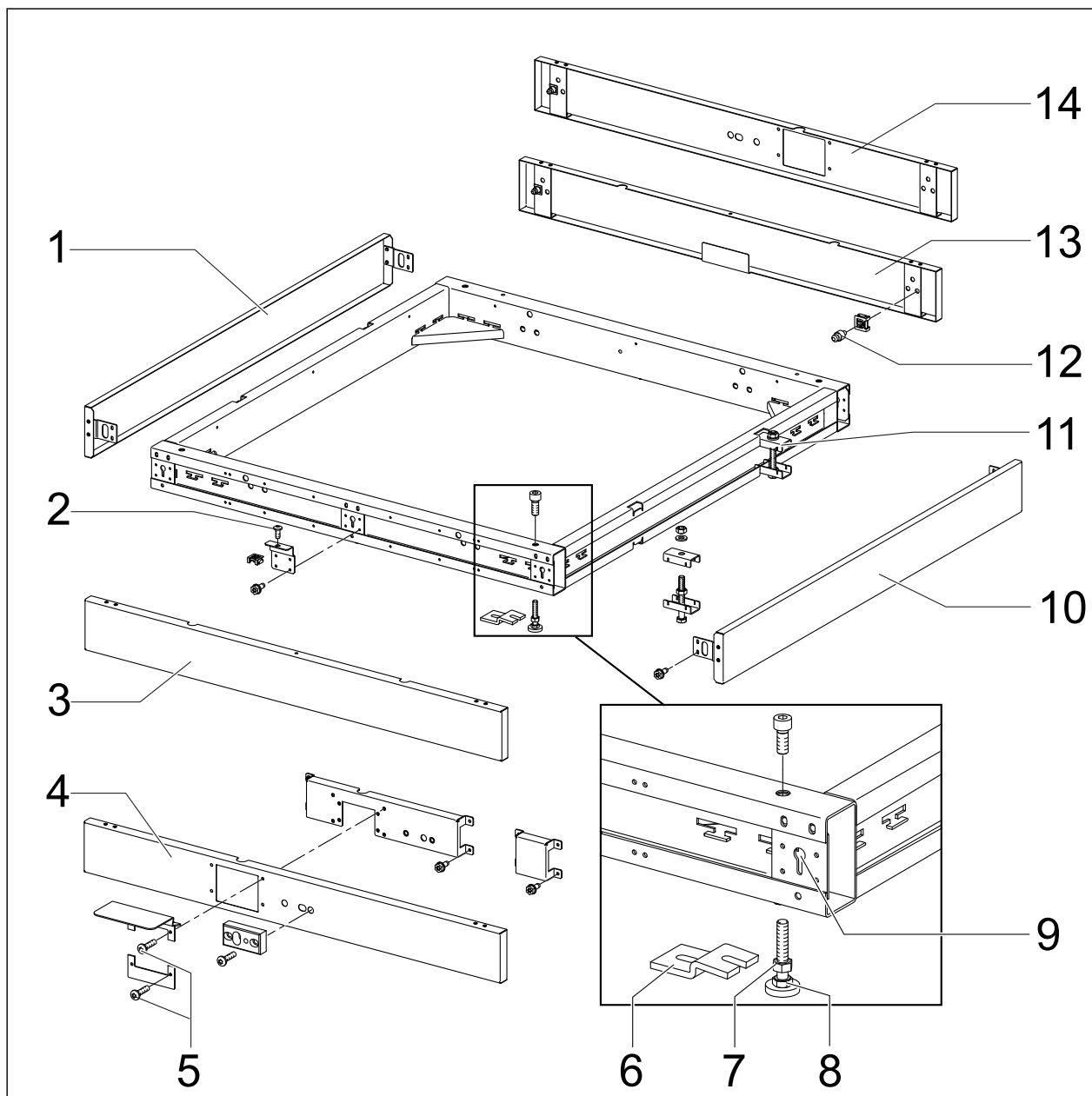## 1

### de

#### Sockel fest (FP)

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## en

### Fixed plinth (FP)

**Reason:** This conversion kit enables the fixed plinth (FP) option to be fitted to PLW 8615, PLW 8616, PLW 8617, PWD 8626 and PWD 8628 machines.

It can only be used with PLW 8615 or PLW 8617 models with just one door if the rear of the appliance can be accessed.

The plinth can be used for drain valve and drain pump drainage solutions.

The fixed plinth (FP) option can only be fitted to machines that can be accessed from both sides. Drain valve and drain pump drainage solutions can be used with the fixed plinth (FP) option.

Please refer to the installation plan for the required connections, drains and infeeds.

- FP900 = suitable for 900 mm wide machines: PLW 8615, PLW 8616, PWD 8626
- FP1150 = suitable for 1150 mm wide machines: PLW 8617, PWD 8628

### Parts required

No.	Mat. no.	Designation
1	10941870	FP900 – Fixed plinth (FP) 900 mm
1	10880750	FP1150 – Fixed plinth (FP) 1150 mm

**In order to carry out this work, the appropriate Technical Service Documentation (TSD) is required.**

The warning and safety instructions given in the applicable model-specific TSD must be complied with.

Work must be carried out in accordance with the instructions given in the applicable model-specific TSD.

### Fitting the fixed plinth (FP)

#### Removing the plinth panel

To level the plinth, remove the plinth panel on both the infeed side and the outfeed side.

- ✚ **Fixed plinth with cutout for attaching a transport conveyor:** Unscrew the 4 fixing screws (5) around the cutout for attaching a conveyor, see Fig. 1.
- ✚ **Fixed plinth without cutout for attaching a transport conveyor:** Unscrew the fixing screws (2) in the centre of the plinth panel from above, see Fig. 1.
- ✚ Lift the plinth panel (4), (5), (13) or (14) up and out of the keyhole brackets (9), see Fig. 1.

#### Levelling the plinth

- ✚ Fit the plinth and level it by adjusting the feet (8) up or down, see Fig. 1.
- ✚ Once the plinth is level, tighten the counter nuts (7) on the feet, see Fig. 1.
- ✚ If you are installing several machines side by side, make sure that the plinth drip trays are aligned absolutely level.
- ✚ Also use the clamp (11) to connect the individual plinths together, see Fig. 1.

#### Fitting optional floor mounts

The floor mounts (6) help to protect the machine in the event of an earthquake, for example, see Fig. 1.

- ✚ Mark out drill holes for the floor mount on the plinth drip tray.
- ✚ Drill holes for the plugs.
- ✚ Remove the drilling dust from the holes.
- ✚ Insert the plugs supplied with the mounts into the holes.

The special plugs supplied with the mounts will subsequently seal the drilled hole to prevent the ingress of moisture.

☞ Secure the floor mounts with the fixing screws.

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**el****en-US****Permanent stand (FP)**

**Reason:** This conversion kit is for installing the optional permanent stand (FP) on models PLW 8615, PLW 8616, PLW 8617, PWD 8626 and PWD 8628.

Use in single-door models PLW 8615 and PLW 8617 is only possible when the rear of the machine is accessible.

The stand can be used with the dump valve or drain pump.

The optional permanent stand (FP) can only be used in machines who have both sides accessible. The optional permanent stand (FP) can be used with the dump valve or drain pump.

See the installation plan for the necessary connections, outputs and inputs.

- FP900 = suitable for 900mm-wide machines: PLW 8615, PLW 8616, PWD 8626
- FP1150 = suitable for 1150mm-wide machines: PLW 8617, PWD 8628

**Parts required**

Quantity	Mat. no.	Designation
1	10941870	FP900 - 900mm permanent stand (FP = "fixed plinth")
1	10880750	FP1150 - 1150mm permanent stand (FP = "fixed plinth")

**The TSD will be needed for all service work**

The warning and safety instructions in the model-specific TSD must be observed.

Service work must be carried out following the instructions in the model-specific TSD.

**Permanent stand (FP) installation****Toekick removal**

In order to align the stand, the toekicks on both the loading and unloading sides must be removed.

- ↪ **For permanent stands with a conveyer-belt cutout:** Remove the 4 retaining screws (5) from the perimeter of the conveyer-belt cutout; see Fig. 1.
- ↪ **For permanent stands without a conveyer-belt cutout:** Remove the retaining screws (2) from the top center of the toekick; see Fig. 1.
- ↪ Lift the toekick ((4), (5), (13) or (14)) up and out of the keyhole retainers (9); see Fig. 1.

### Leveling the stand

- ↪ Line the stand up and level it horizontally by tightening or loosening the feet (8); see Fig. 1.
- ↪ After the stand has been horizontally leveled, tighten the counter nuts (7) on the feet; see Fig. 1.
- ↪ When leveling a row of machines, make sure that the drip pans are absolutely equally level.
- ↪ Additionally, connect the undersides of the stands together using the brackets (11); see Fig. 1.

### Installing the optional floor mounting kit

The floor fasteners (6) prevent the machine from falling over, e.g., during an earthquake; see Fig. 1.

- ↪ Draw the drill holes for the floor fasteners on the drip pan.
- ↪ Drill the holes for the screw anchors.
- ↪ Remove any dust from the holes.
- ↪ Insert the provided screw anchors in the holes.

The provided special screw anchors seal the holes against any incoming moisture.

- ↪ Secure the floor fasteners with the retaining screws.

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