DOS K 85

de	Montageanweisung DOS-Modul
en	Fitting instructions Dispensing module
es	Instrucciones de montaje Módulo DOS
fr	Notice de montage Module DOS
it	Istruzioni di montaggio Modulo DOS
ja	取付説明書 DOSモジュール
ru	Инструкция по монтажу DOS-модуль

de	4
en	14
es	24
fr	34
it	44
ja	54
ru	64

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Product description	14
Caring for the environment	17
Disposal of the packing material	17
Disposing of your old machine	17
Installation	18
Schematic diagram showing the dispensing module	18
Fitting the dispensing module	18
Wallmounting:	18
Surfacemounting:	18
Installation inside an adjacent cabinet:	19
Connection	20
Connecting the dispensing module to the cleaning machine.....	20
Flowmeter.....	20
Strain relief for dispensing hoses	20
Connecting the dispensing hoses.....	20
Colour coding on the dispensing hoses	21
Maintenance	21
Technical data	22

en - Product description

The dispensing module is an external dispensing system that is intended solely for connection to Miele cleaning machines. It dispenses liquid process chemicals (cleaning agent, disinfection solution, neutralising agent or rinsing agent, for example) automatically at the appropriate stage of a programme. The dispensing module does away with the need to fill the dispenser containers in the door of the machine (if there are any). Up to two external dispensing systems can be connected depending on the type of machine.

Suitable process chemicals are available from Miele Professional.

Dispensing modules can be retrofitted and connected if required.

This dispensing module is intended for stationary, indoor use only.

Risk posed by conversions or modifications

Please note that the K 85 dispensing module must only be used for applications described in these instructions.

Any other use, conversion or modification is not permitted and potentially dangerous.

Miele cannot be held liable for damage resulting from incorrect or improper use or operation.

Initial commissioning will be undertaken by the Miele Customer Service Department or authorised service technicians, who will provide you with operating instructions.

Please follow the operating instructions for your cleaning machine and pay particular attention to warning and safety instructions when using and handling chemical agents.


Damage caused by noncompliance with the fitting instructions.

Please read these fitting instructions carefully before fitting the dispensing module.

This prevents both personal injury and damage to the dispensing module.

Keep these fitting instructions in a safe place.


Areas of application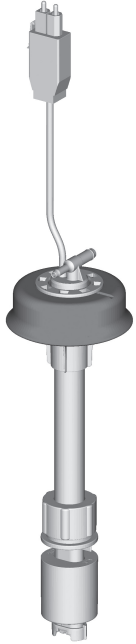

DOS module K 85 flex is suitable for use with the following cleaning machines:

- PG 8562
- PG 8581
- PG 8582
- PG 8583
- PG 8591
- PG 8592
- PG 8593

en - Product description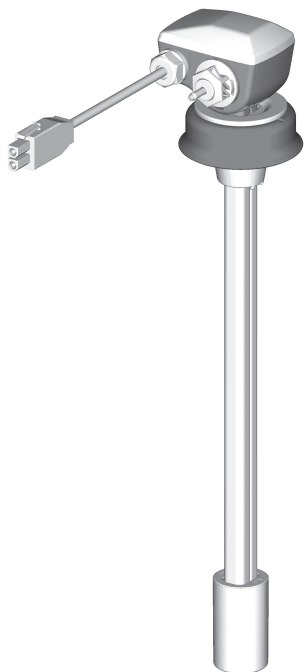

DOS K 85 comfort / DOS K 85/1 comfort


- PWD 8682
- PWD 8692
- PLW 8683
- PLW 8693
- PLW 8683 Safety
- PLW 8693 Safety

DOS K 85 flex suction lances can also be used for the PWD 86xx and PLW 86xx, but with a reduced functional scope which includes level detection only.

Symbols on the DOS module

Flow direction of process chemicals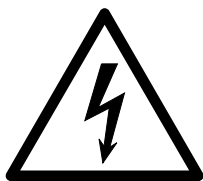

Warning:
Danger of electric shock!

Warning:
Read accompanying documentation!


Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

Disposing of your old machine


Electrical and electronic machines often contain valuable materials. They also contain materials which, if handled or disposed of incorrectly, could be potentially hazardous to human health and to the environment. They are, however, essential for the correct functioning of your machine. Please do not therefore dispose of it with your household waste.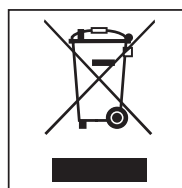

Please dispose of it at your local community waste collection / recycling centre or contact your dealer for advice.

Ensure that it presents no danger to children while being stored for disposal.

Schematic diagram showing the dispensing module

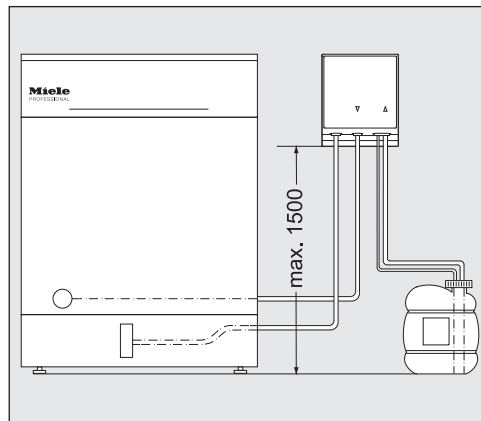
⚠ Dispensing problems caused by fitting the canister incorrectly. The container must not be placed on top of or above the cleaning machine. Make sure that the dispensing hose is not kinked or trapped.

Place the canister on the floor next to the cleaning machine or inside an adjacent cabinet.

Fitting the dispensing module

The dispensing module is supplied with two screws (4.2 x 25 mm) and two plugs (6 mm) for fitting.

- Select a suitably accessible position, e.g. on an adjacent wall or inside a cabinet, for fitting the dispensing module.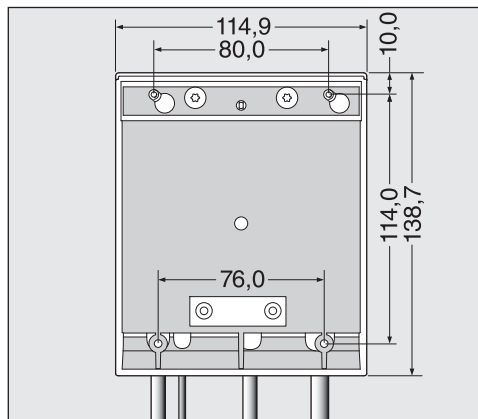

Wall mounting:

- Using the drawing as a guide, fit the screws the correct distance apart (80.0 mm) and suspend the dispensing module from these.

Surface mounting:

Two slots 76.0 mm apart are provided at the bottom of the dispensing module for fixing the module to a thin carrier or plate. Screws 3.9 mm in diameter (not supplied) are required to secure the module.

- Drill two holes through the plate and screw the dispensing module firmly in place from underneath.


**Installation inside
an adjacent cabinet:**

If more than one dispensing module is being used, or if the liquid agent canisters are to be stored with the module, a suitable cabinet is available from Miele.

Make sure hoses are laid without kinks.

Connecting the dispensing module to the cleaning machine


Flowmeter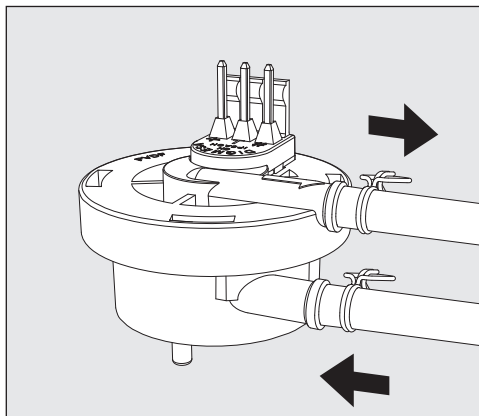

The K 85 dispensing module is supplied with a flowmeter which monitors the dispensing of liquid media.

Highly viscous chemical agents can cause problems with dispenser monitoring. In this case the dispenser monitoring system for laboratory applications should be switched off by the Miele Customer Service Department.

Strain relief for dispensing hoses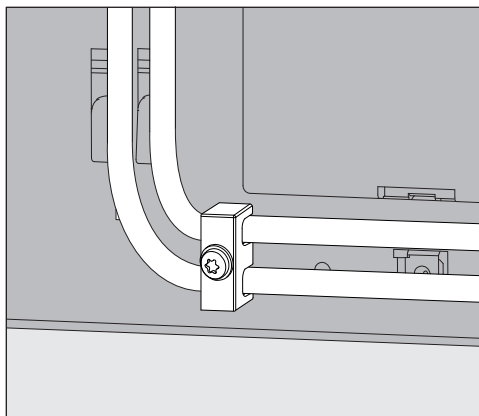


The strain relief provided with the dispensing modules for the dispensing hoses is fitted to the bottom right-hand side of the back of the machine to prevent the hoses from dislodging themselves from the dispenser connections during use.

Connecting the dispensing hoses

The dispensing module is tested at the factory for functionality. Any residues of the harmless test medium are a result of this testing and not evidence of previous use.

- Pull the safety cap off the end of the hose and remove the protective foil from the siphon.
- Connect the hose with the dispenser connection to the cleaning machine.
- Connect the module to the machine's power supply.

Connections for the dispensing hoses are located on the back of the cleaning machine. These are described in detail in the operating instructions for your cleaning machine.

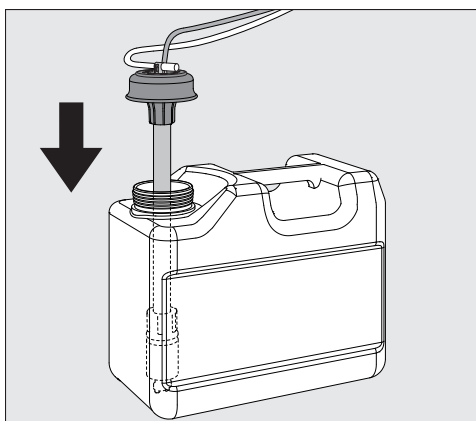- Fill or replace the container, fit the siphon and secure the lid.
- Feed the siphon into the canister until it reaches the bottom.

Follow the instructions for venting the dispensing system and setting the dosage in the operating instructions.

Colour coding on the dispensing hoses

To differentiate between the containers when using more than one DOS module, the hoses must be colour coded using the coloured labels supplied.

- Stick the label near the siphon and one of the same colour close to the free end of the hose to enable clear identification.

Miele use and recommend the following for DOS K 85:

- blue: for cleaning agent
- red: for neutralising agent
- green: for disinfecting agent / additional second cleaning agent
- white: for acidic chemical agent
- yellow: for special agents in laboratory applications

Maintenance

Maintenance of the dispensing system can be performed by Miele Customer Service at the same time as maintenance of the cleaning machine.

en - Technical data

Height	138.7 mm	
Width	114.9 mm	
Depth	114.2 mm	
Weight (net)	1.5 kg (incl. suction lance, hose and cable)	
Connection requirements	Power supply via cleaning machine	
Voltage, rated load, fuse rating via on-site connection	200–230 V 50 Hz 12 W 16 A	200–240 V 60 Hz 10 W 16 A
Flow rate (nominal rating)	55 ml/min	
Ambient temperature	5 °C to 40 °C	
Relative humidity Maximum Linear decreasing to	80 % for temperatures up to 31 °C 50 % for temperatures up to 40 °C	
Installation site height	1.500 m	
Degree of soiling	2	
Safety class (according to DIN EN 60529)	IP21	
Overvoltage category (according to IEC 60664)	II	
Voltage fluctuations	+/- 10 %	
Test certifications	VDE, EMC radio suppressed, CSA	
CE mark	2014/35/EC Low Voltage Directive	
Manufacturer's address	Miele & Cie. KG, Carl-Miele-Straße 29, 33332 Gütersloh, Germany	

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