# A 605

|           |   |
|-----------|---|
| <b>de</b> | Gebrauchsanweisung Injektormodul              |
| <b>cs</b> | Návod k obsluze Injektorový modul             |
| <b>da</b> | Brugsanvisning Injektormodul                  |
| <b>en</b> | Operating instructions Injector module        |
| <b>es</b> | Instrucciones de manejo Módulo inyector       |
| <b>fi</b> | Käyttöohje Suorasuihkumoduuli                 |
| <b>fr</b> | Mode d'emploi Module à injection              |
| <b>it</b> | Istruzioni d'uso Modulo a iniezione           |
| <b>ja</b> | 取扱説明書 インジェクターモジュール                            |
| <b>ko</b> | 사용 설명서 인젝터 모듈                                 |
| <b>nl</b> | Gebruiksaanwijzing injectormodule             |
| <b>no</b> | Bruksanvisning injektormodul                  |
| <b>pl</b> | Instrukcja użytkowania Moduł iniekcyjny       |
| <b>pt</b> | Instruções de utilização Módulo injetor       |
| <b>ru</b> | Инструкция по эксплуатации Инжекторный модуль |
| <b>sk</b> | Návod na použitie Injektorový modul           |
| <b>sv</b> | Bruksanvisning injektormodul                  |
| <b>tr</b> | Kullanım Kılavuzu Enjektör Modülü             |
| <b>uk</b> | Інструкція з експлуатації Інжекторний модуль  |
| <b>zh</b> | 操作说明书 注射器模块                                   |

**de** ..... 3  
**cs** ..... 9  
**da** ..... 15  
**en** ..... 21  
**es** ..... 27  
**fi** ..... 33  
**fr** ..... 39  
**it** ..... 45  
**ja** ..... 51  
**ko** ..... 57  
**nl** ..... 63  
**no** ..... 69  
**pl** ..... 75  
**pt** ..... 81  
**ru** ..... 87  
**sk** ..... 93  
**sv** ..... 99  
**tr** .....105  
**uk** .....111  
**zh** .....117

### Warnungen

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### Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

### Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

### Additional information and comments

Additional information and comments are contained in a simple frame.

### Operating steps

Operating steps are indicated by a black square bullet point.

#### Example:

- Select an option using the arrow buttons and save your choice with *OK*.

### Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

#### Example:

Settings  menu.

## en - Appropriate use

---

This module can be used to reprocess machine-reprocessible laboratory glassware and utensils in a Miele washer-disinfector for laboratory glassware and utensils. Follow the operating instructions for your washer-disinfector as well as the instructions of the glassware and utensil manufacturer on how to reprocess their items by machine.

The A 605 module is designed for reprocessing centrifuge tubes, phials, test tubes for fraction collectors and auto sampler tubes.


The module can be used in an A 503 mobile unit.

The washer-disinfector is generally referred to as “the machine” in these operating instructions. Reprocessible laboratory glassware and utensils are referred to as “items” if they are not more closely defined.

### **Queries and technical problems**

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your washer-disinfector or go to [www.miele-professional.com](http://www.miele-professional.com).

## Load carriers

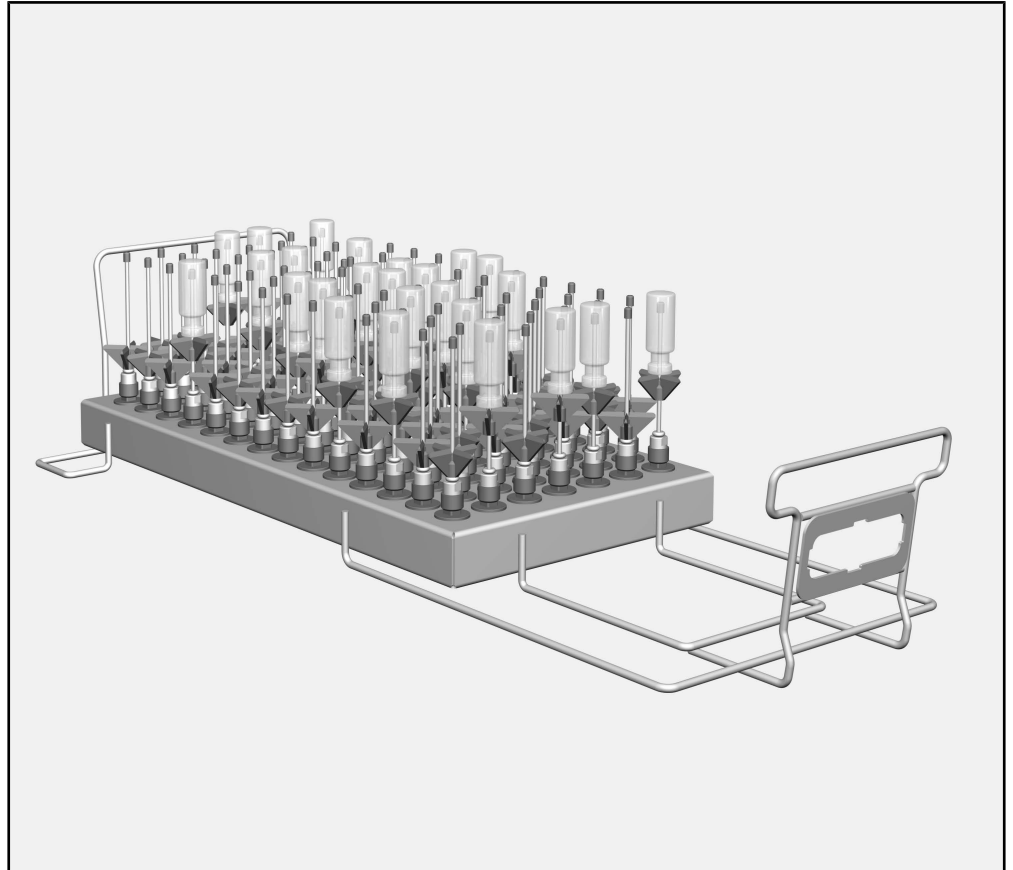- A 605 module, height 160 mm, width 255 mm, depth 611 mm

The module is suitable for items with a maximum diameter of 20 mm and a maximum height of 160 mm.

### Items supplied

- 98 x ID 110 injector nozzles with plastic supports. Length 110 mm,  $\varnothing$  2.5 mm

### Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in landfill sites.

## en - Warning and Safety instructions

---

Read the operating instructions carefully before using this load carrier. This will help protect users from personal injury, and help protect the load carrier from damage.  
Keep these operating instructions in a safe place.

 Please also read the operating instructions for your washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ The module is approved solely for the applications specified in the operating instructions. Components such as nozzles may only be replaced with Miele accessories or genuine Miele original spare parts. Alterations or conversion of the machine, or using it for purposes other than those for which it was intended, are not permitted and could be dangerous.
- ▶ New load carriers must be cleaned in the washer-disinfector without a load prior to first use.
- ▶ Inspect all load carriers as described in the “Maintenance” section in the operating instructions for the cleaning machine.
- ▶ Only items which have been declared by their manufacturer as suitable for machine reprocessing may be processed. The manufacturer's specific reprocessing instructions must be observed.
- ▶ Broken glass can result in serious injury when loading or unloading. Damaged glass items must not be reprocessed in the machine.
- ▶ Only ever place empty, unloaded modules in the mobile unit. Check that they are correctly engaged before loading them. Modules must be completely emptied before removing them. Placing loaded modules in mobile units or removing them can cause damage to the wash load and cause injury should glassware be broken.
- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

Miele cannot be held liable for damage caused by non-compliance with these Warning and Safety Instructions.

The operating instructions for the loading carriage describe how to insert and remove the modules.

### **To check before loading the machine and before starting a programme**

- Check that fittings such as irrigation sleeves and injector nozzles are securely screwed into position.


⚠ Make sure that all screw connectors are fitted with injector nozzles, adapters, irrigation sleeves or blind stoppers to ensure that all fittings in use are supplied with sufficient standardised pressure.

Damaged fittings such as injector nozzles, adapters and irrigation sleeves must not be used.


Fittings not equipped with wash items do not have to be replaced blind stoppers.

- Is the module correctly docked to the water supply in the load carrier?


### Preparing the load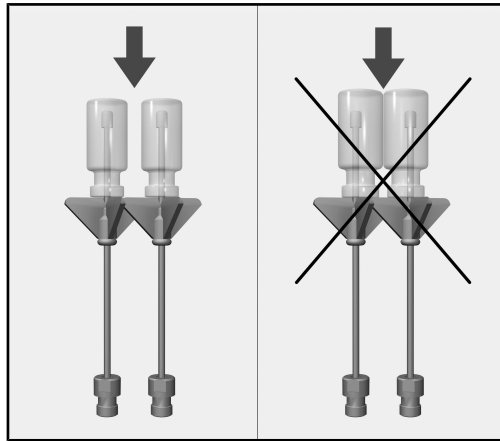

Individual items of glassware must not touch each other.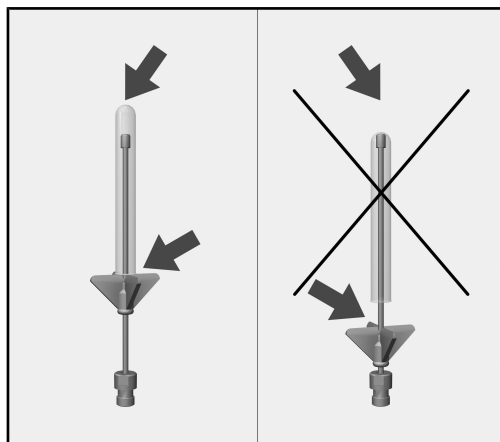

If the tip of an injector nozzle touches the bottom of a piece of glassware, the plastic support has to be pushed upwards.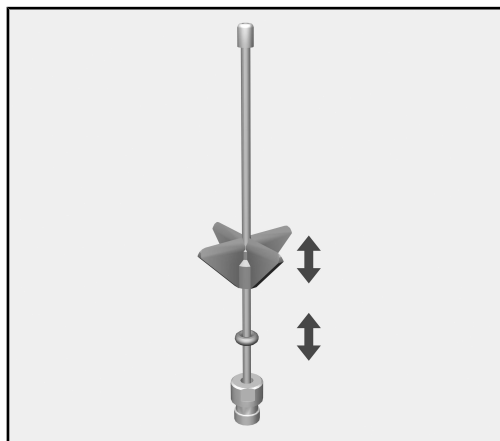

- Push the plastic support upwards to the height required. If necessary use a pair of pointed pliers to do so.
- Push the O-ring up to the plastic support to secure it in position.

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