PG 8595

de	Gebrauchsanweisung Aqua-Purificator
da	Brugsanvisning Demineraliseringsanlæg
en	Operating instructions Aqua Purificator
es	Instrucciones de manejo Aqua-Purificator
fi	Käyttöohje Aqua Purificator
fr	Mode d'emploi Déminéralisateur
hr	Upute za uporabu Uređaj za pročišćavanje vode Aqua-Purificator
it	Istruzioni d'uso Aqua-Purificator
nl	Gebruiksaanwijzing Aqua-Purificator
no	Bruksanvisning Aqua-purificator
pt	Instruções de utilização Aqua-Purificator
sl	Navodila za uporabo Aqua-Purificator
sv	Bruksanvisning Aqua Purificator

de 4
da 18
en 32
es 46
fi 60
fr 74
hr 88
it102
nl116
no130
pt144
sl158
sv172

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

Guide to the machine	32
Product description	33
Guide to the appliance	33
Partial demineralisation cartridges with flow meter.....	33
Full demineralisation cartridges with conductivity meter	33
Warning and Safety instructions	34
Disposal.....	35
Caring for the environment	36
Disposal of the packing material	36
Disposing of your old machine.....	36
FM flow meter and CM conductivity meter	37
FM Flow meter (partial demineralisation)	37
CM Conductivity meter (full demineralisation)	38
Connecting and installing cartridges	39
Connection options	39
Connect the cartridge using the screw connectors.	39
Connecting cartridges with quick release couplings	39
Connecting cartridges to the mains water supply and the washer-disinfector.....	40
Venting the cartridges	40
Inserting a cartridge	40
Removing a cartridge	41
Removing a cartridge	41
LP 2800 empty cartridge	42
Filling an empty LP 2800 cartridge with single use resin	42
Unclamping the lid	42
Filling with resin	42
Wall mounted FlowMeter / ConductivityMeter	44
Maintenance.....	44

en - Guide to the machine


Front view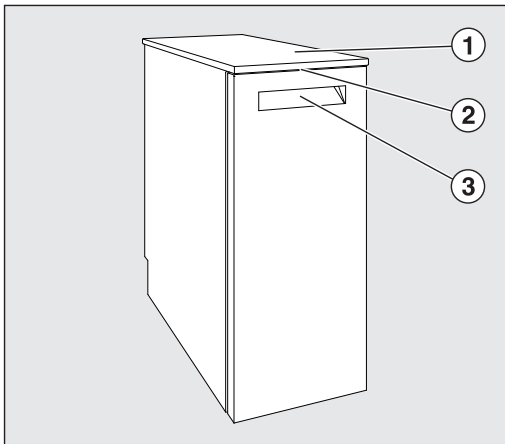

Rear view

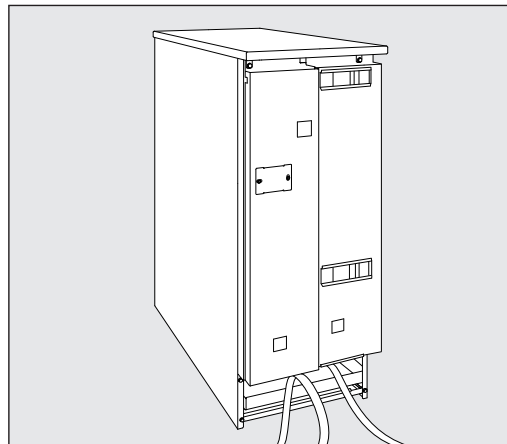- ① Cabinet lid
(remove for building under)
- ② LED for status monitoring when door is closed
- ③ Handle for opening the door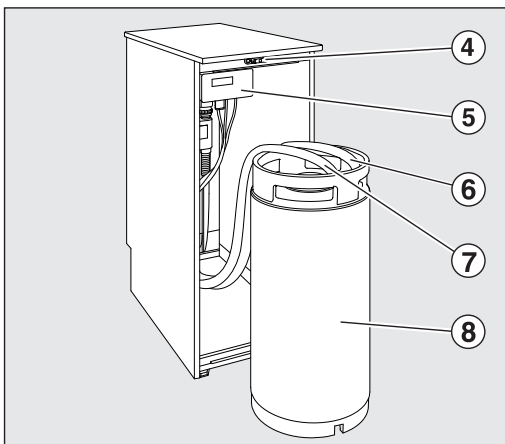

- ④ Door lock
- ⑤ Display unit FM (partially demineralised),
CM (fully demineralised)
- ⑥ “Outlet” connection
- ⑦ “Inlet” connection
- ⑧ Water demineralisation cartridge

Guide to the appliance

The PG Aqua Purificator is a water processing system and can be operated with two different water demineralisation cartridges.

Partial demineralisation cartridges with flow meter

With partial demineralisation, carbonate hardness is reduced by binding calcium and magnesium ions, thereby reducing the mineral content of the water.

The flow meter indicates when the cartridge requires changing.

- **TE P 2000 partial demineralisation cartridge**

Filled with 12.5 l resin.

- **TE P 2800 partial demineralisation cartridge**

Filled with 19 l resin.

Full demineralisation cartridges with conductivity meter

Full demineralisation removes all salts and minerals from water.

The conductivity meter shows the conductivity level of the demineralised water in $\mu\text{S}/\text{cm}$ (microsiemens per centimetre) to enable quality control of the demineralised water.

- **VE P 2000 full demineralisation cartridge**

Filled with 12.5 l resin.

- **VE P 2800 full demineralisation cartridge**

Filled with 19 l resin.

The cartridges are filled with renewable resin which can be regenerated after depletion. The whole cartridge can be sent to a regeneration station for reprocessing.

The PG 8595 Aqua Purificator has to be commissioned before first use and operators trained in its use by a Miele technician or authorised person only.

Please follow the operating instructions for the water demineralisation cartridges, conductivity meter, flow meter and the washer-disinfector.

en - Warning and Safety instructions

The PG 8595 Aqua Purificator complies with current safety requirements. Inappropriate use can, however, lead to personal injury and damage to property.

To avoid the risk of accidents and damage to the unit, please read these instructions carefully before using it for the first time. They contain important information on its installation, safe use and maintenance. Keep these instructions in a safe place and pass them on to any future owner.

Technical safety

- ▶ The Aqua Purificator must only be installed, commissioned, maintained and then repaired by the Miele Service Department or an authorised Miele Service dealer. Installation, repairs and other work by unqualified persons could be dangerous and Miele will not be held liable.
- ▶ The unit must only be operated by appropriately trained personnel. They must be instructed and trained regularly.
- ▶ Check the Aqua Purificator before use for any visible damage. Do not use a damaged unit. It could be dangerous.
- ▶ In the event of a damaged or leaking Aqua Purificator, switch it off immediately and call the Miele Service Department or an authorised Miele Service dealer.
- ▶ National plumbing regulations must be observed.

Correct use

- ▶ The water from demineralisation cartridges is not drinking water.
- ▶ Never clean the Aqua Purificator or near vicinity with a water hose or a pressure washer.
- ▶ Ensure that the outer casing of the machine does not come into contact with solutions or steam containing acid to avoid any corrosion damage.
- ▶ When replacing resin, use only tested and approved single use resin.
- ▶ This unit must only be connected to Miele-approved machines.

Safety with children

- ▶ Supervise children in the vicinity of the Aqua Purificator and do not let them play with it. There is a danger, among other things, of getting locked inside it.
- ▶ Children must not operate the Aqua Purificator.

Disposal

- ▶ For safety and environmental reasons, all residual resin and water must be disposed of in accordance with relevant safety regulations. Disable the door lock so that children cannot become trapped inside before disposing of the Aqua Purificator appropriately.
- ▶ Single use resin or resin which cannot be reactivated must be disposed of together with refuse or other similar industrial waste in accordance with EWC Waste disposal code 190905 for saturated or spent ion exchange resins.

Miele cannot be held liable for damage caused by non-compliance with these Warning and Safety instructions.


Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in landfill sites.

Disposing of your old machine

Electrical and electronic machines often contain valuable materials. They also contain materials which, if handled or disposed of incorrectly, could be potentially hazardous to human health and to the environment. They are, however, essential for the correct functioning of your machine. Please do not therefore dispose of it with your household waste.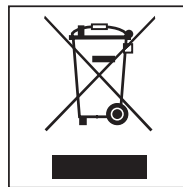

Please dispose of it at your local community waste collection / recycling centre or contact your dealer for advice.


Ensure that it presents no danger to children while being stored for disposal.

FM Flow meter (partial demineralisation)

A throughflow monitor ③ measures the amount of processed water and displays this in the FM flow meter ②.

If the set amount is exceeded, or if there is a fault, an optical LED lights up ① and an acoustic signal is given.

If the demineralisation cartridge is depleted it must be replaced with a new one and the spent resin sent away for reprocessing.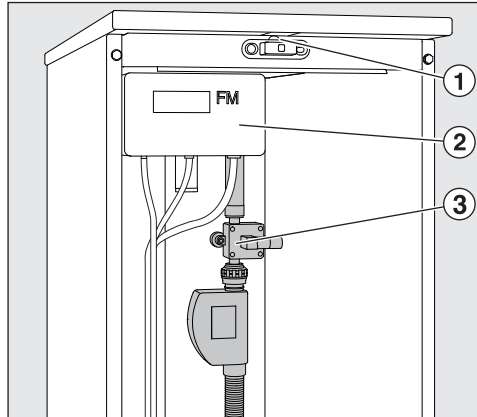

- **TE P 2000 and TE P 2800 cartridges:**

Send cartridges filled with spent resin away for reprocessing.

- **LP 2800 cartridge:**


Replace the resin. The E 316 Resin replacement set is supplied with its own instruction manual.

CM Conductivity meter (full demineralisation)

A conductivity module ③ measures the conductivity level of the demineralised water in a range from 0 - 199.9 $\mu\text{S}/\text{cm}$ (microsiemens per centimetre) and displays this in the conductivity meter ②. This enables quality control of the demineralised water.

If the defined limit value is reached, or if a fault occurs, an optical LED lights up ① and an acoustic signal is given.

If the demineralisation cartridge is depleted it must be replaced with a new one and the spent resin sent away for reprocessing.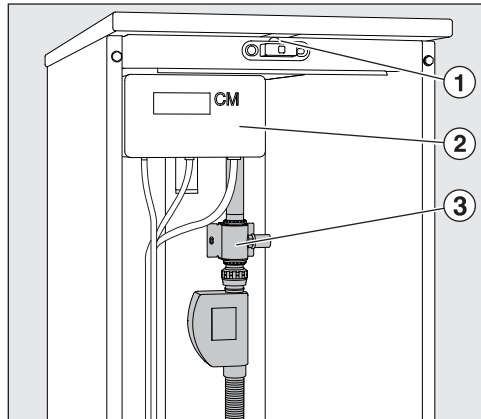

- **VE P 2000 and VE P 2800 cartridges:**

Send cartridges filled with spent resin away for reprocessing.


- **LP 2800 cartridge:**

Replace the resin. The E 316 Resin replacement set is supplied with its own instruction manual.

Connection options

- Open the door by pulling it upwards and disengage it from the hinges to remove it.
- Unscrew the safety caps from the cartridge connectors.

Connect the cartridge using the screw connectors.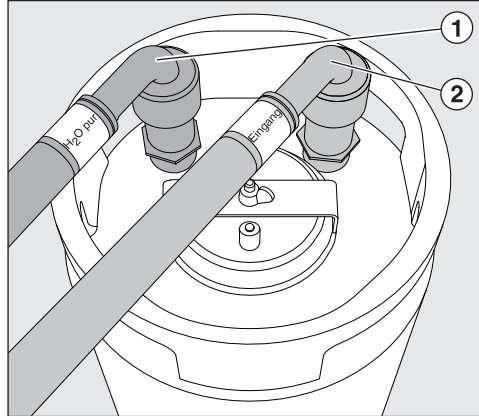

- Connect the hoses according to the colour coding on the cartridge connectors,
“Inlet (Eingang)” (red) to “In” ②,
“H2O pur” (green) to “Out” ①.

Connecting cartridges with quick release couplings

The cartridges and hoses can be fitted with quick release couplings (optional accessory) for rapid changing and reconnection.

When installing for the first time, screw the male part of the quick release couplings to the threaded cartridge connectors (in-out). Check for leaks before using for the first time.

- Connect the hoses with the relevant colour coding to the cartridge couplings and press down until they click firmly into place.
- To release a quick release coupling pull the ring on the end of the hose upwards and pull the hose off.

en - Connecting and installing cartridges

Connecting cartridges to the mains water supply and the washer-disinfector


- Guide the hoses through the openings provided on the back of the Aqua Purificator.
- Ensure that the hoses are not kinked when doing this.
- Connect the “H2O pur” (green) hose to the solenoid valve for AD water or to the AD water inlet hose for the washer-disinfector.
- Connect the “Inlet (Eingang)” (red) hose to the on-site cold water mains supply stopcock.

Important

Water pressure (flow pressure) must be between 250 - 1000 kPa pressure (2.5 - 10 bar).

The water temperature must not exceed 20 °C.

Venting the cartridges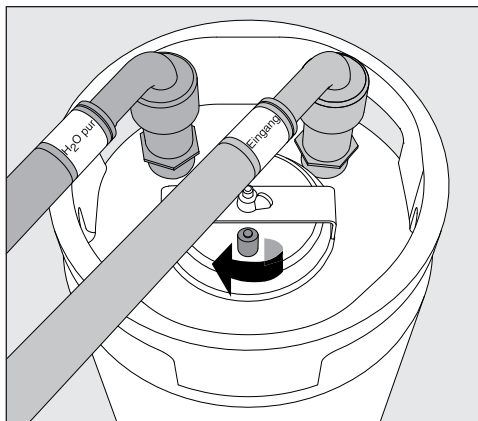

- Open the valve by turning it to the left.
- Open the cold water stopcock until water escapes from the valve. until water escapes from the vent valve.
- Then close the valve.

The demineralisation cartridge is now ready for use.

Inserting a cartridge

- Place the connected cartridges into the back of the Aqua Purificator. The front area can be used to accommodate a spare cartridge.

Removing a cartridge

- Close the cold water stopcock.
- Open the stopcock for drawing manual AD water (if available), or:
- Open the valve on the cartridge to release any pressure.
While doing this, water can escape which should be wiped up with a cloth.
- Unscrew the hoses.
- Upend the cartridge to drain the residual water into a suitable outlet.

en - LP 2800 empty cartridge

The LP 2800 cartridge is supplied unfilled and must be filled on site with 19 l of single use resin prior to use (see “Filling an empty LP 2800 cartridge with single use resin”).

When depleted replace the spent resin with new single use resin.

Single use resin or resin which cannot be reactivated must be disposed of together with refuse or other similar industrial waste in accordance with EWC Waste disposal code 190905 for saturated or spent ion exchange resins.


For easy filling and refilling of an empty LP 2800 cartridge, we recommend the following as basic equipment in addition to the cartridge:

- The original Miele E 316 refill set for single use resin, consisting of a 30 l plastic barrel with special lid and funnel.
- 2 x 10 l bags of single use resin for demineralisation with filter bag for depleted resin, E 315.


Filling an empty LP 2800 cartridge with single use resin

The empty LP 2800 cartridge can be either filled with resin for partial demineralisation or for full demineralisation (E 315).

Unclamping the lid

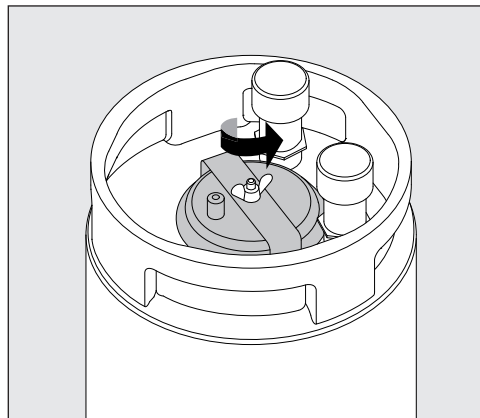
- Loosen the wing nut (hold the lid while doing this so that it does not fall into the cartridge).
- Remove the clamp.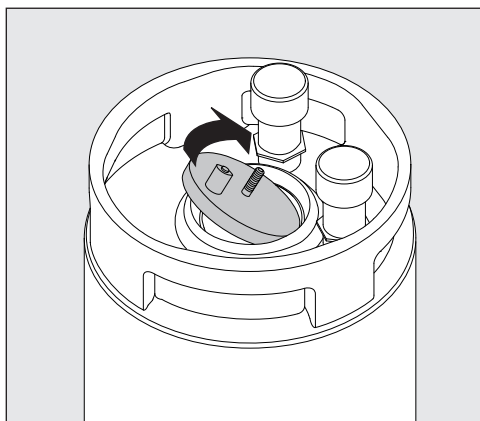

- Open the lid.

Filling with resin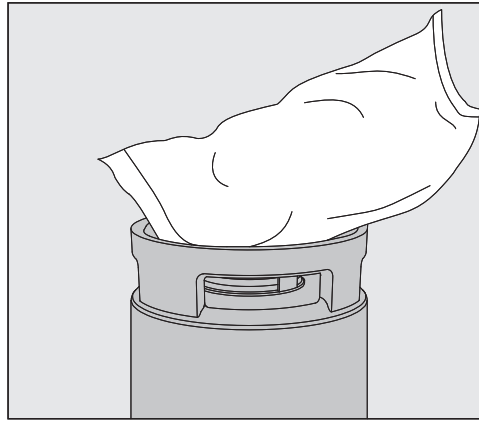


- Place the funnel in the cartridge.
- Cut the corners off both bags of single use resin and empty the contents into the cartridge.
- Remove any residual resin from the mouth of the cartridge.
- Replace the lid, hold it tightly by the valve and secure it with the clamp and wing nut.

en - Wall mounted FlowMeter / ConductivityMeter

The water quality display devices can be mounted outside the Aqua Purificator on the wall.

A suitable wall bracket is supplied with the display devices.

For alternative installation the sequence of individual components must be observed:

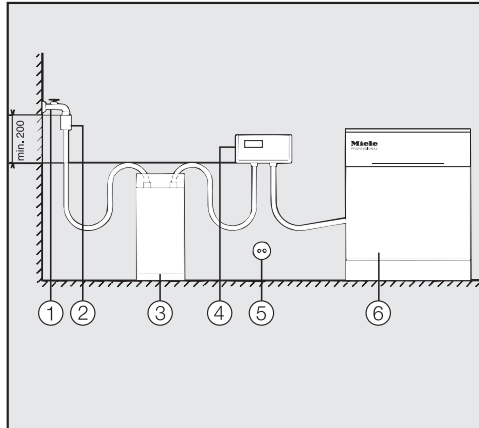The display device is always connected behind the water demineralisation cartridge, i.e. in the following order:

water stopcock ① - non-return valve with pipe aerator according to DIN EN 1717 ② - water demineralisation cartridge ③ - display device FM/CM ④ - washer-disinfector .

Electrical connection ⑤ is made via a switched plug and socket.

Maintenance

The PG 8595 Aqua Purificator should be serviced at the same time as the washer-disinfector by Miele Service.

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