A 613

de	Gebrauchsanweisung Injektormodul
cs	Návod k obsluze Injektorový modul
da	Brugsanvisning Injektormodul
el	Οδηγίες χρήσης μονάδα με εγχυτήρες
en	Operating instructions Injector module
es	Instrucciones de manejo Módulo inyector
fi	Käyttöohje Suorasuihkumoduuli
fr	Mode d'emploi Module à injection
it	Istruzioni d'uso Modulo iniettore
ja	取扱説明書 インジェクターモジュール
ko	사용 설명서 인젝터 모듈
nl	Gebruiksaanwijzing injectormodule
no	Bruksanvisning injektormodul
pl	Instrukcja użytkowania Moduł iniekcyjny
pt	Instruções de utilização Módulo injetor
ru	Инструкция по эксплуатации Инжекторный модуль
sk	Návod na použitie Injektorový modul
sv	Bruksanvisning injektormodul
tr	Kullanım kılavuzu Enjektör modülü
uk	Інструкція з експлуатації Інжекторний модуль
zh	操作说明书 注射器模块

de	4
cs	10
da	16
el	22
en	28
es	34
fi	40
fr	46
it	52
ja	58
ko	64
nl	70
no	76
pl	82
pt	88
ru	94
sk	100
sv	106
tr	112
uk	118
zh	124

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Notes about these instructions	28
Appropriate use	29
Queries and technical problems	29
Items supplied	30
Load carriers.....	30
Disposal of the packing material	30
Warnings and safety notes	31
Areas of application.....	32
To check before loading the machine and before starting a programme	32
Loading pipettes	32
Volumetric pipettes	32

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

- Select an option using the arrow buttons and save your choice with *OK*.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings  menu.

This module can be used to reprocess machine-reprocessable laboratory glassware and utensils in a Miele washer-disinfector for laboratory glassware and utensils. Follow the operating instructions for your washer-disinfector as well as the instructions of the glassware and utensil manufacturer on how to reprocess their items by machine.

The A 613 module is designed for reprocessing volumetric pipettes and measuring pipettes.


The module can be used in an A 503 mobile unit.

The washer-disinfector is generally referred to as “the machine” in these operating instructions. Reprocessable laboratory glassware and utensils are referred to as “items” if they are not more closely defined.

Queries and technical problems

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your dishwasher or the washer-disinfector or go to www.miele-professional.com.

Load carriers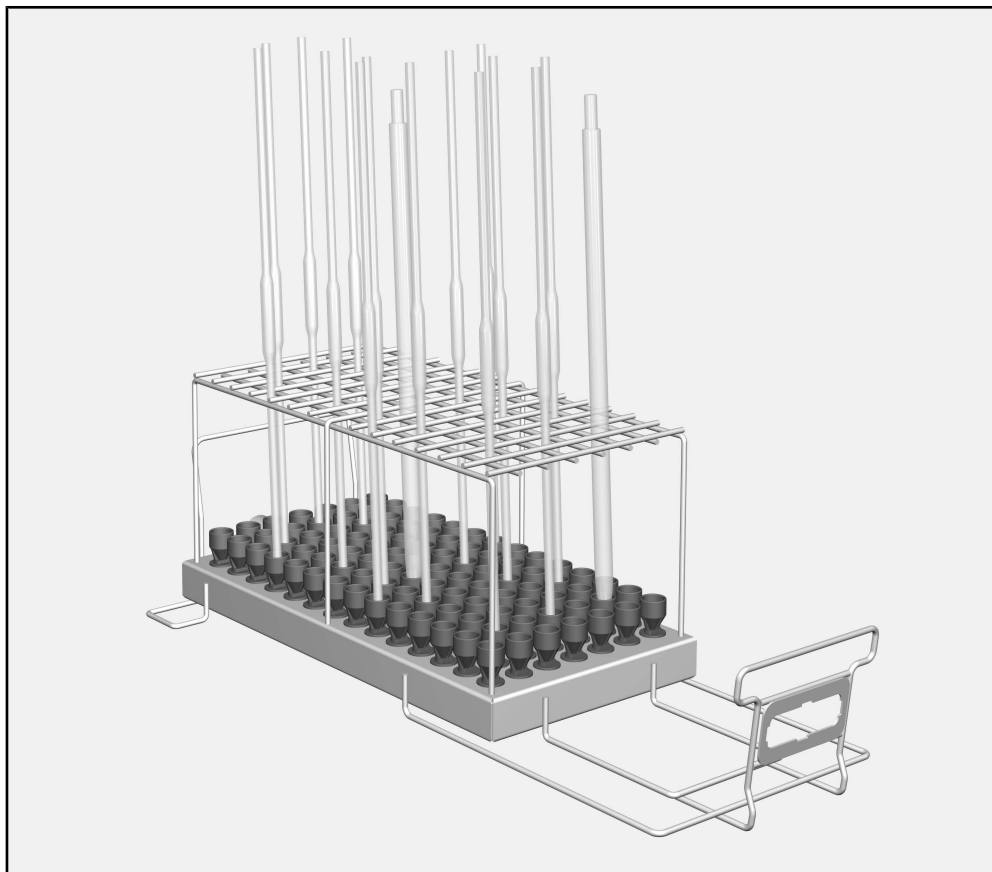

- A 613 module, height 195 mm, width 255 mm, depth 614 mm

The module is suitable for:

- Measuring pipettes with a volume from 1 ml to 50 ml
- Volumetric pipettes with a volume from 1 ml to 100 ml

Capacity depends on the maximum diameter of the pipettes.

Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

Read the operating instructions carefully before using this load carrier. This will help protect users from personal injury, and help protect the load carrier from damage.
Keep these operating instructions in a safe place.

 Please also read the operating instructions for your dishwasher or the washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ The module is approved solely for the applications specified in the operating instructions. Components such as nozzles may only be replaced with Miele accessories or genuine Miele original spare parts. Alterations or conversion of the machine, or using it for purposes other than those for which it was intended, are not permitted and could be dangerous.
- ▶ New load carriers must be cleaned in the washer-disinfector without a load prior to first use.
- ▶ Inspect all load carriers as described in the “Maintenance” section in the operating instructions for the cleaning machine.
- ▶ Only items which have been declared by their manufacturer as suitable for machine reprocessing may be processed. The manufacturer's specific reprocessing instructions must be observed.
- ▶ Broken glass and ceramics can result in serious injury when loading or unloading. Damaged glass or ceramic load items must not be reprocessed in the machine.
- ▶ Only ever place empty, unloaded modules in the mobile unit. Check that they are correctly engaged before loading them. Modules must be completely emptied before removing them. Placing loaded modules in mobile units or removing them can cause damage to the wash load and cause injury should glassware be broken.
- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

Miele cannot be held liable for damage or injury caused by non-compliance with the warning and safety notes.

en - Areas of application


The operating instructions for the loading carriage describe how to insert and remove the modules.

To check before loading the machine and before starting a programme

- Is the module correctly docked to the water supply in the load carrier?

Loading pipettes


The module can take up to 98 pipettes with a maximum length of 585 mm. During reprocessing, the bodies of volumetric pipettes must not touch each other as this can reduce the capacity of the module.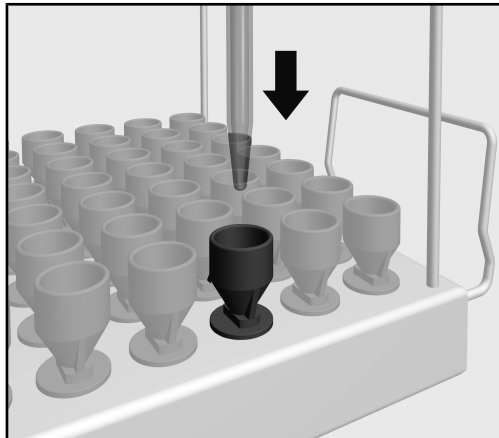

- Insert the pipettes through the frame and into the funnels with the pointed end facing downwards.

Volumetric pipettes

The bodies of volumetric pipettes must not touch each other.

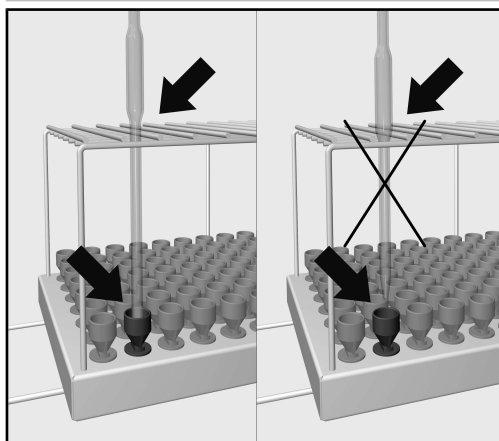- If the body of a pipette does not fit through the frame, the outflow tube will need to be long enough for the pipette to reach the funnel.

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