APFD 210

de	Flaschenkorb Gebrauchsanweisung
da	Flaskekurv Brugsanvisning
en	Bottle basket Operating instructions
fr	Porte-bouteilles Mode d'emploi
nl	Flessenrek Gebruiksaanwijzing
es	Botellero Instrucciones de manejo
pt	Cesto para garrafas Instruções de utilização
it	Cesto per bottiglie Istruzioni d'uso
fi	Pullovaunu Käyttöohje
no	flaskekurv Bruksanvisning
sv	Flaskkorg Bruksanvisning

de	4
da	11
en	18
fr	25
nl	32
es	39
pt	46
it	53
fi	60
no	67
sv	74

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
Scope of delivery	18
APFD 210	18
Overview APFD 210	19
Application technology	21
Removal of lower spray arm from ProfiLine dishwasher	21
Fitting the APFD 211 adapter	21
Removal of lower spray arm from fresh water dishwasher	22
Fitting the APFD 212 adapter	22
Preparing the load items	23

en - Scope of delivery


APFD 210

Bottle basket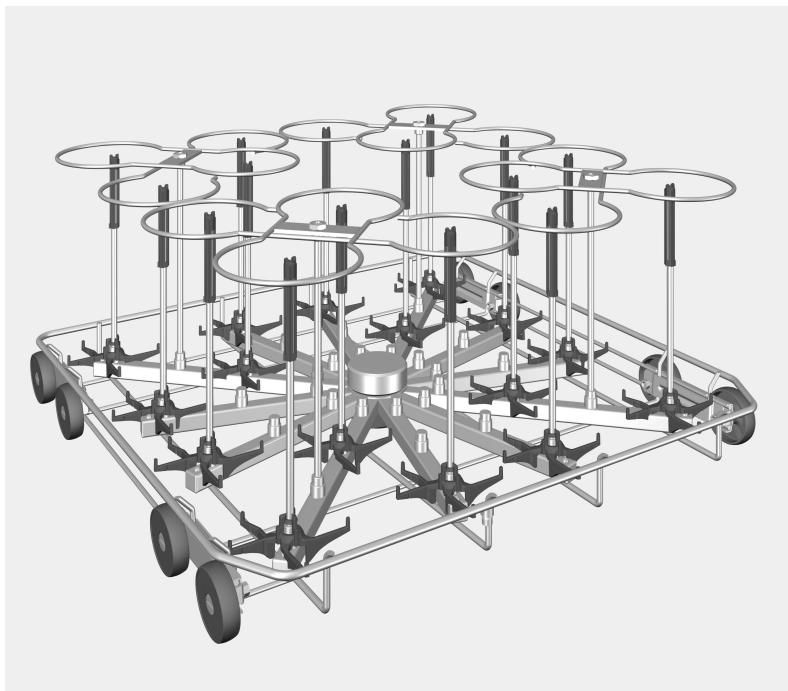

The bottle basket can be used in the following appliances:

ProfiLine dishwasher (APFD 211 adapter)	Fresh water dishwasher (APFD 212 adapter)
PFD 101	PG 8055
PFD 102	PG 8056
PFD 103	PG 8057 TD
PFD 104	PG 8058
PG 8130	PG 8059
PG 8131	PG 8060
PG 8133	PG 8061
PG 8110	PG 8096
PG 8132	PG 8099

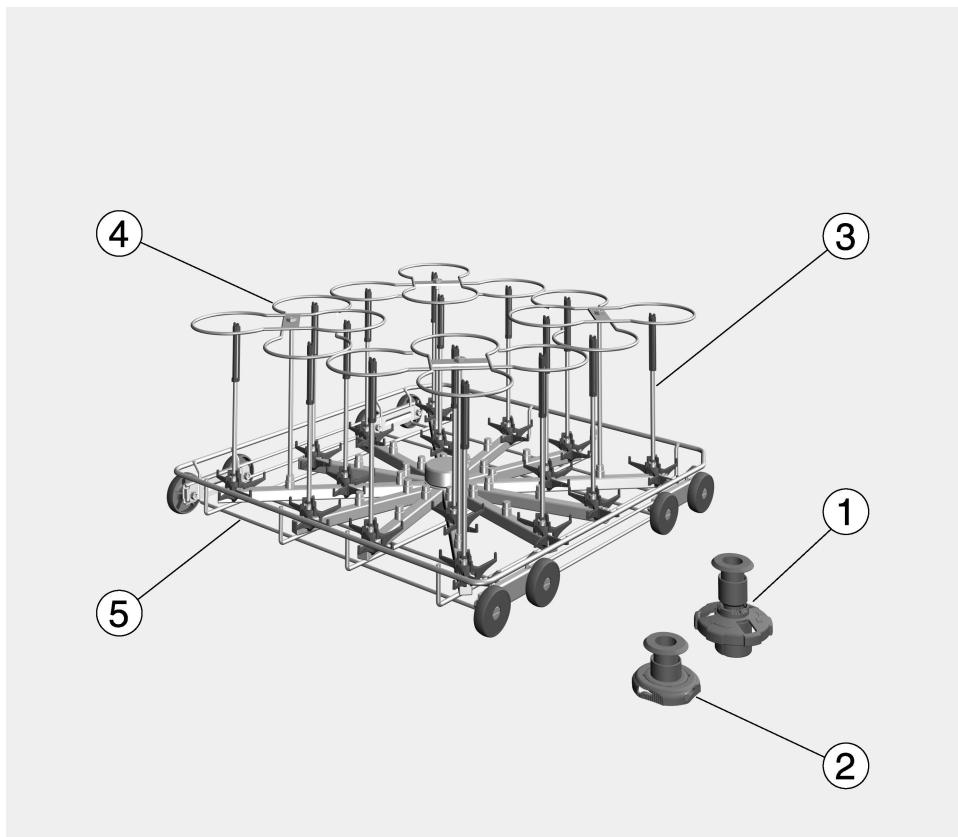- ① APFD 211 for ProfiLine dishwashers
(not included in the scope of delivery)
- ② APFD 212 for fresh water dishwashers
(not included in the scope of delivery)
- ③ Injector nozzle
- ④ Bottle holders
- ⑤ Lower basket

en - Overview APFD 210

Dimensions

Height	235.2 mm
Width	251.2 mm
Depth	515.5 mm
Max. number of nozzles	16
Height of load items, ProfiLine dish-washer with upper basket	255 mm ± 20 mm (position of upper basket bottom/middle/top)
Height of load items, fresh water dish-washer with upper basket	225 mm/275 mm (position of upper basket bottom/top)
Height of load items, fresh water dish-washer without upper basket	500 mm
Bottle diameter	max. 95 mm
Inner diameter of neck of bottle	min. 12 mm


Removal of lower spray arm from ProfiLine dishwasher

The lower spray arm must be removed before the APFD 211 adapter can be fitted in the wash cabinet of the ProfiLine dishwasher.

- Switch off the dishwasher.

Removing the lower spray arm

- Remove the lower basket.
- Rotate the screw thread of the lower spray arm anti-clockwise ① and pull the spray arm and the screw thread up and out of the wash cabinet ②.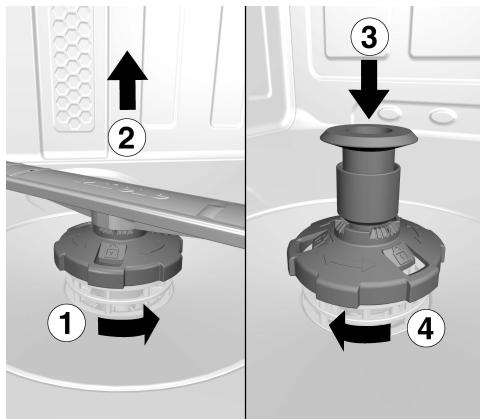

The lower spray arm ② is permanently attached to the screw thread ①. The adapter ③ can only be fitted with a permanently attached screw thread ④.

Fitting the APFD 211 adapter

- Fit the adapter with the screw thread ③ and ensure that the filter combination lies flat in the base of the wash cabinet.

- Rotate the screw thread of the adapter clockwise ④ until the arrow in the viewing window points to the lock symbol .
- Check that the adapter connection is secure.

Damage due to soiling entering the circulation system.

The adapter is used to secure the filter combination. Without filters, coarse soiling can enter the circulation system and cause a blockage. Do not run a wash programme without the adapter or without the filters.


Removal of lower spray arm from fresh water dishwasher

The lower spray arm must be removed before the APFD 212 adapter can be fitted in the wash cabinet of the fresh water dishwasher.

- Switch off the dishwasher.

Removing the lower spray arm

- Remove the lower basket.
- Rotate the screw thread of the lower spray arm anti-clockwise ① and pull the spray arm and the screw thread up and out of the wash cabinet ②.

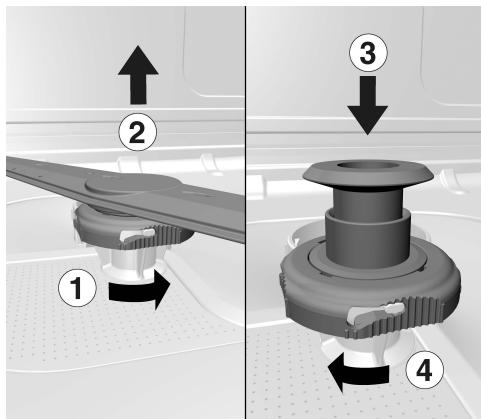The lower spray arm ② is permanently attached to the screw thread ①. The adapter ③ can only be fitted with a permanently attached screw thread ④.

Fitting the APFD 212 adapter

- Fit the adapter with the screw thread ③ and ensure that the filter combination lies flat in the base of the wash cabinet.

- Rotate the screw thread of the adapter clockwise ④ until the arrow in the viewing window points to the lock symbol .
- Check that the adapter connection is secure.

Damage due to soiling entering the circulation system.

The adapter is used to secure the filter combination. Without filters, coarse soiling can enter the circulation system and cause a blockage. Do not run a wash programme without the adapter or without the filters.

Preparing the load items

Remove any coarse soiling from the load items.

Check that the load items are suitable for cleaning in a dishwasher.

Notes

 Risk of damage caused by acids and solvents.

Acid or solvent residues, especially hydrochloric acid or chlorides, must not get inside the wash cabinet.

Do not use acids or solvents in the wash cabinet.

 Damage caused by load items soiled with ash, sand, wax, lubricating grease or paint.

These items will damage ProfiLine and fresh water dishwashers.

Do not wash items soiled with ash, sand, wax, lubricating grease or paint in ProfiLine or fresh water dishwashers.

Position the bottles in the designated locations in the bottle basket.

Please note:

■ Load the bottles so that water can access all surfaces.

The bottle basket has 16 bottle holders.

■ Load the bottles onto the injector nozzles with the openings pointing down.

The bottle basket and holders can only accommodate bottles with a maximum diameter of 95 mm.

The maximum number of bottles that can be loaded for cleaning is 16.

Insufficient cleaning

The height of the bottles loaded on the nozzles at the front ① must not exceed 200 mm.

Liquid dispensing is recommended if all of the bottles loaded for cleaning are taller than 200 mm.

■ Start the wash programme.

Recommended wash programme:
glassware

For more information, see the operating instructions for the dishwasher you are using.

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