A 200

de	Gebrauchsanweisung Injektorwagen
da	Brugsanvisning Injektorvogn
en	Operating instructions Injector unit
es	Instrucciones de manejo Carros inyectores
fi	Käyttöohje Suorasuihkuvaunu
fr	Mode d'emploi Chariot à injection
hr	Upute za uporabu Injektorska kolica
it	Istruzioni d'uso Carrello iniettore
nl	Gebruiksaanwijzing injectorwagen
no	Bruksanvisning injektorvogn
pt	Instruções de utilização Carro injetor
sl	Navodila za uporabo Injektorski voziček
sv	Bruksanvisning Injektorvagn

de	4
da	11
en	18
es	25
fi	32
fr	39
hr	46
it	53
nl	60
no	67
pt	74
sl	81
sv	88

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Notes about these instructions	18
Appropriate use	19
Queries and technical problems	19
Items supplied	20
Disposal of the packing material	20
Warnings and safety notes	21
Areas of application.....	22
To check before loading the machine and before starting a programme	22
Loading pipettes	22

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

- Select an option using the arrow buttons and save your choice with *OK*.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings  menu.

Re-usable laboratory glassware and laboratory utensils can be reprocessed using this mobile unit in a Miele laboratory glassware washer-disinfector. Follow the instructions for your washer-disinfector as well as the instructions of the glassware manufacturer on how to reprocess their items by machine.

The A 200 mobile injector unit is set up for reprocessing pipettes.

Use the Pipettes programme.


The mobile unit can be used in the following washer-disinfectors:

- | | |
|---------------|--------------|
| - PLW 8683 | - PG 8583 |
| - PLW 8683 CD | - PG 8583 CD |
| - PLW 8693 | - PG 8593 |
| - PLW 8636 | |
| - LAB 5010 | |
| - LAB 5011 | |
| - LAB 5020 | |

The washer-disinfector is generally referred to as “the machine” in these operating instructions. Reprocessable laboratory glassware and utensils are referred to as “items” if they are not more closely defined.

Queries and technical problems

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your dishwasher or the washer-disinfector or go to www.miele-professional.com.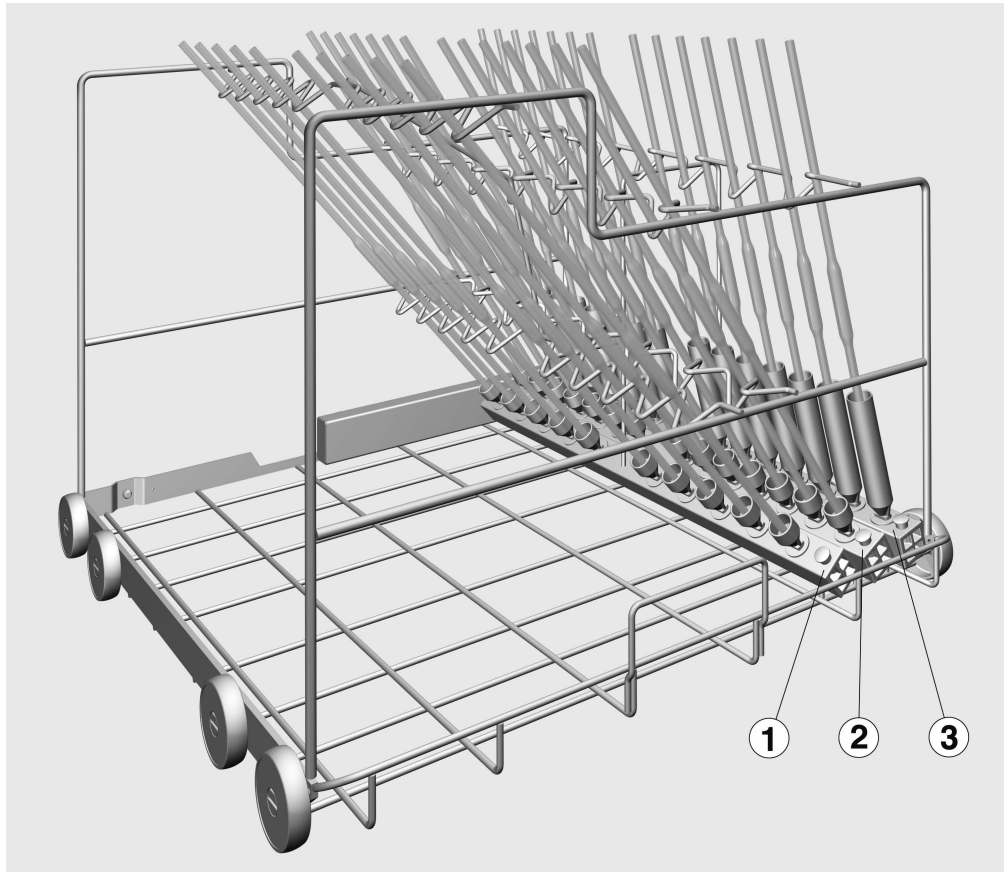

- Mobile injector unit A 200, height 397 mm, width 531 mm, depth 546 mm, with 3 injector manifolds.

Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

To avoid the risk of accidents and damage to this mobile unit please read these instructions carefully before using it for the first time. Keep these instructions in a safe place where they are accessible to users at all times.

 Please also read the operating instructions for your dishwasher or the washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ The mobile unit is approved solely for the applications specified in the “Areas of application” chapter of these operating instructions. Applications that go beyond those listed and use of the mobile unit with additional components from the Miele product range are described in the operating instructions for the components or are to be agreed with Miele.
- ▶ Inspect all mobile units, baskets, modules and inserts daily as described in the “Maintenance” section in the Operating instructions for the washer-disinfector.
- ▶ Only items which have been declared by their manufacturer as suitable for machine reprocessing may be processed. The manufacturer's specific reprocessing instructions must be observed.
- ▶ Broken glass can result in serious injury when loading or unloading. Broken glass items must not be processed in the machine.
- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

Miele cannot be held liable for damage or injury caused by non-compliance with the warning and safety notes.

Before placing pipettes in the injector unit, please make sure that all hollow sections are free of blockages.

To check before loading the machine and before starting a programme


- Check that the mobile unit is correctly docked to the water path in the chamber.

Loading pipettes


The mobile injector unit has three injector manifolds that can be loaded as follows:

Row	Max. length [mm]	Qty	Volume [ml]
1	600	10	10 to 100
2	500	14	10 to 25
3	440	14	1 to 10

The pipettes must not protrude beyond the sides of the mobile unit.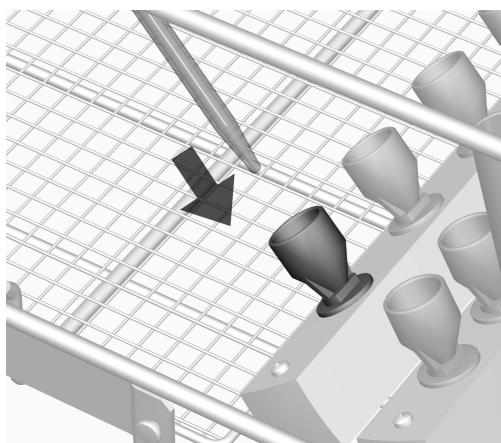

- Insert the pipettes into the funnels with the pointed end facing downwards.
- Place the upper end of the pipettes in the corresponding indentations on the rods.

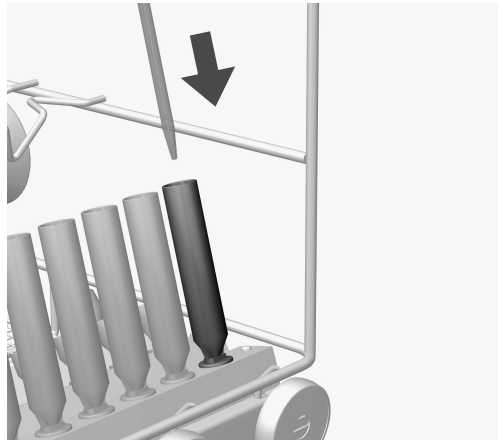- Place pipettes with a volume of 1 to 5 ml in the irrigation sleeves in row ③ of the mobile injector unit.

The wash pressure could push these pipettes out of the funnels in rows ① and ②.

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