A 623

de	Gebrauchsanweisung Injektormodul
cs	Návod k obsluze Injektorový modul
da	Brugsanvisning Injektormodul
en	Operating instructions Injector module
es	Instrucciones de manejo Módulo inyector
fi	Käyttöohje Suorasuihkumoduuli
fr	Mode d'emploi Module à injection
it	Istruzioni d'uso Modulo a iniezione
ja	取扱説明書 インジェクターモジュール
ko	사용 설명서 인젝터 모듈
nl	Gebruiksaanwijzing injectormodule
no	Bruksanvisning injektormodul
pl	Instrukcja użytkowania Moduł iniekcyjny
pt	Instruções de utilização Módulo injetor
ru	Инструкция по эксплуатации Инжекторный модуль
sk	Návod na použitie Injektorový modul
sv	Bruksanvisning injektormodul
tr	Kullanım Kılavuzu Enjektör Modülü
uk	Інструкція з експлуатації Інжекторний модуль
zh	操作说明书 注射器模块

de	4
cs	13
da	22
en	31
es	40
fi	50
fr	59
it	68
ja	77
ko	86
nl	96
no	105
pl	115
pt	124
ru	133
sk	142
sv	152
tr	161
uk	171
zh	181

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en - Contents

Notes about these instructions	31
Intended use	32
Queries and technical problems.....	32
Items supplied	33
Load carriers	33
Disposal of the packing material	33
Warning and Safety instructions	34
Areas of application	35
Before loading the machine and before starting a programme	35
Preparing the load	35
Inserting the measuring cylinder	35
Installation	37
Tools required:	37
Screwing in nozzles.....	37
Screwing the injector nozzle onto the module	37
Mounting the holder frames	38

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property.
Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

■ Select an option using the arrow buttons and save your choice with *OK*.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings  menu.

This module can be used to reprocess machine-reprocessible laboratory glassware and utensils in a Miele washer-disinfector for laboratory glassware and utensils. Follow the operating instructions for your washer-disinfector as well as the instructions of the glassware and utensil manufacturer on how to reprocess their items by machine.

The A 623 injector module is designed for reprocessing measuring cylinders with a volume of 1000 ml and 2000 ml. The module can hold both short and tall construction type measuring cylinders.


The module can be used in an A 503 mobile unit.

The washer-disinfector is generally referred to as “the machine” in these operating instructions. Reprocessible laboratory glassware and utensils are referred to as “items” if they are not more closely defined.

Queries and technical problems

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your washer-disinfector or go to www.miele-professional.com.

Load carriers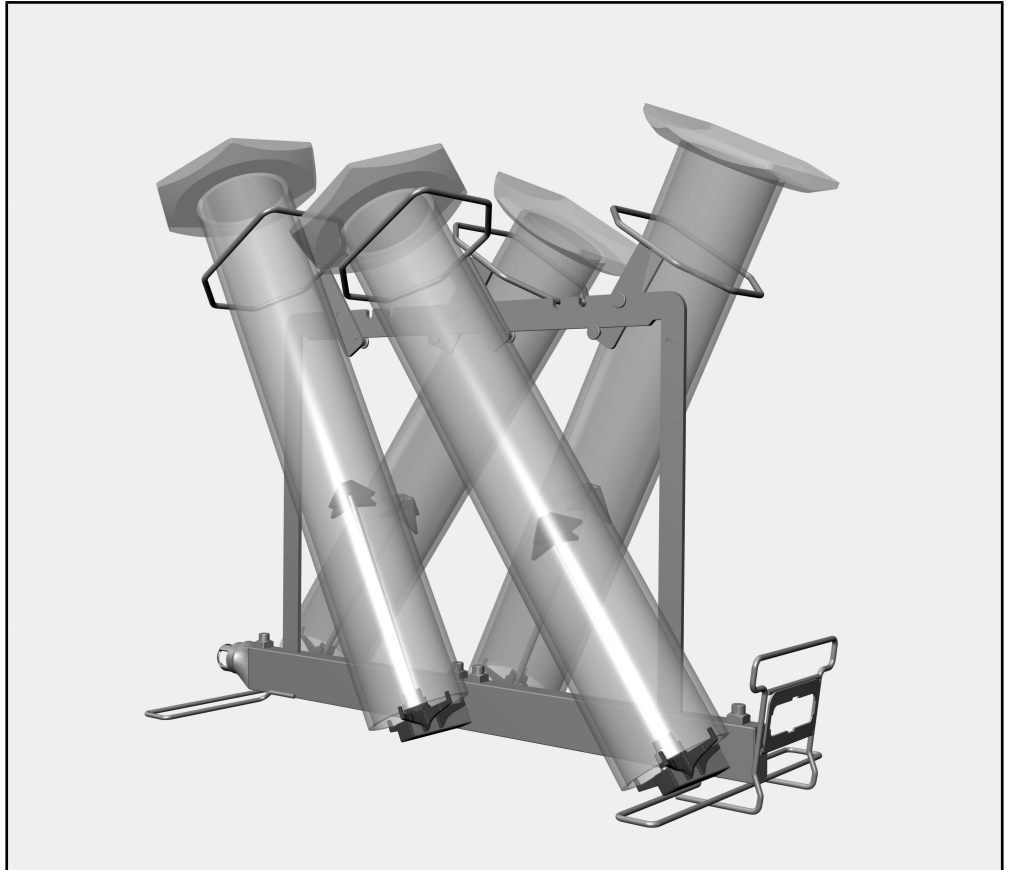

- A 623 module, height 437 mm, width 255 mm, depth 608 mm

Items supplied

- 4 injector nozzles with plastic supports for reprocessing measuring cylinders, length 217 mm, Ø 6 mm
- 4 holder frames with coating; for measuring cylinders up to max. Ø 110 mm

Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in landfill sites.

en - Warning and Safety instructions

Read the operating instructions carefully before using this load carrier. This will help protect users from personal injury, and help protect the load carrier from damage.
Keep these operating instructions in a safe place.

 Please also read the operating instructions for your washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ The module is approved solely for the applications specified in the operating instructions. Components such as nozzles may only be replaced with Miele accessories or genuine Miele original spare parts. Alterations or conversion of the machine, or using it for purposes other than those for which it was intended, are not permitted and could be dangerous.
- ▶ New load carriers must be cleaned in the washer-disinfector without a load prior to first use.
- ▶ Inspect all load carriers as described in the “Maintenance” section in the operating instructions for the cleaning machine.
- ▶ Only items which have been declared by their manufacturer as suitable for machine reprocessing may be processed. The manufacturer's specific reprocessing instructions must be observed.
- ▶ Broken glass can result in serious injury when loading or unloading. Damaged glass items must not be reprocessed in the machine.
- ▶ Only ever place empty, unloaded modules in the mobile unit. Check that they are correctly engaged before loading them. Modules must be completely emptied before removing them. Placing loaded modules in mobile units or removing them can cause damage to the wash load and cause injury should glassware be broken.
- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

The operating instructions for the loading carriage describe how to insert and remove the modules.

Before loading the machine and before starting a programme

- Check that fittings such as irrigation sleeves and injector nozzles are securely screwed into position.

⚠ Make sure that all screw connectors are fitted with injector nozzles, adapters, irrigation sleeves or blind stoppers to ensure that all fittings in use are supplied with sufficient standardised pressure.

Damaged fittings such as injector nozzles, adapters and irrigation sleeves must not be used.

Fittings not equipped with wash items do not have to be replaced blind stoppers.


- Is the module correctly docked to the water supply in the load carrier?

Preparing the load

⚠ Damage to the measuring cylinders.

If the measuring cylinders touch each other during reprocessing or rest against the wall of the wash cabinet or against the door of the cleaning machine, damage can occur at the points of contact. For example, the items may become scratched or the glass may break. Arrange the measuring cylinders so that they are not touching each other or resting against the wall of the wash cabinet or against the door of the cleaning machine.

Inserting the measuring cylinder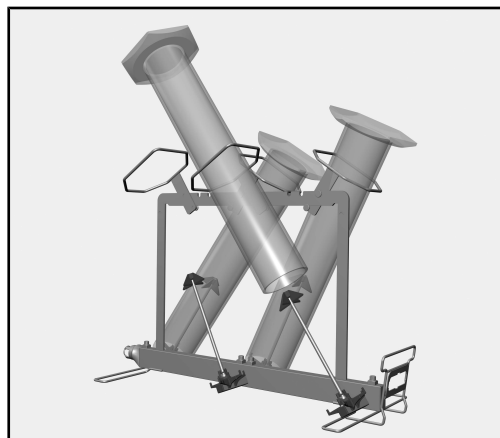


- Guide the measuring cylinder through the holder frame and over the nozzle and place it on the support for the injector nozzle.

en - Areas of application

Depending on the height of the measuring cylinders, the following must be observed:

Tall measuring cylinders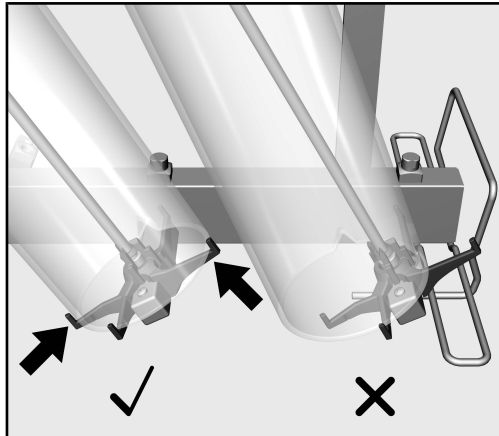


The openings of the measuring cylinders should rest between the support hooks on the injector nozzles to stabilise the cylinders during reprocessing.

If the measuring cylinders cannot be placed between the support hooks, the angle between the measuring cylinders and injector nozzles must be adjusted. To do this:

- Move the holder frame into another recess on the strut

Short measuring cylinders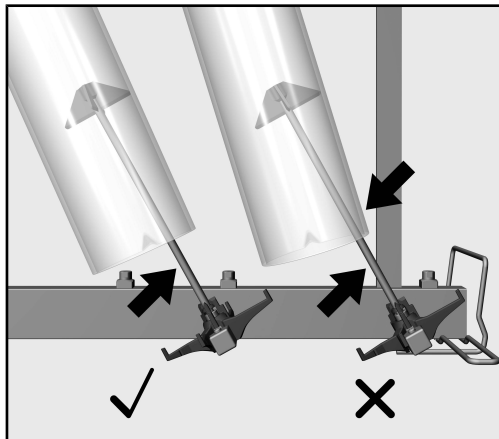

The measuring cylinders must rest on the plastic supports of the injector nozzles. However, they must not be allowed to touch the nozzles themselves.

If the measuring cylinders are touching the nozzles, the angle between the measuring cylinders and the injector nozzles must be altered. To do this:

- Move the holder frame into another recess on the strut

or


- Slide the plastic support on the injector nozzle downwards

Tools required:

- 9 mm Spanner (WAF 9)

Screwing in nozzles

Screwing the injector nozzle onto the module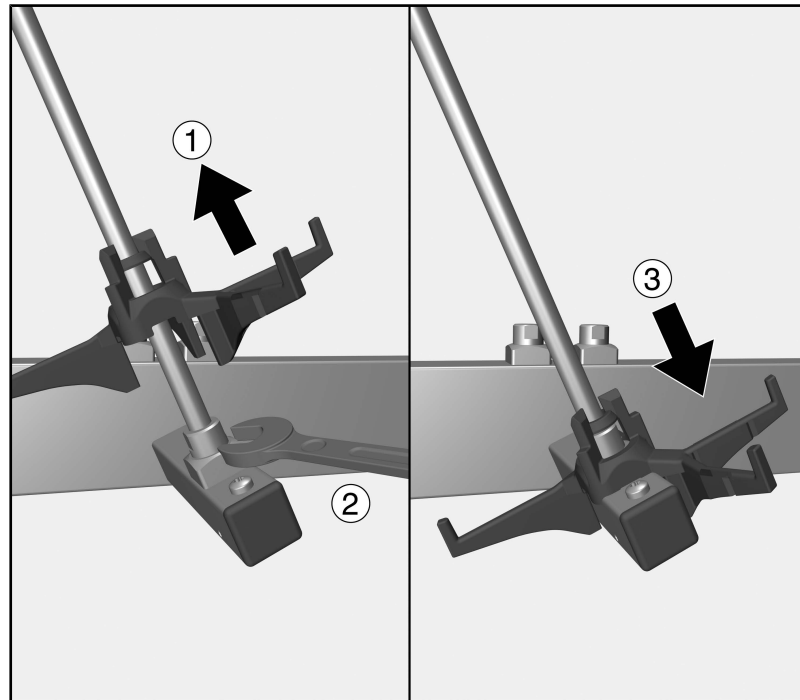


- Lift up the support on the injector nozzle and screw the injector nozzles into the required position ①.
- Use the open spanner to tighten the injector nozzle ②.
- Slide the support on the injector nozzle downwards until it rests on the water supply pipe ③.

Screw the other injector nozzles into the required positions.

Mounting the holder frames

2 holder frames are placed on each side of the upper strut of the module. The slanted corners of the holder frames centre the measuring cylinders during reprocessing. After mounting, they must point downwards.

The centre holder frames must be mounted on both sides first.

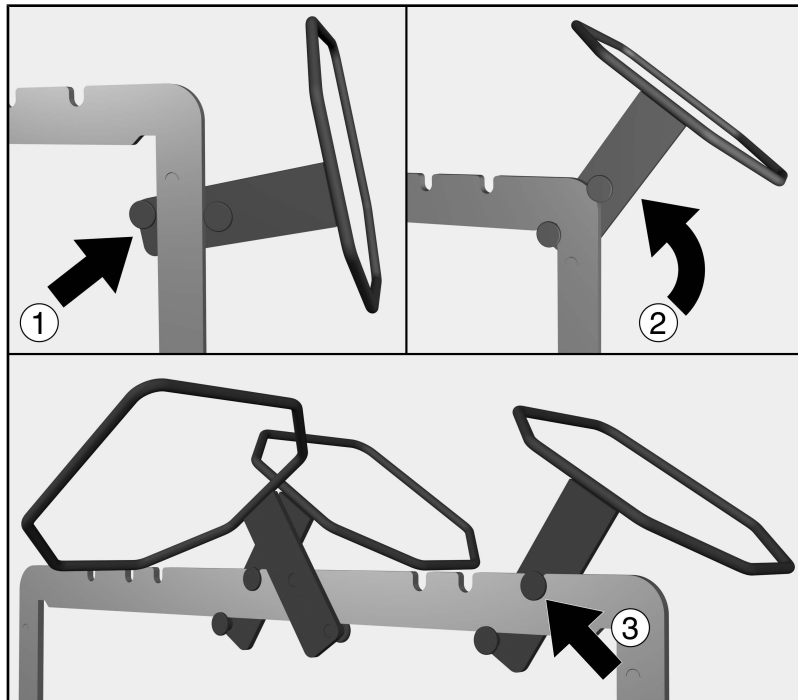- Position the holder frame on one of the side struts of the module so that the pins of the frame reach around the strut ①. The holder frame must be facing outwards and the slanted corners must be facing downwards.
- Slide the holder frame upwards along the strut ②.
- Secure the holder frame in a recess ③.
- Place the other holder frames in the required positions.

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