


de	Zusätzliches Dosiersystem	en	Additional dispenser system	cs	Doplňkový dávkovací systém
da	Ekstra doseringssystem	el	Επιπρόσθετο σύστημα δοσομέτρησης	en-US	Additional dispenser system
es	Sistema de dosificación adicional	fr	Système de dosage supplémentaire	it	Sistema di dosaggio aggiuntivo
nl	Extra doseersysteem	no	Ekstra doseringssystem	pt	Sistema de dosagem adicional
ru	Дополнительная система дозирования	sv	Extra doseringssystem		

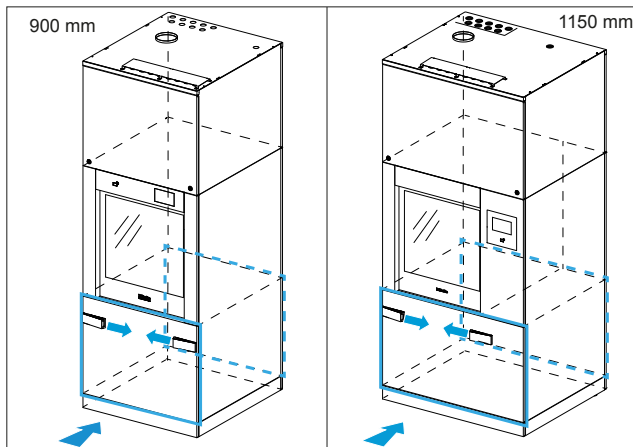
1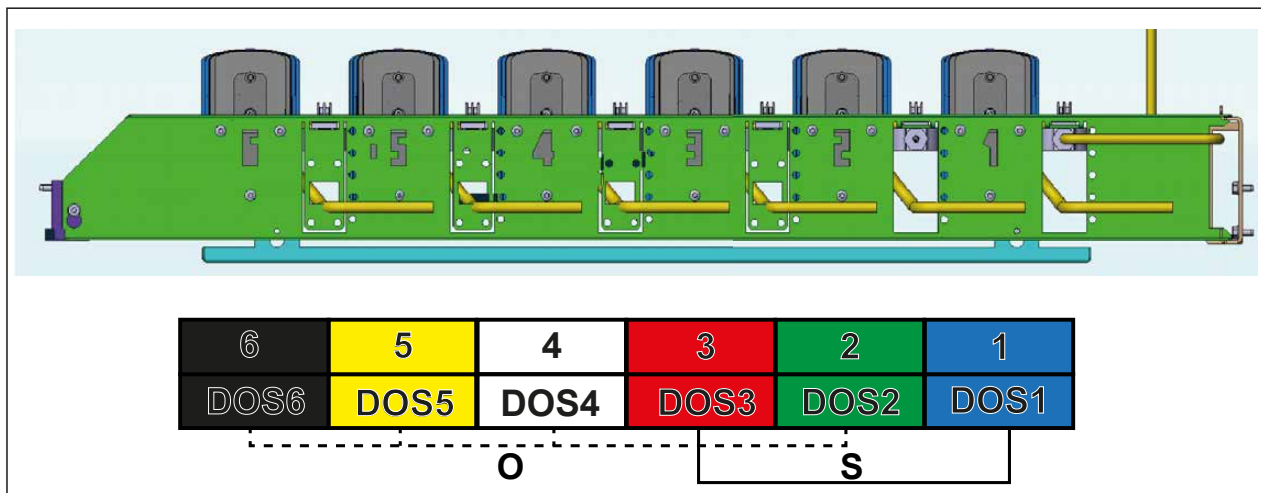


2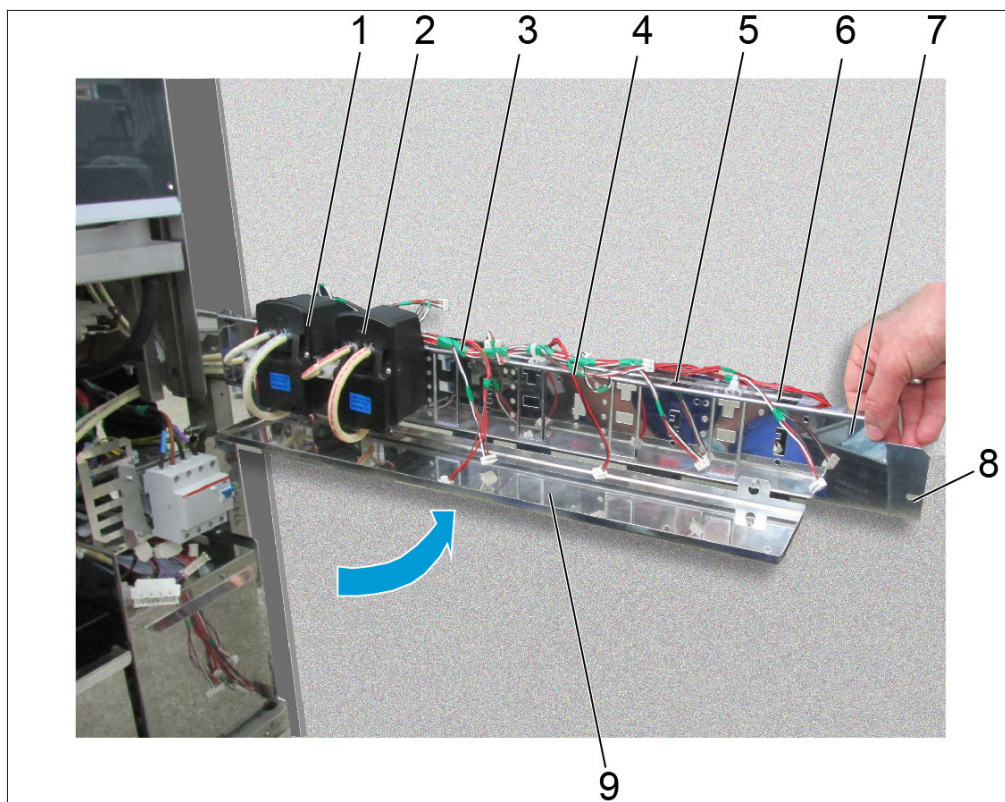


3

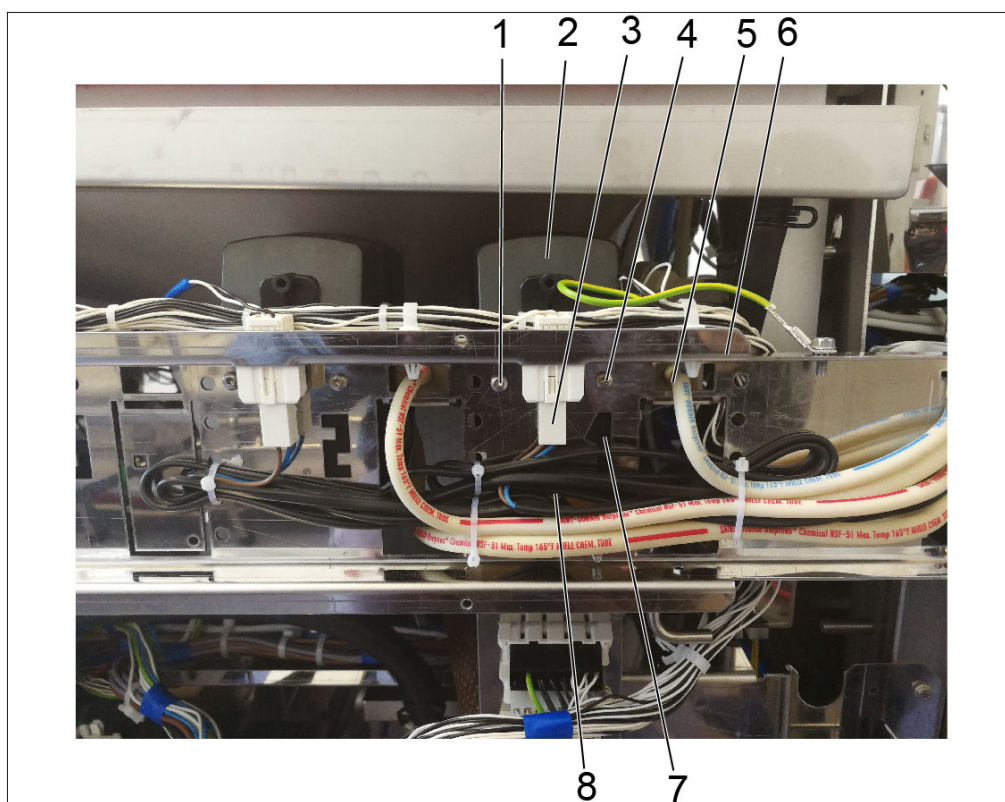
4

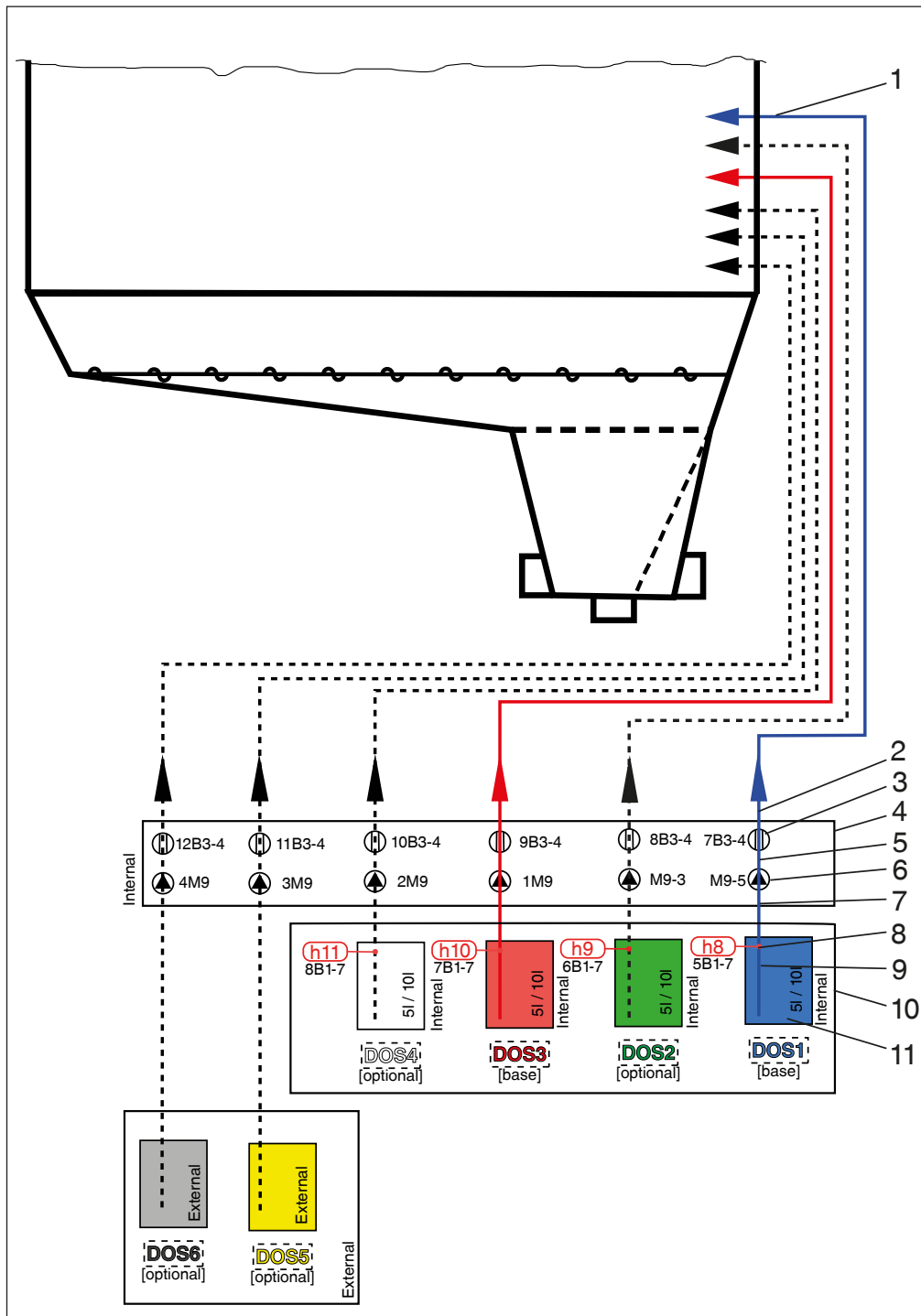
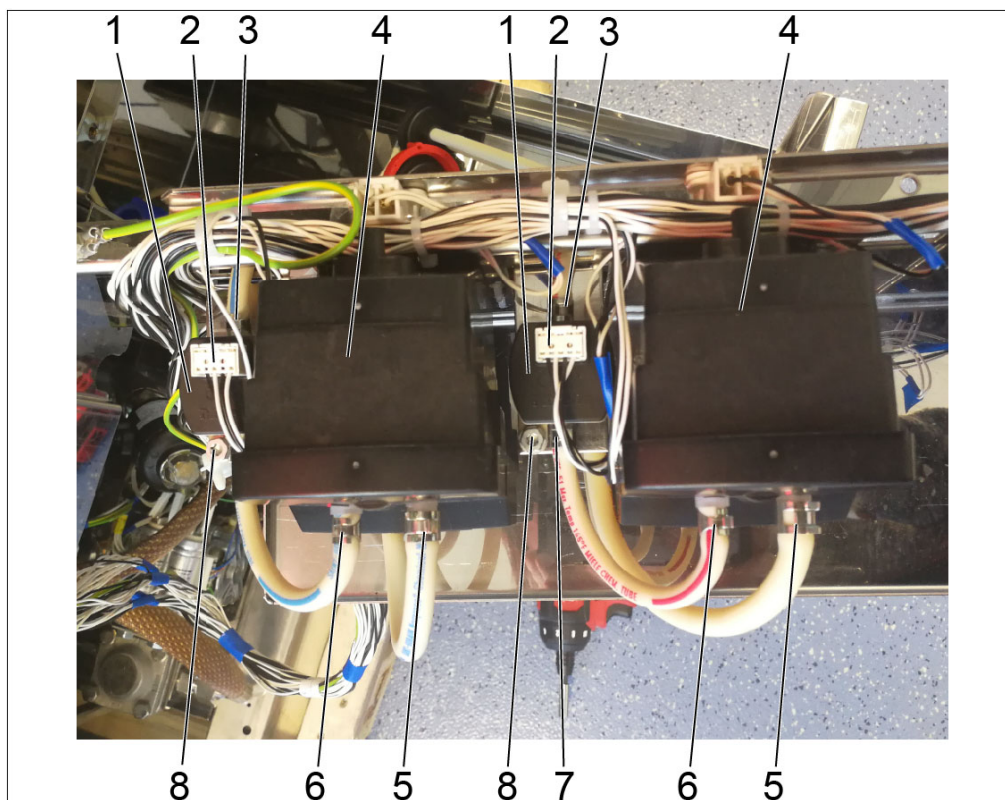
6

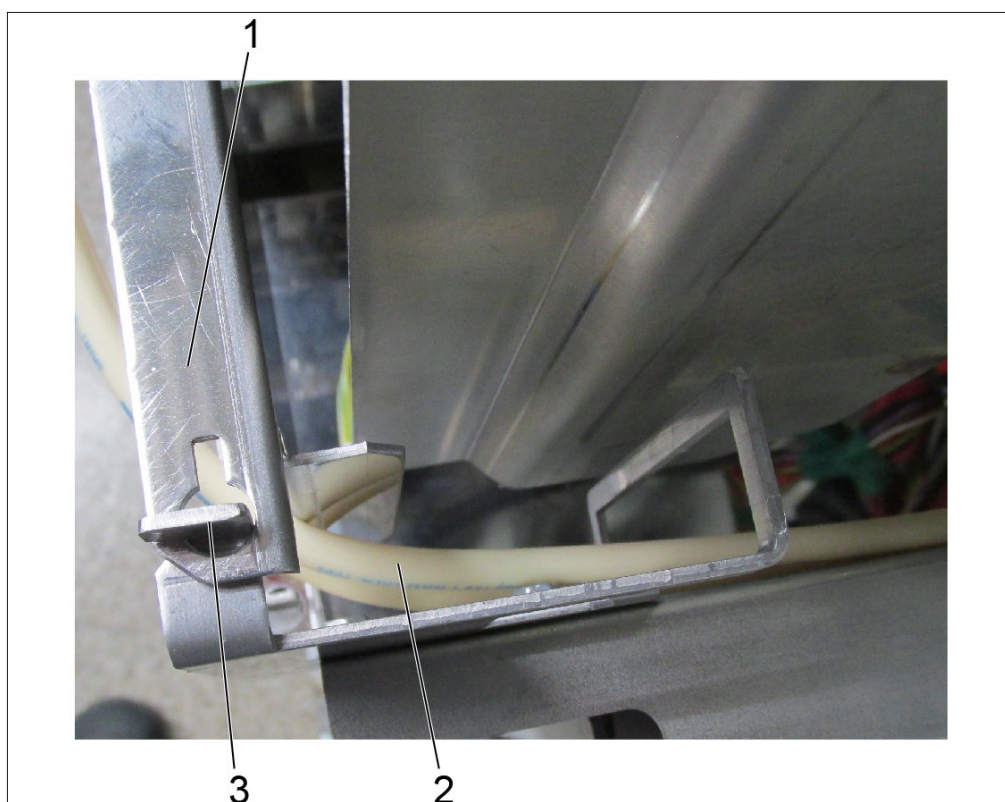7

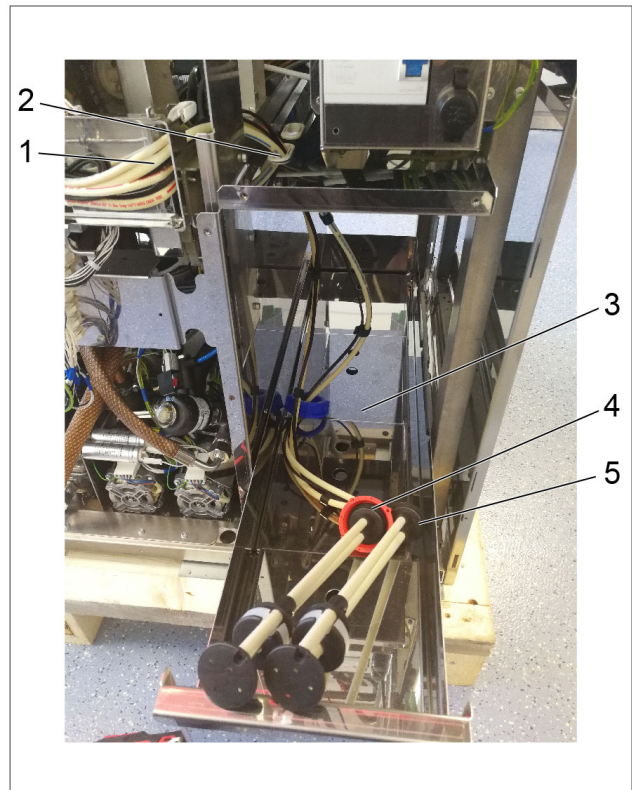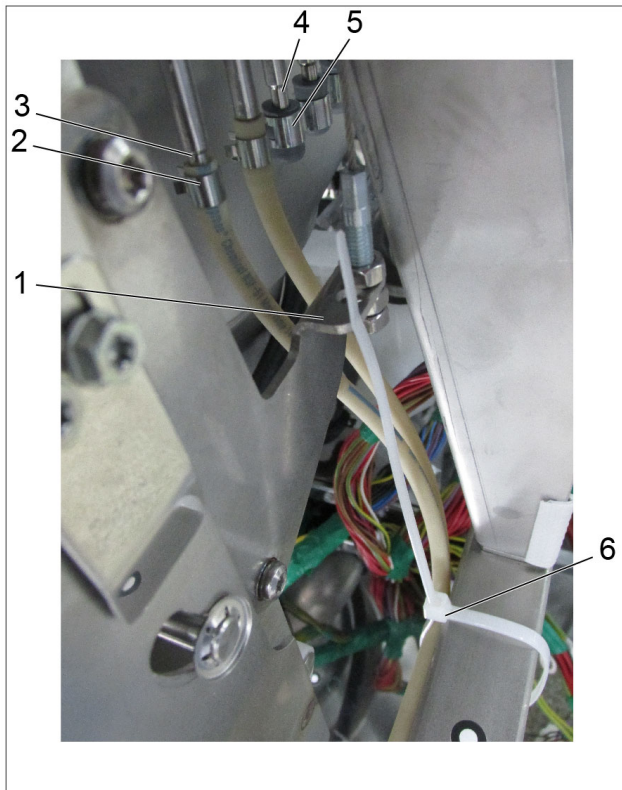
8

10

9

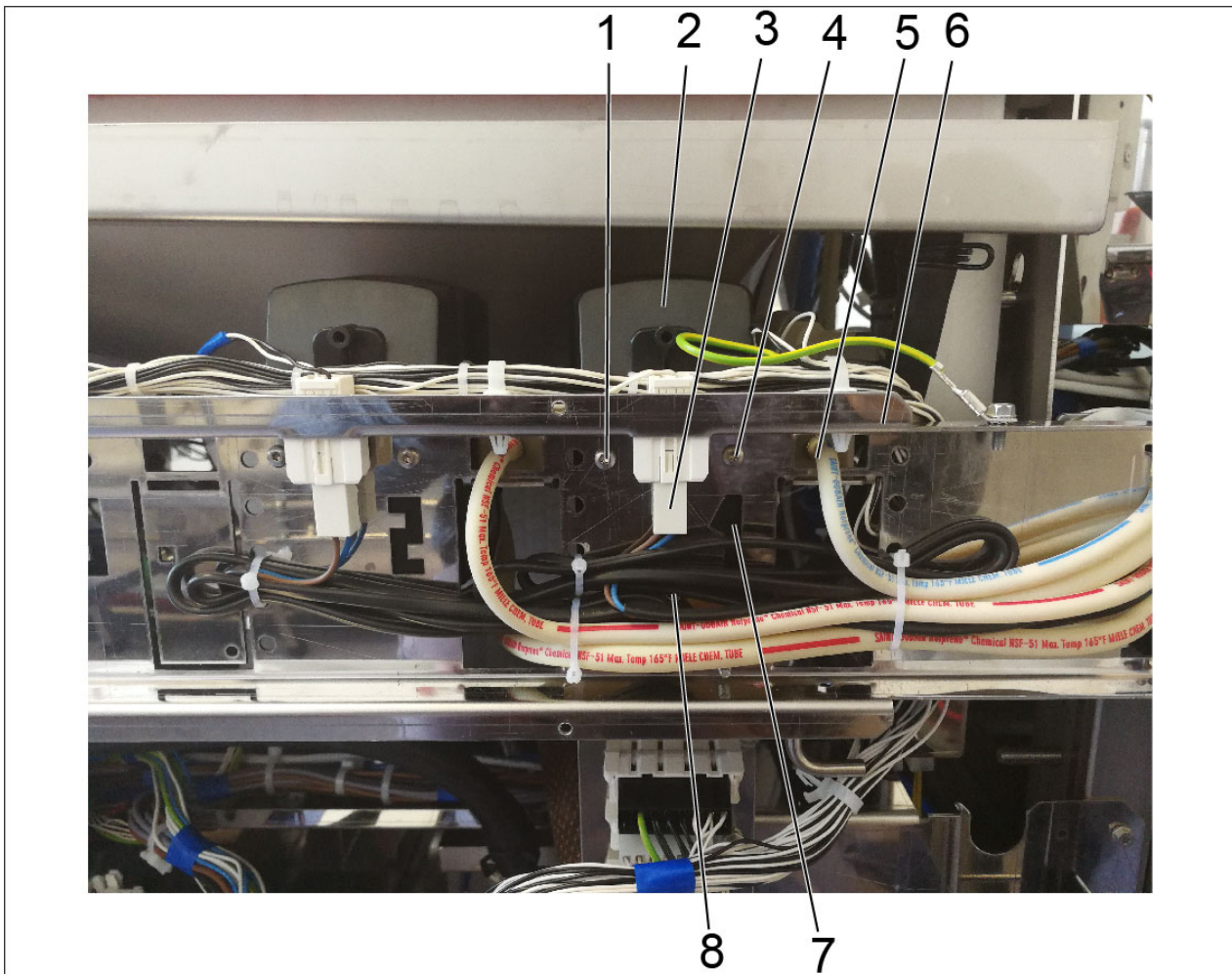11

de

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en

Additional dispenser system

Reason: The conversion kit enables an additional internal dispenser system to be fitted in models PLW 8615, PLW 8616, PLW 8617, PWD 8626 or PWD 8628. There are optionally two different versions for 5 l dispenser container (retrofitting kit - Liquid dispensing, DS 5) or 10 l dispenser container (retrofitting kit - Liquid dispensing, DS 10). The conversion kits differ only in the type of suction tube included.

Retrofitting kit - Liquid dispensing DS 5 - Suction tube

Included parts

No.	Mat. no.	Designation
1	09065773	Suction tube 200 mm

Retrofitting kit - Liquid dispensing DS 10 - Suction tube

Included parts

No.	Mat. no.	Designation
1	09065743	Suction tube 333 mm

Both kits contain the following identical parts:

Included parts

No.	Mat. no.	Designation
1	04012531	Black cap DRM64
1	10873120	Connection cable
1	10895470	Hose, white 5.5 - 1750 mm
1	10895240	Hose, white 3.7 - 1000 mm
1	11043800	Hose clip 8 mm
4	09720740	Hose clip 6 mm
1	09983790	Hose clip 9 mm
1	10303990	Dispenser pump 24 V
3	10719980	Raised-head self-tapping screw 3.5 x 10 EDST
1	10448530	Flow meter

No.	Mat. no.	Designation
1	07246730	Bolt DM4 x 10
2	04658581	Adhesive label, red
2	04658591	Adhesive label, blue
2	06739710	Sticker, yellow
2	04658561	Adhesive label, green
2	04658571	Adhesive label, white
2	05693210	Adhesive label, black
3	03069180	Cable holder
6	04982150	Cable holder

In order to carry out this work, the appropriate Technical Service Documentation (TSD) is required.

The warning and safety instructions given in the applicable model-specific TSD must be complied with.

Work must be carried out in accordance with the instructions given in the applicable model-specific TSD.

Disconnect the machine from the mains and ensure power cannot be switched on again in error.

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Do not lay cables over sharp edges or hot surfaces.

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Bottom part service panel opening

The bottom part service panel is secured with 2 bolts which each clip into the frame at the sides of the machine.

✚ Slide the service panel retaining bolts towards the centre of the machine at the same time, see Fig. 1.

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Gaining access to dispenser systems

Dispenser system colour coding, see Fig. 2.

1. Dispenser position 1 for detergent, colour coding blue (S = Standard basic version)
2. Dispenser position 2, colour coding green (O = Optional)
3. Dispenser position 3 for neutralisation agent, colour coding red (S = Standard basic version)
4. Dispenser position 4, colour coding white (O = Optional)
5. Dispenser position 5, colour coding yellow (O = Optional) without level monitoring
6. Dispenser position 6, colour coding black (O = Optional) without level monitoring

✚ Hold the dispenser cover plate firmly and at the same time remove the dispenser cover plate fixing screws.

✚ Remove the dispenser cover plate.

✚ Remove the fixing screw (8) for the dispenser holding plate (9) on the left, see Fig. 3.

✚ Open the dispenser holding plate, see Fig. 3.

Dispenser pump fitting

✚ Place the dispenser pump (2) on the appropriate dispenser position (7) and secure it with the fixing screws (1), (4) and (8), see Fig. 4.

✚ Unwind the connection plug for the dispenser pump (2) from the machine wiring harness and connect it at the bottom to the dispenser pump (2), see Fig. 4.

✚ Fit the appropriate colour coding on the dispenser pump.

Fitting dispenser hoses

Depending on the machine variant, the dispenser drawer with the dispenser containers is located either under the dispenser system or fitted in a neighbouring unit. As a result the routes for laying the dispenser hoses vary slightly.

The dispenser hose from the dispenser probe to the dispenser pump has a larger diameter than the connection hose between the dispenser pump and flow meter, and the dispenser hose between the flow meter and dispenser connection to the cabinet.

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- ✂ Cut dispenser hoses with a straight cut using a utility knife or similar tool.
- ✂ Do not use scissors, shears or similar to cut the dispenser hoses.

Connection plan legend, see Fig. 5.

1. Dispenser connections to cabinet
2. Narrow dispenser hose: Flow meter - Dispenser connection
3. Flow meter
4. Dispensing system holding plate
5. Narrow dispenser hose (approx. 200 mm long): Dispenser pump - Flow meter
6. Dispenser pump
7. Wide dispenser hose: Dispenser probe - Dispenser pump
8. Fill level monitor on dispenser probe
9. Dispenser probe
10. Dispenser drawer
11. Dispenser container

- ✂ Use the paths taken by the hoses already fitted in the machine as a guide.
- ✂ Always lay the hoses with gentle curves without kinking or straining them.

Laying dispenser hose connections between dispenser pump and flow meter

- ✂ Cut of a 200 mm piece of the narrow dispenser hose.
- ✂ Fit a hose clip (6 mm) on each end of the 200 mm dispenser hose.
- ✂ Connect the dispenser hose to the narrow hose connector (6) on the dispenser pump (4) and secure it with a hose clip, see Fig. 6.
- ✂ Connect the other end of the hose to the hose connector (7) on the flow meter (1) on the side with the fixing screws (8) and secure it with a hose clip, see Fig. 6.
- ✂ Fit a hose clip (6 mm) on one end of the remaining narrow dispenser hose.

Laying dispenser hose between flow meter and dispenser connection on cabinet

- ✂ Lay the remaining narrow dispenser hose to the machine dispenser connection (3) through the dispenser holding plate to the flow meter, see Fig. 6.
- ✂ Connect the dispenser hose to the open hose connector (3) on the flow meter (1) and secure it with a hose clip (6 mm), see Fig. 6.
- ✂ Bend out the metal plate lug to secure the flow meter.
- ✂ Secure the flow meter with the fixing screw (8), see Fig. 6.
- ✂ Lay the dispenser hose (2) along the dispenser holder plate (1) towards the hinged joint (3), see Fig. 7.
- ✂ Pass the dispenser hose at the hinged joint (3) through the opening, see Fig. 7.

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✂ Lay hoses under the door holding bracket.

- ✂ Lay the dispenser hose along the existing dispenser hoses towards the dispenser connections on the cabinet (4) and fit a hose clip (6 mm) on the end, see Fig. 8.
- ✂ Lay the dispenser hose under the door holding bracket (1), see Fig. 8.
- ✂ Remove the dispenser connection cover (5), see Fig. 8.
- ✂ Fit the dispenser hose end fully onto the dispenser connection (4) and secure it with a hose clip (6 mm), see Fig. 8.
- ✂ Fasten the dispenser hose with cable ties at suitable points. At the same time ensure that the dispenser hoses are not closed.
- ✂ Fit the appropriate colour coding on the dispenser hose.

Laying dispenser hose between suction probe and dispenser pump

- ✂ Fit a hose clip (8 mm) on one end of the wide dispenser hose.
- ✂ Connect the dispenser hose to the hose connector (2) on the dispenser probe in the dispenser drawer and secure it with a hose clip (9 mm), see Fig. 9.
- ✂ Lay the dispenser hose from the dispenser probe (4) along the existing dispenser hoses (1) to the dispenser pump, see Fig. 10.
- ✂ Fit a hose clip (8 mm) on the end of the dispenser hose.
- ✂ Connect the dispenser hose to the wide hose connector (9) on the dispenser pump (4) and secure it with a hose clip (8 mm), see Fig. 6.
- ✂ Fit the appropriate colour coding on the dispenser hose and dispenser probe.

Laying cable connection for suction probe fill level monitor

- ✂ Connect the fill level monitor (1) connection plug to the suction probe, see Fig. 9.
- ✂ Lay the fill level monitor connection cable along the appropriate dispenser hose (2), see Fig. 9.
- ✂ Secure the fill level monitor connection cable with cable holders (3) to the appropriate dispenser hose (2), see Fig. 9.
- ✂ Connect the fill level monitor connection plug (3) to the wiring harness plug connector in the machine, see Fig. 11.

Dispenser system activation

- ✂ Log in with a service technician with administrator rights.
- ✂ Select the System menu.
- ✂ Select the Machine configuration menu.
- ✂ Select the Dispensing submenu.
- ✂ Select the appropriate dispenser system.
- ✂ Set the dispenser system parameters.

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el**en-US**

Additional dispenser system

Reason: This conversion kit enables an additional internal dispenser system to be installed in models PLW 8615, PLW 8616, PLW 8617, PWD 8626 or PWD 8628. There are optionally two different versions for 5-liter dispensing containers (retrofit kit - Liquid dispensing, DS 5) or 10-liter dispensing containers (retrofit kit - Liquid dispensing, DS 10). The conversion kits differ only in the type of suction tube included.

Retrofit kit - Liquid dispensing DS 5 - Suction tube**Parts included**

Quantity	Mat. no.	Designation
1	09065773	Suction tube, 200 mm

Retrofit kit - Liquid dispensing DS 10 - Suction tube**Parts included**

Quantity	Mat. no.	Designation
1	09065743	Suction tube, 333 mm

Both kits contain the following identical parts:**Parts included**

Quantity	Mat. no.	Designation
1	04012531	Black cap DRM64
1	10873120	Connection cable
1	10895470	White hose, 5.5 - 1750 mm
1	10895240	White hose, 3.7 - 1000 mm
1	11043800	Hose clip, 8 mm
4	09720740	Hose clip, 6 mm
1	09983790	Hose clip, 9 mm
1	10303990	Dispensing pump, 24V
3	10719980	Raised-head self-tapping screw, 3.5 x 10 EDST
1	10448530	Flow meter
1	07246730	DM 4x10 bolt
2	04658581	Adhesive label, red
2	04658591	Adhesive label, blue
2	06739710	Sticker, yellow
2	04658561	Adhesive label, green
2	04658571	Adhesive label, white
2	05693210	Adhesive label, black
3	03069180	Cable tie
6	04982150	Cable tie

The TSD will be needed for all service work

The warning and safety instructions in the model-specific TSD must be observed.

Service work must be carried out following the instructions in the model-specific TSD.

Disconnect the machine from power and ensure that power cannot be switched on again in error.

Wear protective gloves and goggles.

Chemically resistant protective gloves and tightly fitting protective goggles must be worn when working with process chemicals.

Do not lay cables or wires over sharp edges or hot surfaces.

Lay the hoses without kinking or straining them.

Opening the bottom service panel

The bottom service panel is secured with 2 bolts which each clip into the frame at the sides of the machine.

✚ Slide the service panel retaining bolts towards the center of the machine at the same time; see Fig. 1.

✚ Open the service panel slightly at the top and remove it.

Accessing the dispensing systems

Dispensing-system color coding, see Fig. 2.

1. Dispensing position 1 for detergent, color coding blue (S = Standard basic version)
2. Dispensing position 2, color coding green (O = Optional)
3. Dispensing position 3 for neutralization agent, color coding red (S = Standard basic version)
4. Dispensing position 4, color coding white (O = Optional)
5. Dispensing position 5, color coding yellow (O = Optional), without level monitoring
6. Dispensing position 6, color coding black (O = Optional), without level monitoring

✚ Hold the dispensing cover plate firmly and remove the retaining screws.

✚ Remove the dispensing cover plate.

✚ Remove the retaining screw (8) securing the left side of the dispenser mounting bracket (9); see Fig. 3.

✚ Open the dispenser mounting bracket; see Fig. 3.

Dispensing-pump installation

✚ Place the dispensing pump (2) in its appropriate position (7) and secure it with the retaining screws (1), (4) and (8); see Fig. 4.

✚ Unwind the connection plug for the dispenser pump (2) from the machine wiring harness and connect it at the bottom to the dispenser pump (2); see Fig. 4.

✚ Place the appropriate color-coded sticker on the dispensing pump.

Installing dispenser hoses

Depending on the machine variant, the dispenser drawer with the dispensing containers is located either under the dispensing system or in a neighboring cabinet. As a result, the routes for laying the dispenser hoses will vary slightly.

The dispenser hose from the dispenser probe to the dispensing pump has a larger diameter than the connection hose between the dispensing pump and flow meter, and the dispenser hose between the flow meter and dispenser connection to the wash chamber.

NOTE

Risk of leakage at the hose connectors due to incorrectly cut dispenser hoses.

✚ Cut dispenser hoses with a straight cut using a utility knife or similar tool.

✚ Do not use scissors, shears or similar to cut the dispenser hoses.

Connection plan legend, see Fig. 5.

1. Dispenser connections to the wash chamber
2. Narrow dispenser hose: Flow meter - Dispenser connection
3. Flow meter

4. Dispensing-system mounting bracket
5. Narrow dispenser hose (approx. 200 mm long): Dispensing pump - Flow meter
6. Dispensing pump
7. Wide dispenser hose: Dispenser probe - Dispensing pump
8. Fill-level monitor on dispenser probe
9. Dispenser probe
10. Dispenser drawer
11. Dispensing container

- ✂ Use the paths taken by the hoses already installed in the machine as a guide.
- ✂ Always lay the hoses with gentle curves without kinking or straining them.

Laying dispenser hose connections between the dispensing pump and flow meter

- ✂ Cut off a 200 mm piece of the narrow dispenser hose.
- ✂ Install a hose clip (6 mm) on each end of the 200 mm dispenser hose.
- ✂ Connect the dispenser hose to the narrow hose connector (6) on the dispenser pump (4) and secure it with a hose clip; see Fig. 6.
- ✂ Connect the other end of the hose to the hose connector (7) on the flow meter (1) on the side with the retaining screws (8) and secure it with a hose clip; see Fig. 6.
- ✂ Install a hose clip (6 mm) on one end of the remaining narrow dispenser hose.

Laying dispenser hose connections between the flow meter and the dispenser connection on the wash chamber

- ✂ Lay the remaining narrow dispenser hose to the machine dispenser connection (3) through the dispenser mounting bracket to the flow meter; see Fig. 6.
- ✂ Connect the dispenser hose to the open hose connector (3) on the flow meter (1) and secure it with a hose clip (6 mm); see Fig. 6.
- ✂ Bend out the lug on the metal plate to secure the flow meter.
- ✂ Secure the flow meter with the retaining screw (8); see Fig. 6.
- ✂ Lay the dispenser hose (2) along the dispenser mounting bracket (1) towards the hinged joint (3); see Fig. 7.
- ✂ Feed the dispenser hose at the hinged joint (3) through the opening; see Fig. 7.

NOTE

Risk of hoses being torn off when laid over the door holding bracket.

- ✂ Lay hoses under the door holding bracket.

- ✂ Lay the dispenser hose along the existing dispenser hoses towards the dispenser connections on the cabinet (4) and install a hose clip (6 mm) on the end; see Fig. 8.
- ✂ Lay the dispenser hose under the door holding bracket (1); see Fig. 8.
- ✂ Remove the dispenser connection cover (5); see Fig. 8.
- ✂ Install the dispenser hose end fully onto the dispenser connection (4) and secure it with a hose clip (6 mm); see Fig. 8.
- ✂ Fasten the dispenser hose with cable ties at suitable points. At the same time, ensure that the dispenser hoses are not closed.
- ✂ Place the appropriate color-coded sticker on the dispenser hose.

Laying dispenser hose connections between the suction probe and dispensing pump

- ✂ Install a hose clip (8 mm) on one end of the wide dispenser hose.

- ✚ Connect the dispenser hose to the hose connector (2) on the dispenser probe in the dispenser drawer and secure it with a hose clip (9 mm); see Fig. 9.
- ✚ Lay the dispenser hose from the dispenser probe (4) along the existing dispenser hoses (1) to the dispenser pump; see Fig. 10.
- ✚ Install a hose clip (8 mm) on the end of the dispenser hose.
- ✚ Connect the dispenser hose to the wide hose connector (9) on the dispenser pump (4) and secure it with a hose clip (8 mm); see Fig. 6.
- ✚ Place the appropriate color-coded stickers on the dispenser hose and dispenser probe.

Laying cable connections for the suction probe fill-level monitor

- ✚ Connect the fill-level monitor (1) connection plug to the suction probe; see Fig. 9.
- ✚ Lay the fill-level monitor connection cable along the appropriate dispenser hose (2); see Fig. 9.
- ✚ Secure the fill-level monitor connection cable with cable holders (3) to the appropriate dispenser hose (2); see Fig. 9.
- ✚ Connect the fill-level monitor connection plug (3) to the wiring harness plug connector in the machine; see Fig. 11.

Dispensing-system activation

- ✚ Log in as a service technician with administrator rights.
- ✚ Select the System menu.
- ✚ Select the Machine configuration menu.
- ✚ Select the Dispensing submenu.
- ✚ Select the appropriate dispensing system.
- ✚ Set the dispensing-system parameters.

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