A 503

de	Gebrauchsanweisung Wagen
da	Brugsanvisning Vogn
en	Operating instructions: Mobile unit
es	Instrucciones de manejo Carros
fi	Käyttöohje Vaunu
fr	Mode d'emploi Chariot
it	Istruzioni d'uso Carrelli
nl	Gebruiksaanwijzing Wagen
no	Bruksanvisning vogn
pt	Instruções de utilização Carro
sv	Bruksanvisning Vagn

de 4
da 15
en 26
es 37
fi 48
fr 59
it 70
nl 81
no 92
pt103
sv114

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Notes about these instructions	26
Intended use	27
Queries and technical problems.....	27
Items supplied	28
Load carriers	28
Disposal of the packing material	28
Optional accessories.....	29
Warning and Safety instructions	30
Medical devices.....	30
Laboratory glassware and utensils.....	30
Areas of application	31
To check before loading the machine and before starting a programme	31
Programme selection	31
Modules.....	32
Strip	32
Bar	33
Inserting modules.....	34
Removing modules.....	34
Inserting the A 606 2/3 level.....	35
Removing the A 606 2/3 level	35

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

- Select an option using the arrow buttons and save your choice with *OK*.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings  menu.

This mobile unit can be used to reprocess machine-reprocessible medical devices, laboratory glassware and utensils in a Miele washer-disinfector. Follow the operating instructions for your washer-disinfector as well as the instructions of the medical device or glassware and utensil manufacturer on how to reprocess their items by machine.

The A 503 mobile unit can be used with or without modules that are suitable for the intended purpose as well as with other inserts.

The mobile unit can be used in the following washer-disinfectors:


- PLW 8615
- PLW 8616
- PLW 8617
- PWD 8626
- PWD 8628

The washer-disinfector is generally referred to as “the machine” in these operating instructions. Reprocessible medical devices, laboratory glassware and utensils are referred to as “items” if they are not more closely defined.

Queries and technical problems

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your washer-disinfector or go to www.miele-professional.com.

Load carriers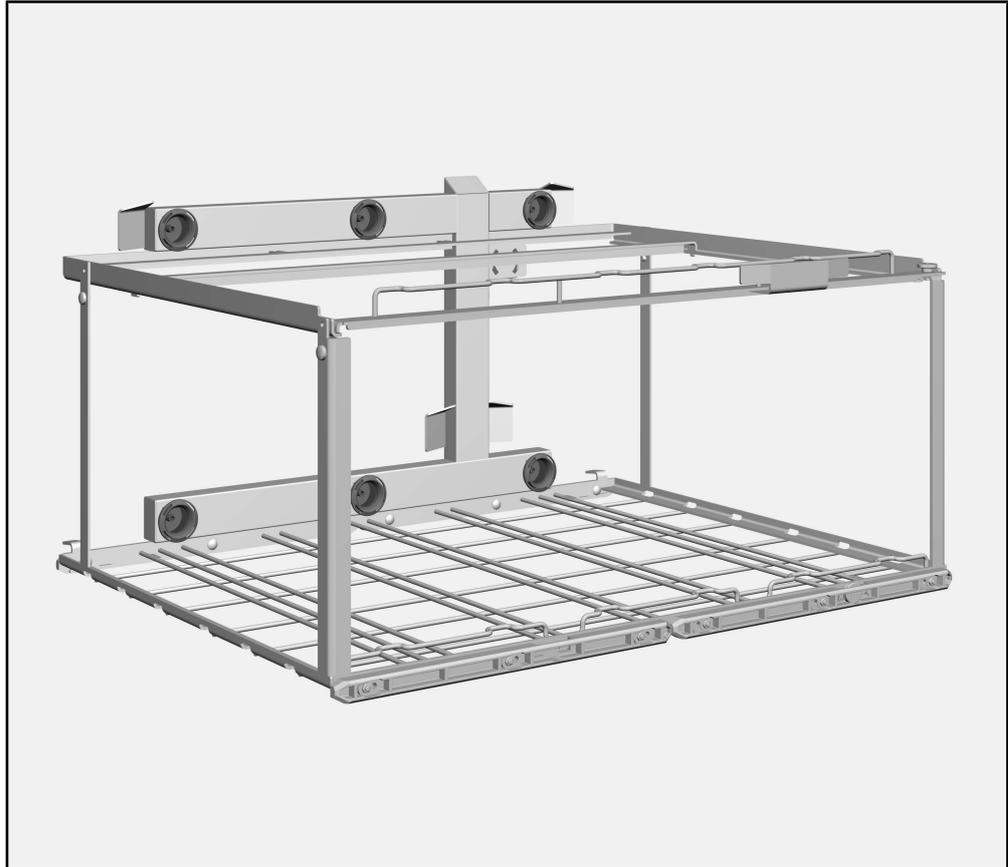

Mobile unit A 503 , height 433 mm, width 650 mm, depth 790 mm

Without inserted modules and bars in the top level, the loading dimensions are height 670 mm, width 565 mm, depth 770 mm.

Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in landfill sites.

Optional accessories

The following and other accessories are available to order from Miele:

Medical devices

- Module A 606, insertable 2/3 level with spray arm, height 97 mm, width 505 mm, depth 606 mm
- Module A 611 for hollow shafted instruments, height 187 mm, width 2454 mm, depth 614 mm
- Module A 618 for a maximum of 10 breathing bags or breathing masks, height 352 mm, width 142 mm, depth 614 mm
- Module A 619 for a maximum of 10 intubation materials, height 138 mm, width 142 mm, depth 614 mm

Laboratory glassware and utensils

- Module A 605, e.g. for vials, small test tubes or centrifuge tubes, height 140 mm, width 255 mm, depth 611 mm
- Module A 612 for large laboratory glassware, height 361 mm, width 255 mm, depth 614 mm
- Module A 613 for pipettes up to 450 mm long, height 195 mm, width 255 mm, depth 614 mm
- Module A 620 for a maximum of 10 items of laboratory glassware, height 113 mm, width 142 mm, depth 614 mm
- Module A 621 for a maximum of 20 items of laboratory glassware, height 113 mm, width 184 mm, depth 614 mm
- Module A 622 for a maximum of 36 items of laboratory glassware, height 113 mm, width 208 mm, depth 614 mm

en - Warning and Safety instructions

To avoid the risk of accidents and damage to this mobile unit please read these instructions carefully before using it for the first time.
Keep these instructions in a safe place where they are accessible to users at all times.

 Please also read the operating instructions for your washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ The mobile unit is approved solely for the applications specified in the “Areas of application” chapter of these operating instructions. Applications that go beyond those listed and use of the mobile unit with additional components from the Miele product range are described in the operating instructions for the components or are to be agreed with Miele.
- ▶ New mobile units and baskets must be washed in the washer-disinfector without a wash load prior to first use.
- ▶ Inspect all mobile units, baskets, modules and inserts daily as described in the “Maintenance” section in the Operating instructions for the washer-disinfector.
- ▶ Broken glass can result in serious injury when loading or unloading. Damaged glass items must not be reprocessed in the machine.

Medical devices

- ▶ The reprocessing programme used should include thermal disinfection, and the chemical agents used must be suitable for the material quality of the wash load.
- ▶ The last rinse, as a minimum, must be carried out with DI water.
- ▶ Visual checks should be undertaken to check the cleaning result. Additional protein analysis checks should also be carried out e.g. weekly, on a random sampling basis.

Laboratory glassware and utensils

- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

Miele cannot be held liable for damage caused by non-compliance with these Warning and Safety Instructions.

For validation purposes please make sure that you follow the loading instructions given on the template.

To check before loading the machine and before starting a programme

- Check that the mobile unit is correctly docked to the water path in the chamber.
- Are the inserted modules correctly connected to the water supply of the mobile unit?

Programme selection

When selecting a programme manually on the machine, the relevant mobile unit must be selected first. Pictograms representing the mobile units are shown on the display.

The pictogram of the mobile unit without inserted modules can be found on a bar on the front of the mobile unit.

Depending on the modules inserted, several pictograms are displayed for the A 503 mobile unit.

Once you have selected a pictogram, the list of assigned reprocessing programmes is displayed.

A 503 + module	Pictogram
A 503	
A 503 + A 605	
A 503 + A 606	
A 503 + A 606 + 4 x A 611	
A 503 + 6 x A 611	
A 503 + A 612	
A 503 + A 613	
A 503 + A 614	
A 503 + A 618 + A 619	
A 503 + A 620, A 621, A 622	

en - Areas of application

Risk of fire due to overheating

When the A 503 mobile unit is used without modules, all water connections are closed. During drying this will cause a build-up of hot air. The drying units can overheat and catch fire.

The A 503 mobile unit must only be used without modules in programmes that do not include a drying programme block.

Modules


Up to 6 modules can be accommodated in the A 503 mobile unit. Among other things, the possible number of modules depends on the height of the load in the mobile unit's bottom level.

Example:


For reprocessing large laboratory glassware, long pipettes or breathing bags, three modules are inserted into the bottom level.

For reprocessing small test tubes, vials or intubation material, three modules are inserted into each of the top and bottom levels.

Strip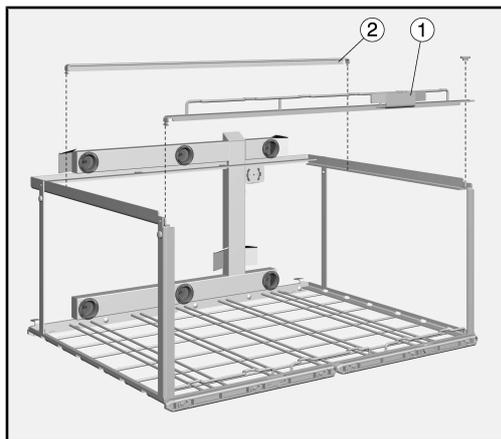

The strip ① on the mobile unit can be removed in order to insert particularly high modules into the bottom level.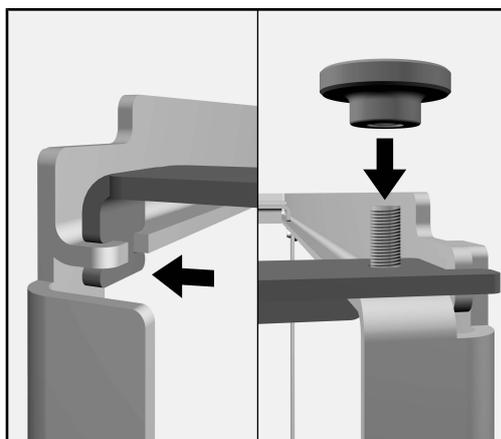

- Unscrew the knurled nuts on the right side of the strip.
- Lift the strip from the screw and pull it out of the guide on the left side.

- After the modules have been inserted into the bottom level, follow the instructions in reverse to reinsert the strip.


⚠ Damage due to bending.

Without the front strip, the mobile unit can bend during transport.

Only use the mobile unit with the strip in place.

Only carry the mobile unit with the strip in place.

Bar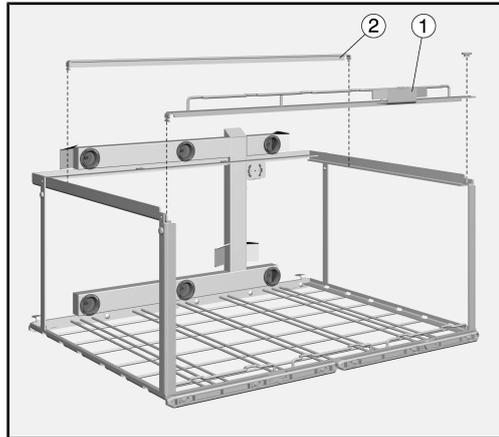

A bar ② is inserted into the top level of the mobile unit. The bar prevents modules from tipping into the bottom level when they are disconnected from the water connection.

The bar must be removed if a higher load is to be reprocessed in the bottom level.


- Pull the bar out of the recess on the right side.
- Pull the left side of the bar out of the guide.

If the load in the bottom level allows, the bar must always be inserted in the top level.


en - Areas of application

Inserting modules

A bar can be inserted into the top level. The bar supports modules that are disconnected from the water connection of the top level. This prevents modules from falling onto the load in the mobile unit's bottom level.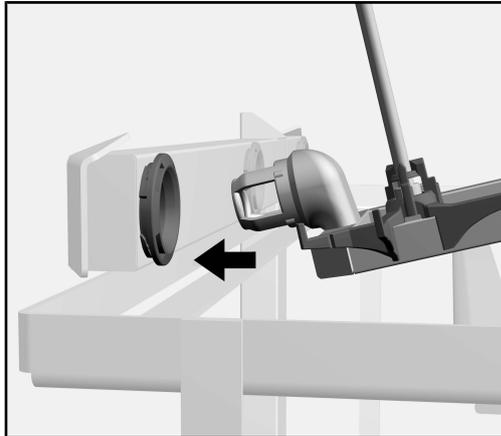

- Plug the module connector into the water connection of the mobile unit at a flat angle.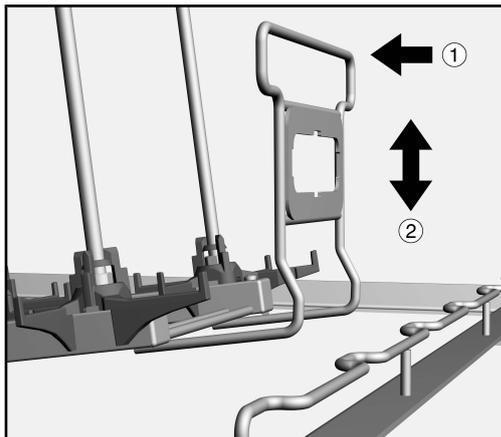

- Then push the module's retainer slightly towards the water inlet ① and lower the module ② to lock it in place behind the struts of the mobile unit.


Modules with holder rack

In the bottom level of the mobile unit, insert the modules in the two outer positions first so that they can be pushed behind the mobile unit's side bars. Then insert the module in the middle position.

Removing modules

- Push the module retainer slightly towards the water inlet ① and pull it out from behind the struts of the mobile unit ②.
- Disconnect the module from the water connection on the mobile unit and remove it.

Inserting the A 606 2/3 level

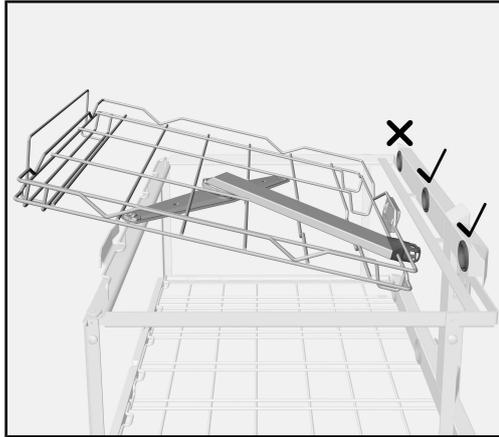The A 606 2/3 level can be inserted into the top level of the A 503 mobile unit.

The level connector can be connected to the middle or right water connection of the mobile unit. Another module can be inserted in addition to the 2/3 level.

To ensure that the spray arm of the 2/3 level is not blocked, the mobile unit must be used without the bar in the top level.

- Remove the bar from the top level of the mobile unit.
- Plug the level connector into the water connection of the mobile unit at a flat angle.
- Then push the level's retainer slightly towards the water inlet and lower the level to lock the module in place behind the struts of the mobile unit.

Removing the A 606 2/3 level

- Push the level's retainer slightly towards the water inlet and pull it out from behind the struts of the mobile unit.
- Disconnect the level from the mobile unit's water connection and remove it.

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