A 505

de	Gebrauchsanweisung Wagen
da	Brugsanvisning Vogn
el	Οδηγίες χρήσης ειδικό συρόμενο βαγονέτο με εγχυτήρες
en	Operating instructions: Mobile unit
es	Instrucciones de manejo Carros
fi	Käyttöohje Vaunu
fr	Mode d'emploi Chariot
it	Istruzioni d'uso Carrelli
nl	Gebruiksaanwijzing Wagen
no	Bruksanvisning vogn
pt	Instruções de utilização Carro
sv	Bruksanvisning Vagn

de	4
da	14
el	24
en	34
es	44
fi	54
fr	64
it	74
nl	84
no	94
pt	104
sv	114

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Notes about these instructions	34
Intended use	35
Queries and technical problems	35
Items supplied	36
Load carriers.....	36
Disposal of the packing material	36
Optional accessories.....	37
Warnings and safety notes	38
Medical devices.....	38
Laboratory glassware and utensils.....	38
Areas of application.....	39
To check before loading the machine and before starting a programme	39
Programme selection.....	39
Medical devices.....	40
Reprocessing hollow items	41
Laboratory glassware and utensils.....	42
Watch glasses	42
Small items	42

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

- Select an option using the arrow buttons and save your choice with *OK*.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings  menu.

This mobile unit can be used to reprocess machine-reprocessible medical devices, laboratory glassware and utensils in a Miele washer-disinfector. Follow the operating instructions for your washer-disinfector as well as the instructions of the medical device or glassware and utensil manufacturer on how to reprocess their items by machine.

The A 505 mobile unit can be loaded with inserts for baby bottles, DIN mesh trays or a variety of mesh trays and inserts for reprocessing medical devices, laboratory glassware and utensils.

The mobile unit can be used in the following washer-disinfectors:


- PLW 8615
- PLW 8616
- PLW 8617
- PWD 8626
- PWD 8628

The washer-disinfector is generally referred to as “the machine” in these operating instructions. Reprocessible medical devices are referred to as “load items” if they are not more closely defined.

Queries and technical problems

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your dishwasher or the washer-disinfector or go to www.miele-professional.com.

Load carriers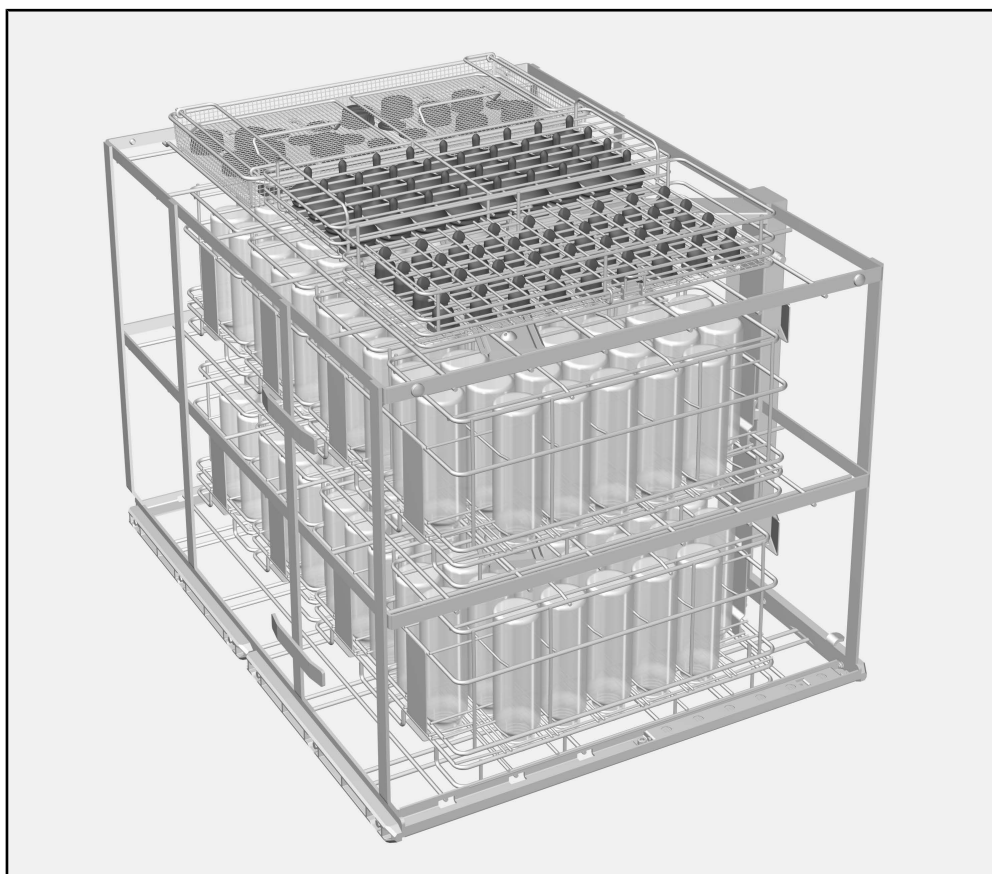

Mobile unit A 505 , height 524 mm, width 650 mm, depth 790 mm

A 505 *	Loading dimensions	
	[height x width x depth in mm]	
	Level 1	202 x 585 x 780
	Level 2	202 x 590 x 780
Level 3	132 x 590 x 780	

* Levels are numbered from the bottom up

Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

Optional accessories

The following and other accessories are available to order from Miele:

- E 142, DIN mesh tray with a 5 mm grid and 2 hinged handles, height 45/55 mm, width, 255 mm, depth 480 mm
- A 810 Cover net, metal frame with replaceable silicone net A 811, 428 x 220 mm

Medical devices

- E 492, insert with 9 holders for kidney dishes, spacing 49 mm, height 120 mm, width 256 mm, depth 474 mm
- E 135, insert for 19 baby bottles up to 250 ml, height 194 mm, width 192 mm, depth 447 mm
- E 364, insert for 36 wide-necked teats, height 77 mm, width 215 mm, depth 445 mm
- E 364, insert for 36 screw-on teats, height 63 mm, width 215 mm, depth 445 mm
- E 936, insert with 15 holders for lamp handles, height 90 mm, width 245 mm, depth 500 mm

Laboratory glassware and utensils

- E 104/1, insert for approx. 200 test tubes up to 12 x 105 mm, including lid A 13, height 132 (152) mm, width 200 mm, depth 320 mm.
- E 144, insert for 18 glass beakers up to 250 ml, 18 x 3 rods, height 131 mm, width 200 mm, depth 445 mm.
- E 402, insert for 44 watch glasses Ø 80 - 125 mm, spacing approx. 15 mm, height 53 mm, width 200 mm, depth 445 mm.
- E 969, insert for funnels, glass beakers, wide-necked glassware, etc., height 67 (122) mm, width 363 mm, depth 533 mm.

en - Warnings and safety notes

To avoid the risk of accidents and damage to this mobile unit please read these instructions carefully before using it for the first time. Keep these instructions in a safe place where they are accessible to users at all times.

 Please also read the operating instructions for your dishwasher or the washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ The mobile unit is approved solely for the applications specified in the “Areas of application” chapter of these operating instructions. Applications that go beyond those listed and use of the mobile unit with additional components from the Miele product range are described in the operating instructions for the components or are to be agreed with Miele.
- ▶ New mobile units and baskets must be washed in the washer-disinfector without a wash load prior to first use.
- ▶ Inspect all mobile units, baskets, modules and inserts daily as described in the “Maintenance” section in the Operating instructions for the washer-disinfector.
- ▶ Broken glass and ceramics can result in serious injury when loading or unloading. Damaged glass or ceramic load items must not be reprocessed in the machine.

Medical devices

- ▶ The reprocessing programme used should include thermal disinfection, and the process chemicals used must be suitable for the material quality of the load items.
- ▶ The last rinse, as a minimum, must be carried out with DI water.
- ▶ Visual checks should be undertaken to check the cleaning result. Additional protein analysis checks should also be carried out e.g. weekly, on a random sampling basis.

Laboratory glassware and utensils

- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

Miele cannot be held liable for damage or injury caused by non-compliance with the warning and safety notes.

For validation purposes please make sure that you follow the loading instructions given on the template.

To check before loading the machine and before starting a programme

- Are the spray arms able to rotate freely?
- Check spray arm jets for blockages.
- Check that the mobile unit is correctly docked to the water path in the chamber.

Programme selection

When selecting a programme manually on the machine, the relevant mobile unit must be selected first. Pictograms representing the mobile units are shown on the display.

The pictogram can be found in the bottom right on the front of the mobile unit, next to the RFID tag.

For the A 505 mobile unit, the pictogram below is displayed.


Once you have selected the pictogram, the list of assigned reprocessing programmes is displayed.

Mobile unit	Pictogram
A 505	


Reprocessing
baby bottles

Medical devices

Maximum load: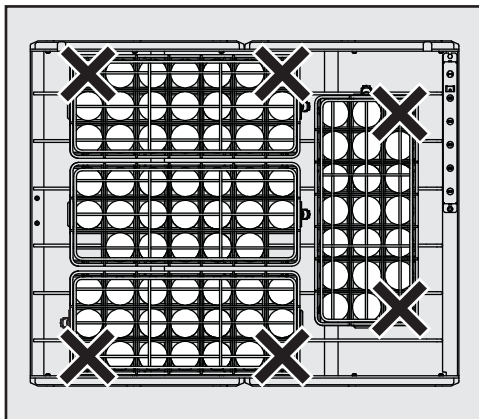


Level 1

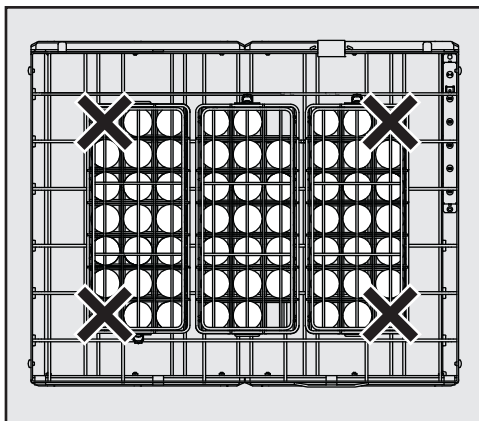Level 2 and 3

The two blocked corners in the inserts must be located over the positions marked with an X on the drawing. The jets of water cannot get into the bottles properly in these positions to clean them.

Reprocessing hollow items

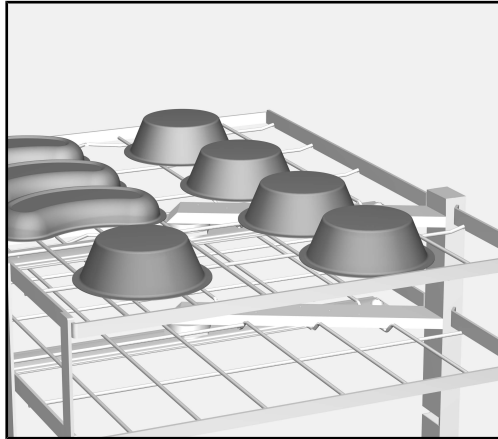- Hollow items such as beakers and kidney dishes should be loaded with the open side facing downwards. Lightweight items should be secured with a cover net, e.g. the A 810.
- Special inserts should be used if necessary for reprocessing hollow items. These inserts can be purchased from Miele.

The hollow items must not hinder the reprocessing of other items and must be secured so that they cannot be flipped over by the force of water during reprocessing.

If hollow items are placed with the open side facing upwards, suds solution can collect in them. This can spill and impair reprocessing results.

Laboratory glassware and utensils

Items with wide necks, e.g. glass beakers, wide-necked Erlenmeyer flasks and petri dishes, or cylindrical items, e.g. test tubes, can be cleaned and rinsed inside and out using rotating spray arms.

- Load wide-necked glassware with the open side facing downwards.

The laboratory glassware must not hinder the reprocessing of other items and must be secured so that it cannot be flipped over by the force of water during reprocessing. Secure light items with a cover net such as the A 810 or reprocess them in mesh trays with lids.

If laboratory glassware is placed with the open side facing upwards, suds solution can collect in it. This can spill and impair reprocessing results.

- Special inserts should be used if necessary for reprocessing laboratory glassware and utensils. Various inserts for use with different items can be purchased from Miele. Some examples are given below.

Watch glasses

Special inserts are available for reprocessing watch glasses, e.g. the E 403 for up to 105 watch glasses.

Small items

Reprocess small items such as stoppers, lids, spatulas, magnetic stirring rods, etc. in mesh trays with lids, e.g. the E 146. Alternatively a mesh tray with a sufficiently large cover net can be used.

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