APLW 868

APLW 869

de	Montageanweisung Ladestation und Akku
da	Monteringsanvisning Ladestation og batteri
en	Installation instructions Charging station and battery
fr	Notice de montage Base de chargement et batterie
it	Istruzioni di montaggio Stazione di ricarica e batteria
no	Monteringsveiledning Ladestasjon og batteri
cs	Montážní návod Nabíjecí stanice a akumulátor
fi	Asennusohje Latausasema ja akku
nl	Montagehandleiding Laadstation en accu
pl	Instrukcja montażu Stacja ładowająca i akumulator
sv	Monteringsanvisning Laddningsstation och batteri

de	4
da	12
en	20
fr	28
it	36
no	44
cs	52
fi	60
nl	68
pl	76
sv	84

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Notes about these instructions.....	20
Queries and technical problems.....	20
Warning and Safety instructions	21
Safety notes and warnings for the rechargeable battery and charger	21
Product description.....	22
APLW 868 charging station.....	22
Caring for the environment	23
Disposal of the packing material	23
Disposing of your old machine.....	23
Installation.....	24
APLW 868/APLW 869 installation	24
Technical data.....	26

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property.

Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

- Select an option using the arrow buttons and save your choice with OK.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings  menu.

Queries and technical problems

In the event of queries or technical problems, please contact Miele. Contact details can be found at the end of the operating instructions for your cleaning machine or at www.miele.com/professional.



Safety notes and warnings for the rechargeable battery and charger

- The rechargeable battery and charger must be used for the variable transfer trolley. For safety reasons, the transfer trolley must not be used with a rechargeable battery or charger produced by another manufacturer.
- If you are using a mains plug adapter, make sure that it is approved for continuous operation.
- Stop using the rechargeable battery immediately if you notice an unusual smell from it, if liquid is leaking from it, if it is significantly overheating or if it is discoloured or deformed. Dispose of the rechargeable battery immediately.
- Batteries will deplete over time. They can then leak corrosive fluids. Avoid contact with the eyes and skin. In the event of contact, rinse thoroughly with water and seek medical advice.
- Disposing of the rechargeable battery: remove the rechargeable battery from the transfer trolley. Insulate the metal contacts with adhesive tape to prevent short circuits. Dispose of the rechargeable battery in a municipal recycling facility. Do not dispose of the rechargeable battery with your household waste.
- Do not allow the rechargeable battery to come into contact with liquids.
- Do not cause a short circuit by bridging the contacts of the rechargeable battery, either unintentionally or deliberately.
- Do not dismantle the rechargeable battery.
- Do not drop or throw the rechargeable battery. For safety reasons, do not continue to use a rechargeable battery which has been damaged.
- Keep the rechargeable battery away from naked flames or other heat sources. Do not heat the rechargeable battery. Do not expose the rechargeable battery to direct sunlight.

APLW 868 charging station

The external charger has been developed specifically for use as a wall-mounted charging station for the APLW 869 battery. It is a switched-mode power supply charger (SMPS) which makes the battery charging process more efficient.

The general operation of the APLW 869 battery is described in the operating instructions for the ATT 86.

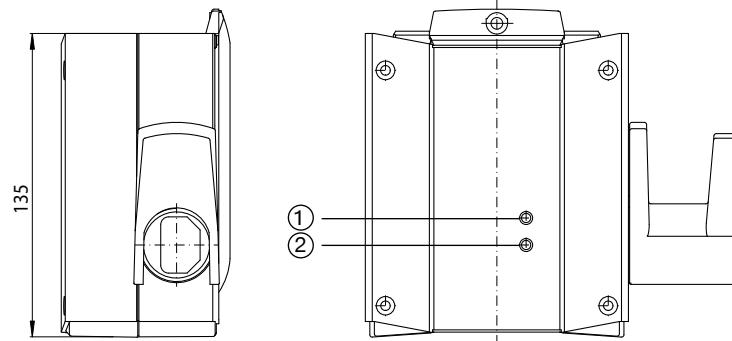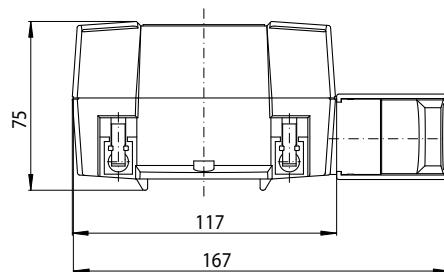① Green LED (lights up when the power supply is connected)

② Yellow LED (lights up when the battery is charging)

The battery module takes less time to charge with the charger.

Risk posed by conversions or modifications

Please note that the charging station and the battery must only be used for applications described in these operating instructions. Alterations or conversion of the charging station, or using it for purposes other than those for which it was intended, are not permitted and could be dangerous.

Miele cannot be held liable for damage caused by improper or incorrect use or operation.

Initial commissioning will be undertaken by the Miele Customer Service Department or an authorised technician, who will provide you with operating instructions.

Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in landfill sites.

Disposing of your old machine

Electrical and electronic machines often contain valuable materials. They also contain materials which, if handled or disposed of incorrectly, could be potentially hazardous to human health and to the environment. They are, however, essential for the correct functioning of your machine. Please do not therefore dispose of it with your household waste.

Please dispose of it at your local community waste collection / recycling centre or contact your dealer for advice.


Ensure that it presents no danger to children while being stored for disposal.

APLW 868/APLW 869 installation

The APLW 868 charging station with wall bracket must be installed vertically upright.

The screws for fastening the wall bracket to the wall are not included in the scope of delivery.

- Select an accessible position on the wall for the charging station.
- Use the diagram to position the screws (not included in the scope of delivery) at the correct intervals on the wall (140 mm), no. ①.
- Use the mounting screw (included in the scope of delivery) to attach the charger to the wall bracket, no. ②.
- Hook the charging station (now attached to the wall bracket) onto the screws on the wall by the lugs, no. ①.

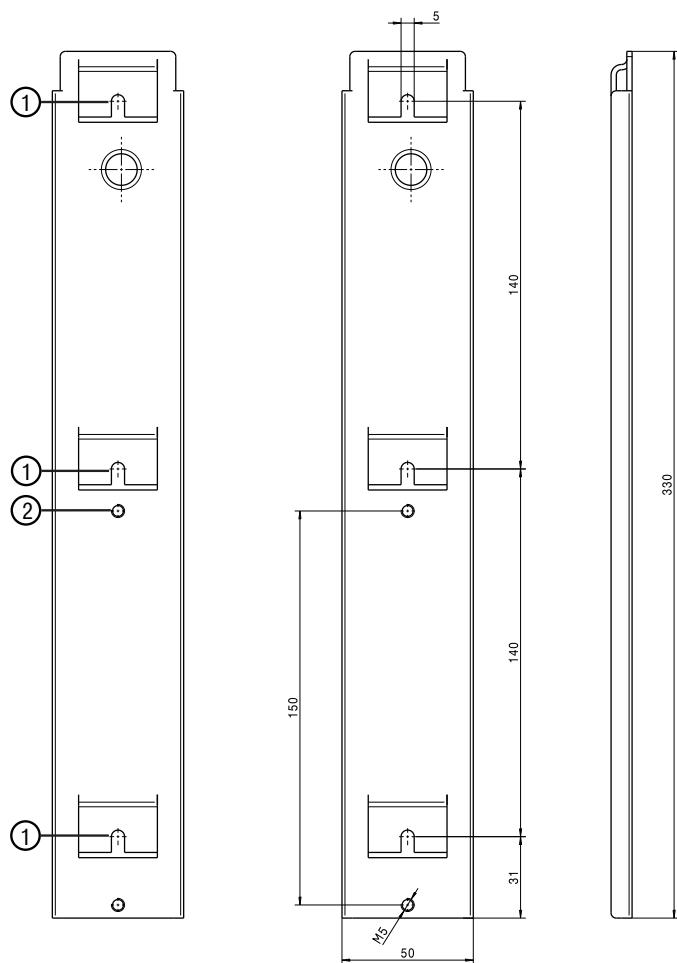Tighten the mounting screws for the charger to a maximum tightening torque of 1 Nm.

- Plug the APLW 869 into the charging station, positioned vertically upright.

Charging is automatic. Once the battery is fully charged, all of the LEDs go out.

- Unplug the battery from the charging station.

en - Technical data

APLW 868

Height	135 mm
Width	167 mm
Depth	75 mm
Charging duration for the battery	Approx. 5 hours
Mains voltage	100–240 V/50–60 Hz
Charging current	Up to 650 mA
Rated current consumption	Max. 400 mA
Electricity consumption (standby)	Max. 2.5 W
Electricity consumption (charging)	Max. 19 W
Ambient temperature	5 °C to 40 °C
Storage temperature	-10 °C to 50 °C
Relative humidity non-condensing	20 % for temperatures up to 80 %
Air pressure	700 to 1060 hPa
Maximum height above sea level	3000 m
Protection rating	IPX5 (vertical installation only)
Approvals	IEC60601-1, ANSI/AAMI ES60601 and CAN/CSA-22.2 No 60601-1

APLW 869

Height	239 mm
Width	117 mm
Depth	75 mm
Switch-on duration	10 % or 2 minutes of continuous operation followed by an 18-minute pause
Ambient temperature	5 °C to 40 °C
Storage temperature	-15 °C to 40 °C
Inflammable materials	UL94-V0
Protection rating	IPX5 (vertical installation only)
Approvals	IEC60601-1, ANSI/AAMI ES60601 and CAN/CSA-22.2 No 60601-1

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Miele Werk Bürmoos GmbH