# A 306/1

|           |                                         |
|-----------|-----------------------------------------|
| <b>da</b> | Brugsanvisning Injektormodul            |
| <b>de</b> | Gebrauchsanweisung Injektormodul        |
| <b>en</b> | Operating instructions Injector module  |
| <b>es</b> | Instrucciones de manejo Módulo inyector |
| <b>fi</b> | Käyttöohje Suorasuihkumoduuli           |
| <b>fr</b> | Mode d'emploi Module à injection        |
| <b>hr</b> | Upute za uporabu Injektorski modul      |
| <b>it</b> | Istruzioni d'uso Modulo iniettore       |
| <b>nl</b> | Gebruiksaanwijzing injectormodule       |
| <b>no</b> | Bruksanvisning injektormodul            |
| <b>pt</b> | Instruções de utilização Módulo injetor |
| <b>sl</b> | Navodila za uporabo Injektorski modul   |
| <b>sv</b> | Bruksanvisning injektormodul            |

|                 |     |
|-----------------|-----|
| <b>da</b> ..... | 4   |
| <b>de</b> ..... | 15  |
| <b>en</b> ..... | 26  |
| <b>es</b> ..... | 37  |
| <b>fi</b> ..... | 48  |
| <b>fr</b> ..... | 59  |
| <b>hr</b> ..... | 70  |
| <b>it</b> ..... | 81  |
| <b>nl</b> ..... | 92  |
| <b>no</b> ..... | 103 |
| <b>pt</b> ..... | 114 |
| <b>sl</b> ..... | 125 |
| <b>sv</b> ..... | 136 |

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|                                                                           |           |
|---------------------------------------------------------------------------|-----------|
| <b>Notes about these instructions .....</b>                               | <b>26</b> |
| Queries and technical problems .....                                      | 26        |
| <b>Appropriate use .....</b>                                              | <b>27</b> |
| <b>Items supplied .....</b>                                               | <b>28</b> |
| Scope of delivery A 306/1 .....                                           | 28        |
| Disposal of the packing material .....                                    | 28        |
| <b>Warnings and safety notes .....</b>                                    | <b>29</b> |
| <b>Installation .....</b>                                                 | <b>30</b> |
| Injector nozzles .....                                                    | 30        |
| Tools required .....                                                      | 30        |
| Screwing in injector nozzles .....                                        | 30        |
| Mounting the holder frames .....                                          | 31        |
| <b>Areas of application .....</b>                                         | <b>32</b> |
| Injector modules .....                                                    | 32        |
| Inserting the module .....                                                | 32        |
| Removing an injector module .....                                         | 32        |
| Preparing the load .....                                                  | 33        |
| Measuring cylinder, tall .....                                            | 34        |
| Measuring cylinder, short .....                                           | 34        |
| To check before loading the machine and before starting a programme ..... | 35        |
| Reprocessing .....                                                        | 35        |

### Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

### Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

### Additional information and comments

Additional information and comments are contained in a simple frame.

### Operating steps

Operating steps are indicated by a black square bullet point.

#### Example:

- Select an option using the arrow buttons and save your choice with *OK*.

### Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

#### Example:

Settings  menu.

### Queries and technical problems

In the event of queries or technical problems, please contact Miele. Contact details can be found at the end of the operating instructions for your cleaning machine or at [www.miele.com/professional](http://www.miele.com/professional).

This module can be used to reprocess machine-reprocessable laboratory glassware and utensils in a Miele washer-disinfector for laboratory glassware and utensils. Follow the operating instructions for your washer-disinfector as well as the instructions of the glassware and utensil manufacturer on how to reprocess their items by machine.

The A 306/1 module is designed for reprocessing measuring cylinders with a volume of 1000 ml and 2000 ml. The module can hold both short and tall construction type measuring cylinders up to a max. length of 500 mm.

Because of its height, this module can only be used in a lower basket. It is not possible to use an upper basket at the same time.

This module is suitable for use in the following:


- Lower basket A 150
- APLW 106 basket

Special reprocessing programmes are required for reprocessing laboratory glassware and utensils by machine. These programmes are installed ex-works on the following washer-disinfectors or can be installed, created or activated by the Miele Customer Service Department where necessary:

- |               |              |
|---------------|--------------|
| - PLW 8604    | - PG 8504    |
| - PLW 8636    | - PG 8583    |
| - PLW 8683    | - PG 8583 CD |
| - PLW 8683 CD | - PG 8593    |
| - PLW 8693    |              |
| - LAB 5010    |              |
| - LAB 5011    |              |
| - LAB 5020    |              |

The washer-disinfector is generally referred to as “the machine” in these operating instructions. Reprocessable laboratory glassware and utensils are referred to as “items” if they are not more closely defined.

### Scope of delivery A 306/1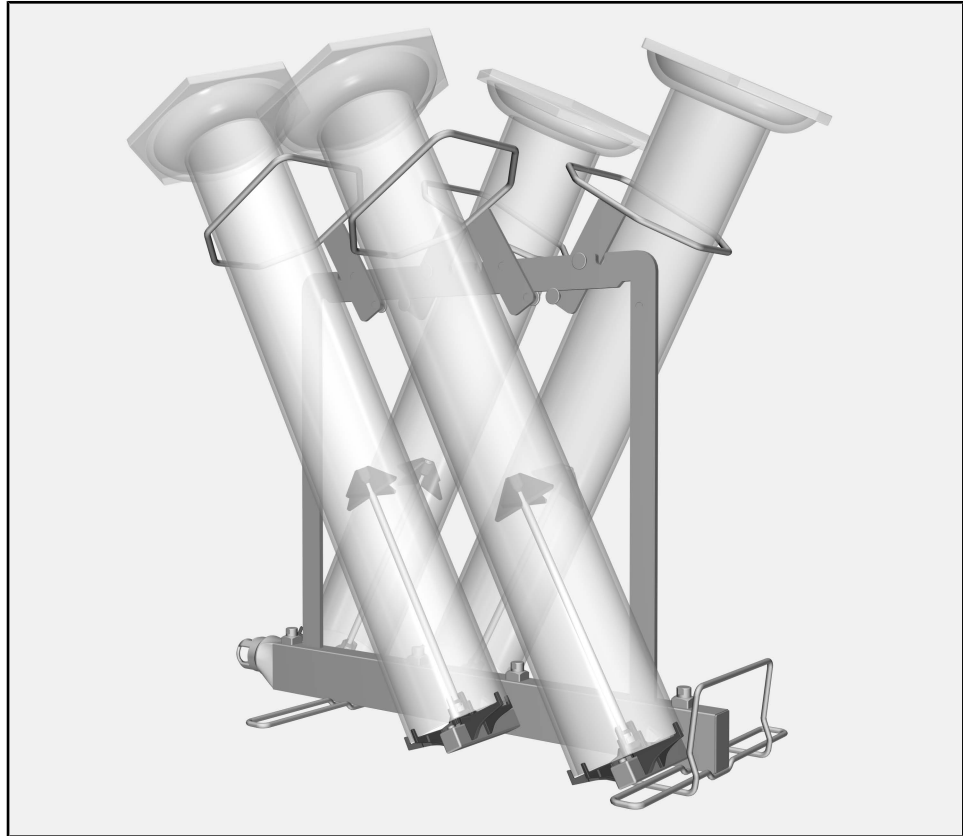

- Injector module A 306/1
  - Height 418 mm, width 235 mm, depth 471 mm
  - Max. vertical clearance (from injector nozzle supports) 500 mm

### Items supplied

- 4 injector nozzles with plastic supports for reprocessing measuring cylinders, length 217 mm, Ø 6 mm
- 4 holder frames with coating; for measuring cylinders up to max. Ø 110 mm

### Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

To avoid the risk of accidents and damage to this module please read these instructions carefully before using it for the first time. Keep these instructions in a safe place where they are accessible to users at all times.

 Please also read the operating instructions for your dishwasher or the washer-disinfector and pay particular attention to the Warning and Safety instructions.


- ▶ The module is approved solely for the applications specified in the “Areas of application” chapter of these operating instructions. Components such as nozzles may only be replaced with Miele components or genuine original spare parts.
- ▶ New load carriers must be cleaned in the washer-disinfector without a load prior to first use.
- ▶ Inspect all mobile units, baskets, modules and inserts as described in the “Maintenance” section of the operating instructions for your machine.
- ▶ Only items which have been declared by their manufacturer as suitable for machine reprocessing may be processed. The manufacturer's specific reprocessing instructions must be observed.
- ▶ Broken glass can result in serious injury when loading or unloading. Broken glass items must not be processed in the machine.
- ▶ Only place empty, unloaded modules in baskets. Check that they are correctly engaged before loading them. Modules must be completely emptied before taking them out. Placing loaded modules in baskets or removing them can cause damage to the wash load and cause injury should glassware be broken.
- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

### Injector nozzles

#### Tools required

- 9 mm Spanner (WAF 9)


#### Screwing in injector nozzles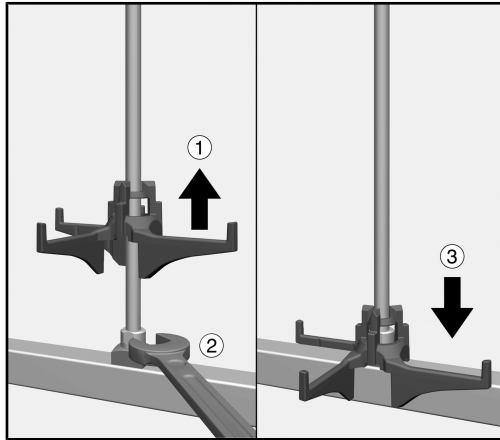

- Lift up the injector nozzle support and screw the nozzle into the required position ①.
- Use the spanner to tighten the injector nozzle ②.
- Slide the support downwards until it rests on the water supply pipe ③.

## Mounting the holder frames

The holder frames are mounted above the nozzles on the upper cross strut of the module, with 2 frames mounted on each side. The slanted corners of the holder frames centre the measuring cylinders during re-processing. After mounting they must point downwards.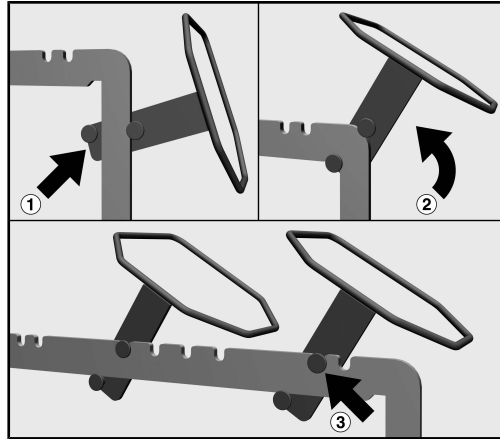


- Position the holder frame on one of the side struts of the module. The pins of the holder frame must reach around the strut of the module. The frame must be facing outwards and the slanted corners must be facing downwards ①.
- Move the holder frame upwards onto the upper cross strut of the module ②.
- Hook the holder frame into one of the notches ③.
- The frames must be mounted on each side in turn. Repeat the process on the other side.

### Injector modules

⚠ Only place empty, unloaded modules in baskets. Check that they are correctly engaged before loading them. Modules must be completely emptied before taking them out. Placing loaded modules in baskets or removing them can cause damage to the wash load and cause injury should glassware be broken.

#### Inserting the module

Ensure that the module is correctly orientated when inserting it. For example, the nozzles must point upwards.

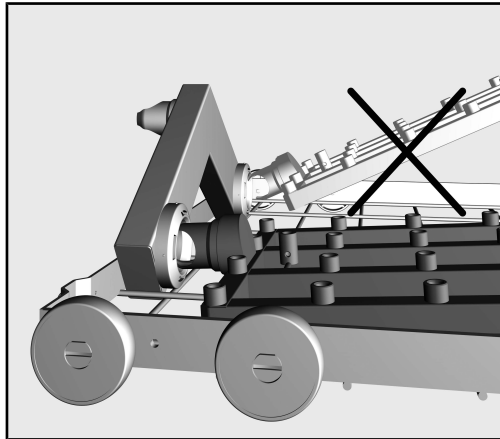Example of a flat insertion angle

- Insert the module into the water connection with the connector at a flat angle.
- Then engage the module by pushing it down onto the struts of the basket.

#### Removing an injector module

⚠ Do not hold the basket by the injector module when removing it from the cleaning machine. The module could come loose or the water connection could be damaged.

- Release the module from the locking mechanism and pull it out of the water connection at a flat angle. If the module is removed at too steep an angle, the lever action may damage the connection point and the connector.

### Preparing the load

⚠ Damage to the measuring cylinders.


If the measuring cylinders touch each other during reprocessing or rest against the wash cabinet walls, damage can occur at the points of contact. For example, the items may become scratched or the glass may break.

Arrange the measuring cylinders so that they are not touching each other or resting against the wash cabinet walls.

To prevent damage, adjust the position of the measuring cylinders as required by moving the holder frames or the plastic supports on the nozzles or change the load arrangement of the module.

Plastic supporting feet can be removed as required and reprocessed separately.

Long measuring cylinders or measuring cylinders with extra-wide supporting feet must not be reprocessed in the cleaning machine.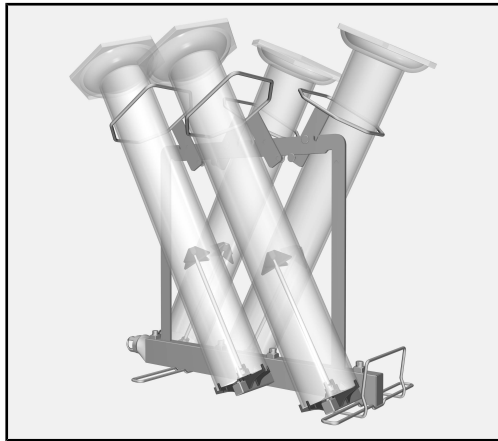


- Guide the measuring cylinders through the holder frames and position the cylinders on the injector nozzles.

Depending on the construction type of the measuring cylinder, the following points must be observed.

## en - Areas of application


---

### Measuring cylinder, tall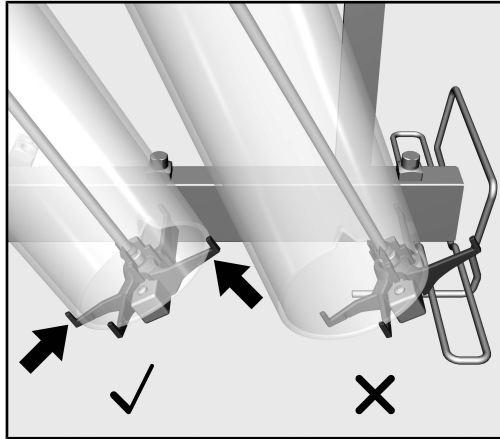

The openings of the measuring cylinders must rest completely on the injector nozzle supports. The cylinders must not touch the module or the injector nozzles in order to prevent damage such as glass breakage during reprocessing. Move the holder frames as required so that the openings rest fully on the supports.

### Measuring cylinder, short

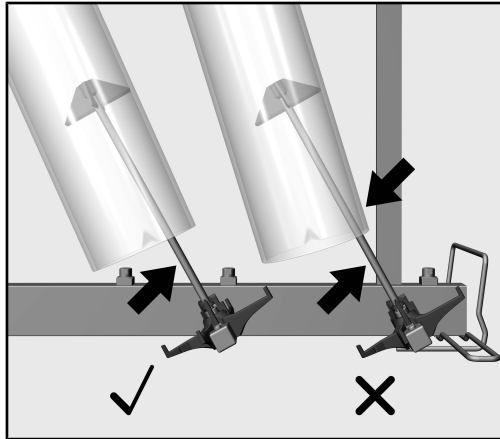The measuring cylinders must rest on the plastic supports of the injector nozzles. However, they must not be allowed to touch the nozzles themselves. If the measuring cylinders rest on the nozzles during reprocessing, damage such as glass breakage may occur. Move the holder frames or the plastic supports on the nozzles as required in order to adjust the positions of the measuring cylinders.

### To check before loading the machine and before starting a programme

- Check that fittings such as irrigation sleeves and injector nozzles are securely screwed into position.

⚠ Make sure that all screw connectors are fitted with injector nozzles, adapters, irrigation sleeves or blind stoppers to ensure that all fittings in use are supplied with sufficient standardised pressure. Damaged fittings such as injector nozzles, adapters and irrigation sleeves must not be used.

Fittings not equipped with wash items do not have to be replaced blind stoppers.

- Check that the baskets and modules are correctly connected to the water supply.
- Make sure that the measuring cylinders are not touching each other or resting against the wash cabinet walls.

### Reprocessing

- Slowly and carefully insert the lower basket with the loaded module into the cleaning machine.  
Make sure that the supporting feet of the measuring cylinders do not stick out into the door opening or the door seal or collide with the wash cabinet walls.
- Carefully close the cabinet door.  
Protruding measuring cylinders which extend into the door area may be damaged when the door is closed or during reprocessing.

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