APLW 511

de	Gebrauchsanweisung Wagen
da	Brugsanvisning Vogn
el	Οδηγίες χρήσης ειδικό συρόμενο βαγονέτο με εγχυτήρες
en	Operating instructions: Mobile unit
es	Instrucciones de manejo Carros
fi	Käyttöohje Vaunu
fr	Mode d'emploi Chariot
it	Istruzioni d'uso Carrelli
nl	Gebruiksaanwijzing Wagen
no	Bruksanvisning vogn
pt	Instruções de utilização Carro
sv	Bruksanvisning Vagn

de	4
da	16
el	28
en	40
es	52
fi	64
fr	76
it	88
nl	100
no	112
pt	124
sv	136

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Notes about these instructions	40
Appropriate use	41
Queries and technical problems	41
Items supplied	42
Load carriers.....	42
Disposal of the packing material	42
Optional accessories.....	43
Warnings and safety notes	44
Areas of application.....	45
To check before loading the machine and before starting a programme	45
Reprocessing programmes	45
Modules	45
Removable bars	46
Supports	47
Inserting modules.....	48
Removing modules.....	48
Modules with holder rack.....	49
Inserting the A 606 2/3 level	50
Removing the A 606 2/3 level.....	50

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

- Select an option using the arrow buttons and save your choice with *OK*.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings  menu.

Re-usable laboratory glassware and laboratory utensils can be reprocessed using this mobile unit in a Miele laboratory glassware washer-disinfector. Follow the instructions for your washer-disinfector as well as the instructions of the glassware manufacturer on how to reprocess their items by machine.

The APLW 511 mobile unit can be used with or without modules that are suitable for the intended purpose as well as with other inserts.

The mobile unit can be used in the following washer-disinfectors:

- PLW 8615
- PLW 8616
- PLW 8617

The washer-disinfector is generally referred to as “the machine” in these operating instructions. Reprocessable laboratory glassware and utensils are referred to as “items” if they are not more closely defined.

Queries and technical problems

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your dishwasher or the washer-disinfector or go to www.miele-professional.com.

Load carriers

APLW 511 mobile unit, height 534 mm, width 650 mm, depth 790 mm

Without modules inserted and without supports on the middle and top levels, the loading dimensions are height 670 mm, width 565 mm, depth 770 mm.

Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

Optional accessories

The following and other accessories are available to order from Miele:

- Module A 605, e.g. for vials, small test tubes or centrifuge tubes, height 140 mm, width 255 mm, depth 611 mm
- Module A 606, insertable 2/3 level with spray arm, height 97 mm, width 505 mm, depth 606 mm
- Module A 613 for pipettes up to 450 mm long, height 195 mm, width 255 mm, depth 614 mm
- Module A 620 for a maximum of 10 items of laboratory glassware, height 113 mm, width 142 mm, depth 614 mm
- Module A 621 for a maximum of 20 items of laboratory glassware, height 113 mm, width 184 mm, depth 614 mm
- Module A 622 for a maximum of 36 items of laboratory glassware, height 113 mm, width 208 mm, depth 614 mm

en - Warnings and safety notes

To avoid the risk of accidents and damage to this mobile unit please read these instructions carefully before using it for the first time. Keep these instructions in a safe place where they are accessible to users at all times.

 Please also read the operating instructions for your dishwasher or the washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ The mobile unit is approved solely for the applications specified in the “Areas of application” chapter of these operating instructions. Applications that go beyond those listed and use of the mobile unit with additional components from the Miele product range are described in the operating instructions for the components or are to be agreed with Miele.
- ▶ New mobile units and baskets must be washed in the washer-disinfector without a wash load prior to first use.
- ▶ Inspect all mobile units, baskets, modules and inserts daily as described in the “Maintenance” section in the Operating instructions for the washer-disinfector.
- ▶ Broken glass and ceramics can result in serious injury when loading or unloading. Damaged glass or ceramic load items must not be reprocessed in the machine.
- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

Miele cannot be held liable for damage or injury caused by non-compliance with the warning and safety notes.

To check before loading the machine and before starting a programme

- Check that the mobile unit is correctly docked to the water path in the chamber.
- Are the inserted modules correctly connected to the water supply of the mobile unit?

Risk of fire due to overheating

When the APLW 511 mobile unit is used without modules, all water connections are closed. During drying this will cause a build-up of hot air. The drying units can overheat and catch fire.

The APLW 511 mobile unit must only be used without modules in programmes that do not include a drying programme block.

Reprocessing programmes

New reprocessing programmes adapted for the mobile unit's loading templates are available for the APLW 511.

The new reprocessing programmes are available from the manufacturer of the cleaning machine.

The factory default setting assigns the standard programmes for the PLW 8615, PLW 8616 and PLW 8617 cleaning machines to the A 503 mobile unit. The standard programs must be copied and adapted for the APLW 511 mobile unit.

Modules

Up to 9 modules on 3 levels can be accommodated in the APLW 511 mobile unit. Among other things, the possible number of modules depends on the height of the load items on the bottom and middle levels of the mobile unit.

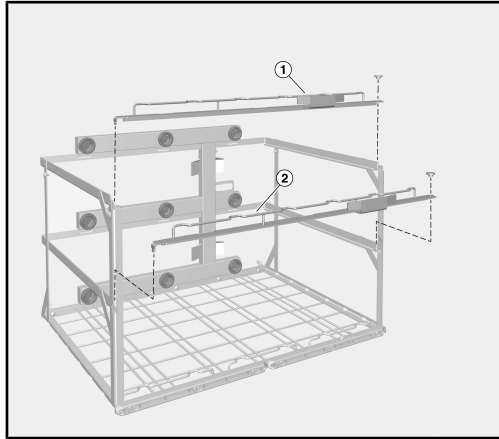
Examples:

For reprocessing long pipettes, 3 modules are placed on the bottom level.

For reprocessing small items of laboratory glassware, 3 modules can be placed on each level.


For reprocessing phials, one A 605 module can be placed on each level. The three A 605 modules must be arranged diagonally, not one on top of the other. A 620, A 621 or A 622 modules can be added in the available space on each level.

Removable bars

The bars ① and ② on the mobile unit can be removed to accommodate modules with tall load items on the middle or bottom level.

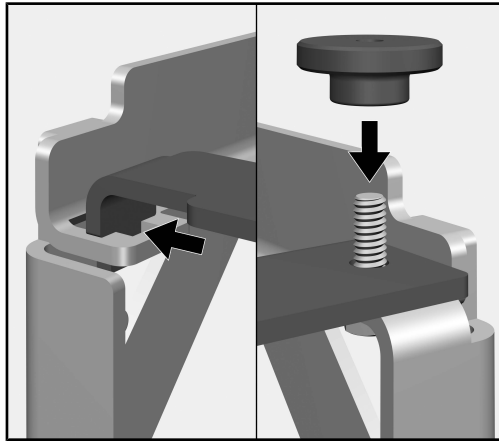To do this:

- Unscrew the knurled nuts on the right side of the bar.
- Lift the bar from the screw, slide it to the left and pull it out of the guide on the left side.
- After the modules have been removed, follow the instructions in reverse to refit the bar.

The bar ② must be fitted to the mobile unit in order to insert modules on the middle level. Therefore, the modules on the bottom level must be loaded and the bar ② must be fitted first.

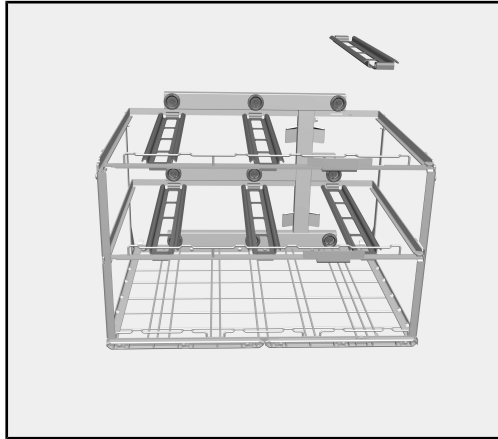
⚠ Damage due to bending.

Without the front bars, the mobile unit can bend during transport.


Only use the mobile unit with the bars in place.

Only carry the mobile unit with the bars in place.

Supports

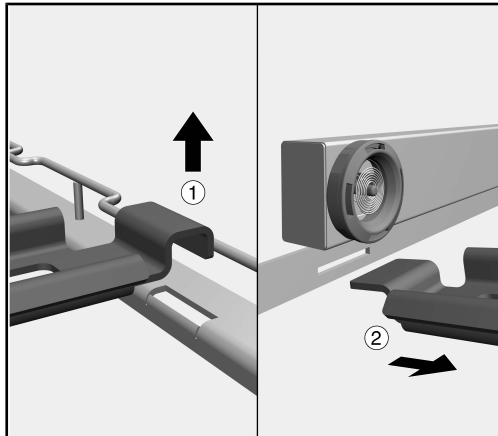
Supports are fitted on the middle and top levels of the mobile unit. The supports prevent modules from tipping down when they are disconnected from the water connection. The supports must be removed if taller load items need be reprocessed on the bottom or middle level. To do this:


- Pull the support up and out of the recess in the bar ①.
- Pull the support out of the bar underneath the water inlet ②.

The supports should always be fitted whenever the load items on the lower levels allow.

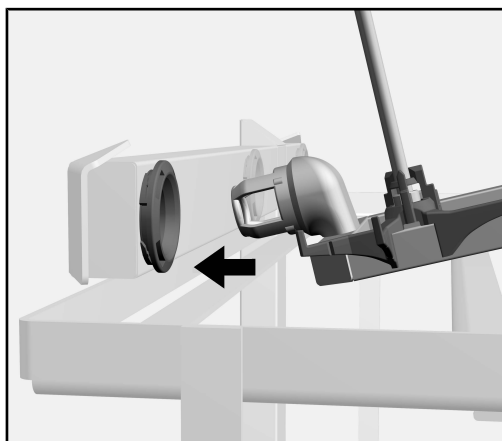Inserting modules

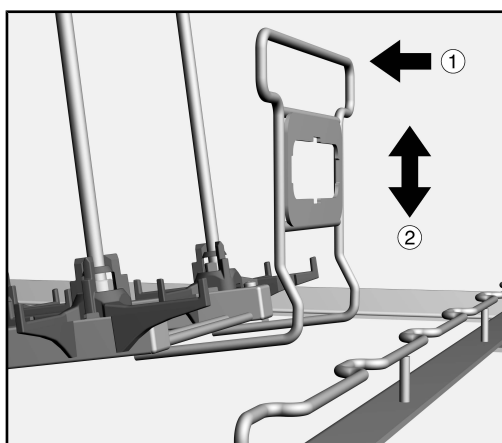
- Attach the back end of the module to the support and push towards the water connection.
- Plug the module connector into the water connection of the mobile unit at a flat angle.

- Then push the module retainer slightly towards the water inlet ① and lower the module ② to lock it in place behind the struts of the mobile unit.
- Push the module retainer slightly towards the water inlet ① and pull it out from behind the struts of the mobile unit ②.
- Disconnect the module from the water connection on the mobile unit and remove it.

Removing modules

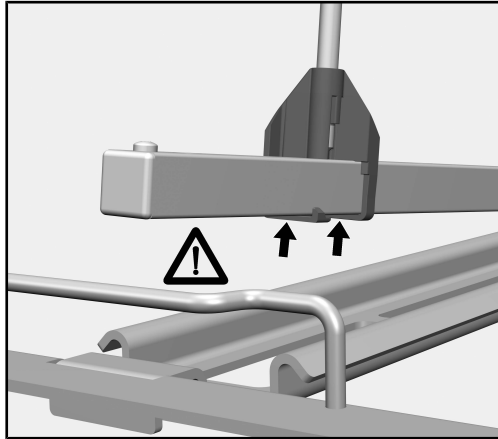Modules with holder rack

 Risk of damage.

The holders on the holder rack can break off if they bang against the struts on the bar of the mobile unit.

Inserting modules:

Do not lower the module until the holders are located behind the struts on the mobile unit.

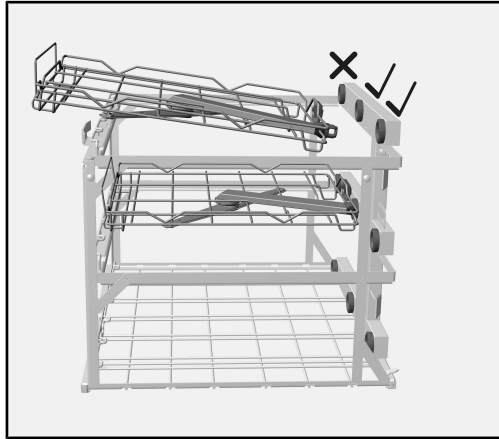
Removing modules:

Lift the module until the holders are located above the struts on the mobile unit.

In the bottom and middle levels of the mobile unit, insert the modules in the two outer positions first so that they can be pushed behind the bars on the side of the mobile unit. Then insert the module in the middle position.

en - Areas of application

Inserting the A 606 2/3 level

The A 606 2/3 level can be inserted in the middle and top levels of the APLW 511 mobile unit.

The 2/3 level connector can be connected to the middle or right water connection on the mobile unit.

To avoid blocking the spray arm of the 2/3 level, the supports must be removed at the relevant positions.

- Remove the relevant supports from the middle or top level of the mobile unit.
- Plug the 2/3 level connector into the water connection on the mobile unit at a flat angle.
- Then push the retainer on the 2/3 level slightly towards the water inlet and lower the 2/3 level to lock the module in place behind the struts of the mobile unit.

Another module with injector nozzles, e.g. the A 620, can be inserted in addition to the 2/3 level. Alternatively, a module for tall load items can be inserted in the bottom level, e.g. the A 613 for pipettes.

Removing the A 606 2/3 level

- Push the level's retainer slightly towards the water inlet and pull it out from behind the struts of the mobile unit.
- Disconnect the level from the mobile unit's water connection and remove it.

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