


Installation plan:

PG 8592 - PG 8593

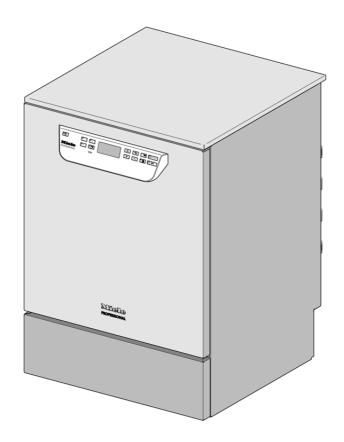

Mat. no. Version 10 344 170 00 **US**

Technical data sheet



PG 8592 - 8593

Legend:



Circled, bold-type abbreviations: Connection required



Abbreviations surrounded by broken circle: Connection optional or required, depending on model

PG 8592:

EL	Electrical connection PG 8592	Voltage (default) Rated load	2 AC 208V 60 Hz 6.2 kW	2 AC 208V 60 Hz 6.2 kW
		Fuse rating	2 pole 30 Amp	2 pole 30 Amp
		Minimum cross-section	3 × AWG 10	3 × 5.2 mm ²
		Length of electrical cord (UL 921) with plug	approx. 5' 9"	approx. 1.8 m
		Electrical connection	NEMA L6-30	NEMA L6-30
F	Electrical connection PG 8592	Voltage (convertible)	3 AC 208V 60 Hz	3 AC 208V 60 Hz
		Rated load	6.2 kW	6.2 kW
		Fuse rating	3 pole 20 Amp	3 pole 20 Amp
		Minimum cross-section	4 × AWG 12	4 × 3.3 mm ²
		Length of electrical cord (UL 921) with plug	approx. 5' 9"	approx. 1.8 m
		Electrical connection	NEMA L15-20	NEMA L15-20

PG 8593:

EL	Electrical connection PG 8593	Voltage (default) Rated load Fuse rating Minimum cross-section Length of electrical cord (UL 921) with plug Electrical connection	3 AC 208V 60 Hz 6.2 kW 3 pole 20 Amp 4 × AWG 12 approx. 5' 9" NEMA L15-20	3 AC 208V 60 Hz 6.2 kW 3 pole 20 Amp 4 × 3.3 mm ² approx. 1.8 m NEMA L15-20
EL	Electrical connection PG 8593	Voltage (convertible) Rated load Fuse rating Minimum cross-section Length of electrical cord (UL 921) with plug Electrical connection	2 AC 208V 60 Hz 6.2 kW 2 pole 30 Amp 3 x AWG 10 approx. 5' 9" NEMA L6-30	2 AC 208V 60 Hz 6.2 kW 2 pole 30 Amp 3 × 5.2 mm ² approx. 1.8 m NEMA L6-30

Connection to the electrical supply must be carried out in accordance with local and national safety regulations. The power cord must be protected from the risk of thermal damage.

It is recommended to make an electrical connection via a plug and socket so that electrical safety checks can be carried out easily. The socket must be accessible after machine installation.

For hard-wired machines, a main switch must be provided on site, which must completely isolate the machine from the power supply with a contact gap of at least 1/8" (3 mm).

The plug and socket as well as the main switch must be easily accessible after the machine has been installed.

Installation plan PG 8592 – PG 8593 M.-Nr. 10 344 170 / 00 Date created: 06.11.2015 Page 2 / 6

(A)	Equipotential bonding and grounding Communication modules (optional)	For added safety the machine should be protected with a residual current device with a trip current of 30 mA. If necessary, an equipotential bond with good contact connection must be provided in accordance with all appropriate national and local regulations. Equipotential bonding and grounding must be completed before the machine is commissioned. There is a screw connection point at the back of the washer-disinfector, to which equipotential bonding can be connected. Optional ethernet and RS232 communication modules are available. They are not supplied with the machine. The fuse box for the modules should be installed near the machine for transferring and printing process data. They must be installed and connected in accordance with local codes and standards.			
	Cold Water				
Cold Water Connections Use supplied Y-adapter if only one cold water connection is					
	(machine comes with two cold	Max. temperature	68 °F	20 °C	
	water connections)	Max. water hardness	73 gpg	70 °dH	
	,	Recommended flow pressure	29 psi	200 kPa	
		Min. flow pressure with extended water intake	14.5 psi	100 kPa	
		Max. pressure	145 psi	1,000 kPa	
		Throughput	2 gal/min	7.5 l/min	
		Connection thread - on site (flat sealing)	3/4" Male Garden Hose Thread	3/4" Male Garden Hose Thread	
		Length of connection hose (supplied)	approx. 5' 7"	approx. 1.7 m	
		Length of the steam condenser hose	approx. 5' 7"	approx. 1.7 m	
6000	Hot Water	Max. temperature	150 °F	65 °C	
WW	Connection	Max. water hardness	73 gpg	70 °dH	
		Recommended flow pressure	29 psi	200 kPa	
		Min. flow pressure with extended water intake	5.8 psi	40 kPa	
		Max. pressure	145 psi	1,000 kPa	
		Throughput	2 gal/min	7.5 l/min	
		Connection thread - on site (flat sealing)	3/4" Male Garden Hose Thread	3/4" Male Garden Hose Thread	
		Length of connection hose (supplied)	approx. 5' 7"	approx. 1.7 m	
(VE)	Demineralized	Max. temperature	150 °F	65 °C	
	water	Recommended flow pressure	29 psi	200 kPa	
		Min. flow pressure with extended water intake	4.4 psi	30 kPa	
		Max. pressure	145 psi	1,000 kPa	
		Connection thread - on site (flat sealing)	3/4" Male Garden Hose Thread	3/4" Male Garden Hose Thread	
		Throughput	2 gal/min	7.5 l/min	
		Length of connection hose (supplied)	approx. 5' 7"	approx. 1.7 m	
Requirements for unpressurized demineralized water system:					
		Minimum flow pressure (DI unpressurized)	1.3 psi	8.5 kPa	
		Maximum pressure (DI unpressurized)	8.7 psi	60 kPa	
		Machine connection nipple (outer dia. x I)	1/4" x 3/16"	6 mm x 30 mm	
		Hose from supply container to machine to be installed on site.			
i					

Installation plan PG 8592 – PG 8593 Date created: 06.11.2015

(AVA)	Drain Water	Drainage temperature	200 °F	93 °C
W		Length of drain hose	approx. 4' 6"	approx. 1.4 m
		Drain hose, max. drainage length	approx. 13' 1"	approx. 4.0 m
		Drain pump head height from floor level, max.	approx. 3' 3"	approx. 1 m
		Drain hose delivery, max.	4.2 gal/min	16 l/min
		On-site hose connector (external diameter x length)	7/8" × 1 3/16"	22 mm × 30 mm
		It is advisable to lay the drain hose in a curve min. 23 the drainage point.	5/8" (600 mm) above floor level to	
DS	External dispensing (optional)	Connection for up to two dispensing units for liquid agents on the back panel.		
	(note machine model)			
	Machine feet	Height-adjustable	0-2 3/8"	0-60 mm
		Foot diameter	1 3/8"	35 mm
		Entry point into housing	8 M	8 M
	Heat Dissipation	Heat transfer by emission	1,194 btu/h	0.35 kWh/h
	to Room	Load on removal	1,365 btu/h	0.40 kWh/h
	Machine data	Height	32 7/8"	835 mm
		Height without lid	32 5/16"	820 mm
		Width	23 5/8"	600 mm
		Width without lid	23 9/16"	598 mm
		Depth	23 5/8"	600 mm
		Depth without lid	23 9/16"	598 mm
		Door height	24 1/2"	622 mm
		Net weight	172 lbs	78 kg
		Dynamic floor load	1,200 N	1,200 N
		Width including transport pallet	26 3/8"	670 mm
		Depth including transport pallet	29 3/8"	740 mm
		Height including transport pallet	36 1/4"	920 mm
		Sound emissions in dB (A),	< 70 dB	< 70 dB
		Sound pressure LpA, washing and drying	< 10 ub	< 10 UB
	Set-up Conditions	Ambient temperature	40-105 °F	5-40 °C
		Max. humidity	80%	80%
		Rel. humidity, declining proportionally up to 104°F (40°C)	50%	50%
		Max. altitude above sea level	4,921'	1,500 m

Please read and observe the instructions in the operating instruction manual and the service documents regarding installation and setting up of this medical washer/disinfector.


Installation should only be carried out by authorized installers in accordance with valid regulations, relevant standards and health and safety codes.

The machine must be commissioned and operators trained in its use by Miele Service or by an approved Miele Service Technician only.

Condensate can build up in the area surrounding the machine. Any cabinetry and fixtures in the room must be suitable.


If the machine is installed under a countertop (built-under model), a stainless steel panel must be installed above the door opening to protect the countertop from moisture damage. A suitable stainless steel cover plate is available to order from Miele.

Installation plan PG 8592 – PG 8593 M.-Nr. 10 344 170 / 00 Date created: 06.11.2015 Page 4 / 6

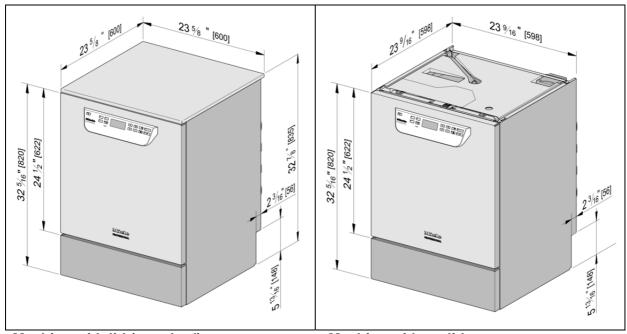


Machine with lid (standard)

Machine without lid

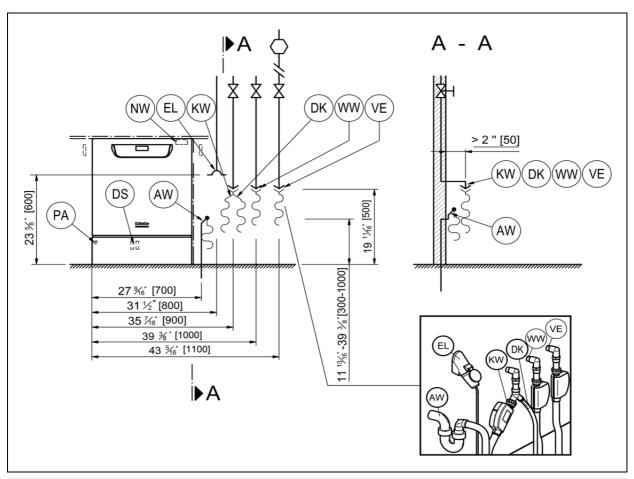

AW	Drain water	NW	Network and printer connection
DK	Cold water connection, steam condenser	PA	Equipotential bonding
DS	External dispensing system connection	VE	Demineralized water connection
EL	Electrical connection	WW	Hot water connection
KW	Cold water connection		

Installation plan PG 8592 – PG 8593 Date created: 06.11.2015

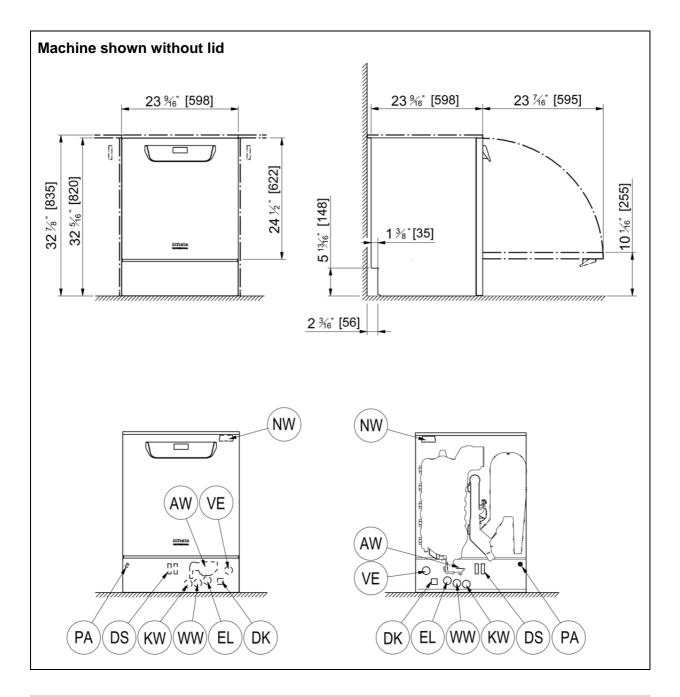

AW	Drain water	NW	Network and printer connection
DK	Cold water connection, steam condenser	PA	Equipotential bonding
DS	External dispensing system connection	VE	Demineralized water connection
EL	Electrical connection	WW	Hot water connection
KW	Cold water connection		