# APFD 200

|           |                               |
|-----------|-------------------------------|
| <b>de</b> | Gebrauchsanweisung Einsatz    |
| <b>da</b> | Brugsanvisning Indsats        |
| <b>el</b> | Οδηγίες χρήσης Ένθετο         |
| <b>en</b> | Operating instructions Insert |
| <b>fi</b> | Käyttöohje Teline             |
| <b>fr</b> | Mode d'emploi Complément      |
| <b>it</b> | Istruzioni d'uso Inserto      |
| <b>nl</b> | Gebruiksaanwijzing inzet      |
| <b>no</b> | Bruksanvisning innsats        |
| <b>sv</b> | Bruksanvisning Insats         |

**de** ..... 4  
**da** ..... 16  
**el** ..... 28  
**en** ..... 41  
**fi** ..... 53  
**fr** ..... 65  
**it** ..... 77  
**nl** ..... 89  
**no** ..... 101  
**sv** ..... 113

---

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## en - Contents

---

|  |           |
|--|-----------|
| <b>Notes about these instructions</b> .....        | <b>41</b> |
| Symbols and conventions used in this document..... | 41        |
| Definition of terms .....                          | 41        |
| <b>Appropriate use</b> .....                       | <b>42</b> |
| Queries and technical problems .....               | 42        |
| <b>Items supplied</b> .....                        | <b>43</b> |
| Load carriers.....                                 | 43        |
| Disposal of the packing material .....             | 43        |
| Optional accessories.....                          | 43        |
| <b>Warnings and safety notes</b> .....             | <b>44</b> |
| <b>Areas of application</b> .....                  | <b>45</b> |
| Colour coding.....                                 | 45        |
| Preparing the load .....                           | 46        |
| Preparing the insert.....                          | 46        |
| Pre-cleaning air regulators.....                   | 46        |
| Dismantling air regulators.....                    | 46        |
| Inserting air regulators .....                     | 46        |
| Connecting the hose .....                          | 47        |
| Completely loading the insert.....                 | 49        |
| Filling the pressure chamber .....                 | 49        |
| Placing the insert in the load carrier .....       | 50        |
| After reprocessing .....                           | 50        |
| <b>Maintenance</b> .....                           | <b>51</b> |
| Routine check .....                                | 51        |

### Symbols and conventions used in this document

#### Warnings

 Warnings contain information which is important for safety. This alerts you to the potential danger of injury to people or damage to property.

Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

#### Notes

Notes provide information of particular importance that must be observed.

#### Additional information and comments

Additional information and comments are contained in a simple frame.

#### Operating steps

Operating steps are indicated by a black square bullet point.

##### Example:

■ Select an option using the arrow buttons and save your choice with *OK*.

##### Example:

■ Select an option.

#### Display

Display text can be identified from the special font.

##### Example:

Save.

### Definition of terms

#### Cleaning machine

In these operating instructions, the dishwasher or the washer-disinfector is referred to as “the machine”.

#### Load items

The term “load items” is used wherever the items to be processed are not defined in any further detail.

#### Wash water

The term “wash water” is used for the mixture of water and process chemicals.

## en - Appropriate use

---

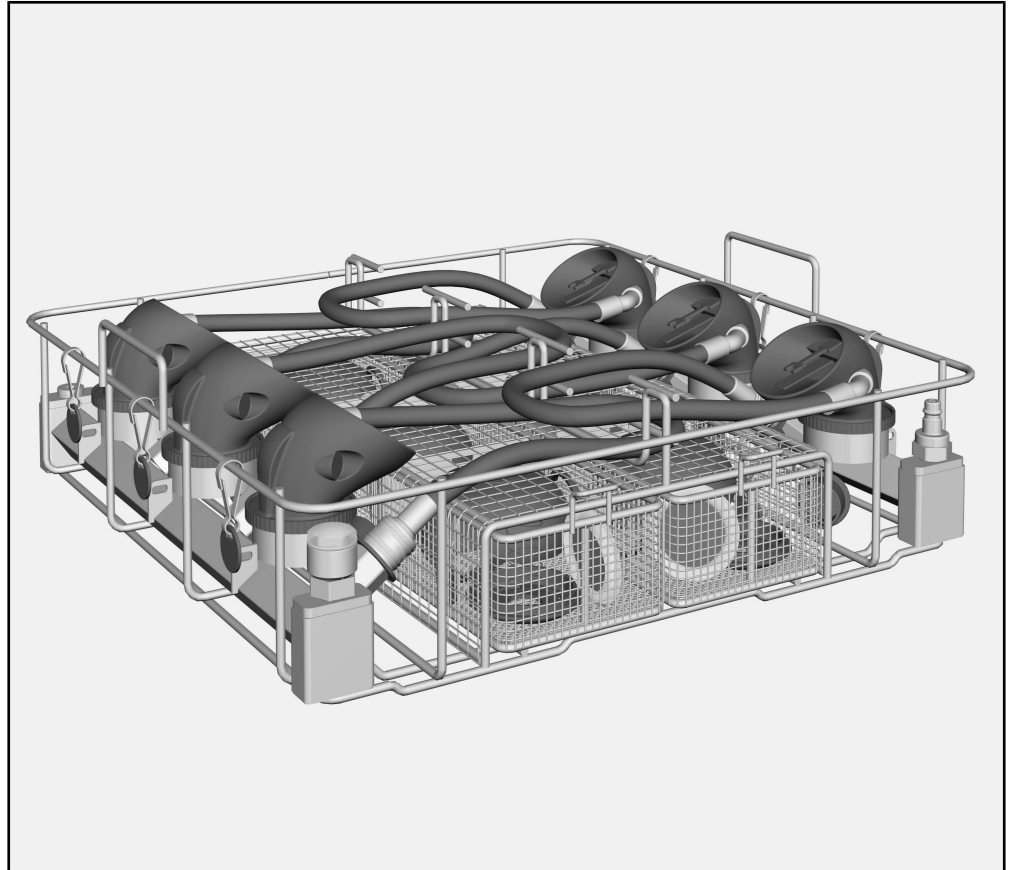
The APFD 200 insert can be used to reprocess machine-reprocessable air regulators in a Miele dishwasher or a machine. Follow the operating instructions for your dishwasher or the machine as well as the instructions issued by the manufacturer of the air regulators.

The APFD 200 insert can be used together with the A 101 upper basket or the A 151 lower basket.

### **Queries and technical problems**

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your dishwasher or the washer-disinfector or go to [www.miele-professional.com](http://www.miele-professional.com).

## Load carriers

- APFD 200 insert, height 153 mm, width 484 mm, depth 452 mm

### Items supplied

- 2 x compartment for small parts, divided into three parts
- 6 x clip labels in 6 colours

### Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

### Optional accessories

The following and other accessories are available to order from Miele:

- APFD 202, 6 adapters for loading specially shaped air regulators, e.g. from Interspiro.

## en - Warnings and safety notes

---

Read the operating instructions carefully before using this insert.  
Keep these operating instructions in a safe place.

 Please also read the operating instructions for your dishwasher or the washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ The inserts are only approved for use in the applications stated in the operating instructions. Applications that go beyond those listed and use of the inserts with additional components from the Miele range of components are described in the operating instructions of the components or are to be agreed with Miele. Conversions, modifications and any other use are not permitted.
- ▶ Prior to being used for the first time, new inserts must undergo thorough cleaning in the dishwasher or the washer-disinfector without load items.
- ▶ Inspect all mobile units, baskets, modules and inserts as described in the “Maintenance” section of the operating instructions for your machine.
- ▶ Only items which have been declared by their manufacturer as suitable for machine reprocessing may be processed. The manufacturer's specific reprocessing instructions must be observed.
- ▶ Follow the warnings and instructions provided by the breathing apparatus manufacturer.
- ▶ Any seals on the load items should be protected against prolonged contact with wash water. Therefore, always close unused connections with protective caps or blind covers.
- ▶ The last rinse, as a minimum, must be carried out with DI water.
- ▶ Disposable items must not be reprocessed.

The APFD 200 insert for reprocessing air regulators must be supplied with pressure for reprocessing. The pressure ensures that wash water cannot enter the air regulators during reprocessing, assuming the correct adapters are used.

**⚠ Risk of infection**

The compressed air must correspond to the quality of breathing air as a minimum.

For this, observe the national and regional regulations concerning compressed air.

**⚠ Risk of injury caused by compressed air**

A flexible compressed air connection with defined quick-release couplings must be provided on site.

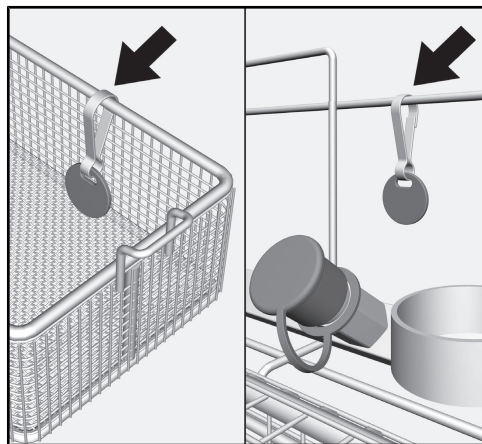
Use a compressed air connection with one of the following quick-release couplings:

- CEJN series 341
- Parker Rectus series 25S


### Colour coding

The small parts of a dismantled air regulator are reprocessed in a compartment of the small parts basket. Colour-code the compartments and the ring holders or the load item carriers to ensure that the small parts can be assigned to the right air regulators after reprocessing.

Version with carabiner:

Version with clip label:

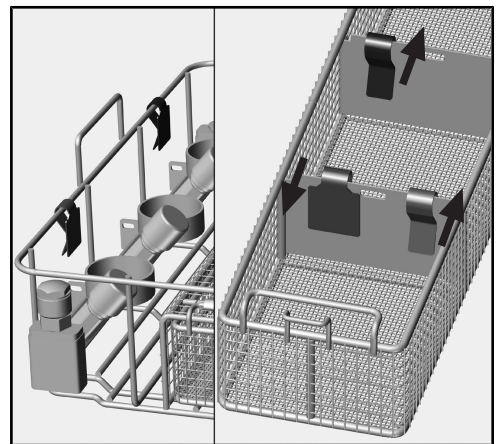For each colour, attach one tag to the ring holders or the load item carriers and one tag to the small parts basket compartments.


If you are using clip labels, they must be attached to the dividers inside the small parts compartments. We recommend that the colour-coded tag is attached to the side facing the surface (see arrows in image above).

When using other systems for identification, ensure that these are suitably chemical-resistant and temperature-resistant.

### Preparing the load
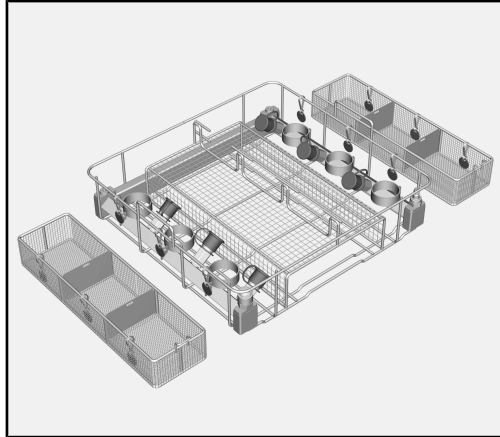
#### Preparing the insert

- Place the APFD 200 insert on the worktop. The insert must be stable.

- Remove the small-part baskets and place these on the left and right next to the APFD 200 insert.


The quick-release couplings on the insert should have the same colour coding as the accompanying compartments of the small-part baskets so that you can assign the air regulator small parts clearly.

#### Pre-cleaning air regulators

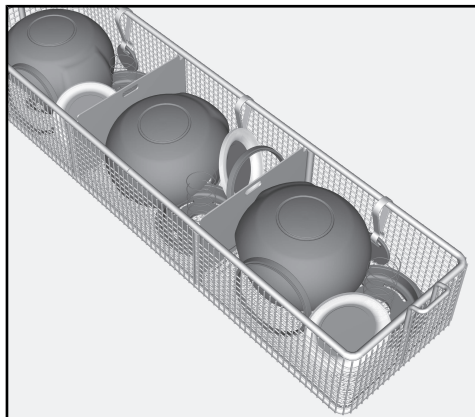- Check whether the air regulator requires manual pre-cleaning, e.g. to remove soot.
- If required, carry this pre-cleaning out in line with the manufacturer's instructions.

#### Dismantling air regulators

- Dismantle the air regulator in line with the manufacturer's instructions.


- Place the small parts in a compartment in a small-part basket.
- Do not position small parts in such a way that they will collect water. For example, insert rubber covers with the curved hollow side facing down.

Note the colour-coding for the small-part basket compartment to place the air regulator in the same colour in the insert.


#### Inserting air regulators

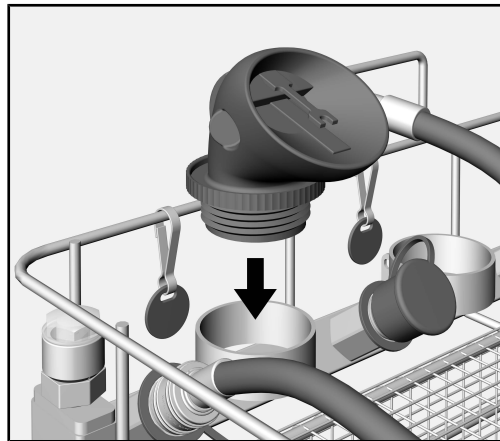- Select the ring holder with the same colour-coding as on the small-part basket.

- Insert the coupling of the air regulator with the opening facing downwards in the ring holder.

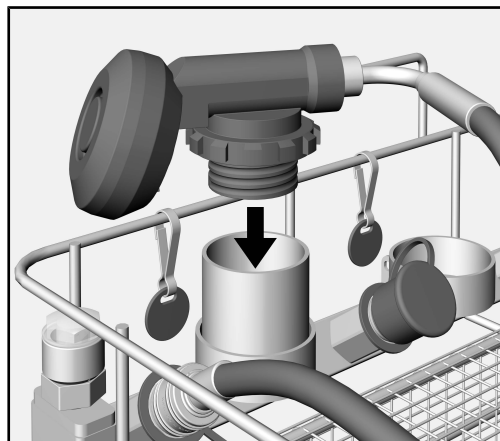For specially shaped air regulators, the ring holders must be equipped with APFD 202 adapters. The adapters extend the ring holders and ensure a better grip.

APFD 202 adapters are available as an optional accessory.

### Connecting the hose

Quick-release couplings

To supply the air regulator with compressed air, the hose must be connected to the insert.

The quick-release couplings on the insert are fitted with protective caps. The protective caps prevent wash water from collecting in the couplings.

#### ⚠ Danger of toxic fumes

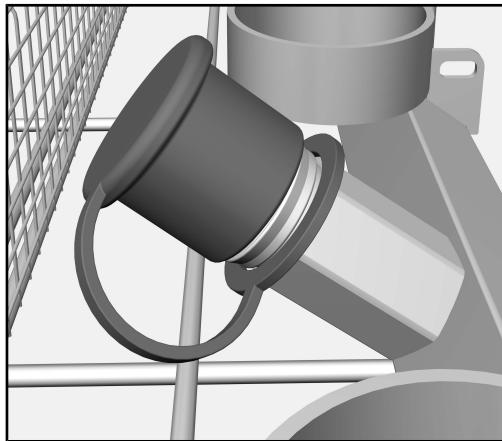
If wash water collects in the quick-release couplings, the pressure can force this into the air regulator when it is connected.

The seals in the quick-release couplings swell up when they are in contact with wash water for an extended period, causing them to become damaged.

For this reason, always seal unused quick-release couplings with protective caps.


- Select the adapter with the same colour-coding as on the ring holder.

## en - Areas of application

- Remove the protective cap from the quick-release coupling.

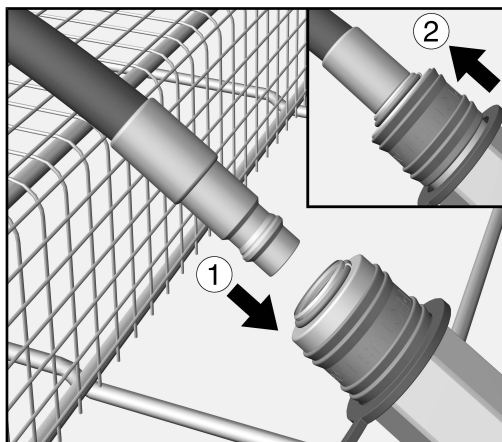Connecting the ad-  
apter

- Insert the hose adapter into the quick-release coupling on the insert ① until it engages with an audible click.


The collar on the coupling is pushed downwards in the process. ②. As soon as the collar has snapped back, the hose is pressure-tight.

**⚠ Danger to health**

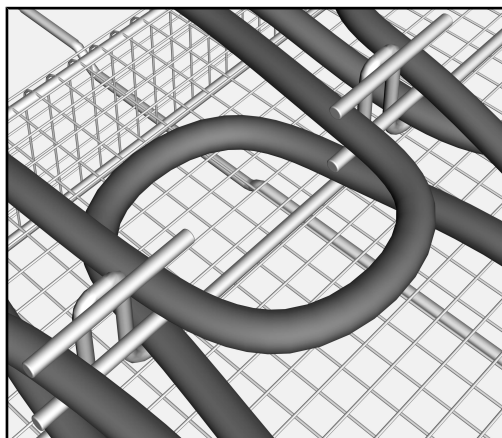Damaged quick-release couplings can force wash water into the air regulator.

Have the quick-release coupling replaced with an original spare part if it is no longer possible to hear the click clearly or the collar does not snap back fully.

Securing the hose

- Lay the hose in an arc shape in the hose guide for the insert.

The locking pins on the hose guide hold the hose in position.

If hoses are not secured, they may block the spray arms during re-processing.

This will affect the reprocessing result.

Secure all hoses using the locking pins on the hose guide.

### Completely loading the insert

- Equip the remaining positions of the insert in the same way.
- Then put the small-parts baskets back in the APFD 200 insert.

### Filling the pressure chamber

⚠ Risk of injury due to compressed air.

The jet of compressed air could be dangerous for the user.

Users must be trained on how to use the compressed air system on the customer's premises.

The compressed air must correspond to the quality of breathing air as a minimum. For this, observe the national and regional regulations.

The pressure chamber is designed for pressure ranging up to 400–600 kPa.

- Connect the compressed air hose on the customer's premises to the quick-release coupling on the pressure chamber of the insert.
- Add pressure to the pressure chamber until the red compressed air indicator pokes out.

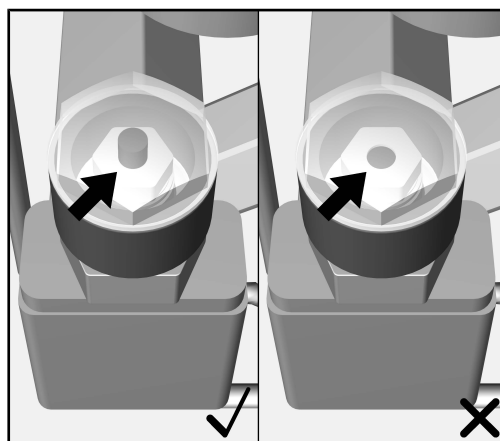
⚠ Damage to the air regulators caused by pressure.

Do not provide more than 600 kPa of pressure to the pressure chamber. Higher pressure can damage the air regulators.

Observe the air regulator manufacturer's instructions.

### Compressed air indicator

The compressed air indicator on the pressure chamber shows whether a sufficient amount of pressure has been provided.

- ✓ If the red indicator of the compressed air indicator pokes out, the pressure is sufficient.
- ✗ If the indicator does not poke out, the pressure is not sufficient.

 Danger to health

If the connection between the quick-release couplings on the pressure chamber and the hoses for the air regulators is not absolutely tight, the pressure will drop over time.

If the pressure drops too much, wash water and particles may get into the air regulators and impair their function.

Therefore, supply the pressure chamber with pressure as close to the programme start as possible.

### **Placing the insert in the load carrier**

- Lift the insert at the side handles.
- Place the insert in the load carrier of the dishwasher or the machine.

### **After reprocessing**

It is essential that you pay attention to the “Application technology” chapter in the operating instructions for the dishwasher or the machine and the instructions specified by the manufacturer of the air regulators.

### Routine check

 Danger to health

Damaged seals in the quick-release couplings can force wash water and particles into the air regulator. This prevents the air regulator from working properly.

Have the quick-release coupling with damaged seals replaced with an original spare part.

The seals and airtightness of the pressure chamber must be checked regularly to make sure the insert is functioning correctly.

#### Daily

- During the daily routine, visually check the seals in the quick-release couplings to ensure they are intact.

#### Weekly

- During the weekly routine, check that the pressure chamber is airtight.

To do this:

- Fit the quick-release couplings on the insert with blind covers.
- Apply compressed air to the pressure chamber of the insert.

The compressed air indicator must remain switched after an hour.

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