# A 200

<b>de</b>	Gebrauchsanweisung Injektorwagen
<b>da</b>	Brugsanvisning Injektorvogn
<b>en</b>	Operating instructions Injector unit
<b>es</b>	Instrucciones de manejo Carros inyectoros
<b>fi</b>	Käyttöohje Suorasuihkuvaunu
<b>fr</b>	Mode d'emploi Chariot à injection
<b>hr</b>	Upute za uporabu Injektorska kolica
<b>it</b>	Istruzioni d'uso Carrello a iniezione
<b>nl</b>	Gebruiksaanwijzing injectorwagen
<b>no</b>	Bruksanvisning injektorvogn
<b>pt</b>	Instruções de utilização Carro injetor
<b>sl</b>	Navodila za uporabo Injektorski voziček
<b>sv</b>	Bruksanvisning Injektorvagn

**de** ..... 4  
**da** ..... 11  
**en** ..... 18  
**es** ..... 25  
**fi** ..... 32  
**fr** ..... 39  
**hr** ..... 46  
**it** ..... 53  
**nl** ..... 60  
**no** ..... 67  
**pt** ..... 74  
**sl** ..... 81  
**sv** ..... 88

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<b>Notes about these instructions</b> .....	18
<b>Correct application</b> .....	19
Queries and technical problems.....	19
<b>Items supplied</b> .....	20
Disposal of the packing material .....	20
<b>Warning and Safety instructions</b> .....	21
<b>Areas of application</b> .....	22
Before loading the machine and before starting a programme .....	22
Loading pipettes.....	22

### Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

### Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

### Additional information and comments

Additional information and comments are contained in a simple frame.

### Operating steps

Operating steps are indicated by a black square bullet point.

#### Example:

- Select an option using the arrow button and save your choice with *OK*.

### Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

#### Example:

Settings  menu.

Re-usable laboratory glassware and laboratory utensils can be re-processed using this mobile unit in a Miele laboratory glassware washer-disinfector. Follow the instructions for your washer-disinfector as well as the instructions of the glassware manufacturer on how to reprocess their items by machine.

The A 200 mobile injector unit is set up for reprocessing pipettes.

Use the **Pipettes** programme.


The mobile unit can be used in the following washer-disinfectors:

- PG 8583
- PG 8583 CD
- PG 8593

Laboratory glassware and laboratory utensils for reprocessing are referred to as the wash load if they are not more closely defined.

### **Queries and technical problems**

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your Washer-Disinfector or go to [www.miele-professional.com](http://www.miele-professional.com).

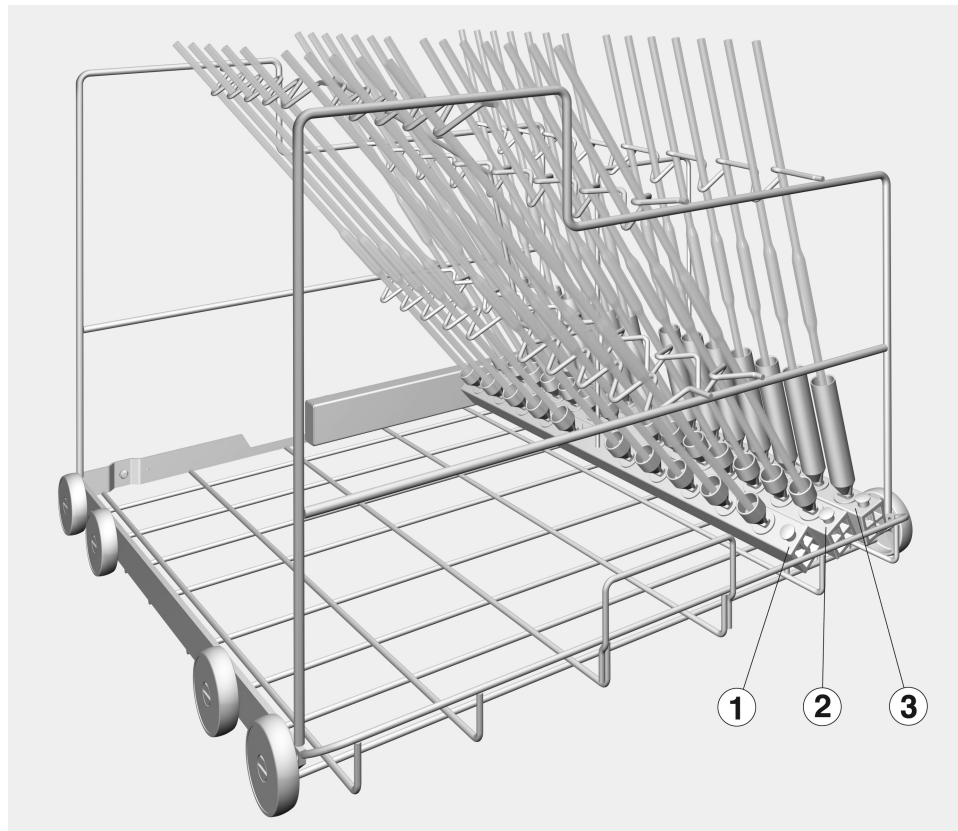– Mobile injector unit A 200, height 397 mm, width 531 mm, depth 546 mm, with 3 injector manifolds.

### **Disposal of the packing material**

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in landfill sites.

## en - Warning and Safety instructions

---

To avoid the risk of accidents and damage to this mobile unit please read these instructions carefully before using it for the first time.  
Keep these instructions in a safe place where they are accessible to users at all times.

 Please also read the Operating instructions for your washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ This mobile injector unit is only for use in the areas of application as defined in the Operating instructions. Components such as injector nozzles must only be replaced with Miele accessories or original parts. Any other application, modification or change is unauthorised and could be dangerous.  
Miele cannot be held liable for damage caused by unauthorised use or incorrect operation.
- ▶ Inspect all mobile units, baskets, modules and inserts daily as described in the "Maintenance" section in the Operating instructions for the washer-disinfector.
- ▶ Only items which have been declared by their manufacturer as suitable for machine reprocessing may be processed. The manufacturer's specific reprocessing instructions must be observed.
- ▶ Broken glass can result in serious injury when loading or unloading. Broken glass items must not be processed in the washer-disinfector.
- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

Miele cannot be held liable for damage caused by non-compliance with these Warning and Safety Instructions.

## en - Areas of application

Before placing pipettes in the injector unit, please make sure that all hollow sections are free of blockages.

### Before loading the machine and before starting a programme


- Check that the mobile unit is correctly connected to the water supply in the washer-disinfector.

### Loading pipettes


This mobile injector unit has three injector manifolds which must be loaded as follows:

Row	Max. length [mm]	Qty	Volume [ml]
1	550	10	10 to 100
2	500	14	10 to 25
3	440	14	1 to 10

Pipettes must not extend beyond the sides of the mobile injector unit.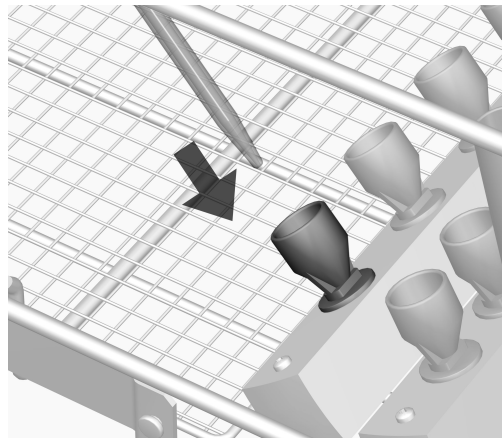

- Insert the pipettes into the funnels with the pointed end facing downwards.
- Fit the top ends of the pipettes into the relevant holders on the support rail.

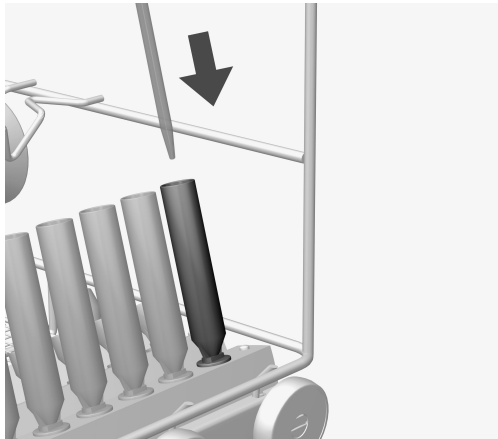- Pipettes with a volume of between 1 and 5 ml should be fitted into irrigation sleeves in row ③ in the mobile injector unit.

The spray pressure could eject them from their funnels if placed in rows ① and ②.

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