A 304

de	Gebrauchsanweisung Injektormodul
da	Brugsanvisning Injektormodul
en	Operating instructions Injector module
es	Instrucciones de manejo Módulo inyector
fi	Käyttöohje Suorasuihkumoduuli
fr	Mode d'emploi Module à injection
hr	Upute za uporabu Injektorski modul
it	Istruzioni d'uso Modulo iniettore
nl	Gebruiksaanwijzing injectormodule
no	Bruksanvisning injektormodul
pt	Instruções de utilização Módulo injetor
sl	Navodila za uporabo Injektorski modul
sv	Bruksanvisning injektormodul

de	4
da	12
en	20
es	28
fi	36
fr	44
hr	52
it	60
nl	68
no	76
pt	84
sl	92
sv	100

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Notes about these instructions	20
Appropriate use	21
Queries and technical problems	21
Items supplied	22
Disposal of the packing material	22
Warnings and safety notes	23
Areas of application.....	24
Injector modules.....	24
Inserting the module	24
Removing an injector module	24
To check before loading the machine and before starting a programme	25
Preparing the load	26

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

- Select an option using the arrow buttons and save your choice with *OK*.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings  menu.

This module can be used to reprocess machine-reprocessible laboratory glassware and utensils in a Miele washer-disinfector for laboratory glassware and utensils. Follow the operating instructions for your washer-disinfector as well as the instructions of the glassware and utensil manufacturer on how to reprocess their items by machine.

The A 304 injector module is designed for reprocessing centrifuge tubes, phials, test tubes for fraction collectors and auto-sampler tubes.

To ensure sufficient water pressure, only one module can be used at a time in each upper and lower basket. The module must always be placed in the right hand position. The remaining connections can be used for other modules.

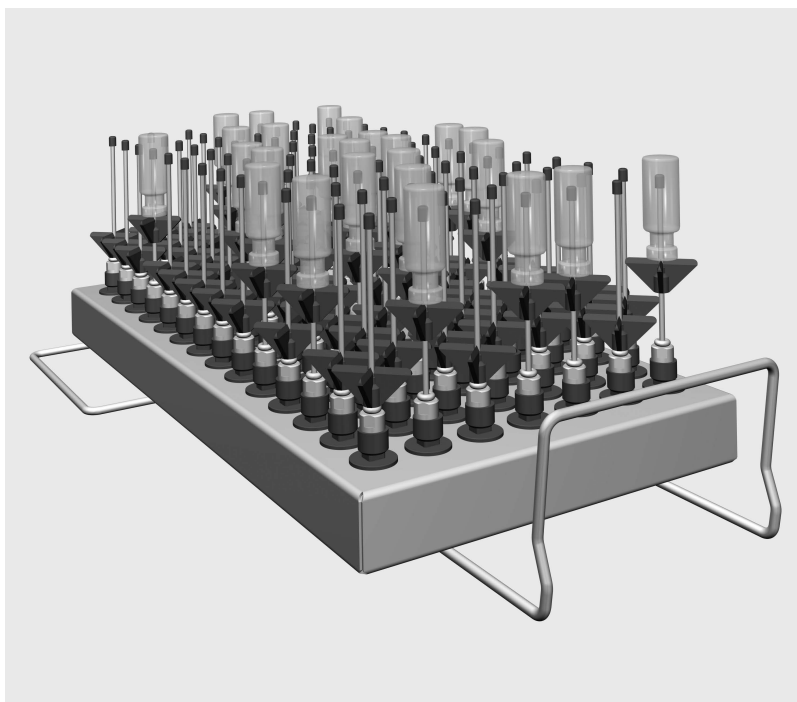
This module is suitable for use in the following:

- A 100 upper basket
- Lower basket A 150
- APLW 106 basket

The washer-disinfector is generally referred to as “the machine” in these operating instructions. Reprocessible laboratory glassware and utensils are referred to as “items” if they are not more closely defined.

Queries and technical problems

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your dishwasher or the washer-disinfector or go to www.miele-professional.com.

- Injector module A 304, height 130 mm, width 222 mm, depth 471 mm.

Items supplied

- 98 x ID 90 injector nozzles with plastic supports. Length 90 mm, Ø 2.5 mm

Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

To avoid the risk of accidents and damage to this module please read these instructions carefully before using it for the first time. Keep these instructions in a safe place where they are accessible to users at all times.

 Please also read the operating instructions for your dishwasher or the washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ The module is approved solely for the applications specified in the “Application technology” chapter of these operating instructions. Components such as nozzles may only be replaced with Miele accessories or genuine Miele original spare parts. Conversions, modifications and any other use are not permitted.
- ▶ Inspect all mobile units, baskets, modules and inserts daily as described in the “Maintenance” section in the Operating instructions for the washer-disinfector.
- ▶ Only items which have been declared by their manufacturer as suitable for machine reprocessing may be processed. The manufacturer's specific reprocessing instructions must be observed.
- ▶ Broken glass can result in serious injury when loading or unloading. Broken glass items must not be processed in the machine.
- ▶ Only place empty, unloaded modules in baskets. Check that they are correctly engaged before loading them. Modules must be completely emptied before taking them out. Placing loaded modules in baskets or removing them can cause damage to the wash load and cause injury should glassware be broken.
- ▶ The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

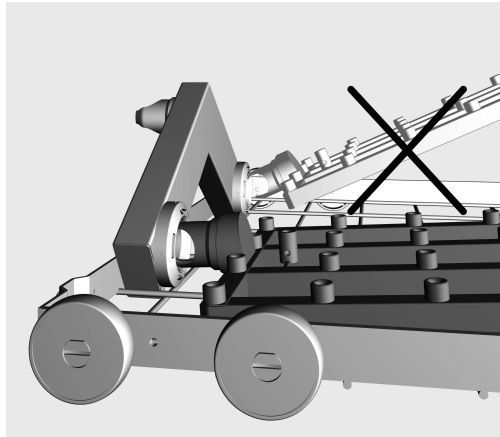
Miele cannot be held liable for damage or injury caused by non-compliance with the warning and safety notes.

Injector modules

⚠ Only place empty, unloaded modules in baskets. Check that they are correctly engaged before loading them. Modules must be completely emptied before taking them out. Placing loaded modules in baskets or removing them can cause damage to the wash load and cause injury should glassware be broken.

Inserting the module

Ensure that the module is correctly orientated when inserting it. For example, the nozzles must point upwards.

Example of a flat insertion angle

- Insert the module into the water connection with the connector at a flat angle.
- Then engage the module by pushing it down onto the struts of the basket.

Removing an injector module

⚠ Do not hold the basket by the injector module when removing it from the cleaning machine. The module could come loose or the water connection could be damaged.

- Release the module from the locking mechanism and pull it out of the water connection at a flat angle. If the module is removed at too steep an angle, the lever action may damage the connection point and the connector.

To check before loading the machine and before starting a programme

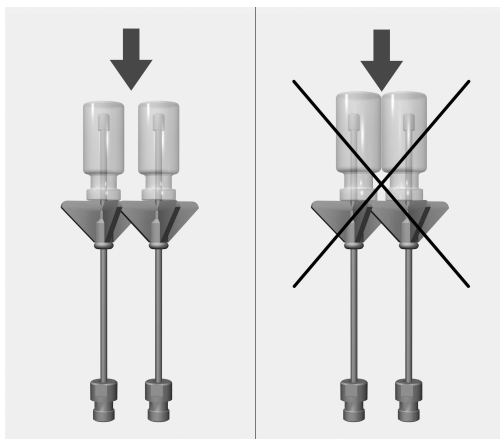
- Check that fittings such as irrigation sleeves and injector nozzles are securely screwed into position.

 Make sure that all screw connectors are fitted with injector nozzles, adapters, irrigation sleeves or blind stoppers to ensure that all fittings in use are supplied with sufficient standardised pressure. Damaged fittings such as injector nozzles, adapters and irrigation sleeves must not be used.

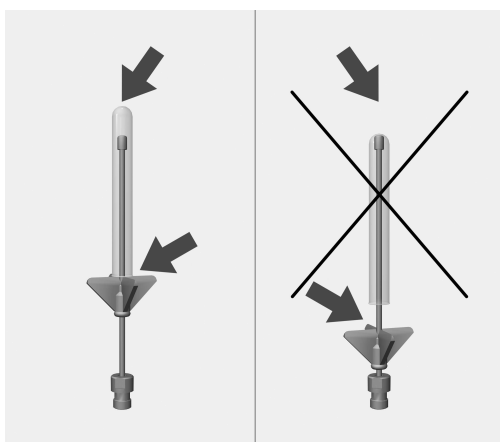
Fittings not equipped with wash items do not have to be replaced blind stoppers.

- Check that the baskets and modules are correctly connected to the water supply.

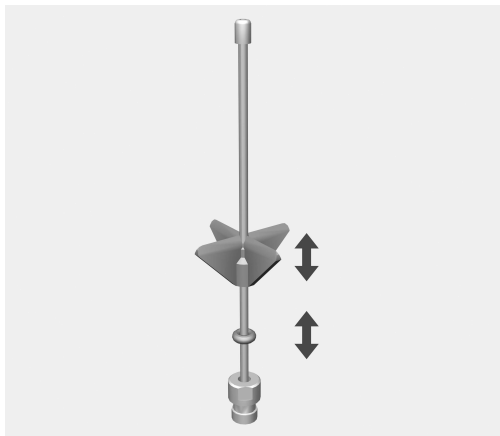
Preparing the load

Individual items of glassware must not touch each other.

If the tip of an injector nozzle touches the bottom of a piece of glassware, the plastic support has to be pushed upwards.

- Push the plastic support upwards to the height required. If necessary use a pair of pointed pliers to do so.
- Push the O-ring up to the plastic support to secure it in position.

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