XKM 3200 WL LMD86

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Notes about these instructions	46
Queries and technical problems	46
Definition of terms	46
Module	46
Intended use	47
Items supplied	48
Scope of delivery	48
Technical data	48
EU Conformity declaration.....	48
Guide to the appliance	49
Module overview	49
Network connection	49
Switching on and off	50
LED indicator.....	50
Reset button	50
Data protection and data security	51
Factory default settings for network configuration	51
Copyright and licences	51
Warnings and safety notes	52
Fitting and installation	53
Installation	53
Module slot.....	53
Inserting the module	54
Laying data cables	54
Network connection	55
Configuring the interface.....	55
Switching the cleaning machine on	55
Web interface	56
Password	56
Caring for the environment	58
Disposal of the packing material	58

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property. Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

■ Select an option using the arrow buttons and save your choice with *OK*.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

menu.

Queries and technical problems

In the event of queries or technical problems, please contact Miele. Contact details can be found at the end of the operating instructions for your cleaning machine or at www.miele.com/professional.

Definition of terms

Machine

In these operating instructions, the laboratory washer is referred to as “the machine”.

Module

The communication module is generally referred to as the module in these operating instructions.

The XKM 3200 WL LMD86 communication module supplied enables an interface to be set up between a Miele washer-disinfector and a terminal device or local network. The data connection is a standard Ethernet connection. The interface serves a number of purposes, including process documentation. Further information on process documentation can be obtained from Miele.

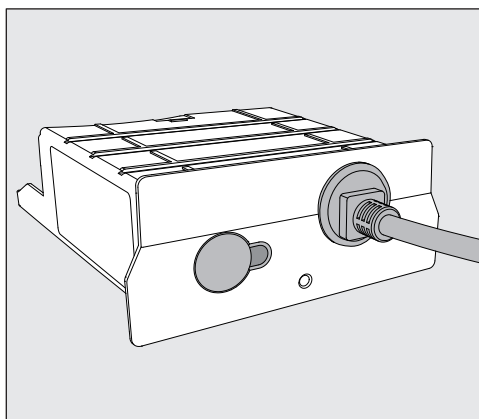
Only terminal devices (PCs, tablets, printers, etc.) compliant with IEC/EN 62368 should be used.

Only specialists are permitted to configure the interface.

The communication module can be used in the following washer-disinfectors:

- PLW 8683
- PLW 8683 CD
- PLW 8693
- PWD 8682
- PWD 8682 C
- PWD 8682 CD
- PWD 8692

Scope of delivery

- XKM 3200 WL LMD86 communication module

Technical data

Voltage, rated load, fuse rating	See data plate
Mains cable	Approx. 1,5 m
Ethernet baud rate	10/100 Mbit/s
Temperature during operation	Up to max. 80 °C
CE marking	2014/53/EU Radio Equipment Directive
Manufacturer address	Miele & Cie. KG, Carl-Miele-Straße 29, 33332 Gütersloh, Germany

EU Conformity declaration

Miele hereby declares that the XKM 3200 WL LMD86 communication module complies with Directive 2014/53/EU.

The complete text of the EU declaration of conformity can be downloaded from the following website by quoting the product name or serial number:

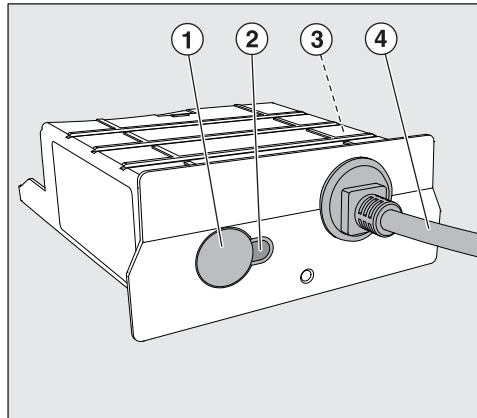
- www.miele.com/professional

> Products > Download

or

> Service & Support > Operating instructions

Module overview

- ① Reset button
- ② LED indicator
Indicates operational readiness after installation and active communication when in operation
- ③ Data plate (located on the side)
- ④ Data cable

Network connection

Integrating the cleaning machine into the on-site network requires the following:

- Ethernet network
- DHCP server or specifications for static IPv4 settings
- Network printer (compatible with Internet Printing Protocol and PostScript 3): printer URL and port

Switching on and off

The module does not have a separate On/Off switch. It is switched on and off together with the cleaning machine.

The module requires approx. 90 s for initialisation after switch-on. Until this has happened data cannot be transferred and interface settings cannot be changed.

LED indicator

The LED indicator (light emitting diode) provides status information about the communication module.

LED	Status
Lit up GREEN	Connection to the network was successful. The module is ready for operation.
Lit up YELLOW	Network is configured but there is no connection to the network.
Lit up RED	The module is not configured.
Off	Power is not being supplied to the module. Check that the module is in the correct position in the module slot.

Reset button

The button on the module can be used to reset the module to the factory default settings.

- Press the button for approx. 5–9 seconds to restore the factory default settings.

Data protection and data security

When you activate the networking function and connect your machine to the Internet, your machine sends the following data to the Miele Cloud:

- Machine serial number
- Machine model and technical features
- Machine status
- Information about the software version of your machine

Initially, this data cannot be assigned to a specific user and is not saved permanently. Data cannot be saved permanently or assigned to a specific user until after you have linked your machine to a user. Data transmission and processing are governed by Miele's strict security standards.

Factory default settings for network configuration

You can reset all of the settings on the communication module or your integrated WiFi module to the factory default settings. The network configuration should be reset whenever an appliance is being disposed of or sold, or if a used appliance is being put into operation. This is the only way to ensure that all personal data has been removed and the previous owner will no longer be able to access the appliance.

Copyright and licences

For the purpose of operating and controlling the communication module, Miele uses proprietary or third-party software that is not covered by open source licensing terms. This software/these software components are protected by copyright. The copyrights held by Miele and third parties must be respected.

Furthermore, this communication module contains software components which are distributed under open source licensing terms. The open source components contained in the machine along with the corresponding copyright notices, copies of the licensing terms valid at the time and any further information can be accessed locally by IP using a web browser (<https://<IP address>/Licenses>). The liability and warranty arrangements for the open source licences displayed in this location only apply in relation to the respective rights holders.

en - Warnings and safety notes

To avoid the risk of accidents and damage to this communication module please read these instructions carefully before using it for the first time.

Keep these instructions in a safe place.

 Please also read the operating instructions for your dishwasher or the washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ This communication module must only be used for applications specified in the operating instructions. Changes or modifications to it and incorrect use of it are not permitted as this can be dangerous.
- ▶ The communication module must only be installed, commissioned and serviced by Miele service or by a suitably qualified person. Repairs are only to be undertaken by Miele Service. Repair work by unqualified persons could be dangerous.
- ▶ Check the communication module carefully before installation for any visible damage. A damaged module could be dangerous and must not be installed.
- ▶ Before installing the communication module compare the connection data (voltage and frequency) on the data plate with that on the data plate of your washer-disinfector. If the data does not match, the module could sustain damage. If in any doubt, consult an electrician.
- ▶ The communication module housing must not be opened on any account.
Touching live electrical components or altering the electrical or mechanical components is dangerous and could cause the module to malfunction.
- ▶ Make sure that the data cable for the communication module is assembled and routed correctly. The cable must not be kinked or trapped, for example. It must not be routed around sharp edges or be placed under tension.

Installation

⚠ Before fitting the communication module compare the connection data (voltage and frequency) on the data plate with that on the data plate of your cleaning machine. If the data does not match, the module could sustain damage. If in any doubt, consult a qualified electrician.

Module slot

- Disconnect the cleaning machine from the mains electricity supply.

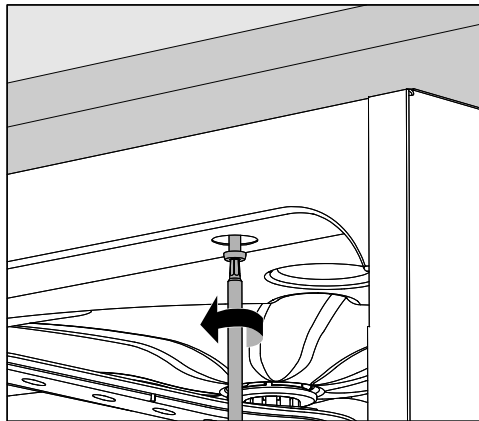
The module is inserted into the module slot on the cleaning machine. The module slot is located on the back of the cleaning machine.

- If the cleaning machine is installed in a niche:

- Pull the cleaning machine forwards out of the niche.

- If the cleaning machine is installed under a worktop:

- Open the door.


- Unscrew the fixing screws.
- Pull the cleaning machine out from under the worktop. Take care not to damage the hoses and cables when moving the machine.

en - Fitting and installation

Inserting the module

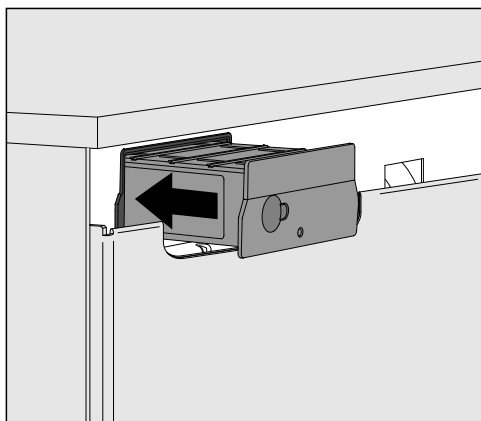- Remove any stickers or protective caps as necessary from the slot aperture.

- Push the module into the module slot. Gently push the module until it clicks into place.

Laying data cables

- Take care to ensure that the cable is laid correctly; it should not be kinked or trapped, for example. Keep the cable away from sharp edges and do not place it under tension.

Only terminal devices (PCs, printers, etc.) which comply with EN/IEC 62368 should be used.

Configuring the interface

⚠ Unauthorized access poses a risk.

Settings in the cleaning machine, e.g. parameters for dispensing process chemicals, may be changed as a result of unauthorised access via the network.

Under no circumstances should it be possible to access the machine via the Internet or other public or unsecured networks, either directly or indirectly (e.g. using port forwarding).

⚠ Please also read the operating instructions for your dishwasher or the washer-disinfector and pay particular attention to the Warning and Safety instructions.

Switching the cleaning machine on

- Connect the cleaning machine to the electricity supply.
The module is supplied with mains voltage by the cleaning machine.
No additional power connection is required.

After around 90 s, the module is ready for operation and the interface can be configured.

⚠ Only specialists are permitted to configure the interface.

⚙ Extended settings

Networking

COM module selection

WiFi

LAN

► COM module selection ► LAN

WiFi / LAN

Connection status

IP configuration

Network printer

Web interface

The XKM 3200 WL LMD86 module has a dedicated web interface via which the network settings can be configured, printers can be set up or the password can be changed.

The interface can be accessed from any workstation on your local network. A configured interface in the controller of the cleaning machine, an up-to-date web browser and an existing network connection to the cleaning machine are required.

The output language is English.

- Enter one of the following addresses in the address bar of your browser (where *<IP-address>* corresponds to the IP address of your cleaning machine).

https://<IP-address>/index.html

Password

When the web interface is opened for the first time, the “Admin” user must set up a password for their profile (once only). A module reset will reset the module settings to the factory default settings; this will include resetting the password for the “Admin” user.

- Change Password

The “Admin” user and the password must always be set at the factory as a first step before further settings can be made. When setting the password for the first time, the field for the old password must be left blank.

Login

User

Password

Login Logout

▼ Change Password

Old Password

New Password

New Password

Apply

Remote Update

Type

Version

New Version

Release Notes

No Update Available

Restart Module

Reset to Factory Defaults

Apply

Network Configuration

Security

SSID

SSID

DHCP ☒

IP Address

Subnet Mask

Gateway

DNS Auto ☒

DNS Server 1

DNS Server 2

Network Printer

Auto Print ☐

Security

Port

URI

Customer Name

Language ▾

Device Language

Format

Duplex

Protocoll Data

Batch Number

Format

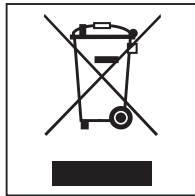
Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

Disposing of your old appliance

Electronic and electrical appliances contain many valuable materials. They also contain certain materials, compounds and components which were essential for their correct functioning and safety. These could be hazardous to your health and to the environment if disposed of with general waste or if handled incorrectly. Please do not, therefore, dispose of your old appliance with general waste.

Instead, please make use of officially designated collection and disposal points to dispose of and recycle electrical and electronic appliances in your local community, with your dealer, with Miele or with the manufacturer. By law, you are solely responsible for deleting any personal data from the appliance prior to disposal. You are legally obliged to remove any old batteries which are not securely enclosed by the appliance and to remove any lamps without destroying them, where this is possible. These must be taken to a suitable collection point where they can be handed in free of charge. Please ensure that your old appliance poses no risk to children while being stored prior to disposal.

Napomene uz upute	60
Pitanja i tehnički problemi	60
Definicija pojmova	60
Modul	60
Opis namjene	61
Sadržaj isporuke	62
Sadržaj isporuke	62
Tehnički podaci	62
Izjava o sukladnosti	62
Opis uređaja	63
Pregled uređaja	63
Integracija u mrežu	63
Uključivanje i isključivanje	64
Kontrolni LED	64
Tipka za resetiranje	64
Zaštita podataka i sigurnost podataka	65
Konfiguracija mreže Tvorničke postavke	65
Autorska prava i licence	65
Sigurnosne napomene i upozorenja	66
Ugradnja i instalacija	67
Ugradnja	67
Otvor za modul	67
Postavljanje modula	68
Polaganje podatkovnog kabela	68
Umrežavanje	69
Konfiguracija sučelja	69
Uključivanje uređaja za pranje	69
Internetsko sučelje	70
Zaporka	70
Vaš doprinos zaštiti okoliša	72
Zbrinjavanje transportne ambalaže	72

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