
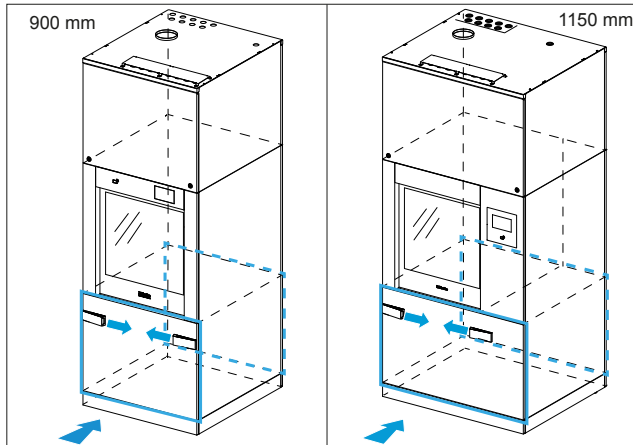
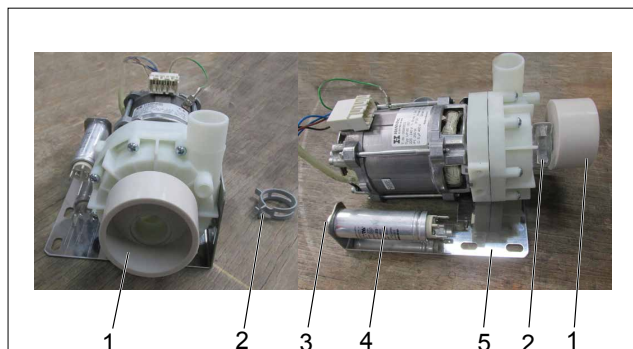
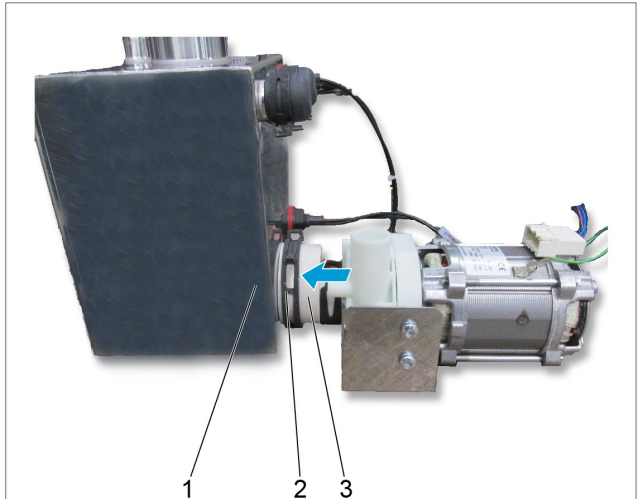
de	Ablaufpumpe (DP)	en	Drain pump (DP)	cs	Vypouštěcí čerpadlo (DP)
da	Afløbspumpe (DP)	el	Αντλία αποχέτευσης (DP)	en-US	Drain pump (DP)
es	Bomba de desagüe (DP)	fr	Pompe de vidange (DP)	it	Pompa di scarico (DP)
nl	Afvoerpomp (DP)	no	Avløpspumpe (DP)	pt	Bomba de esgoto (DP)
sv	Avloppspump (DP)				

1

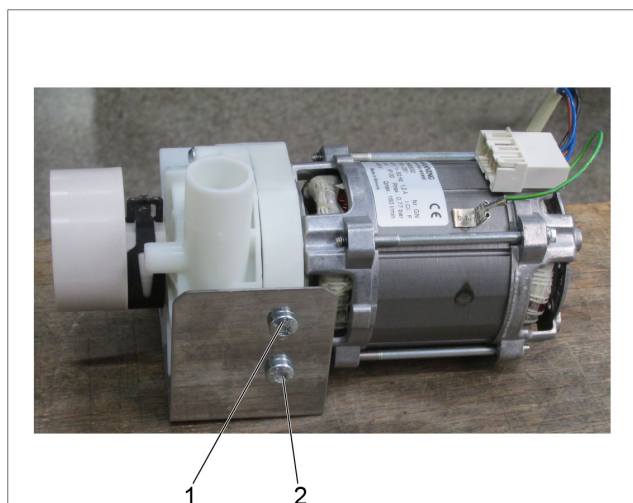
2


3

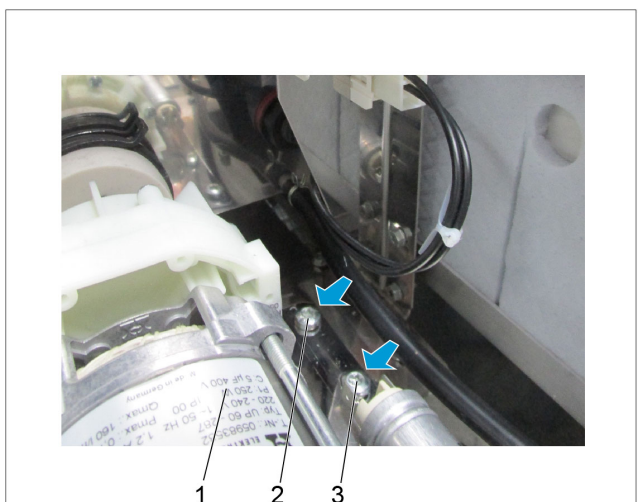
4

5

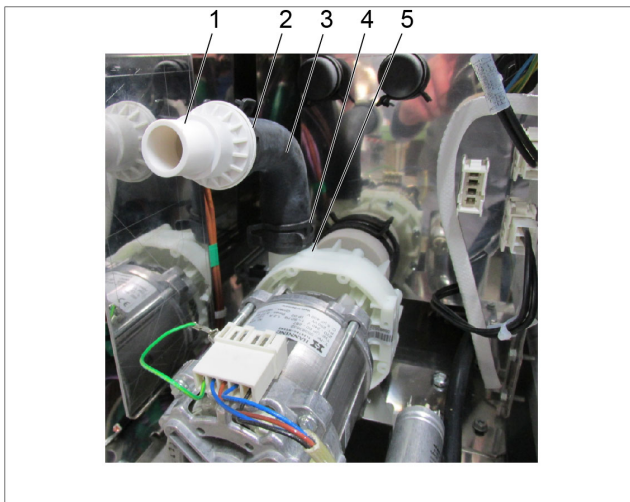6

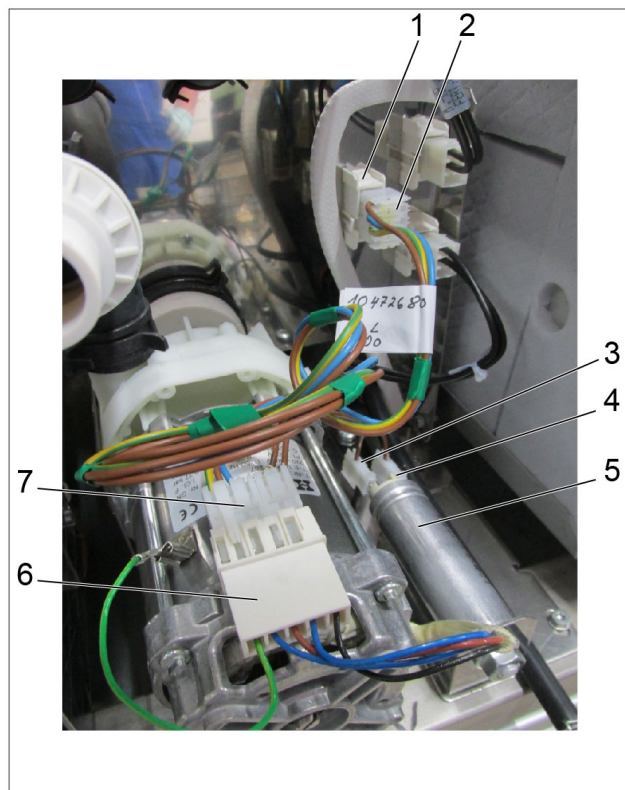7

de

Ablaufpumpe (DP)

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In order to carry out this work, the appropriate Technical Service Documentation (TSD) is required.

The warning and safety instructions given in the applicable model-specific TSD must be complied with.

Work must be carried out in accordance with the instructions given in the applicable model-specific TSD.

Disconnect the machine from the mains and ensure power cannot be switched on again in error.

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Do not lay cables over sharp edges or hot surfaces.

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Bottom part service panel opening

The bottom part service panel is secured with 2 bolts which each clip into the frame at the sides of the machine.

✚ Slide the service panel retaining bolts towards the centre of the machine at the same time, see Fig. 1.

✚ Open the service panel slightly at the top and remove it.

Drain pump fitting

Drain pump fitting preparatory work

✚ Fit the spring clip (2) on the small side of the adapter (1), see Fig. 2.

✚ Slide the adapter (1) onto the drain pump connection and secure it with the spring clip (2), see Fig. 2.

✚ Fit the interference suppression capacitor (4) with the M8 nut (3) to the holding bracket (5), see Fig. 2.

✚ Fit the drain pump on the holding bracket (5), see Fig. 2.

✚ Fit the holding bracket with two fixing screws (1) and (2) onto the drain pump, see Fig. 3.

Fitting drain pump in machine

✚ Fit the large side of the adapter (3) onto the drain connection of the intermediate container (1) and secure it with a hose clip (2), see Fig. 4.

✚ Fit the drain pump (1) with the fixing screws (2) and (3), see Fig. 5.

✚ Insert the non-return valve (1) in the connection hose (3) to the drain pump and secure it with a spring clip (2), see Fig. 6.

✚ Fit the connection hose (3) on the drain pump drain connection (5) and secure it with a hose clip (4), see Fig. 6.

Drain pump minor wiring harness fitting

- ✚ Connect the connection plug (7) in the connection socket (6) of the drain pump, see Fig. 7.
- ✚ Connect the connection plugs (3) and (4) to the connections of the interference suppression capacitor (5), see Fig. 7.
- ✚ Connect the connection plug (2) in the connection socket (1) of the machine wiring harness, see Fig. 7.

Splash protection plate fitting

- ✚ Fit the splash protection plate (3) with the fixing screws (1) and (2) to the machine frame, see Fig. 8.

Drain hose fitting

- ✚ Fit the drain hose onto the non-return valve (4) and secure it with a hose clip, see Fig. 8.
- ✚ Lay the drain hose to the on-site drain.

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en-US

Drain pump (DP)

Reason: This conversion kit enables the drain pump (DP) option to be installed in models PLW 8615, PLW 8616, PLW 8617, PWD 8626 or PWD 8628.

Parts required

Quantity	Mat. no.	Designation
1	10931330	Drain pump (DP - Drain pump) PLW/PWD 86xx

Parts included

Quantity	Mat. no.	Designation
1	10469320	Adapter
1	07651460	Hose clip FBS 60/12 (suds container adapter)
1	05983532	Pressure pump UP60 220 - 240 V 50Hz
1	09973120	Mounting bracket (for drain pump)
4	06499370	Raised-head screw M6 x 12 (for drain pump and mounting bracket)
1	10714411	Protective panel
2	07246730	DM 4x10 bolt
1	01022310	Edge protection (0.150 m)
1	06847870	Connection hose
2	05146920	Hose clip DN34 15 mm
1	05433341	Capacitor, 5µF 400V
1	07638120	M8 nut (interference suppression filter)
1	04127480	Hose clip 29 (pump connection hose)
1	06496094	Non-return valve
1	09706850	Drain hose, 1.5 m
2	01627501	Hose clip, 19-31
1	10472680	Minor wiring harness, water drainage, 8615-28

The TSD will be needed for all service work

The warning and safety instructions in the model-specific TSD must be observed.

Service work must be carried out following the instructions in the model-specific TSD.

Disconnect the machine from power and ensure that power cannot be switched on again in error.

Wear protective gloves and goggles.

Do not lay cables or wires over sharp edges or hot surfaces.

Lay the hoses without kinking or straining them.

Opening the bottom service panel

The bottom service panel is secured with 2 bolts which each clip into the frame at the sides of the machine.

- ✚ Slide the service panel retaining bolts towards the center of the machine at the same time; see Fig. 1.
- ✚ Open the service panel slightly at the top and remove it.

Drain pump installation**Preparing the drain pump for installation**

- ✚ Install the spring clip (2) on the small side of the adapter (1); see Fig. 2.
- ✚ Slide the adapter (1) onto the drain pump connection and secure it with the spring clip (2); see Fig. 2.
- ✚ Secure the interference suppression capacitor (4) to the mounting bracket (5) with the M8 nut (3); see Fig. 2.
- ✚ Install the drain pump on the mounting bracket (5); see Fig. 2.
- ✚ Install the mounting bracket with two retaining screws (1) and (2) on the drain pump; see Fig. 3.

Installing the drain pump in the machine

- ✚ Install the large side of the adapter (3) onto the drain connection of the intermediate container (1) and secure it with a hose clip (2); see Fig. 4.
- ✚ Install the drain pump (1) with the retaining screws (2) and (3); see Fig. 5.
- ✚ Insert the non-return valve (1) in the connection hose (3) to the drain pump and secure it with a spring clip (2); see Fig. 6.
- ✚ Install the connection hose (3) on the drain pump drain connection (5) and secure it with a hose clip (4); see Fig. 6.

Installing the drain pump minor wiring harness

- ✚ Connect the connection plug (7) in the connection socket (6) of the drain pump; see Fig. 7.
- ✚ Connect the connection plugs (3) and (4) to the connections of the interference suppression capacitor (5); see Fig. 7.
- ✚ Connect the connection plug (2) in the connection socket (1) of the machine wiring harness; see Fig. 7.

Splash protection plate installation

- ✚ Install the splash protection plate (3) with the retaining screws (1) and (2) to the machine frame; see Fig. 8.

Drain hose installation

- ✚ Install the drain hose on the non-return valve (4) and secure it with a hose clip; see Fig. 8.
- ✚ Lay the drain hose to the on-site drain.

Drain pump activation

- ✚ Connect the service computer to the machine.
- ✚ Log in to the Web User Interface (WebUI) with a service technician with administrator rights.
- ✚ Request write rights in the WebUI.
- ✚ Select the System menu.
- ✚ Select the Machine configuration menu.
- ✚ Select the Drainage submenu.
- ✚ Select the Drain pump option parameter.

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