# PLW 8683

**en** Installation Plan

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**en** ..... 4  
**fr** ..... 17

<b>Information on the installation plan</b> .....	<b>4</b>
Definition of terms .....	4
<b>Appliance dimensions</b> .....	<b>5</b>
<b>Connections</b> .....	<b>6</b>
<b>Installation notes</b> .....	<b>9</b>
Surrounding area .....	9
Vapor barrier for built-under appliances .....	9
Electrical connection .....	9
Equipotential bonding and grounding .....	9
Electromagnetic compatibility (EMC) .....	10
Network connection .....	11
Water connection .....	12
Water inlet .....	12
Drainage .....	12
<b>Technical details</b> .....	<b>13</b>
Electrical connection .....	13
Water inlet .....	13
Drainage .....	14
Dimensions .....	14
Weight .....	14
Emission values .....	15
Heat dissipation .....	15
Sound pressure level .....	15
Operating conditions .....	15
Storage and transportation conditions .....	15

## en - Information on the installation plan

---

For safe installation and commissioning of the laboratory glassware, please read the installation plan, the service documentation, and the operating instructions.

This installation plan includes the dimensions of the laboratory glassware, the technical data, and the requirements to be met by the customer for installation.

The dimensions in the figures are given in mm. Dimensions that deviate from this are labeled accordingly.

The machine features vary depending on the model and can be extended with additional options. The installation plan describes the maximum amount of possible features. Individual connections are not required for each item of optional equipment.

### Installation requirements

Installation and commissioning work should be performed by Miele Customer Service or a suitably authorized technician in accordance with local and national safety regulations.

Installation must comply with valid regulations, legal requirements, relevant standards, and accident prevention regulations.

### Definition of terms

#### The machine

In this installation plan, the laboratory glassware is referred to as “the machine”.

#### Wash water

The term “wash water” refers to water or to a mixture of water and process chemicals.

#### Process chemicals

All media dispensed during a program sequence are generally referred to as process chemicals, e.g., cleaning agents.

#### Device variants



SD stainless steel door  
GD glass door

# en - Appliance dimensions



SD

GD

Front view

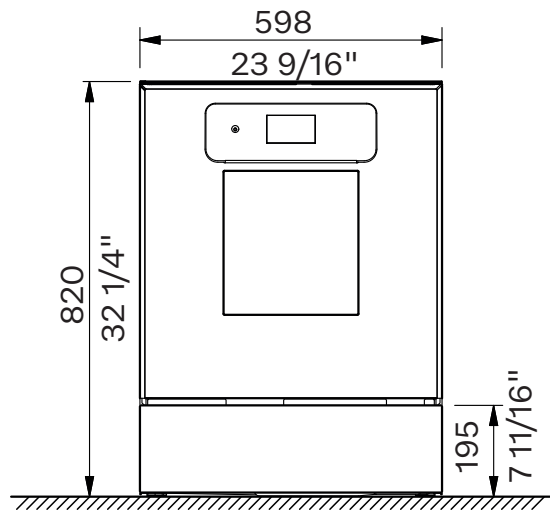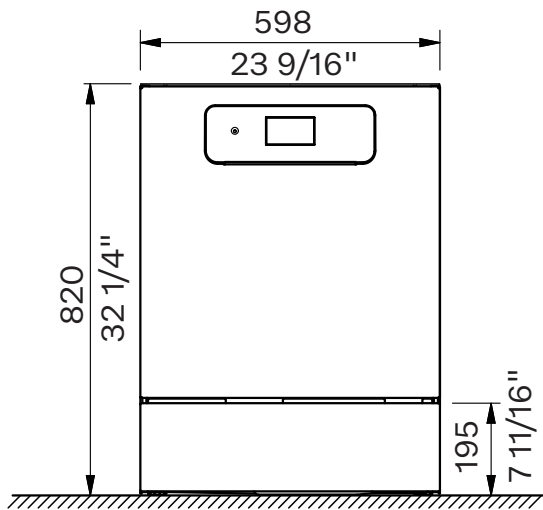
Side view
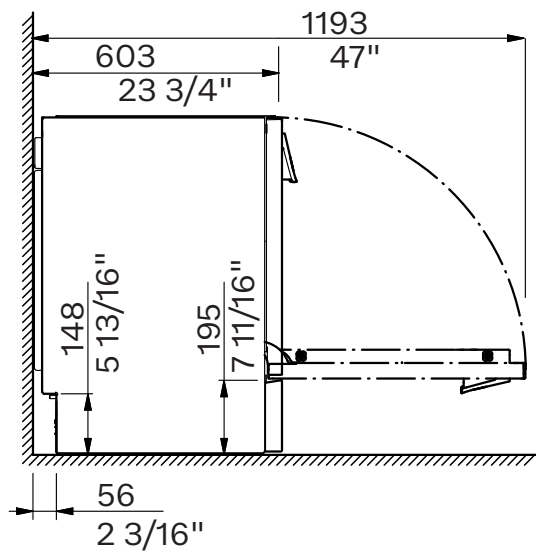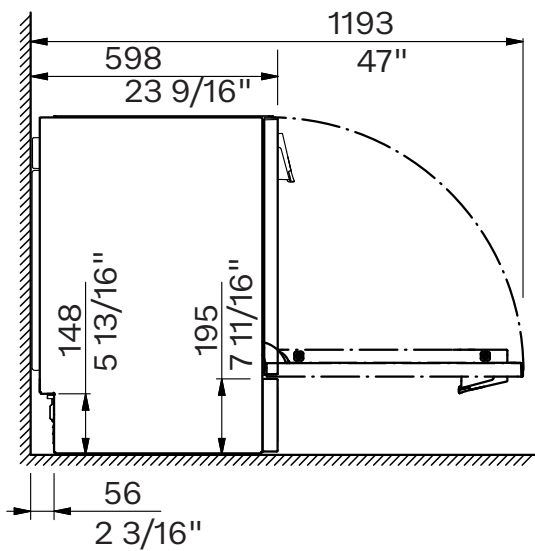
## en - Connections


---

### Rear view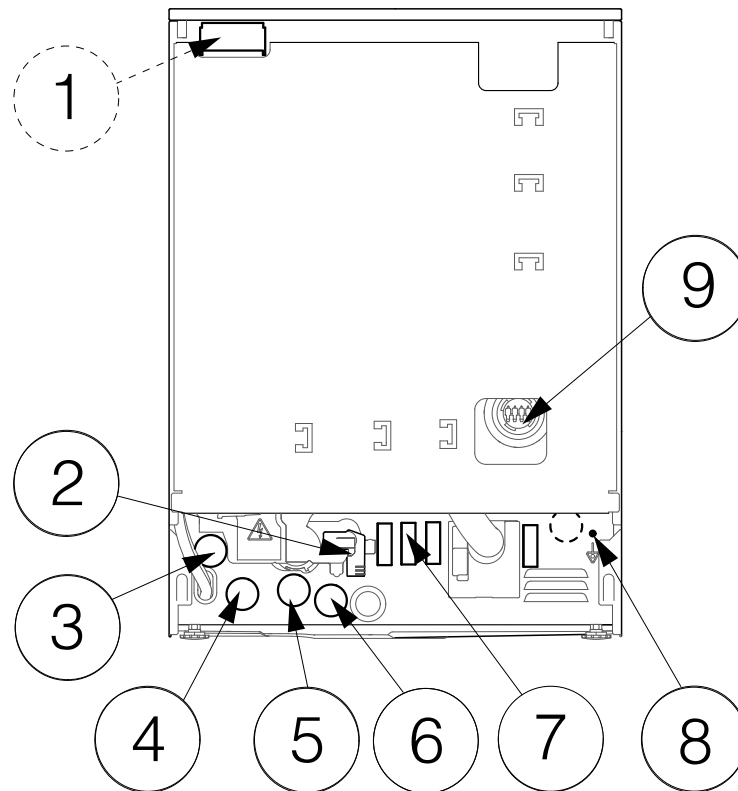


- ① Network connection (optional)
- ② Waste water
- ③ DI water
- ④ Cold water
- ⑤ Power supply cord
- ⑥ Hot water
- ⑦ External dispensing, power supply connection
- ⑧ Equipotential bonding and grounding
- ⑨ External dispensing, connection for dispensing hose

On-site connections

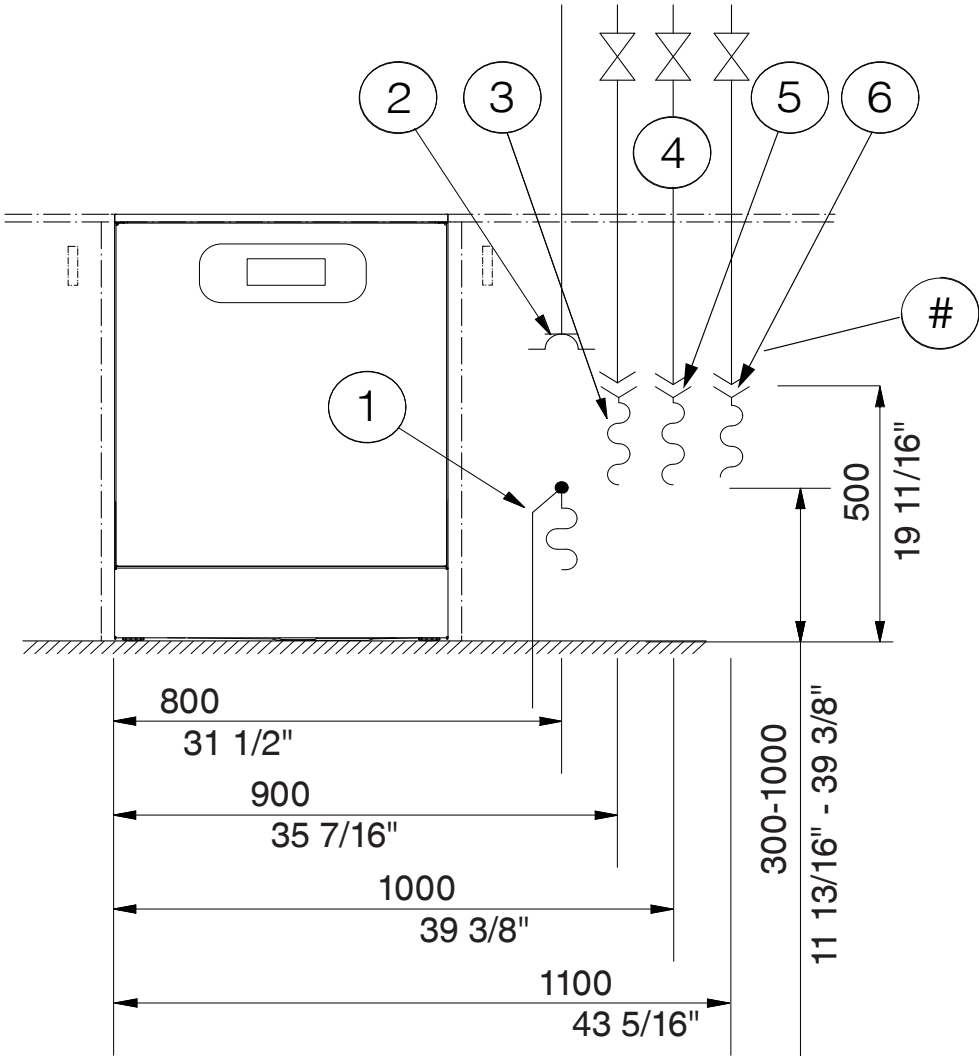
- ① Waste water
- ② Power supply cord
- ③ DI water
- ④ Shut-off valves
- ⑤ Cold water
- ⑥ Hot water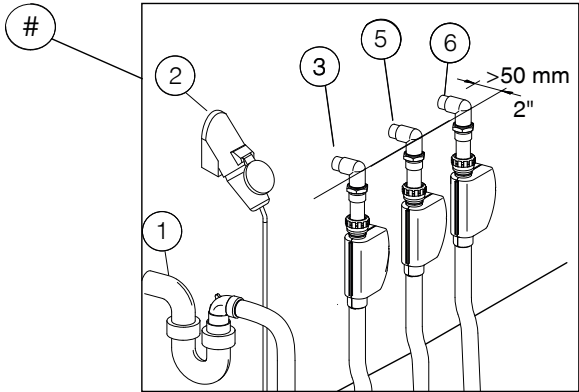

## en - Connections

### External dispensing

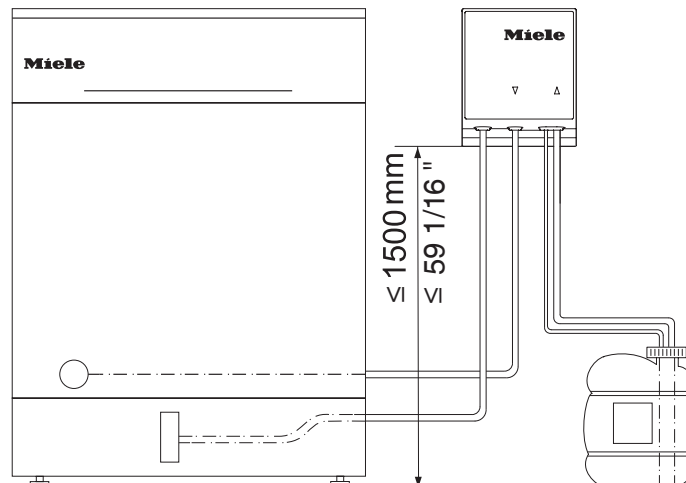Max. delivery head	4' 11" (1.5 m)
Length of dispenser hose, DOS module to suction wand	5' 11" (1.8 m)
Length of dispenser hose, back of appliance to DOS module	9' 2" (2.8 m)
Length of power cord, back of appliance to DOS module	9' 2" (2.8 m)

Place the container on the floor next to the machine or in an adjacent cabinet.  
The container must not be placed on top of or above the machine.

### Surrounding area

Condensate can build up in the area surrounding the machine. Any cabinetry and fittings in the room must, therefore, be suitable for purpose.

### Vapor barrier for built-under appliances

The enclosed vapor barrier protects the countertop from damage caused by water vapor that may escape when the door is opened. Place the vapor barrier above the door under the countertop.

### Electrical connection

All work on the electrical connection must be carried out by Miele Service or a qualified electrician.

The customer's installation of the machine must comply with over-voltage category CAT II in accordance with IEC 61010-1.

The voltage may deviate from the rated voltage by max.  $\pm 10\%$ .

### Power switch

Install a power switch capable of disconnecting the machine at all poles at the customer's site. The power switch must:

- Have a contact gap of at least 1/8" (3 mm).
- Be designed to operate at the rated current of the machine.
- Be able to be locked in the zero position.
- Be accessible after the machine has been installed.

The power cord must be laid protected from the risk of thermal damage.

### Plug connection

The machine should be connected to the electricity supply via a electrical outlet.

### Hard-wired

Alternatively, the machine can be connected via a hard-wired connection.

### Residual current device (RCD)

For added safety, each machine should be protected by a type A residual current device with a trip current of 30 mA.

The residual current device must be installed so that it is easily accessible.

### Equipotential bonding and grounding

The machine has been set up for equipotential bonding and grounding. The connection screw (M 8) is located on the rear of the machine. Equipotential bonding and grounding should be carried out if possible on site.

### **Electromagnetic compatibility (EMC)**

The machine has been tested for electromagnetic compatibility (EMC) in accordance with EN 61326-1 and is suitable for operation in commercial environments, such as hospitals, medical practices, and laboratories plus other similar environments which are connected to the domestic electrical supply.

The machine's high frequency (HF) energy emissions are very low and are therefore unlikely to interfere with other electronic machines in the vicinity.

Flooring in the installation area must be wood, concrete, or tiled. Synthetic flooring must be able to withstand a relative humidity level of at least 30% to minimize the risk of electrostatic discharges.

The quality of the power supply should comply with that found in a typical commercial or hospital environment. Check that the power supply voltage is within a range of +/-10% of its nominal value.

### Network connection

The machine can be integrated into a customer's network via an Ethernet interface.

The operator is responsible for the secure configuration and operation of the network. Creating a secure configuration reduces the scope for the machine to be attacked. Non-secure configurations can make it easier for potential attackers to gain unauthorized access.

 Unauthorized access poses a risk.

Settings in the machine may be changed as a result of unauthorized access via the network.

Under no circumstances should it be possible to access the machine via public or unsecured networks, either directly or indirectly (e.g., using port forwarding).

The IP addresses from 192.168.10.1 to 192.168.10.255 are reserved for Miele Service. They must never be set in the internal network.

Please also note the following recommendations for the network configuration of the machine:

1. Configure the network in which the machine is located with as many restrictions as possible:
  - Only allow machines or persons access to the network if absolutely necessary.
  - Use a VLAN for network segmentation, for example.
2. Make sure to create a secure configuration for all machines connected to the network, e.g., with the help of:
  - The information on network security in the operating instructions for the connected machines.
  - The recommendations of USA: Cybersecurity & Infrastructure Security Agency (CISA - [www.cisa.org](http://www.cisa.org)), Canada: Communications Security Establishment Canada (<https://www.cse-cst.gc.ca>).
3. Pay particular attention to the following areas:
  - user administration
  - staff entering and exiting
  - authorization
  - authentication
  - update management, etc.

### Water connection

#### Water inlet

The machine must only be connected to fully vented pipework. A brief increase in the water pressure can damage components of the machine.

The quality of the incoming water must at least correspond to the drinking water regulations of the country in which the machine is being operated.

The machine must be connected to cold and hot water. If hot water is not available, the inlet hose for hot water must be connected to the cold water supply. For this purpose, a Y-piece can be used on the cold water inlet on the customer side.

To be able to use the programs supplied by the factory, a DI water supply is required.

The connections with shut-off valves at the customer's site must have 3/4" (MGHT) outer threads that seal the surface.

The shut-off valves must be accessible after the machine has been installed.

#### DI water ring line

The machine can be connected to a ring line system for DI water. For this purpose, the machine must be technically adapted and the controls reset by Miele Service or an authorized technician.

#### Drainage

The machine should be connected to a separate on-site drainage system for the machine only. If a separate connection is not available, it is recommended to connect the hose to a dual-chamber siphon. The drain lines must be heat resistant to at least 200°F (93°C).

If the hose is to be directly connected to the drainage system on site, use the hose clip supplied in the accessory pack.

The on-site connector for the drain hose can be used for various hose diameters. If the connector extends more than 1 1/4" (30 mm) into the drain hose, it must be shortened. Otherwise, the drain hose can become clogged.

Lay the drain hose so that it does not kink and is not being subjected to pressure or tugged.

If the on-site drain connection is situated lower than the guide path for the lower basket rollers in the open door, a siphoning effect during a program can cause the appliance to empty itself of water. In this case, lay the drain hose with a bend in it so that its highest point is at least level with the guide path for the lower basket rollers.

## Electrical connection

	Standard	Convertible *
Voltage	3 AC 208 V	2 AC 208 V
Frequency	60 Hz	60 Hz
Fuse rating	3 x 20 A	2 x 30 A
Connection	NEMA L15-20P	NEMA L6-30P
Power cord length	1.9 m/6.2 ft	1.9 m/6.2 ft
Power cord cross-section	4 x AWG12	3 x AWG10
Heater rating	5.4 kW	5.4 kW
Total rated load	6.2 kW	6.2 kW

\* conversion by Miele Service

## Water inlet

Water connections		
- Cold water	1	1
- Hot water	1	1
- DI water	1	1
Water temperature		
- Cold water	5–20 °C	41–68 °F
- Hot water	5–65 °C	41–149 °F
- DI water	5–20 °C	41–68 °F
Connection hose length	1.7 m	51 3/16"
Connection at customer's site with shut-off valve, outer thread that seals the surface	3/4" Male Garden Hose Thread MGHT	3/4" Male Garden Hose Thread MGHT
Water connection pressure	200–1000 kPa	29–145 psi
Flow rate	2.0-7.5 l/min	0.5–1.98 gpm
Water hardness	0 – 10.7 mmol/l	0 – 62.56 gpg
Chloride content	<100 mg/l	<5.85 gpg sodium chloride
DI water		
- Recommended conductivity	<15 µS/cm	<0.0015 Ohms
- pH value	5–8	5–8

## en - Technical details

### Drainage

Drain hose length	1.4 m	55 1/8"
Max. drain hose length	4 m	157 1/2"
Max. delivery head from floor	1 m	39 3/8"
Max. water temperature	93 °C	200 °F
Max. transient flow rate	16 l/min	4.23 gpm
Hose inner diameter (Ø)	22 mm	7/8"
Drain line	≥ DN 50	≥ DN 50

### Dimensions

SD		
Height	820 mm	32 1/4"
Height adjustment	+60 mm	+2 3/8"
Width	598 mm	23 9/16"
Depth	598 mm	23 9/16"
Depth with control panel	639 mm	25 3/16"
Depth with open door	1193 mm	47"
GD		
Height	820 mm	32 1/4"
Height adjustment	+60 mm	+2 3/8"
Width	598 mm	23 9/16"
Depth	603 mm	23 3/4"
Depth with control panel	644 mm	25 3/8"
Depth with open door	1193 mm	47"

### Weight

SD		
Net weight	72 kg	159 lbs
Max. operating weight	128.5 kg	283 lbs
Maximum load capacity of open door	37 kg	82 lbs
Max. individual load (foot Max. individual load (foot Ø 38 mm/1.5"))	320 N	320 N
Max. distributed load (load-bearing capacity)	3.59 kN/m <sup>2</sup>	3.59 kN/m <sup>2</sup>
GD		
Net weight	78 kg	172 lbs
Max. operating weight	137 kg	302 lbs
Maximum load capacity of open door	37 kg	82 lbs
Max. individual load (foot Max. individual load (foot Ø 38 mm/1.5"))	340 N	340 N
Max. distributed load (load-bearing capacity)	3.83 kN/m <sup>2</sup>	3.83 kN/m <sup>2</sup>

## Emission values

### Heat dissipation

Heat output into the room	5.4 MJ/h	5118.2 BTU
---------------------------	----------	------------

### Sound pressure level

Sound pressure level LpA during cleaning and drying	<65 dB(A) re 20 µPa	
---	---------------------	--

## Operating conditions

Ambient temperature	5–40°C	40–105°F
Relative humidity:		
Minimum	10%	10%
Maximum for temperatures up to 201°F (31°C)	80%	80%
Decreasing linearly for temperatures up to 104°F (40°C)	50%	50%
Max. altitude above sea level	2,000 m *	6,561 ft *

If installed above 4,921.26 ft/1,500 m, the boiling point of the wash water will be lower. In this case, Miele Service must adjust the disinfection temperature and the holding time.

## Storage and transportation conditions

Ambient temperature	-20 – +60 °C	-4 – +140 °F
Relative humidity	10–85 %	10–85 %
Air pressure	500–1060 hPa	500–1060 hPa
Weight including packaging max.	91.5 kg	202 lbs
Maximum floor load, transport route	935 kN	935 kN

## Site-access dimensions including transport pallet

Height	920 mm	36 1/4"
Width	670 mm	26 3/8"
Depth	740 mm	29 1/8"

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**U.S.A.**
**Miele, Inc.**
**Miele Limited | limitée**