ADP 86

de	Montageanweisung Pumpenmodul
en	Installation instructions Pump module
es	Instrucciones de montaje Módulo de bombeo
fr	Notice de montage Module de pompe
it	Istruzioni di montaggio Modulo pompa

de	4
en	19
es	34
fr	49
it	64

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en - Contents

Notes about these instructions	19
Symbols and conventions used in this document.....	19
Queries and technical problems	19
Definition of terms	19
Appropriate use	20
General description.....	20
How it works	20
Purpose.....	20
Contraindications	20
Intended user group.....	20
Foreseeable misuse	20
Installation requirements	20
Items supplied	22
Scope of delivery	22
Technical data	23
Product description	24
Appliance overview	24
Warnings and safety notes	25
Installation	27
Installation site	27
Initial commissioning.....	27
Fitting demineralised water hoses.....	27
Fitting the connecting hose to the demineralised water container.....	27
Fitting the demineralised water inlet hose to the cleaning machine	28
Connection to the control unit of the cleaning machine	28
Electrical connection	28
Unplugging the demineralised water hoses from the module.....	29
Disconnecting the connection hose from the demineralised water container	29
Disconnecting the water inlet hose for the cleaning machine.....	29
Operation	30
Switching on and off	30
Fault messages during the pumping process.....	30
Decommissioning	31
Defect.....	31
Permanent decommissioning.....	31
Caring for the environment	32
Disposal of the packing material	32

Symbols and conventions used in this document

Warnings

 Warnings contain information which is important for safety. This alerts you to the potential danger of injury to people or damage to property.
Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Notes provide information of particular importance that must be observed.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

■ Select an option using the arrow buttons and save your choice with *OK*.

Example:

■ Select an option.

Display

Display text can be identified from the special font.

Example:

Save.

Queries and technical problems

In the event of queries or technical problems, please contact Miele. Contact details can be found at the end of the operating instructions for your cleaning machine or at www.miele.com/professional.

Definition of terms

Cleaning machine

In these installation instructions, the washer-disinfector is referred to as “the cleaning machine”.

General description

The pump module is an unfinished machine designed for pumping demineralised water and ultrapure water.

How it works

The pump module is intended exclusively for operation with washer-disinfectors manufactured by Miele. The module is controlled via the washer-disinfector control system.

Purpose

This module is designed exclusively for commercial use.

The pump module can only be operated in combination with Miele washer-disinfectors that can be connected to a ring line system for demineralised water.

Contraindications

The pump module is not suitable for pumping water contaminated with process chemicals or dirt particles or water contaminated in any other way.

Intended user group

All users without exception must be instructed in how to use the module by Miele Customer Service, by a customer service team specially trained and authorised by the manufacturer or by a person who has already received instruction.

Foreseeable misuse

Improper use can be caused by unsuitable water quality, the addition or pumping of (process) chemicals or operation with contaminated water.

Failure of the operator to comply with routine checks and regular service intervals.

Failure to observe the specified installation requirements.

Installation requirements

Installation site


The pump module is intended for installation in laboratories and laboratory-like rooms.

The installation must be carried out in rooms in which ambient conditions meet the following requirements:

- Draught-free and dry
- Equipped with suitable room ventilation
- Solid and even surface, observe floor load-bearing capacity
- No direct sunlight

Conditions of use	Operation (according to IEC/EN 61010-1):	5 °C to 40 °C
	Ambient temperature	80 % for temperatures up to 31 °C
	Max. relative humidity	50 % for temperatures up to 40 °C
	linear decrease to	10 %
	Min. relative humidity	
	Altitude above sea level (according to IEC/EN 61010-1)	Up to 2000 m
	Background noise level	N/A

Scope of delivery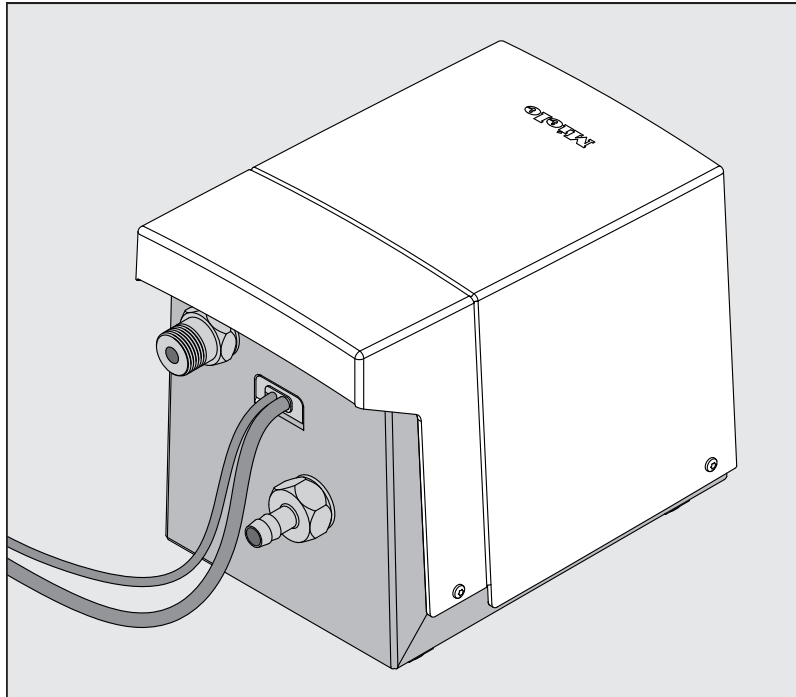


- ADP 86 pump module
- Retrofitting kit for potential-free contact

Height	240 mm
Width	190 mm
Depth	315 mm
Weight	6 kg
Voltage, rated load, fuse rating	See data plate
Mains connection cable	Approx. 2,5 m
Hose sleeve	Ø 13 mm
Operation (according to IEC/EN 61010-1): Ambient temperature Max. relative humidity linear decrease to Min. relative humidity	5 °C to 40 °C 80 % for temperatures up to 31 °C 50 % for temperatures up to 40 °C 10 %
Storage and transportation conditions: Ambient temperature Relative humidity Air pressure	-20 °C to 60 °C 10 % to 85 % 500 hPa to 1060 hPa
Altitude above sea level (according to IEC/EN 61010-1)	Up to 2000 m
Protection category (according to IEC 60529)	IP21
Degree of soiling (according to IEC/EN 61010-1)	2
Overvoltage category (according to IEC 60664)	II
VDE radio suppression, EMC equipment class (according to EN 61236-1)	B
VDE electrical safety	IEC/EN 61010-1, IEC 61010-2-040
CE marking	Machinery Directive 2006/42/EC *)
Manufacturer address	Miele & Cie. KG, Carl-Miele-Straße 29, 33332 Gütersloh, Germany

*) Unfinished machine

Appliance overview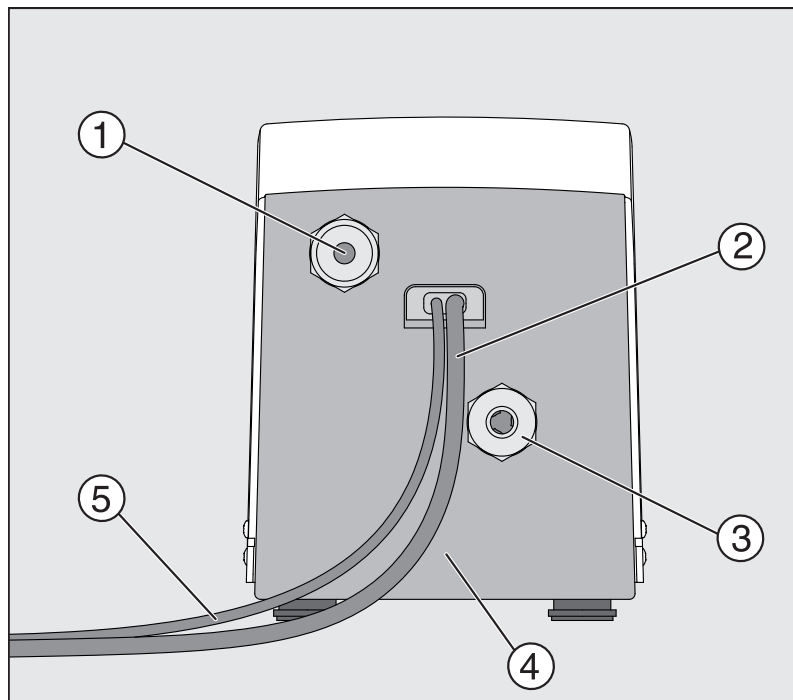

- ① Connection thread for the cleaning machine's demineralised water inlet hose
- ② Mains connection cable
- ③ Hose sleeve for the connecting hose for the demineralised water container
- ④ Data plate
- ⑤ Control lines for connection to the cleaning machine's control system

Read the installation instructions carefully before using this module. Keep these installation instructions in a safe place.

 Please also read the operating instructions for your dishwasher or the washer-disinfector and pay particular attention to the Warning and Safety instructions.

- ▶ The pump module is only approved for use in the applications stated in these installation instructions. Alterations or conversion of the dispensing module, or using it for purposes other than those for which it was intended, are not permitted and could be dangerous.
- ▶ The pump module is only designed for operation with demineralised water or ultrapure water.
The use of organic solvents or inflammable liquids is not permitted. This could cause an explosion or damage rubber or plastic components in the machine and cause liquids to leak out.
- ▶ The pump module is intended for stationary indoor use only.
- ▶ The pump module may only be installed, commissioned and maintained by Miele Customer Service or by a qualified specialist authorised by the module manufacturer. Repairs may only be carried out by Miele Customer Service. Incorrect repairs can cause considerable danger to users.
- ▶ Do not install the pump module in an area where there is any risk of explosion or of freezing conditions.
- ▶ Check the pump module and its electrical cables for any visible damage before installation. A damaged module can jeopardise your safety and must not be installed under any circumstances.
- ▶ Before installing the pump module, compare the connection data (voltage and frequency) on the data plate with that on the data plate of your cleaning machine. If the data does not match, the module could sustain damage. If in any doubt, consult a qualified electrician.
- ▶ The electrical safety of the pump module can only be guaranteed when it is correctly earthed. It is essential that this standard safety requirement is observed and regularly tested. If in any doubt, please have the electrical installation inspected by a qualified specialist.
- ▶ For operation, the pump module must be installed horizontally on a stable surface with the machine feet facing downwards. Operation in an inclined position, lying on its side/upside down or with hose connections pointing upwards is not permitted.
- ▶ During installation, ensure that the mains connection cable and control cables running to the cleaning machine are laid correctly. The cables must not be kinked or trapped, for example. They must not be routed around sharp edges or be placed under tension.
- ▶ After installing the module, make sure that there is no risk of tripping posed by the module, the electrical cables or the connected hoses.
- ▶ To ensure that the pump is ventilated, the pump module must not be covered or placed in an airtight cabinet or box after commissioning.
- ▶ A damaged or leaking pump module is a safety risk. Take the pump module out of operation immediately and inform Miele Customer Service.

en - Warnings and safety notes

- ▶ Label pump modules that have been taken out of operation and secure them against unauthorised restarting. The device may only be put back into operation once it has been successfully repaired by Miele Customer Service or by appropriately qualified specialists.
- ▶ Never open the casing of the pump module. Any contact with live connections or tampering with the electrical or mechanical components can endanger your safety and may lead to module malfunctions.
- ▶ Disconnect the pump module from the power supply before any maintenance work is carried out.
- ▶ Never clean the pump module or near vicinity with a water hose or a pressure washer.

Water damage due to demineralised water.

Uncontrolled leakage of demineralised water can lead to water damage.

Use a connection hose with a suitable diameter for the connection to the demineralised water container (see  "Technical data") and secure it to the pump module with a hose clip.

Screw the water inlet hose for the cleaning machine firmly onto the thread of the pump module.

After initial commissioning and at regular intervals thereafter, check for water leaking from the hose connections.

Installation site

The pump module must be installed in the immediate vicinity of the cleaning machine so that the control lines and the water inlet hose to the cleaning machine can be laid without tension.

The installation site of the pump module must be lower than the outlet of the demineralised water container.

The surface must be stable and level.

A free electrical connection for the mains connection cable of the pump module must be available.

Danger of tripping.

The installation site must be selected so that neither the module nor the hoses or line pose a tripping hazard after the pump module has been installed.

Initial commissioning

Initial commissioning must be carried out by Miele Customer Service or by a qualified specialist authorised by the manufacturer of the pump module.

Customer Service must fit a potential-free contact to the cleaning machine and activate the pump module in the control unit of the cleaning machine.

Fitting demineralised water hoses

Fitting the connecting hose to the demineralised water container

The following items must be kept ready for fitting the connecting hose between the pump module and the demineralised water container:

- Connecting hose to demineralised water container; internal diameter max. 13 mm
- Hose clip for attaching the connecting hose to the pump module
- Suitable tool for fitting the hose clip
- Attach the hose clip to the connecting hose.
- Attach the connecting hose to the hose sleeve of the pump module.
- Use the hose clip to secure the hose to the hose sleeve.

The hose clip is designed to prevent the connecting hose from coming loose or being disconnected accidentally, e.g. when the module is moved during cleaning or maintenance work.

After installation, the connecting hose must not sag or be laid in loops; this is to prevent air collecting in the hose and disrupting the pumping process.

Fitting the demineralised water inlet hose to the cleaning machine

- Check that the seal in the screw thread of the water inlet hose is in place and that there is no visible damage, e.g. cracks.
- Screw the water inlet hose firmly onto the screw thread of the pump module.

Connection to the control unit of the cleaning machine

- Insert the control cable plugs into the designated sockets on the potential-free contact of the cleaning machine.

The sockets on the potential-free contact must be retrofitted by Miele Customer Service. The potential-free contact is located on the back of the cleaning machine. The exact position varies depending on the machine model. You can ask Customer Service about the position if necessary. In order for the pump module to be used, Customer Service must activate the pump module in the control unit of the cleaning machine.

Electrical connection

⚠ Compare electrical connection data.

Before installing the pump module, you must compare the connection data (voltage and frequency) on its data plate with the connection data of your on-site mains connection cable. If the data does not match, the module could sustain damage. If in any doubt, consult a qualified electrician.

⚠ All electrical connections must be carried out by a suitably qualified electrician in accordance with local and national safety regulations.

- The electrical installation must be carried out in accordance with IEC 60364-4-41 or the local regulations.
- The connection to the power supply must be via a suitably rated plug and socket and must comply with national regulations. The socket must be accessible after the machine has been installed. An electrical safety test must be carried out after any maintenance or repair work, for example.
- If the machine is hard-wired to the power supply or connected via a socket, a power switch with all-pole isolation must be installed. The power switch must be designed to operate at the rated current for the machine, must ensure a 3 mm gap between all open contacts and must be able to be locked in the off position. The power switch must be accessible after the machine has been installed.
- Equipotential bonding should be carried out if required.
- The rated loads are indicated on the data plate.

- For added safety, every module should be protected by a type A residual current device with a trip current of 30 mA (DIN VDE 0664). The installation of the residual current device must be carried out on site by the operator.
- The mains connection cable may only be replaced by an original spare part from the manufacturer.
- Connect the pump module to a mains electrical connection on site.

Unplugging the demineralised water hoses from the module

Disconnecting the connection hose from the demineralised water container

- Secure the container against uncontrolled leakage of demineralised water, e.g. empty the container, close shut-off valves (if there are any) or clamp the connecting hose.

There may still be residual water in the connecting hose and in the pump module.

- Place a suitable container underneath the hose sleeve of the pump module to catch any demineralised water that may escape.
- Loosen the hose clip and pull off the hose.

Disconnecting the water inlet hose for the cleaning machine

The water inlet hose for the cleaning machine is equipped with an automatic shut-off valve that prevents uncontrolled water leakage from the hose. Nevertheless, there may still be residual water upstream of the valve and in the pump module.

- Place a suitable container underneath the screw thread of the pump module to catch any demineralised water that may escape.
- Unscrew the water inlet hose.

 Do not tilt or lay the pump module on its side during operation. For operation, the pump module must be installed horizontally on a stable surface with the machine feet facing downwards. Operation in an inclined position, lying on its side/upside down or with hose connections pointing upwards is not permitted.

Switching on and off

The module does not have a separate On/Off switch. It is activated by the cleaning machine when required.

Fault messages during the pumping process

If the flow rate is too low during the pumping process or if air is sucked in, fault messages will appear on the cleaning machine display. The cleaning machine repeats the pumping process several times despite the fault messages.

- If the flow rate is sufficient after the repetitions, the current programme is continued by the cleaning machine without interruption.
- If the flow rate is still too low after the repetitions, the cleaning machine will switch to fault mode and abort the current programme with a fault message.

If the programme aborts, you can try to remedy the fault by taking the following measures:

- Check the fill level in the demineralised water container and top up with demineralised water if necessary.
- Open any shut-off valves or hose clips completely.
- Remove kinks from the hoses.
- Check the hoses and hose connections for leaks.
- Vent the system by starting a DI water rinse (standalone programme; available depending on the equipment variant of the cleaning machine) or start a programme with DI water rinse.
- If the fault occurs again, contact Miele Customer Service.

Defect

In the event of a technical defect, you must take the pump module out of operation.

- Make a note of the fault message or error code shown on the cleaning machine display.
- Disconnect the pump module from the mains.
- Contact Miele Customer Service.
- Label the module as decommissioned.
- Label the cleaning machine so that only programmes without the addition of demineralised water can be used until the defect in the pump module has been remedied.

Permanent decommissioning

- Disconnect the pump module from the mains.
- Disconnect the connection to the cleaning machine.
- Detach the water hoses, see  "Detaching demineralised water hoses from the module".
- If the cleaning machine is to continue to be used, contact Miele Customer Service so that they can deactivate the pump module in the cleaning machine's control unit.


Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in land-fill sites.

Disposing of your old appliance

Electronic and electrical appliances contain many valuable materials. They also contain certain materials, compounds and components which were essential for their correct functioning and safety. These could be hazardous to your health and to the environment if disposed of with general waste or if handled incorrectly. Please do not, therefore, dispose of your old appliance with general waste.

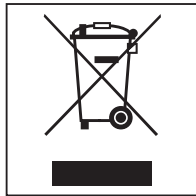Instead, please make use of officially designated collection and disposal points to dispose of and recycle electrical and electronic appliances in your local community, with your dealer, with Miele or with the manufacturer. By law, you are solely responsible for deleting any personal data from the appliance prior to disposal. You are legally obliged to remove any old batteries which are not securely enclosed by the appliance and to remove any lamps without destroying them, where this is possible. These must be taken to a suitable collection point where they can be handed in free of charge. Please ensure that your old appliance poses no risk to children while being stored prior to disposal.

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