E 941/1 mobile unit

Items supplied:

Mobile unit E 941/1, H 421, W 619, D 790 mm
with a magnetic strip for automatic mobile unit recognition

⚠ Please observe the Warning and Safety instructions for washer-disinfector G 7827/28 or PG 8527/28.
If using automatic mobile unit recognition, please make sure the coding is correctly set.


Areas of application:

Modules for laboratory glassware or for orthopaedic surgical instruments can be placed in the E 941/1 mobile unit.
The modules have their own operating instructions.


Up to four modules can be placed in the mobile unit at once, depending on the size of the items being processed.

Before loading the machine and before starting a programme, please check that:


- the sealing caps in the module connections are in good working order,
- the modules are correctly connected to the water intake in the mobile unit,
- the mobile unit is correctly connected to the water intake in the washer-disinfector.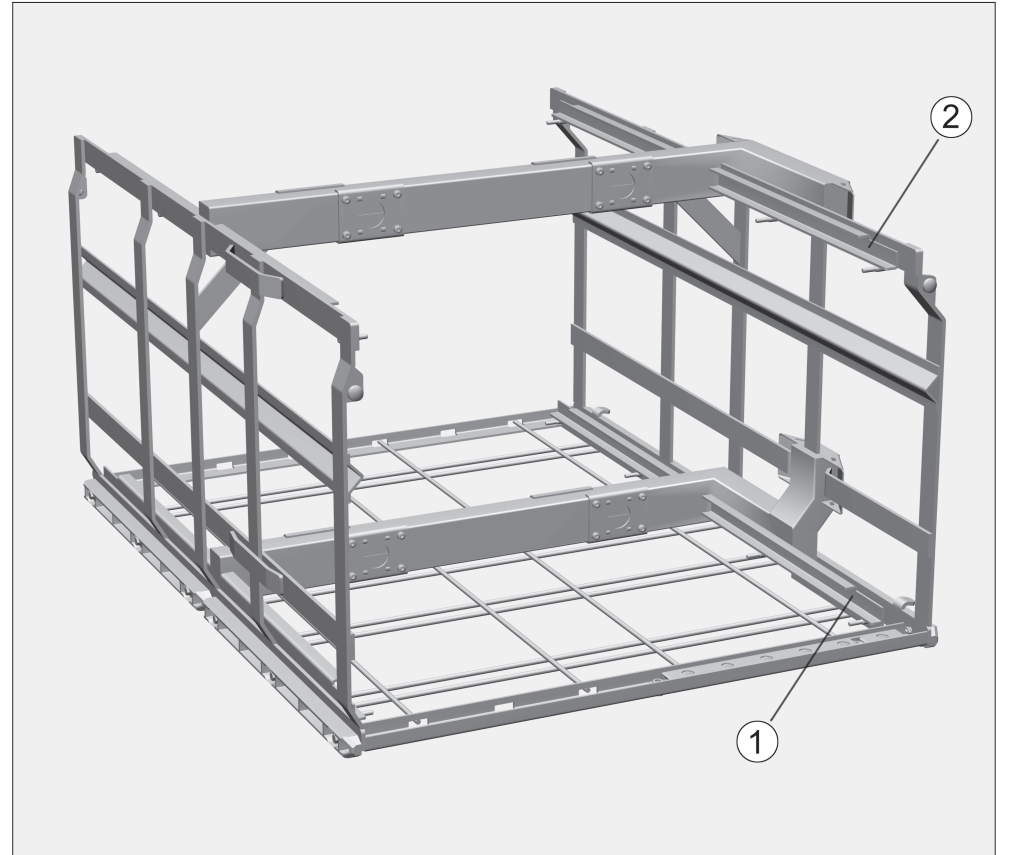


Placing modules into the mobile unit

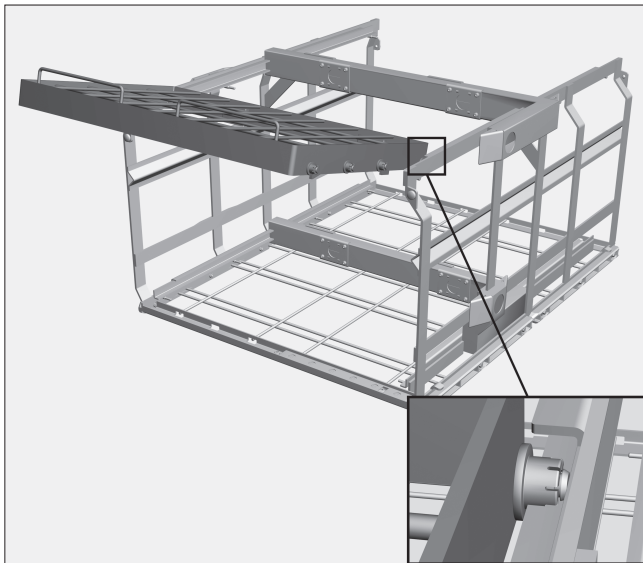- Locate the module rollers on the runners, with the spring-mounted water intakes facing the middle of the mobile unit.
- Push the module completely into the mobile unit.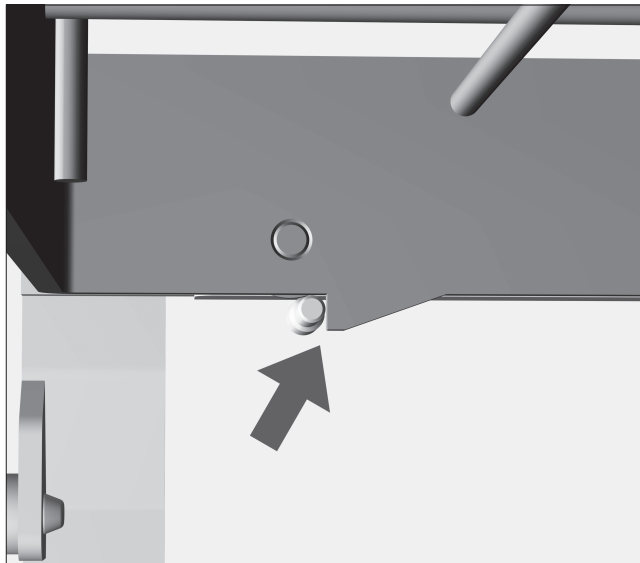

Engage the hooks on the underside of the module frame behind the pins on the runners on either side of the mobile unit.
Make sure it clicks into position correctly.

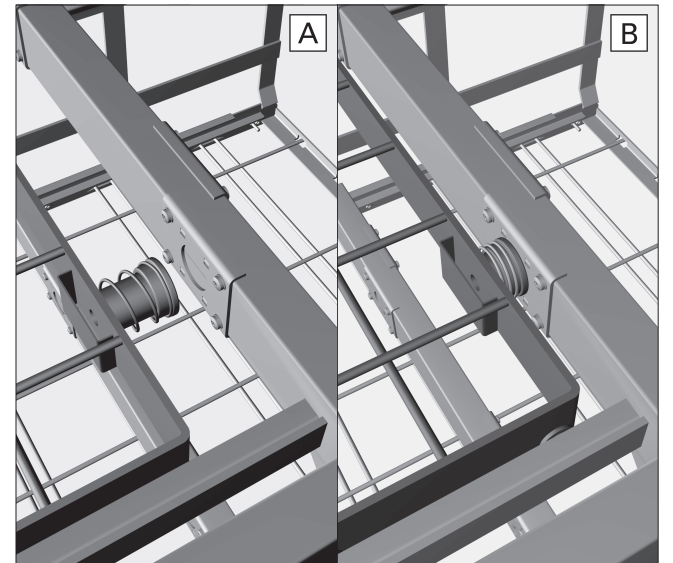When the module is correctly positioned, the water intake on modules with integrated water connection will connect correctly.

Removing modules from the mobile unit

- Lift the module up by the handles until the hooks clear the pins.
- Then pull the module forwards and off the runners.

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