E 741/1 mobile unit

Items supplied:

Mobile unit E 741/1, H 680, W 530, D 600 mm
with a magnetic strip for automatic mobile unit recognition

 Please observe the Warning and Safety instructions for washer-disinfector G 7825/26.
If using automatic mobile unit recognition, please make sure the coding is correctly set.


Areas of application:

Modules for laboratory glassware can be placed in the E 741/1 mobile unit.
The modules have their own operating instructions.



Depending on the size of the items to be processed, up to three modules can be placed in levels ① to ④ at the same time.

Before loading the machine and before starting a programme, please check that:

- the sealing caps in the module connections are in good working order,
- the modules are correctly connected to the water intake in the mobile unit,
- the mobile unit is correctly connected to the water intake in the washer-disinfector.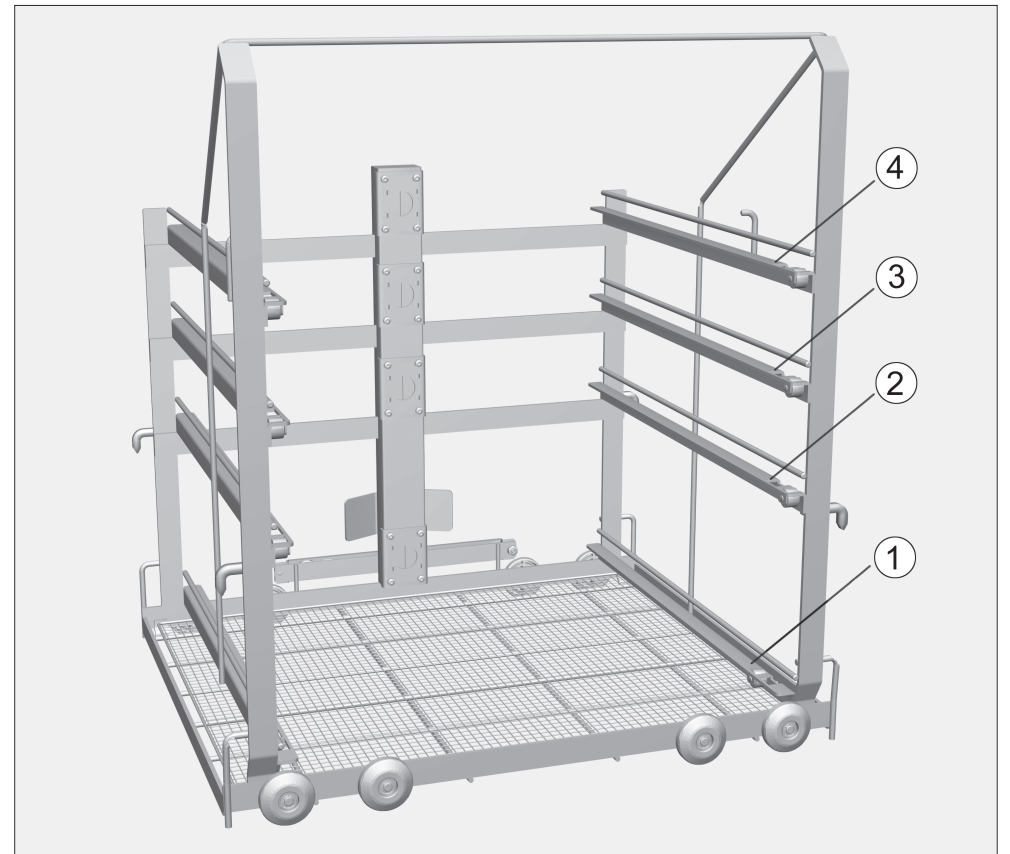

Placing modules into the mobile unit

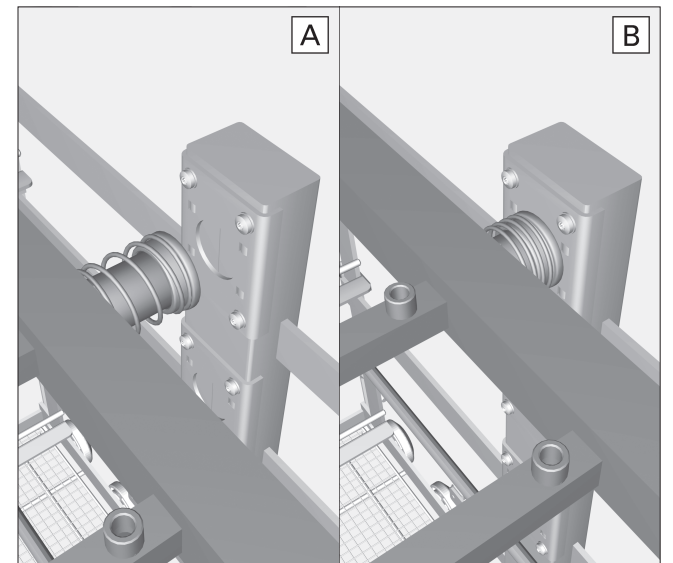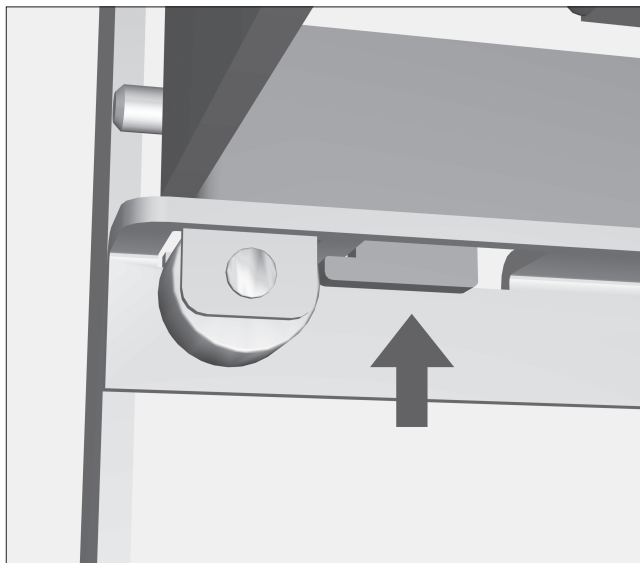- Locate the module rollers on the runners, with the spring-mounted water intake facing towards the mobile unit.
- Push the module completely into the mobile unit.

Engage the catches on the underside of the module frame into the recess in the runners.
Make sure it clicks into position correctly.

If the module is accurately positioned, the water intake on modules with integrated water connection will connect correctly.

Removing modules from the mobile unit

- Push the module gently in the direction of the water intake, and then lift it off the guide rail.

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