DOS-Modul G 10

de	Montageanweisung
en	Fitting instructions
fr	Notice de montage
it	Istruzione di montaggio
es	Instrucciones de montaje
pt	Instruções de montagem
nl	Montage-instructies
da	Monteringsanvisning
sv	Monteringsanvisning
no	Monteringsveiledning
fi	Asennusohje

de	3
en	6
fr	9
it	12
es	15
pt	18
nl	21
da	24
sv	27
no	30
fi	33

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Caring for the environment

Disposal of packing


The transport and protective packing has been selected from materials which are environmentally friendly for disposal and can normally be recycled.

Ensure that any plastic wrappings, bags etc. are disposed safely and kept out of the reach of babies and young children. Danger of suffocation!

Rather than just throwing these materials away, please ensure they are offered for recycling.

Disposal of your old machine

Electrical and electronic appliances / machines often contain materials which, if handled or disposed of incorrectly, could be potentially hazardous to human health and to the environment. They are, however, essential for the correct functioning of your appliance or machine. Please do not therefore dispose of your old machine or appliance with your household waste.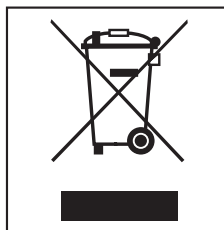

Please dispose of it at your local community waste collection / recycling centre and ensure that it presents no danger to children while being stored for disposal.

It should be unplugged or disconnected from the mains electricity supply by a competent person.

Description of the machine


The DOS-Module G10 is an external dispensing unit. It automatically dispenses liquid cleaning agents at the correct programme stage.

DOS-Modules can also be fitted and connected up after the machine has been installed.

DOS modules must only be set up and connected for the first time by a Miele authorised and trained service technician. Personnel using the DOS module must be trained in its use by a Miele authorised and trained service technician.


Read the warning and safety instructions for the machine to which the DOS-Module is connected.

G 7804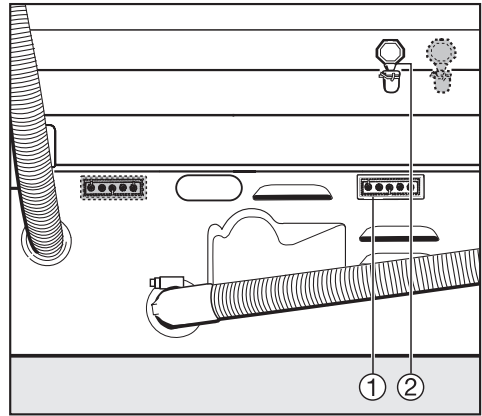

- ① = electrical supply connection neutralising agent
- ② = dispenser hose connection neutralising agent


The connections for the mains cable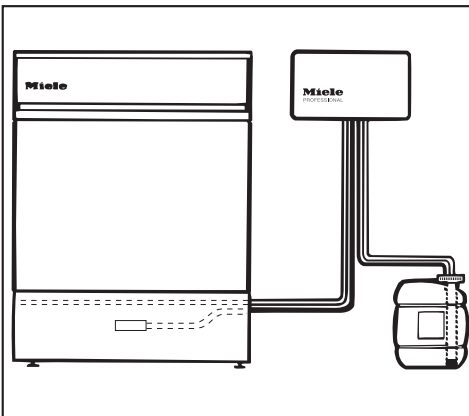

and dispenser hose are on the back of the machine.

Description of the machine


90 cm wide machines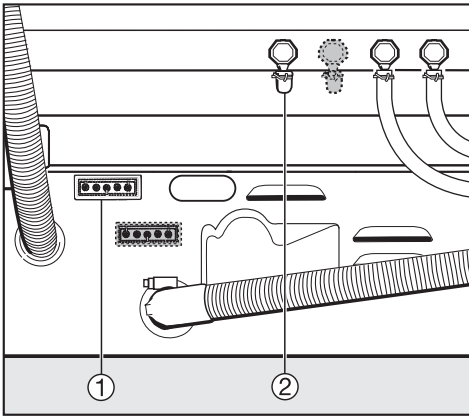

- ① = electrical supply connection
rinsing agent / neutralising agent
- ② = dispenser hose connection
rinsing agent / neutralising agent

For wall mounting the DOS module two screws (4.8 x 32) and two plugs (8 mm) are included with the unit.

Dimensions:

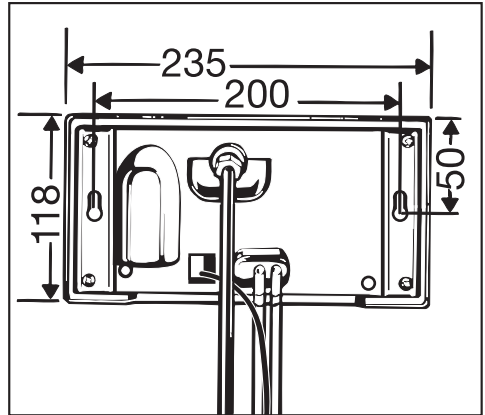- Mount the module in a suitable location on the wall.
- Connect the dispenser pump and electricity supply.
- Adjust the position of the cover caps along the siphon tube so that the tube will reach the bottom of the container.
- Fill the storage container with the relevant agent, feed the siphon tube into the container and tighten the cover cap securely.

DOS-Module venting and dispensing instructions are given in the Operating Instruction booklet for the machine.

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