
E 747 injector module for laboratory glassware

Miele
PROFESSIONAL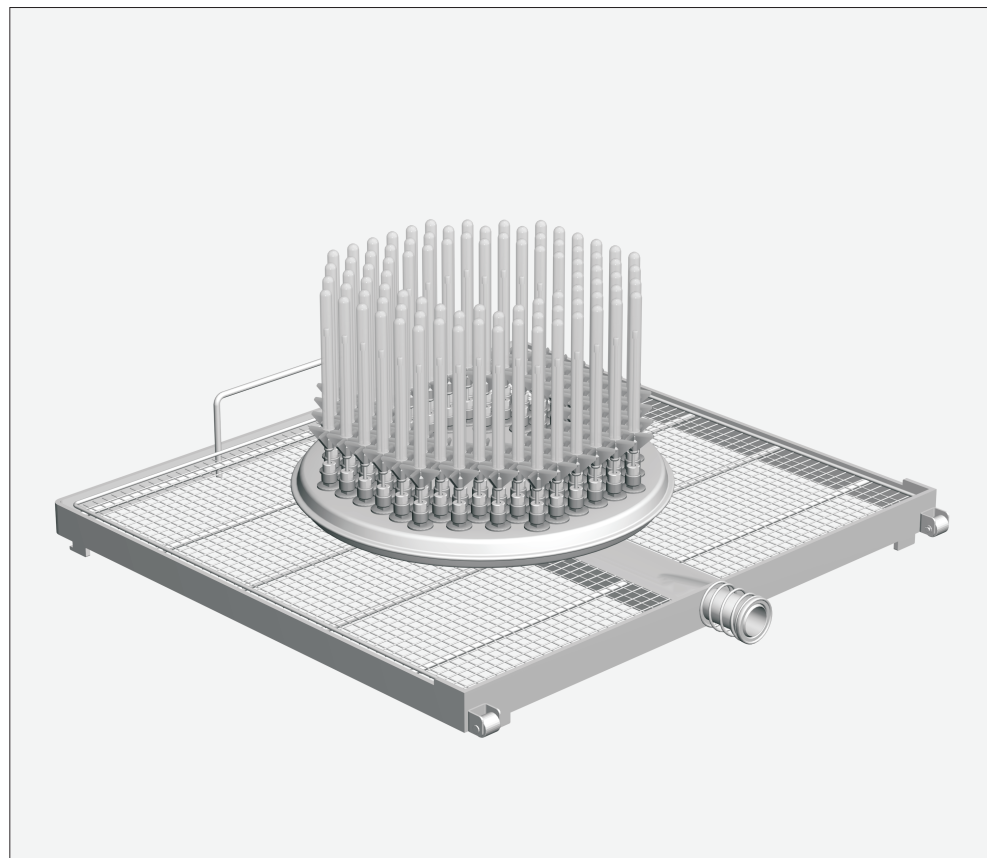

Items supplied:

1 x E 747 injector module, H 168, W 492, D 496 mm

104 x ID 110 jets, 2.5 x 110 mm

Areas of application:

The injector module can be placed on any level in the E 741/1 mobile unit depending on which other modules are being loaded at the same time.


Up to 104 centrifugal tubes, phials and test tubes can be processed in this injector module.

Laboratory glassware, which because of its design cannot be flushed through properly, will not be sufficiently processed in this injector module.

This injector module should only be used for processing laboratory glassware which has been declared as suitable for machine processing by the glassware manufacturer.

⚠ Please observe the Warning and Safety instructions in the Operating Instructions for your washer-disinfector as well as any instructions provided by the glassware manufacturer.
Any broken glassware can cause serious injury.
All glassware should be properly inspected and tested for cleanliness. Do not rely on a visual check alone.


Arranging the laboratory glassware: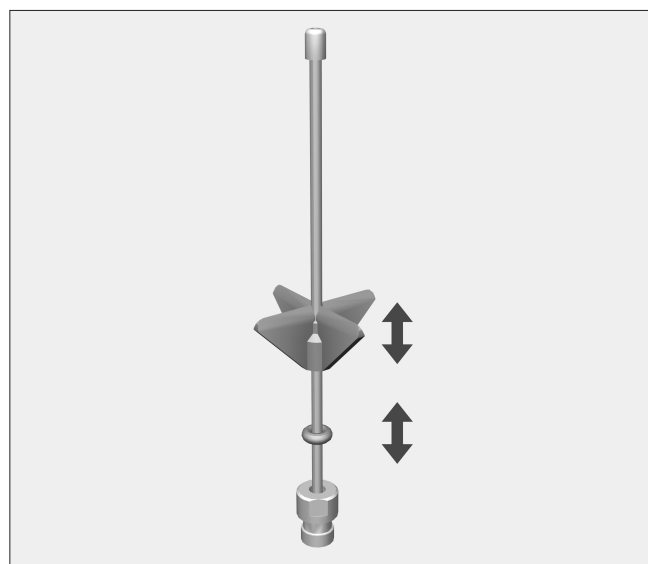

To avoid having to keep adjusting the height of the plastic supports, the injector module should be used for glassware of the same type and size.

The plastic support can be moved in order to adjust the height of the jet to the height of the item being processed. Pliers can be used to do this if necessary.

- Slide the plastic support on the jet to the required height.
- Slide the O ring up under the plastic support to secure its position.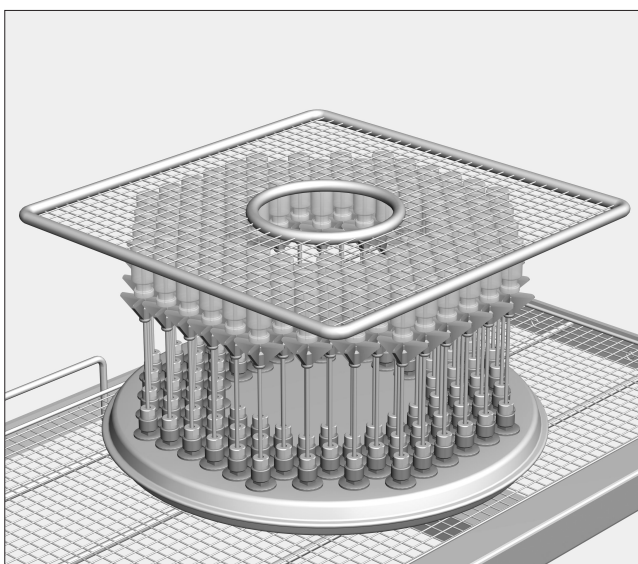


Light items can be secured with an A5 cover.


To avoid the A5 cover slipping during processing, only glassware of a similar type and size should be processed together.

Remove coarse soiling from centrifugal tubes before processing.


⚠ The water intake level must be increased for the relevant reprocessing programmes.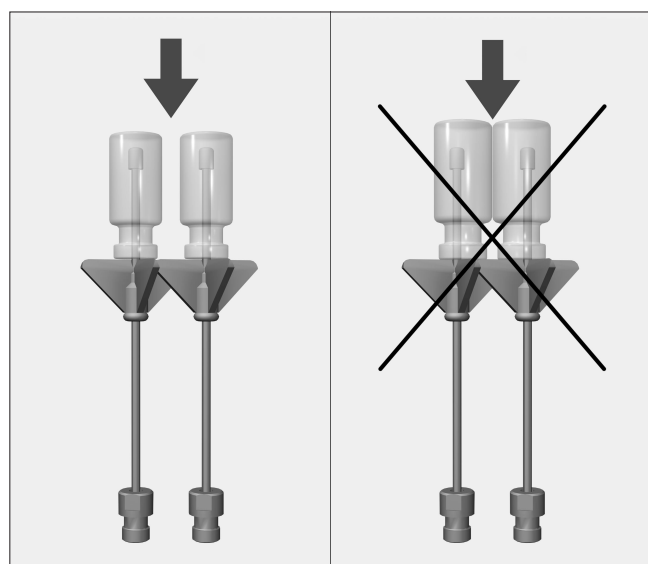

Do not allow individual pieces of glassware to touch each other.

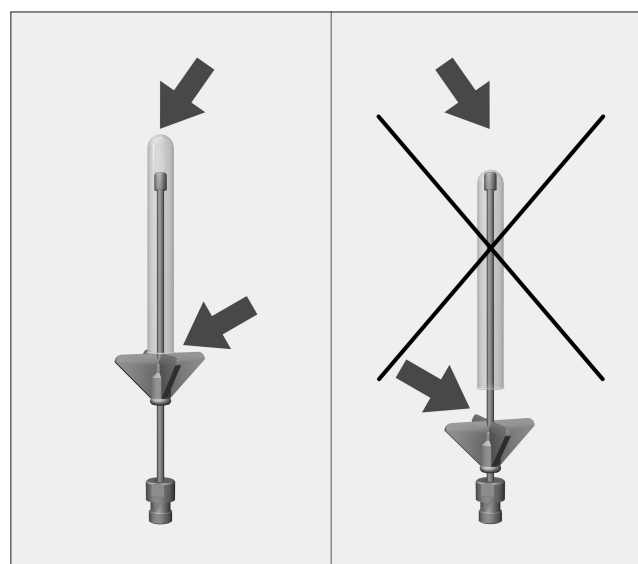If the tip of the jet touches the base of the glassware, slide the plastic support up and secure with the O ring.

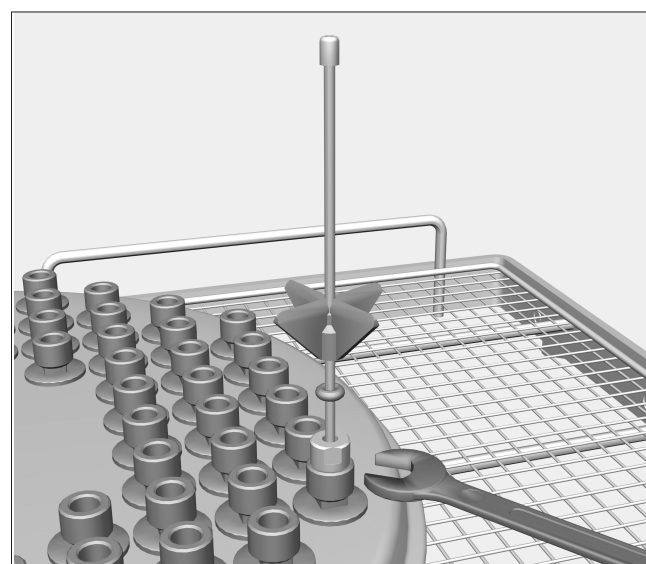Tighten loose jets with an SW 9 spanner.

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