
E 755 injector module for narrow-necked glassware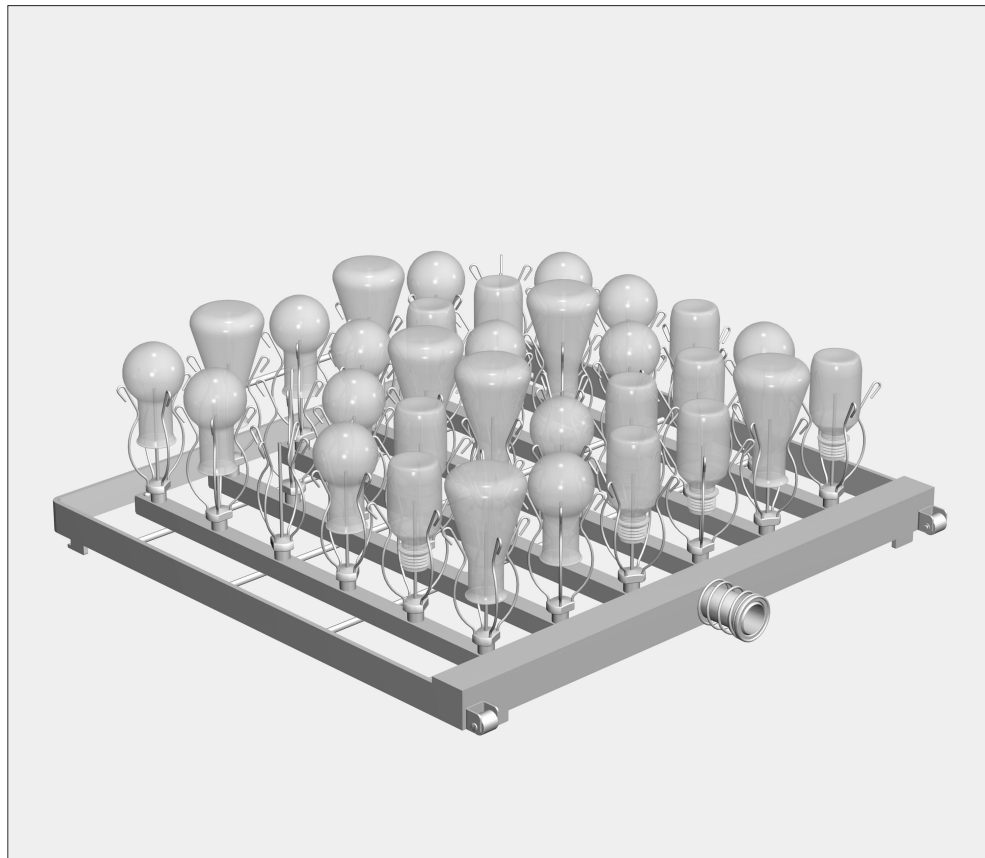

Items supplied:

1 x E 755 injector module, H 130, W 492, D 496 mm

36 jets with E 470 holders, 2.5 x 90 mm

Areas of application:

The injector module can be placed on any level in the E 741/1 mobile unit depending on which other modules are being loaded at the same time.



36 pieces of narrow-necked glassware with a volume of 25-100 ml can be processed in this injector module.

Laboratory glassware, which because of its design cannot be flushed through properly, will not be sufficiently processed in this injector module.

This injector module should only be used for processing laboratory glassware which has been declared as suitable for machine processing by the glassware manufacturer.

⚠ Please observe the Warning and Safety instructions in the Operating Instructions for your washer-disinfector as well as any instructions provided by the glassware manufacturer.
Any broken glassware can cause serious injury.
All glassware should be properly inspected and tested for cleanliness. Do not rely on a visual check alone.

Installation: Tools required: SW 15 spanner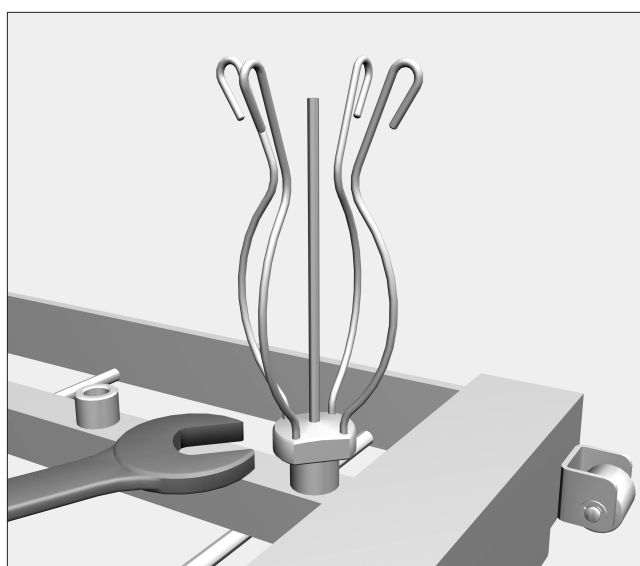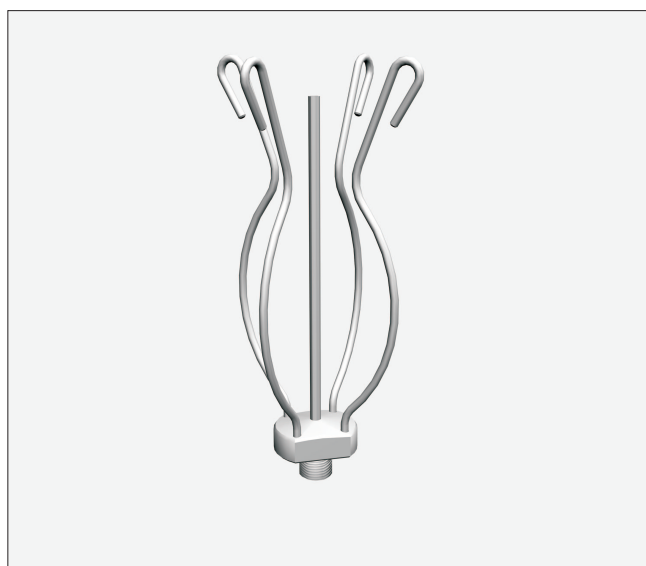


The injector module is supplied with jets and holders.


- Screw the jets into position, and tighten with an SW 15 spanner.

To ensure adequate water pressure, all positions need to be equipped with jets.


Arranging the laboratory glassware: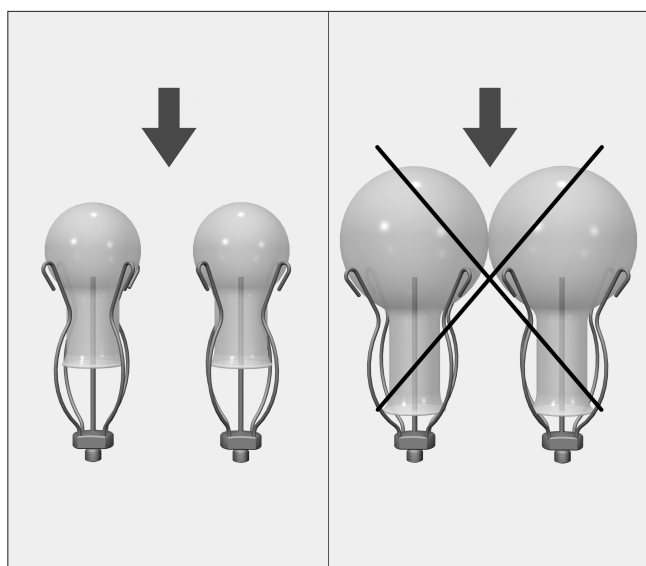

Do not allow individual pieces of glassware to touch each other.

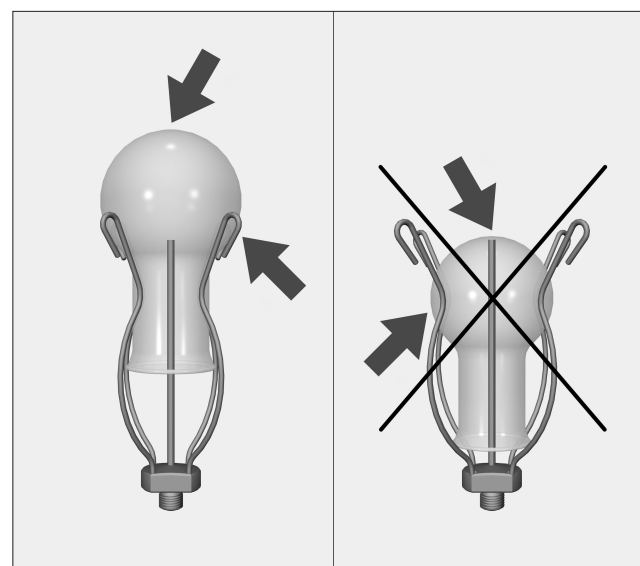The tip of the jet must not touch the bottom of the glassware.

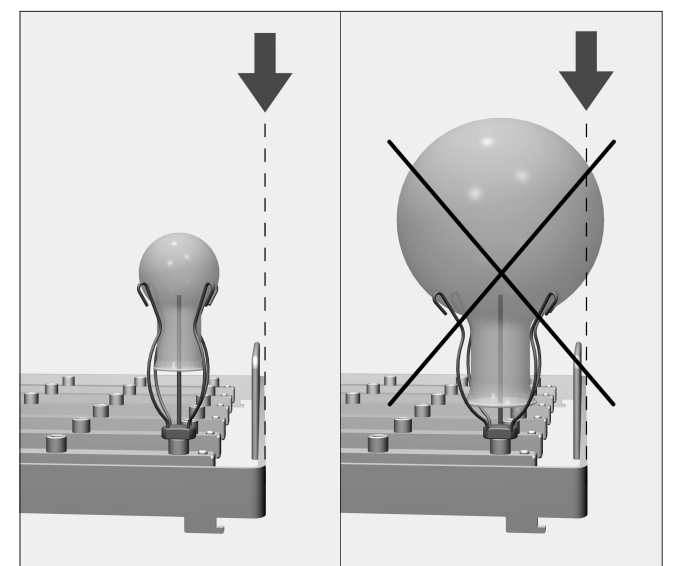Glassware must not extend beyond the sides of the injector module.

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