E 944/2 injector module for narrow-necked laboratory glassware **Miele** PROFESSIONAL

Items supplied:

- 1 x E 944/2 injector module, H 250, W 558, D 352 mm
- 15 x E 352 jets, 6 x 220 mm
- 15 x E 354 clips for 6 x 220 mm jets, consisting of holder + clamp

Areas of application:

The injector module can be placed on either level of the E 941 mobile unit.


The module can be used to process 15 pieces of narrow-necked glassware with a volume of 500-1000 ml.

Laboratory glassware, which because of its design cannot be flushed through properly, will not be sufficiently processed in this injector module.

This injector module should only be used for processing laboratory glassware which has been declared as suitable for machine processing by the glassware manufacturer.


⚠ Please observe the Warning and Safety instructions in the Operating Instructions for your washer-disinfector as well as any instructions provided by the glassware manufacturer.
Any broken glassware can cause serious injury.
All glassware should be properly inspected and tested for cleanliness. Do not rely on a visual check alone.

Installation: Tools required: SW 9 spanner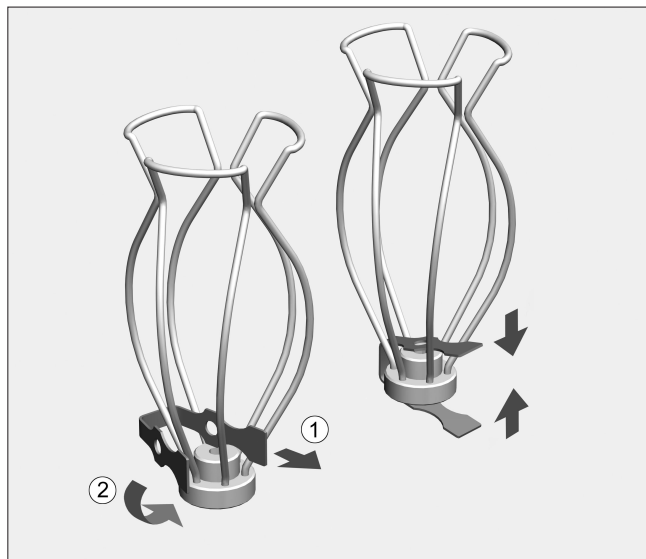

Clamps, holders and jets are supplied with the injector module.


- Push the clamp over the holder as shown above.
- Then squeeze the clamp together and feed it, together with the holder, over the jet.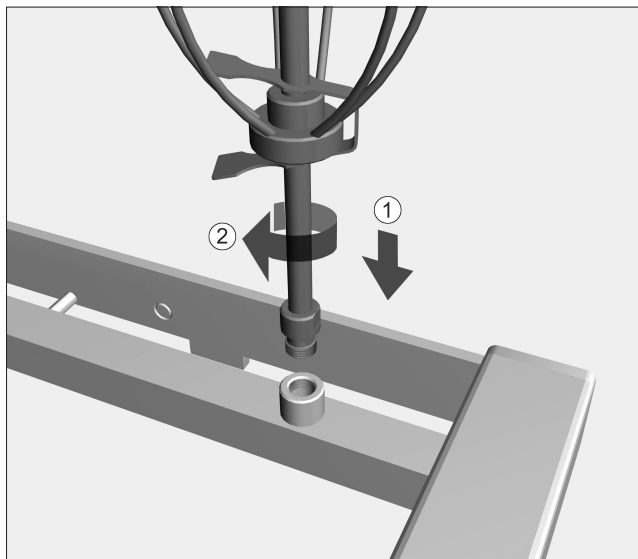


- Screw the jets into position, and tighten with an SW 9 spanner.

To ensure adequate water pressure all positions need to be equipped with jets.


Arranging the laboratory glassware: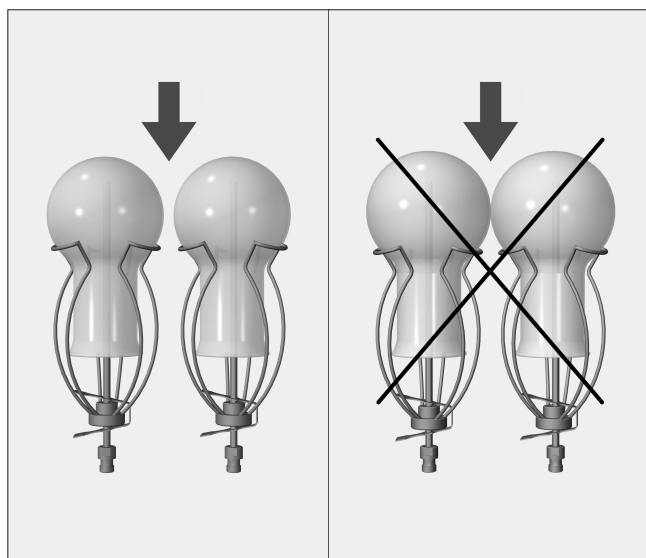

Do not allow individual pieces of glassware to touch each other.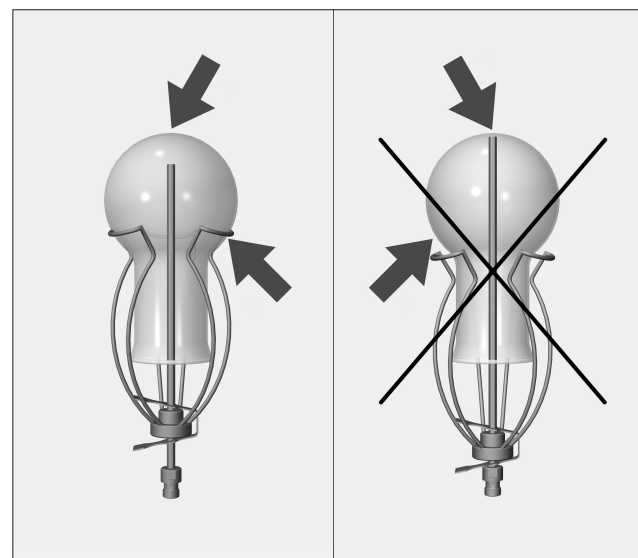

If the tip of the jet touches the base of the glassware, push the holder and clamp upwards.

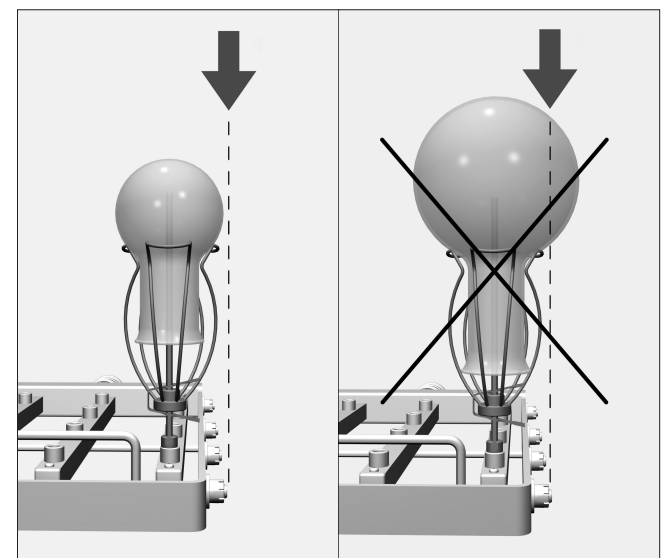Glassware must not extend beyond the sides of the injector module.

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