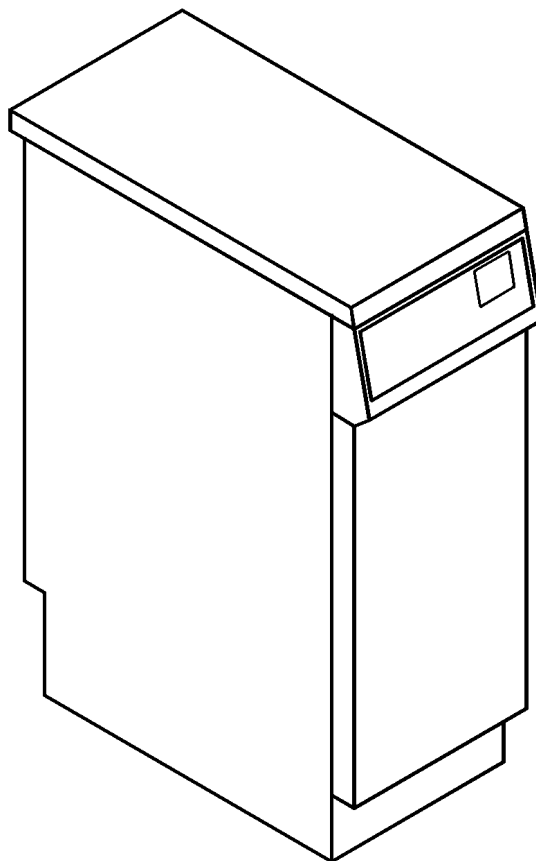
Installationsplan / Installation plan

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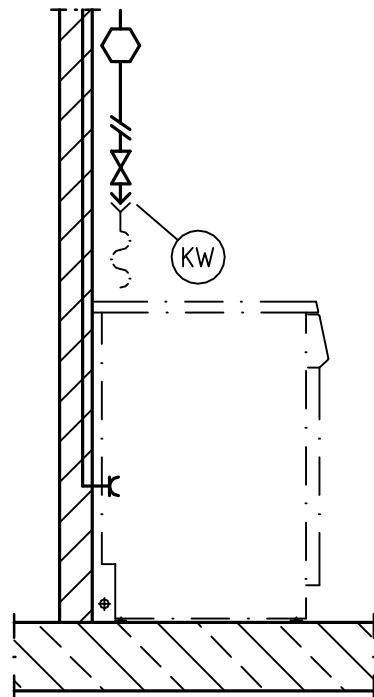
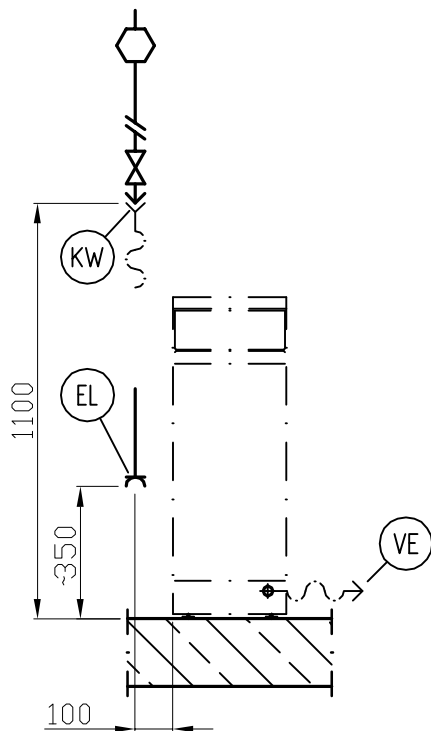
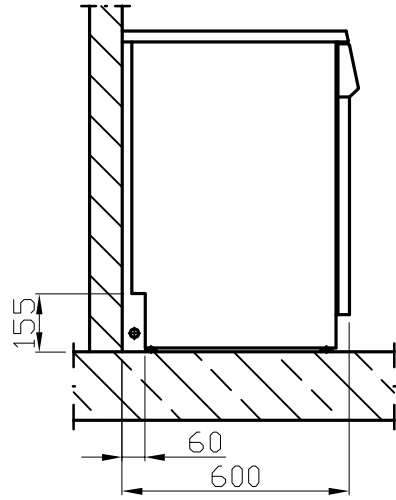
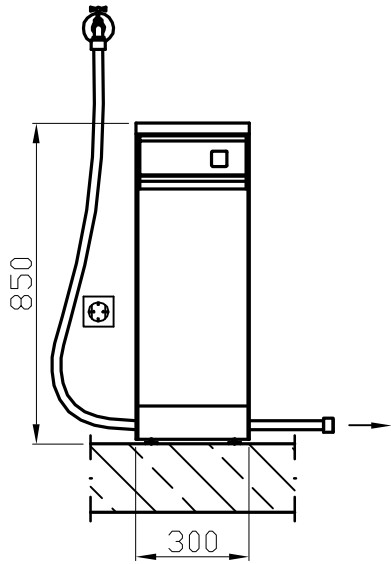
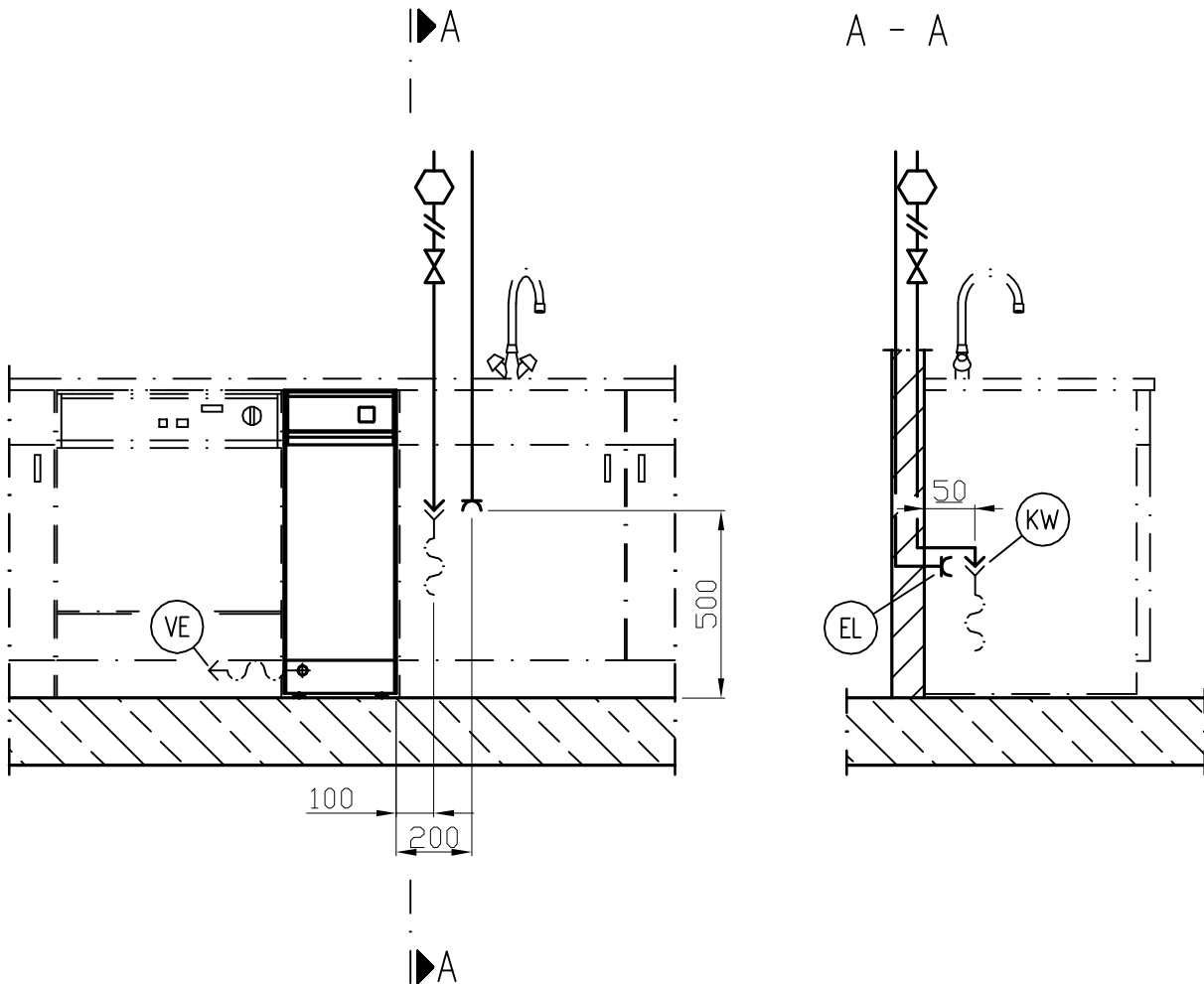
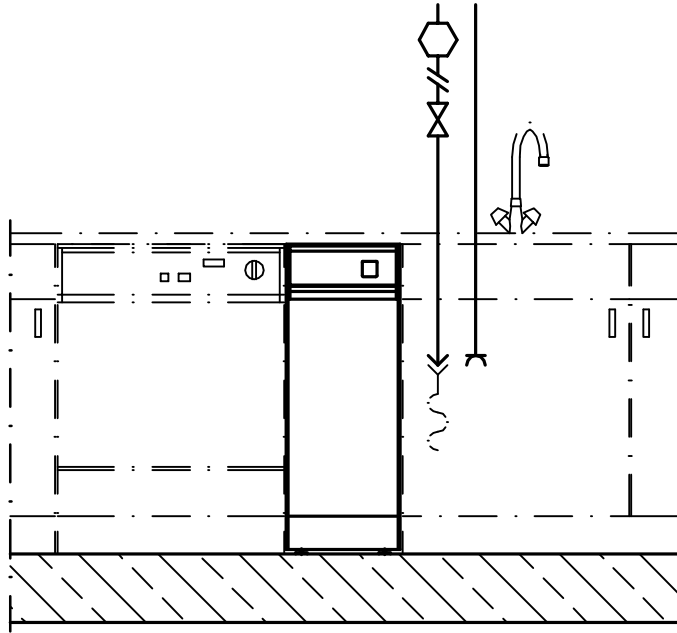
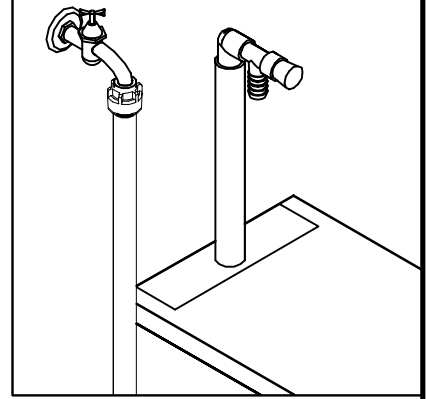
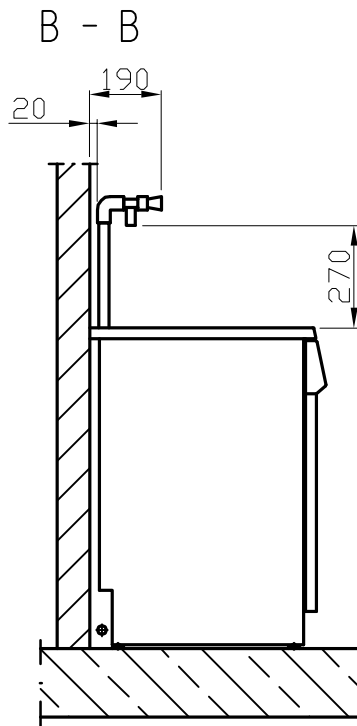
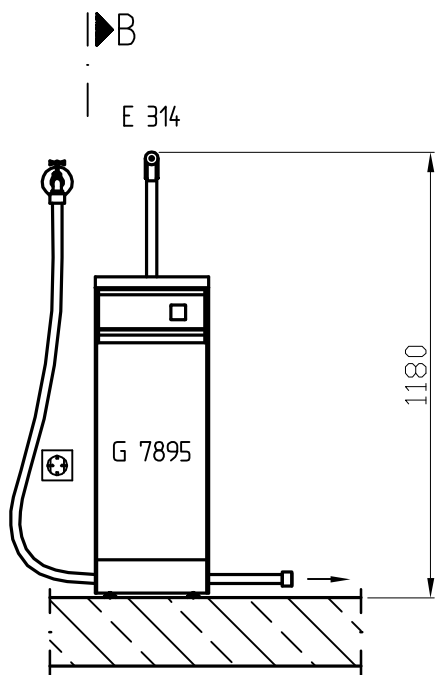
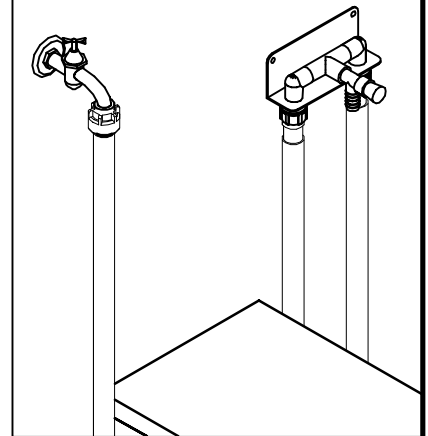
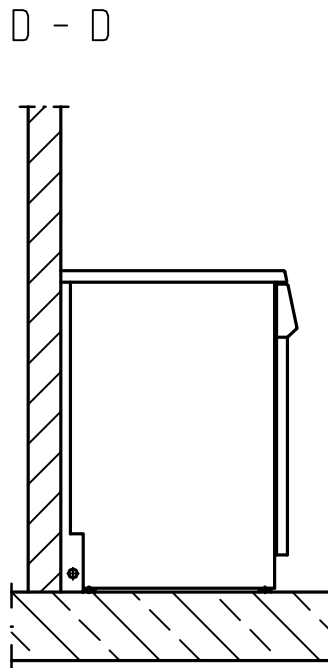
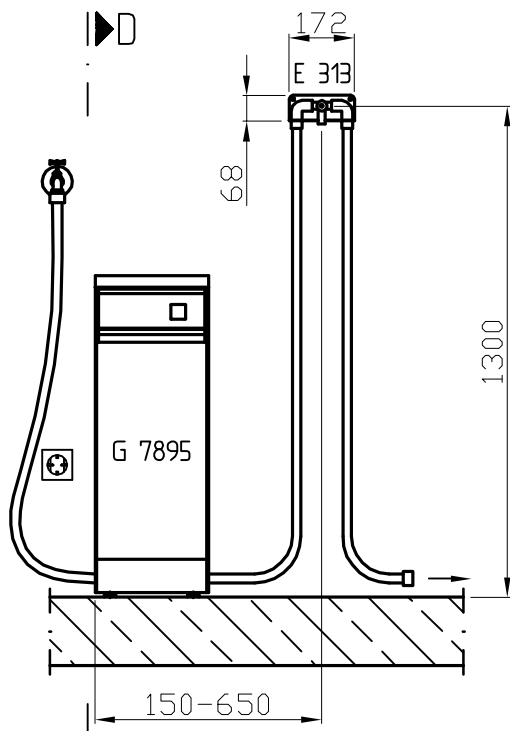
G 7895

Legend:

Highlighted circled abbreviations:
Connection required

(EL)	Electrical connection (German standard models only)	Voltage	V/Hz	AC 208-240/50
		Connected load	W	1.2
		Fuse rating	A	1 × 16
		Cross-sectional area of supply lead	mm ²	3 × 1.5
		Connection cable length (cable supplied with earthed plug)	m	1.7
		<p>It is recommended to make electrical connection via a plug and socket so that electrical safety checks can be carried out easily. The socket must be accessible after machine installation. With a fixed installation, this must be via a main switch to be provided on site which must totally isolate the machine from the mains with a contact gap of at least 3 mm. The fitting of an earth leakage circuit breaker is strongly recommended. If necessary, an equipotential bond with good contact connection must be provided in accordance with all appropriate national and local regulations. The connection cable must be protected against the effects of heat.</p>		
(KW)	Cold water connection (input)	Max. temperature	°C	20
		Max. pressure at demineralising cartridge (cartridge not supplied as standard)	kPa	1000
		Pressure loss at each demineralising cartridge with flow pressure 8 l/Min	kPa	100
		On-site threaded union in accordance with DIN 44 991 (flat seating)	inch	¾ male thread
		Connection hose length, marked blue, (supplied)	mm	1200
		Water connection optional left or right next to machine. Water filter provided must be fitted!		
		The Aqua Purificator must be connected on site via a water safety device as suitable for the on-site conditions.		
(VE)	Demineralised water (output)	Threaded union for connection hose to following machine in accordance with DIN 44991 (flat seating)	inch	¾ male thread
		Connection hose length, marked green, (supplied)	mm	1200
	General data	Height with lid	mm	850
		Building-under height	mm	820
		Width	mm	300
		Depth	mm	600
		Net weight	kg	26.5
		Floor load with one cartridge (with two cartridges)	N	530 (795)
		Min. installation location access width including packaging	mm	400
		Min. installation location access height including packaging	mm	900
	Demineralising cartridge holding capacity, max.	qty.	2	

Installation should only be carried out by suitably qualified persons in accordance with all appropriate local and national safety regulations, accident prevention regulations and applicable standards!
During installation it is essential to follow the installation instructions provided! Modification rights reserved!
Dimensions in mm

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