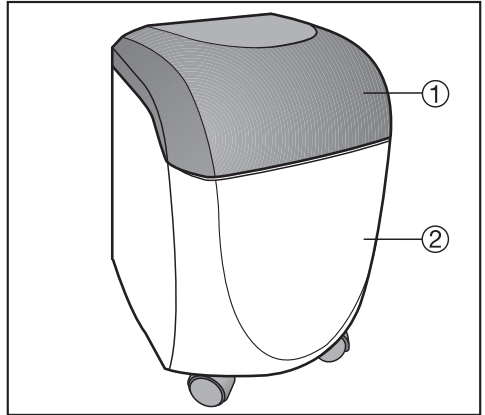
Aqua-Soft-System PG 8597

de	Gebrauchsanweisung
en	Operating instructions
fr	Mode d'emploi
it	Istruzione d'uso
es	Instrucciones de manejo
pt	Instruções de utilização
nl	Gebruiksaanwijzing
da	Brugsanvisning
no	Bruksanvisning

de	3
en	10
fr	17
it	24
es	31
pt	38
nl	45
da	53
no	60

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

en - Description of the appliance

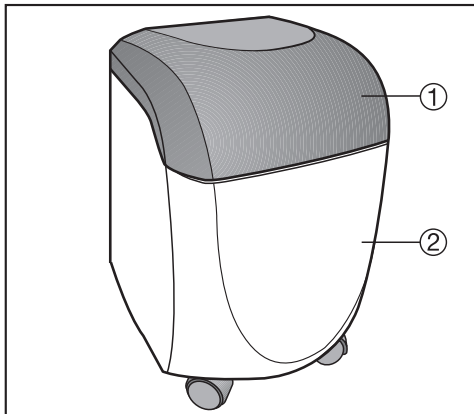- 1 Lid
- 2 Cabinet

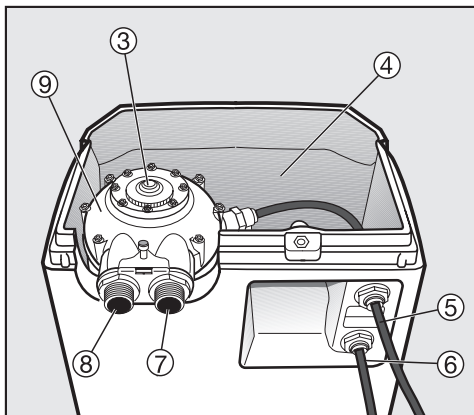- 3 Water hardness range Meter Disk (under lid)
- 4 Salt inlet
- 5 Reactivated water outlet (free outlet)
- 6 Overflow (free outlet)
- 7 Water inlet
- 8 Water outlet to machine
- 9 Valve head with resin cartridge

en - Description of the appliance

How the Aqua-Soft-System works

The Aqua-Soft-System PG 8597 is a "two-bed" system which supplies a constant supply of softened water.

It is a dual-tank system. One tank supplies softened water whilst the other reactivates.

Reactivation is controlled by quantity and is automatically set in motion by the requirement for softened water.

The use of salt is kept to a minimum because reactivation is on demand.

This appliance does not require electricity, solenoid valve or timer to operate.

Moving components only come into contact with softened water so limescale does not develop.

The Aqua-Soft-System PG 8597 must be fitted with the correct water hardness range disk (Meter Disk) for the water supply. A total of 8 different disks are supplied.

The Miele service technician will install the correct hardness range disk (Meter Disk) to set the degree of hardness.

Minimum flow pressure

The Aqua-Soft-System operates with a minimum flow pressure of approx. 250 kPa (2.5 bar) when used in conjunction with either a fresh water machine or a tank machine which does not have a built-in water softener.

Water pressure must not drop below 250 kPa (2.5 bar) even when at full flow.

en - Caring for the environment

Disposing of the packing materials

The packaging is designed to protect the appliance during transportation. It has been selected from materials that are environmentally friendly for disposal and that can be recycled.

Recycling saves on the use of raw materials and reduces the amount of waste going to landfill.

Ensure that any plastic wrappings, bags etc. are disposed of safely and kept out of the reach of babies and young children. Danger of suffocation. Depending on country, your dealer may also be able to take packaging back from you.

Disposal of your old appliance

Old appliances contain materials which can be reclaimed or recycled. Please contact your dealer, local authority or scrap merchant about recycling schemes. Ensure that the appliance presents no danger to children while being stored for disposal.

This Aqua-Soft-System PG 8597 complies with all relevant safety requirements. Incorrect use can, however, lead to personal injury and damage to property.

To avoid the risk of accidents and damage to the appliance please read these instructions carefully before starting to use it. They contain important notes on the installation, safety, operation and care of the appliance.

Keep these instructions in a safe place for reference, and pass them on to any future user.

Correct application

► The Aqua-Soft-System PG 8597 must only be used as described in these instructions. Using it for purposes other than those for which it is designed is not permitted and could be dangerous.

► The manufacturer cannot be held liable for damage caused by non-compliance with these Warning and Safety instructions or resulting from improper or incorrect use or operation.

► Untrained personnel should not be allowed access to the Aqua-Soft-System PG 8597.

► This appliance is not intended for use by persons (including children) with reduced physical, sensory or mental capabilities, or lack of experience and knowledge, unless they are supervised whilst using it or have been shown how to use it correctly by a person responsible for their safety.

Technical safety

► The Aqua-Soft-System PG 8597 must be checked for any visible damage before it is installed. Do not install and use a damaged appliance. A damaged appliance is dangerous.

► The Aqua-Soft-System PG 8597 should be commissioned and then maintained only by a Miele authorised and trained service technician or other suitably qualified and competent approved person.

► Installation, maintenance and repairs should only be carried out by a Miele approved service technician in accordance with local and national safety regulations.

Installation, repairs and other work by unqualified persons could be dangerous and the manufacturer will not be held liable.

► Faulty components must only be replaced by genuine Miele original parts. The manufacturer can only guarantee the safety of the appliance when Miele replacement parts are used.

en - Warning and Safety instructions

▶ To ensure safety, national and local Water Authority Regulations (NWC in the UK/DIN EN 1717 in Germany) must be complied with when installing this equipment. When connection to the water supply is made, an approved backflow prevention device (Type C) must be incorporated between the mains water supply and the water softener cartridge.

The device must be installed at least 250 mm above the upper edge of the Aqua-Soft-System PG 8597. Waste water must be discharged in accordance with national and local water regulations.

▶ A damaged Aqua-Soft-System is dangerous. Stop using it immediately and call the Miele Service Department or an authorised Miele Service Dealer.

Correct use

▶ The water in the appliance must not be used as drinking water.

▶ Never clean the PG 8597 or near vicinity with a water hose or a pressure washer when used in conjunction with an electrically operated appliance.

▶ Ensure that solutions or steam containing hydrochloric acid do not come into contact with the steel outer casing, to avoid any corrosion damage.

▶ While the appliance is under guarantee, repairs should only be undertaken by a service technician authorised by the manufacturer. Otherwise the guarantee will be invalidated.

▶ This appliance must only be used in conjunction with machines specified by Miele as being suitable for such use.

Installation

▶ Do not install the appliance in conjunction with an electrical appliance in areas where there is a risk of explosion.

▶ This appliance may only be used in mobile installations such as ships, if a risk assessment of the installation has been carried out by a suitably qualified engineer.

Safety with children

▶ To avoid the risk of injury, keep children away from the appliance at all times. Do not allow them to play with it or use the controls. They will not understand the potential dangers posed by it.

▶ Older children may only use this appliance when its operation has been clearly explained to them and they are able to use it safely, recognising the dangers of misuse.

Before use

Before the appliance is used for the first time, the Miele approved service technician will need to install the correct hardness range disk (Meter Disk) for the water hardness level in your area.

- Find out the mains water hardness level in your area.
- Please make a note of the water hardness level here.

Hardness level:

The service technician will connect the Aqua-Soft-System to the water supply, fill the resin cartridges and start the system. The built-in salt valve limits the amount of water going into and out of the unit by means of a float switch to ensure there is sufficient salt available for reactivation.

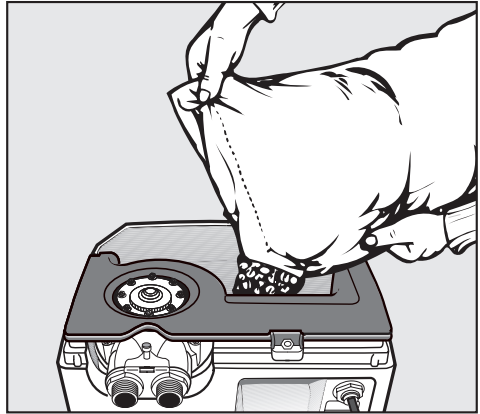
It is important to use good quality reactivation salt. Only use salt tablets or coarse grained reactivation salt (8 - 15 mm grain).

Only use reactivation salt that is suitable for water softeners.

Never use cooking salt, as it contains particles which do not dissolve in water. These particles can cause a blockage in the Aqua-Soft-System.

Filling the salt container

- Remove the lid.

- Fill the salt container with reactivation salt or salt tablets.

Capacity: 20 kg
First fill: 20 kg

- Replace the lid.

To keep evaporation to a minimum make sure that the lid is always replaced securely in position.

The Aqua-Soft-System is now ready for operation and will deliver softened water.

Please contact the Miele Service Department if in any doubt.


en - Technical data

External dimensions:	H 570 / W 360 / D 360 mm
Hard / Soft water connection:	R 3/4"
Resin cartridges:	2 cartridges
Resin content:	2 x 4.5 l +/- 10 % Sybron Chemicals - Dow Chemical Co.
Capacity:	19 l/min
Minimum flow pressure:	
without/with softener in Fresh water machines:	250 kPa (2.5 bar) / 350 kPa (3.5 bar)
without/with softener in Tank system machines:	250 kPa (2.5 bar) / 350 kPa (3.5 bar)
Maximum static pressure:	800 kPa (8.0 bar)
Pressure difference:	maximum 100 kPa (1 bar) at 10 l/min
Temperature:	4 °C - 70 °C
Salt consumption:	0.45 kg NaCl/Reactivate +/- 10 %
Water consumption:	19 l/Reactivate +/- 10 %
Reactivation time:	11 minutes +/- 10 %
Capacity/Diff.:	+/- 10 %
Salt container:	20 kg
Tablet shape:	round, approx. Ø = 25 mm, h = 14 mm

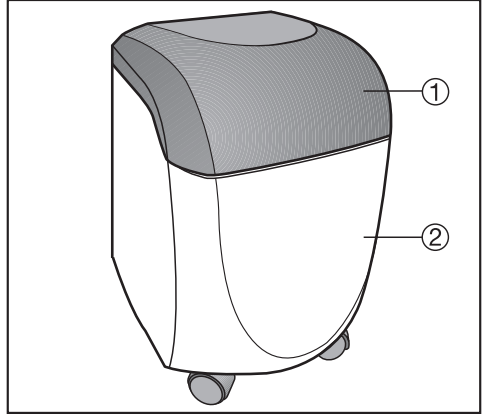Meter Disk	1	2	3	4	5	6	7	8
°d	1-5	6-10	11-15	16-20	21-25	26-30	31-35	36-40
Capacity (l/Reactivate)	1442	721	480	360	288	240	206	180

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