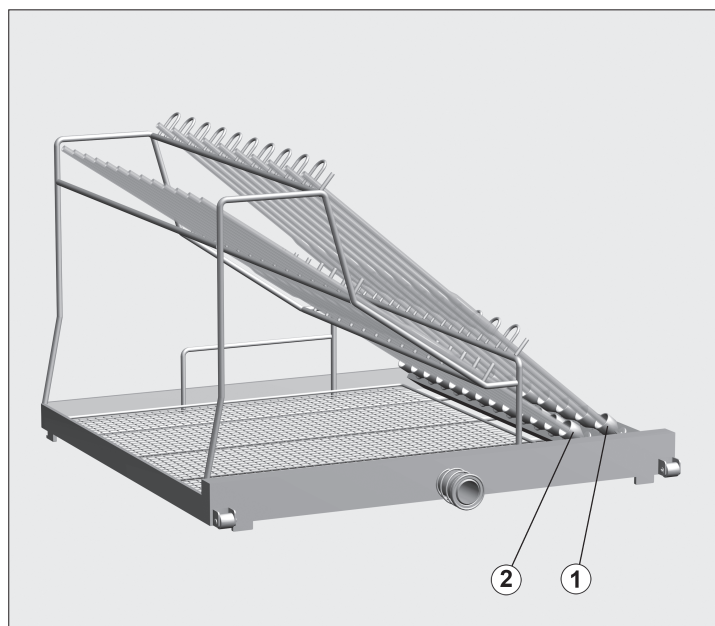
E 746 injector module for pipettes

Items supplied:

1 x E 746 injector module, H 288, W 492, D 496 mm

Intended use:

This injector module is for use in the 1st or 2nd level of the E 741/1 mobile unit.

Capacity:

Row	max. length [mm] *	Quantity	Volume [ml] *	
			Measuring pipettes	Bulb pipettes
1	560	10	10 to 50	10 to 100
2	490	13		

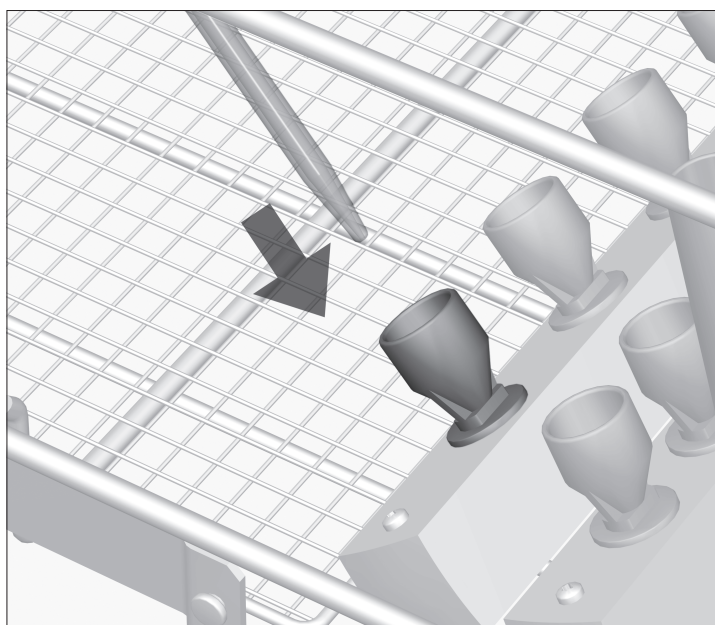
* The pipettes must not extend beyond the sides of the injector module.

This mobile injector unit should only be used for processing laboratory glassware which has been declared as suitable for machine processing by the glassware manufacturer.

⚠ Please observe the Warning and Safety instructions in the Operating Instructions for your washer-disinfector as well as any instructions provided by the glassware manufacturer.

Broken glassware can cause serious injury. All glassware should be properly inspected and tested for cleanliness. Do not rely on a visual check alone.

⚠ The water intake level must be increased for the relevant reprocessing programmes.

Placing pipettes into the injector unit:

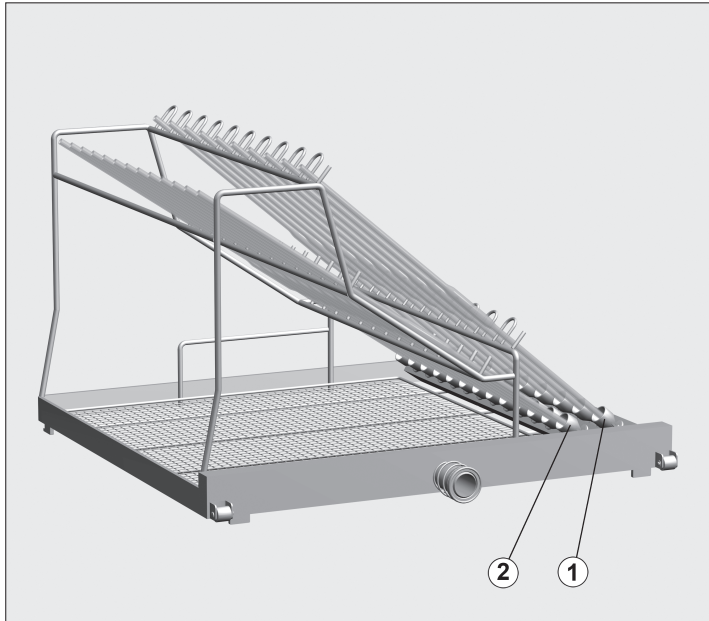
- With the pointed end facing downwards, place each pipette into the appropriate holder in the base.
- The upper ends of each pipette should be supported in the relevant holders on the support rail.

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