
Items supplied:

Mobile injector unit, H 502, W 535, D 515 mm, E 405/1 with connection for hot air drying unit and holder for ML/2 magnetic strip (optional extra) for automatic mobile unit recognition (G 7835, G 7836, PG 8535 and PG 8536)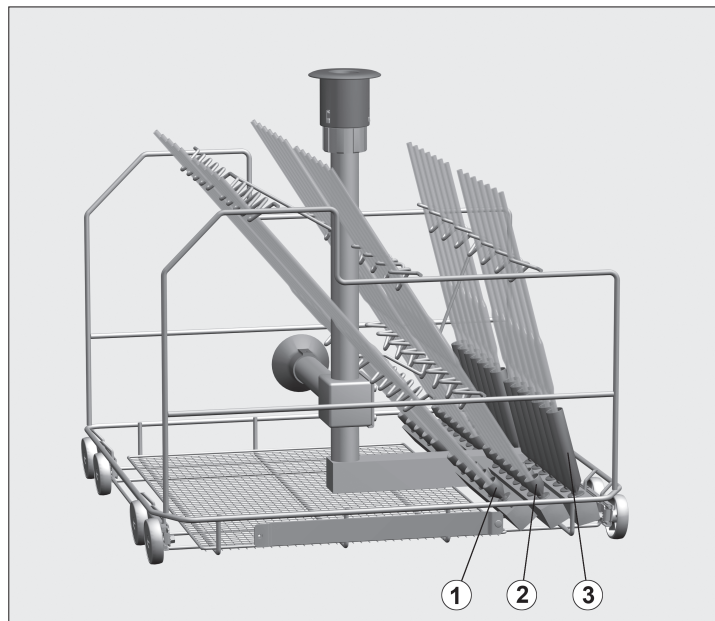

Intended use:

This mobile injector unit can be used for reprocessing a wide range of pipettes. Capacity:

Row	max. length [mm] *	Quantity	Volume [ml]
1	550	10	10 to 100
2	500	14	10 to 25
3	440	14	1 to 10

* The pipettes must not extend above the water intake pipe or behind the sides of the mobile injector unit.

This mobile injector unit should only be used for processing laboratory glassware which has been declared as suitable for machine processing by the glassware manufacturer.

 Please observe the Warning and Safety instructions in the Operating Instructions for your washer-disinfector as well as any instructions provided by the glassware manufacturer.

Broken glassware can cause serious injury.


All glassware should be properly inspected and tested for cleanliness. Do not rely on a visual check alone.

 Use the LAB PIPETTES programme or increase the water intake for the relevant reprocessing programmes as follows:


G 7836, PG 8536 + 5 litres

G 7835, G 7883, G 7883 CD, G 7893, PG 8535 + 6 litres

Placing pipettes into the injector unit: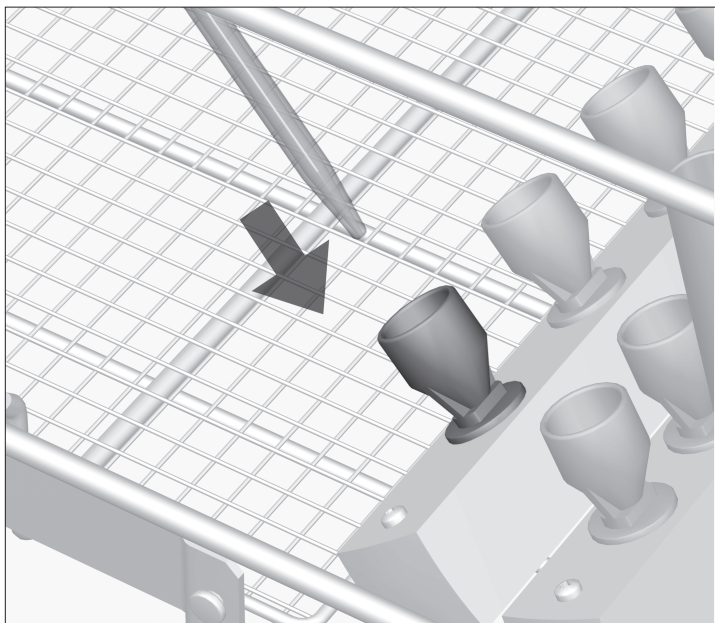

- With the pointed end facing downwards, place each pipette into the appropriate holder in the base.
- The upper ends of each pipette should be supported in the relevant holders on the support rail.

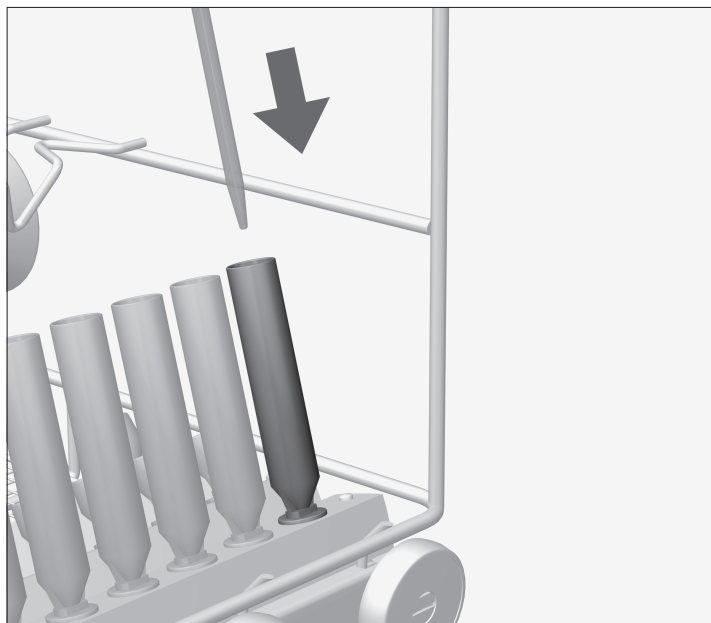- Arrange pipettes with a volume of between 1 and 5 ml in the injector sleeves in row ③ in the mobile injector unit.

The water pressure could push these pipettes out of the funnel shaped holders in rows ① and ②.

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