A 304

de	Gebrauchsanweisung Injektormodul
da	Brugsanvisning Injektormodul
en	Operating instructions Injector module
es	Instrucciones de manejo Módulo inyector
fi	Käyttöohje Suorasuihkumoduuli
fr	Mode d'emploi Module à injection
hr	Upute za uporabu Injektorski modul
it	Istruzioni d'uso Modulo a iniezione
nl	Gebruiksaanwijzing injectormodule
no	Bruksanvisning injektormodul
pt	Instruções de utilização Módulo injetor
sl	Navodila za uporabo Injektorski modul
sv	Bruksanvisning injektormodul

de	4
da	12
en	20
es	28
fi	36
fr	44
hr	52
it	60
nl	68
no	76
pt	84
sl	92
sv	100

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Notes about these instructions	20
Correct application	21
Queries and technical problems	21
Items supplied	22
Disposal of the packing material	22
Warning and Safety instructions	23
Areas of application	24
Injector modules	24
Inserting an injector module	24
Removing an injector module	24
Before loading the machine and before starting a programme	25
Preparing the load	26

Important warnings

 Information which is important for safety is highlighted in a thick framed box with a warning symbol. This alerts you to the potential danger of injury to people or damage to property.

Read these warning notes carefully and observe the procedural instructions and codes of practice they describe.

Notes

Information of particular importance that must be observed is highlighted in a thick framed box.

Additional information and comments

Additional information and comments are contained in a simple frame.

Operating steps

Operating steps are indicated by a black square bullet point.

Example:

- Select an option using the arrow button and save your choice with OK.

Display

Certain functions are shown in display messages using the same font as that used for the function itself in the display.

Example:

Settings  menu.

Re-usable laboratory utensils can be reprocessed using this module in a Miele laboratory washer-disinfector. Follow the instructions for your washer-disinfector as well as the instructions of the laboratory utensil manufacturer on how to reprocess their items by machine.

The A 304 injector module is designed for reprocessing centrifuge tubes, phials, test tubes for fraction collectors and auto-sampler tubes.

To ensure sufficient water pressure, only one module can be used at a time in each upper and lower basket. The module must always be placed in the right hand position. The remaining connections can be used for other modules.


This module is suitable for use in the following:

- Upper basket A 100
- Lower basket A 150

Laboratory glassware and laboratory utensils for reprocessing are referred to as the wash load if they are not more closely defined.

Queries and technical problems

If you have any queries or technical problems that you would like to discuss please contact Miele. Contact details can be found at the back of the Operating Instructions for your Washer-Disinfector or go to www.miele-professional.com.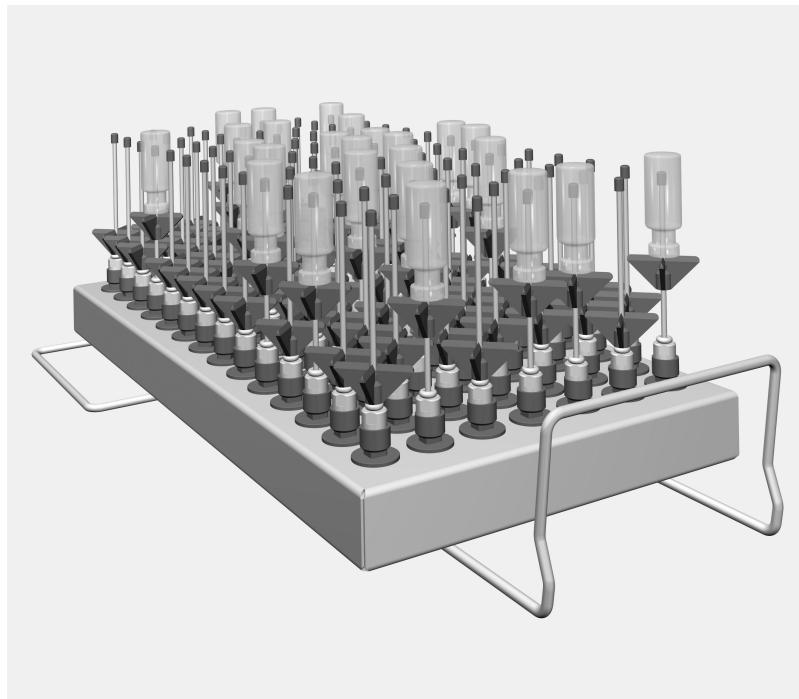

- Injector module A 304, height 130 mm, width 222 mm, depth 471 mm.

Items supplied

- 98 x ID 90 injector nozzles with plastic supports. Length 90 mm, Ø 2.5 mm

Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in landfill sites.

To avoid the risk of accidents and damage to this module please read these instructions carefully before using it for the first time. Keep these instructions in a safe place where they are accessible to users at all times.

⚠ Please also read the Operating instructions for your washer-disinfector and pay particular attention to the Warning and Safety instructions.

- This module is only for use in the areas of application as defined in the Operating instructions. Components such as injector nozzles must only be replaced with Miele accessories or original parts. Any other application, modification or change is unauthorised and could be dangerous.
Miele cannot be held liable for damage caused by unauthorised use or incorrect operation.
- Inspect all mobile units, baskets, modules and inserts daily as described in the "Maintenance" section in the Operating instructions for the washer-disinfector.
- Only items which have been declared by their manufacturer as suitable for machine reprocessing may be processed. The manufacturer's specific reprocessing instructions must be observed.
- Broken glass can result in serious injury when loading or unloading. Broken glass items must not be processed in the washer-disinfector.
- Only place empty, unloaded modules in baskets. Check that they are correctly engaged before loading them.
Modules must be completely emptied before taking them out.
Placing loaded modules in baskets or removing them can cause damage to the wash load and cause injury should glassware be broken.
- The reprocessing result must be checked as necessary using a suitable test method and not just by a visual check.

Miele cannot be held liable for damage caused by non-compliance with these Warning and Safety Instructions.

Injector modules


⚠ Only place empty, unloaded modules in baskets. Check that they are correctly engaged before loading them.

Modules must be completely emptied before taking them out.

Placing loaded modules in baskets or removing them can cause damage to the wash load and cause injury should glassware be broken.

Inserting an injector module

The injector module is only suitable for use in an A 100 upper basket or an A 150 lower basket. The injector nozzles must point upwards.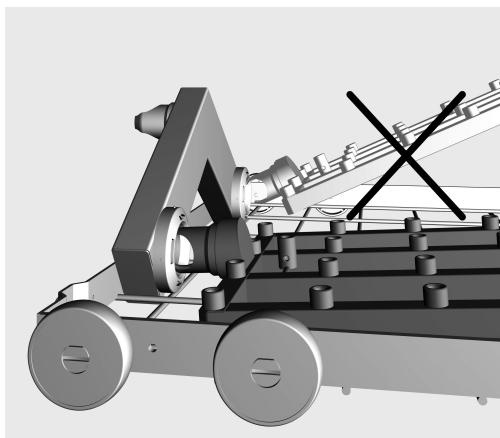

- Carefully slide the injector module in and connect it to the water connection point.
- Then press down on the module to secure it on the rails in the basket.

Removing an injector module

Do not attempt to pull the basket out of the washer-disinfector by the injector module.

The module could come loose or the water connection point could suffer damage.

- Release the module from its position, then carefully pull it off the water connection point. Do not pull it off at too steep an angle as any levering action could damage both the connection point and the coupling.

Before loading the machine and before starting a programme

- Check that fittings such as irrigation sleeves and injector nozzles are securely screwed into position.


 Make sure that all screw connectors are fitted with injector nozzles, adapters, irrigation sleeves or blind stoppers to ensure that all rinsing fittings in use are supplied with sufficient standardised pressure.

Damaged fittings such as injector nozzles, adapters and irrigation sleeves must not be used.


Rinsing positions not in use do not have to be fitted with blind stoppers.

- Check that the baskets and modules are correctly connected to the water supply.


Preparing the load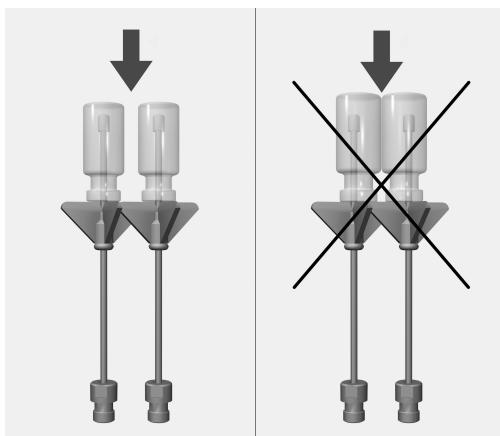

Individual items of glassware must not touch each other.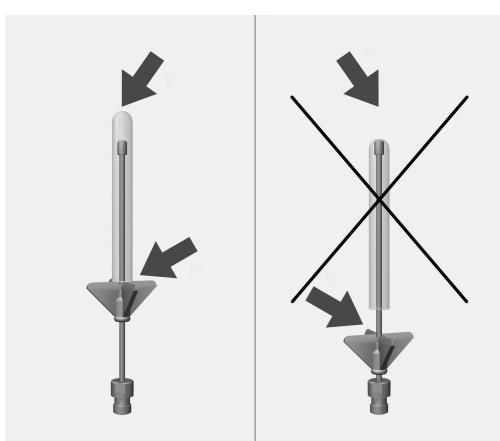

If the tip of an injector nozzle touches the bottom of a piece of glassware the plastic support has to be pushed upwards.

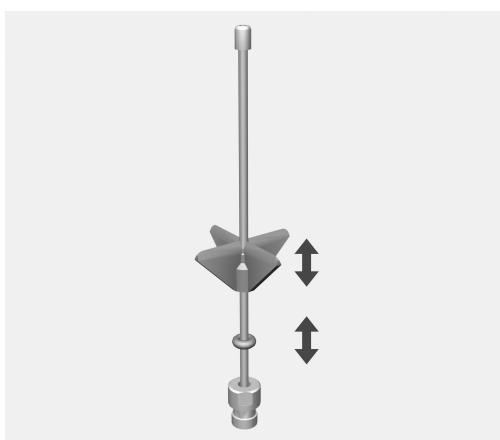- Push the plastic support upwards to the height required. If necessary use a pair of pointed pliers to do so.
- Push the O-ring up to the plastic support to secure it in position.

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